# IKA®

20000005499

Vortex\_1\_012016

## IKA® Vortex 1

Betriebsanleitung	DE	4
Operating instructions	EN	7
Mode d'emploi	FR	10
Indicaciones de seguridad	ES	13
Veiligheidsinstructies	NL	14
Normi di sicurezza	IT	15
Säkerhetsanvisningar	SV	16
Sikkerhedshenvisninger	DA	17
Sikkerhedsanvisninger	NO	18
Turvallisuusohjeet	FI	19
Normas de segurança	PT	20
Wskazówki bezpieczeństwa	PL	21
Bepečnostní upozornění	CS	22
Biztonsági utasítások	HU	23
Varnostna navodila	SL	24
Bezpečnostné pokyny	SK	24
Ohutusjuhised	ET	25
Drošības norādes	LV	26
Saugos reikalavima	LT	27

IKA®-Werke, Germany  
Reg. No. 004343

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**DE**

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## Contents

	Page
Safety instructions	7
Correct use	8
Unpacking	8
Commissioning	8
Maintenance and cleaning	8
Warranty	9
Technical data	9

## Safety instructions

### Read the operation instructions in full before starting up and follow the safety instructions.

Keep the operating instructions in a place where they can be accessed by everyone.

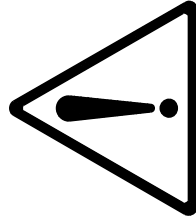
The appliance complies with international laboratory instrument safety regulations.

**Caution:** overfilling with liquid may cause the medium to splash out.

For this reason, only closed test tubes should be used.

Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.

Set up the appliance in a spacious area on an even, stable, clean, non-slip, dry and fireproof surface.

Wear your personal protective equipment in accordance with the hazard category of the medium to be processed.

Otherwise there is a risk of:

- splashing liquids,
- projectile parts,
- body parts, hair, clothing and jewellery getting caught.

Beware of the risk of:

- flammable materials.
- glass breakage as a result of mechanical shaking power.

Only process media that will not react dangerously to the extra energy produced through processing. This also applies to any extra energy produced in other ways, e.g. through light irradiation.

The appliance is maintenance-free and must not be opened.

**Do not** operate the appliance in explosive atmospheres, with hazardous substances or under water.

The feet of the appliance must be clean and undamaged.

Sharp edged vessels cause wear debris on the attachments. This can lead to the destruction of the attachments and the appliance.

## Correct use

The **IKA® Vortex 1** is designed for shaking test tubes. Gently push the test tube into the soft attachment to start shaking. Shaking stops as soon as you lift out the test tube. The shaking rate is set using the front knob (left stop = approx. 1000 rpm - right stop = approx. 2800 rpm).

## Unpacking

Please unpack the device carefully and check it for damage. It is important that any transport damage is detected when the device is unpacked. In the case of any damage a report must be sent immediately (post, rail or forwarder).

As supplied, the package contains:

The **IKA® Vortex 1**, a plug-in power supply unit and an instruction manual.

## Commissioning

Connect the appliance to a socket (please refer to the information in the "Technical Data" section).

## Maintenance and cleaning

The **IKA® Vortex 1** is maintenance-free. It is subject only to the natural wear and tear of components and their statistical failure rate. Clean the devices only with the following cleaning agents.

To remove use:

Dyes	isopropyl alcohol
Constructions materials	water containing tenside / isopropyl alcohol
Cosmetics	water containing tenside / isopropyl alcohol
Foodstuffs	water containing tenside
Fuels	water containing tenside

Wear the proper protective gloves during cleaning of the devices. Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.

Before using another then the recommended method for cleaning or decontamination, the user must ascertain with the manufacturer that this method does not destroy the instrument.

### Repair:

**Please send in instrument for repair only after it has been cleaned and is free from any materials which may constitute a health hazard.**

For repair, please request the "Decontamination Certificate" form **IKA®** or use the download printout of it from **IKA®** website: [www.ika.com](http://www.ika.com).

If you require servicing, return the instrument in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

## Warranty

You have purchased an laboratory machine which meets the highest engineering and quality standards.

In accordance with warranty conditions, the warranty period is 24 months. For claims under the warranty please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

## Technical data

### Plug-in power supply unit

Rated voltage	<b>VAC</b>	100 - 240
Rated frequency	<b>Hz</b>	50 - 60
Output voltage	<b>VDC</b>	12
Absorbed power	<b>W</b>	max. 30
Output power	<b>W</b>	7,2
Power input	<b>W</b>	1,5
<u>Drive</u>		
Speed adjustable	<b>rpm</b>	1000 .... 2800 ±25%
Operating time	<b>%</b>	100
Protection class		II
Pollution degree		2
Overvoltage category		II
Protection class acc. to DIN EN 60529		IP 40
Ambient humidity (relative)	<b>%</b>	80
Ambient temperature	<b>°C</b>	5 to 40 (in operation)
Operation at the terrestrial altitude	<b>m</b>	max. 2000
Transport temperature	<b>°C</b>	above sea level
Storage temperature	<b>°C</b>	-25 to +70
Weight	<b>kg</b>	-10 to +70
Installed weight (max.)	<b>kg</b>	0,55
Test tube size	<b>mm</b>	0,1
Dimensions (B x H x T)	<b>mm</b>	ø 30
Type of movement		95 x 70 x 110
Orbital diameter	<b>mm</b>	orbital
Touchfunction		4,5
		yes

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**IKA® -Werke GmbH & Co.KG**