Biosigma SpA

a Dominique Dutscher Company

Via Valletta, 6 | 30010 Cantarana di Cona (VE), Italy |
Tel. ++39 0426 302224 (r.a.) | Fax ++39 0426 302228 |

E-mail info@biosigmaeu.com | <http://www.biosigma.com> | www.biosigma.it

M521D

24/10/2024

Pag. 1 di 3

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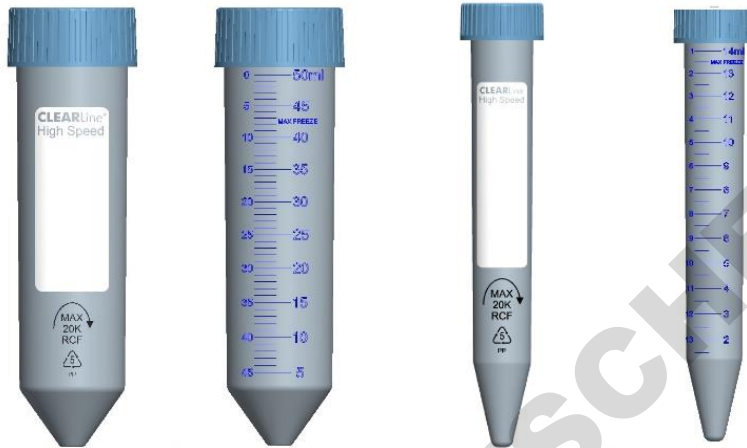
STERILE CENTRIFUGE TUBES with SCREW CAP

Human DNA, DNase, RNase, and bacterial Endotoxin free

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DESCRIPTION

- *Centrifuge tubes are made from transparent polypropylene and designed to withstand high centrifugal forces, making them suitable for a wide range of laboratory applications. These tubes are engineered to endure rigorous conditions, ensuring reliable and high-quality results. They are particularly effective for the sedimentation of cells in cell cultures, the isolation and purification of bacteria in microbiology, as well as the centrifugation, purification, and precipitation of nucleic acids and proteins in the field of molecular biology.*
- *Material: centrifuge tube made of Polypropylene (PP) and cap made of Polyethylene (PE) with inner ring for a perfect seal.*
- *The translucent wall facilitates clear observation of the contents.*
- *Two volumes are available: 15 ml and 50 ml.*
- *The tubes have been tested for centrifugation up to 20,000g (Sorvall X Pro Series Centrifuge).*
- *Indication of the maximum volume for storage at low temperatures.*
- *Clearly visible ascending and descending graduation lines printed on the tube body, with the graduation engraved directly into the plastic of the conical part of the tube.*
- *White writing area for better sample identification.*
- *Serigraphy is resistant to contact with acids, alcohols, and solvents.*
- *The recommended working temperature range is from -80°C to +120°C with appropriate precautions. The suitability of these tubes for storage below 0°C depends on both the solution and the storage conditions. It is strongly recommended to conduct a real-test to assess the suitability of the tubes for frozen storage. Due to their low temperature resistance, tube caps typically made of polyethylene cannot be recommended for autoclave sterilization. If autoclave sterilization is desired, it is advisable to unscrew the caps almost completely or remove them entirely before starting the sterilization process. Pressure differences that occur in the autoclave could cause tube collapse if the cap is screwed tightly. It is important to ensure that the sterilization method used is suitable for all components: tubes and caps. Otherwise, it is advisable to sterilize them separately using the most appropriate method for each. It should be noted that sterilization may result in a reduction in the mechanical, functional, or aesthetic properties of the product.*
- *Available either packaged loose in zip-lock bags or arranged in recyclable polystyrene racks. The racks are not autoclavable.*
- *The internal packaging includes all product-related information such as Catalog Code, Lot Number, and Expiry Date.*
- *Sterilized by electron beam (E-Beam).*
- *Free from human DNA, DNase, RNase, and bacterial Endotoxins. Non-cytotoxic.*
- *Metal free.*

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RECOMMENDATION USE

- *For generic laboratory use.*
- *In case of content spillage, use gloves: risk of contamination.*
- *Do not bring the device near flames or sources of heat that could damage it.*
- *Do not use after expiry date.*
- *Sterile product: Do not use if package damaged.*
- *Do not vary the intended purpose of the product.*
- *Product not suitable for children.*
- *Store in dry place, temperature range: min. -10°C max +50°C.*
- *Disposal: Use appropriate P.P.E. and dispose of according to current regulations.*

DOMINIQUE DUTSCHER SAS