

Protein A/G SpinTrap Buffer Kit

Instructions

Introduction

Important

Read these instructions carefully before using the products.

Intended use

The products are intended for research use only, and shall not be used in any clinical or *in vitro* procedures for diagnostic purposes.

Background

Protein A/G SpinTrap™ Buffer Kit is designed for protein enrichment using Protein A HP SpinTrap or Protein G HP SpinTrap columns.

- The kit is sufficient for 16 reactions and contains reagents for both the Classic and Crosslink protocols (with or without elution of antibody together with target protein). Some of the reagents are delivered in two bottles for increased stability.
- The Classic and Crosslink protocols are available in the user instruction delivered with Protein A HP SpinTrap, Protein G HP SpinTrap and Ab SpinTrap - see [Ordering information, on page 2](#).
- The kit consists of both stock solutions and reagents ready for use. Working solutions are prepared by adding distilled water directly into the stock solution bottle according to the Dilution Table.
- The kit eliminates time-consuming buffer preparation and thus promotes fast, reproducible and convenient enrichment of target protein from a complex protein mix.

Note: For stability reasons urea and DMP (Dimethyl pimelimidate dihydrochloride) are not included in the kit. Urea and DMP are used in the Crosslink protocol.

Kit content

Buffer	Buffer content	Formulation	No. of bottles	Volume
Binding/ Washing Buffer (TBS)	0.5 M Tris	10×	2	2 × 5 ml
	1.5 M NaCl pH 7.5			
Elution Buffer (Classic)	2.5% HAc	Ready to use	1	20 ml
Elution Buffer (Crosslink)	1 M Glycine-HCl pH 2.9	10×	1	3 ml
Crosslink Solution A	2 M Triethanolamine pH 8.9	10×	1	4 ml
Crosslink Solution B	1 M Ethanolamine	10×	2	2 × 1 ml

Buffer	Buffer content	Formulation	No. of bottles	Volume
	pH 8.9			

Dilution Table

Buffer	Stock solution volume	Water added	Final concentration
Binding/ Washing Buffer (TBS)	5 ml	45 ml	50 mM Tris 150 mM NaCl pH 7.5 ¹
Elution Buffer (Crosslink)	3 ml	27 ml	0.1 M Glycine- HCl pH 2.9 ²
Crosslink Solution A	4 ml	36 ml	200 mM Triethanolamine pH 8.9 ³
Crosslink Solution B	1 ml	9 ml	100 mM Ethanolamine pH 8.9

Preparations for Crosslink protocol

1. Washing Buffer is TBS (50 mM Tris; 150 mM NaCl) pH 7.5 with 2 M urea. Add urea to a final concentration of 2 M to a portion of the working solution (2 ml required/column)
2. Elution Buffer is 0.1 M Glycine-HCl pH 2.9 with 2 M urea. Add urea to a final concentration of 2 M to a portion of the working solution (1 ml required/column)
3. Triethanolamine is used for buffer change, cross-linking and washing. For cross-linking, DMP (Dimethyl pimelimidate dihydrochloride) is added to a final concentration of 50 mM. Add DMP to a portion of the working solution (400 µl required/column)

Ordering information

Products

Description	Quantity	Product code
Protein A/G SpinTrap Buffer Kit	1	28913567

Related products

Description	Quantity	Product code
Protein A HP SpinTrap	16 columns	28903132
Protein G HP SpinTrap	16 columns	28903134
Ab SpinTrap	50 columns	28408397

Literature

Title	Product code
Data File Ab SpinTrap	28902030
Data File Protein A HP SpinTrap	28906789
Data File Protein G HP SpinTrap	28906790
Antibody Purification Handbook	18103746
Affinity Chromatography Handbook	18102229

cytiva.com

Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.

SpinTrap is a trademark of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

All other third-party trademarks are the property of their respective owners.

© 2020–2021 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request. Contact your local Cytiva representative for the most current information.

For local office contact information, visit cytiva.com/contact

28923247 AD V:5 10/2021

