RESOLUCIÓN DE PROBLEMAS

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MALFUNCTION AND SOLUTION

No.	Phenomenon cause	Cause	Solution
1	No flame	A.gas have been used out or not enough	A.refilling
		B.no flame after refilling	B.wait for 10minute or adjust the switch if necessary
		C.the adjustable leve is too near the "YELLOW"	C.turn adjustable lever to "BLUE"
		D.the pipe or eoretric leak out	D.replace the accessory
2	The flame is too large	The gas is in large quantity	Turn the button to "-" till normal position
3	The flame is too samll	A.the gas is too little	A.turn the button to "+", till normal position
		B.air pressure is too low or the gas is volatilized	B.use hand to warm it till normal
		C.gas in the tank is not enough	C.refilling
4	Flame is in tampon shape	A.accessory is blocked	A.replace the accessory
		B.turn adjustable lever to "YELLOW"	B.the adjustable lever is too near to "BLUE"

Notice:No one can repair the lighter except honest engineer

TECHNICAL PARAMETER

Product size	base:141.5x67wx69h
Gross weight:	323.6g
Apply to gas pressure:	0.36x0.42Mpa
Burning time:	40-50mins
Flame tempature	1300 °C
Flame chan ging	The minimum of the flame is 20mm,the maxinmun of the55mm under the 22 degrees of the room temperature
Gas-tank capacity	8.5g

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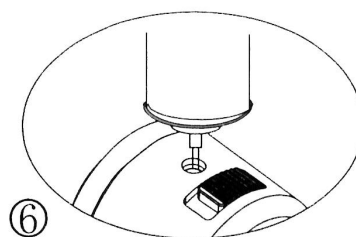
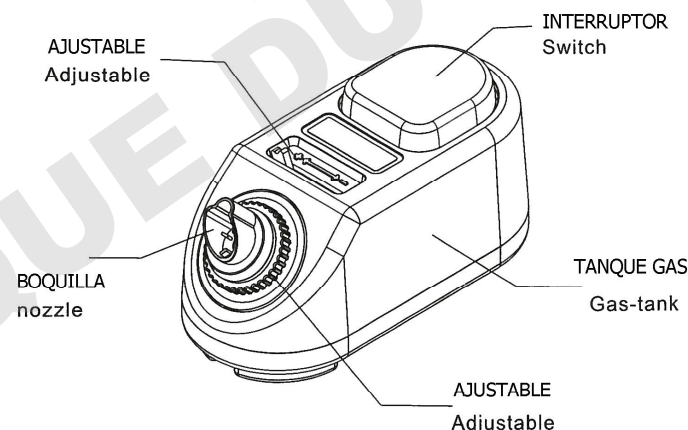
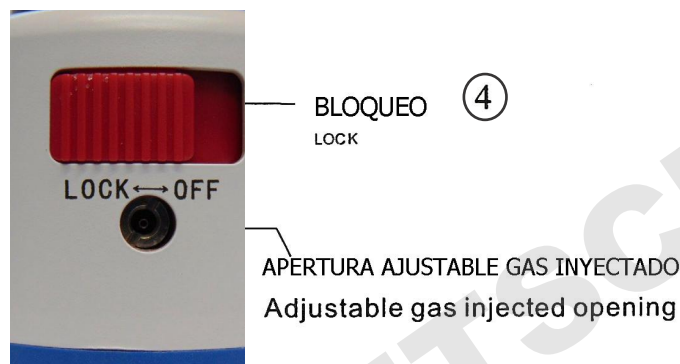
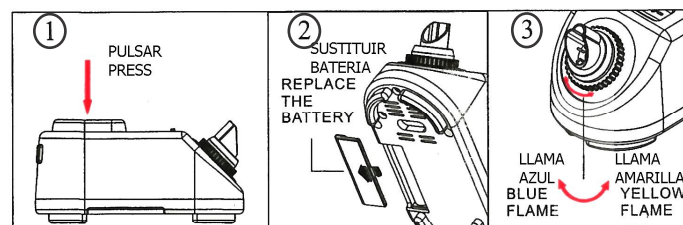
**OPERATION:****HOW TO IGNITE (Fig.1)**

The adjustable lever must be raised. Press the switch to turn on. When you stop pressing, it will turn off.

REPLACING THE BATTERY (Fig.2)

Battery Specifications LR6: 5 1.5V

As shown in the picture, open the cover, replace the battery according to the original position, and turn on the burner.

Note: Be careful not to reverse the polarity.

ADJUSTING THE FLAME (Fig.3)

When the flame is on, you will need to adjust it: BLUE when you need a longer flame and YELLOW with a shorter flame.

ADJUSTMENT SIZE

This lighter can be adjusted according to the requirements of different work situations or fire.

"+" or "-" denote the size of the flame. If the flame is too large, it can be adjusted without being less than 15mm or the flame will turn off automatically.

HOW TO LOCK (Fig.4)

Unlock the burner (OFF) for use. Once its daily use is over, lock the ignition (LOCK) again. For safety check that the lock is locked.

WARNINGS:

- The length of the flame should not be too large.
- The burner may not be on for a long time and should not be turned upside down.
- Wait 10 minutes after refilling and do not shake it vigorously.
- Keep away from flames, sunlight or temperatures above 40°C.
- Keep out of the reach of children.
- Keep away from face and clothing.

FILLING THE GAS (Fig.6)

Put the burner upside down and place the recharger valve in the gas inlet. Press for a few seconds. Keep away from fire during filling and do not light the burner.

WARNINGS:

This product has a self-filling function. The gas could be spilled out when reaching the safe quantity. Please stop refilling it or the burner could not work properly and could even cause damages.