POSITIVE CONTROL (PG + OT)

Ref. ZE/PGOT5

Descripción

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Ref. ZE/PGOT5

Description

Positive control vials with penicilin G and oxytetracycline for use with Duplex BT Scan, BT Scan Strip and Beta Scan.

Kit components

	ZE/PGOT5
Vials	5
Kit instructions	✓
Product certificate	✓

Stability and Storage

This reagent should be stored under refrigeration (4-12°C) until use. Once the powder has been dissolved, the preparation should be kept in freezer for a maximum of one week. To avoid freeze-thaw cycles it is recommended to divide the volume in aliquots adapted to the use (for example 0.5 ml).

Safety

Good laboratory practices are recommended when using this kit. A Safety Data Sheet (SDS) is available from your local distributor or ZEULAB by request.

Preparation

Add 5 ml of a negative milk sample to the vial and stir gently. Leave the solution for 15 min at room temperature and mix the vial few times during that time for full dissolution.

Note. The negative milk sample must not contain antibiotic residues. These samples should have been previously tested as negative. It is recommended that the negative sample belong to the same matrix and specie as the samples to be analysed.

ZEULAB, S.L.

C/ Bari, 25 dpdo. • 50197 Zaragoza (SPAIN)

Tel.: +34 976 731 533

info@zeulab.com • www.zeulab.com