EN

ES

FR

DE


Defender[®] Series Base User Guide

IT

PT

SE

NL

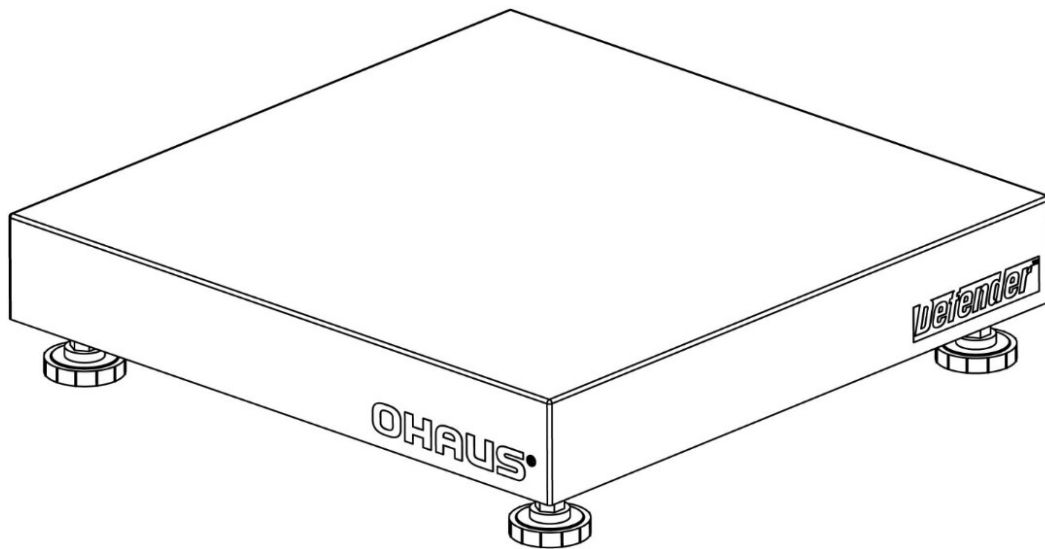DK

PL

CZ

HU

INTRODUCTION

This manual covers installation, and maintenance instructions for the Ohaus Defender® Series Base. Please read this manual completely before installation and operation.

SAFETY PRECAUTIONS

For safe and dependable operation of this product, please comply with the following precautions:

- Operate the base only under ambient conditions specified in these instructions.
- Ensure that the load cell cable does not pose an obstruction or tripping hazard.
- Do not operate the base in hazardous environments or unstable locations.
- Do not drop loads on the base.
- Do not lift the base by the top frame; always lift from the bottom frame when moving the base.
- Service should only be performed by authorized personnel.

INSTALLATION

Unpacking

Unpack and inspect the product to make sure that all components have been included. The package includes the following:

- Defender™ Series Base
- Weighing Pan
- User Guide


When purchased as a complete Defender Series Scale, the package will also include:

- Indicator
- Column Assembly

ASSEMBLY

Weighing Pan

Remove the shipping spacers, and place the weighing pan securely over the rubber load pads on the top frame of the base.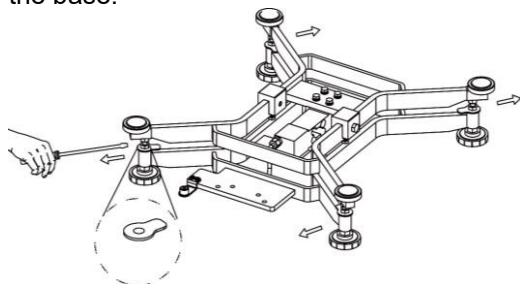

Note: Not included with some models.

Wiring Connections

When connecting the base to an Ohaus indicator equipped with the matching circular connector, push the base connector onto the indicator connector and turn the locking ring clockwise to lock it in place.

When connecting the base to an indicator that does not have the matching connector, install the optional load cell cable adapter (P/N 30429931 sold separately) to the indicator using the wiring codes in Table 1. Then attach the load cell cable to the cable adapter.

Note: As an alternate wiring method, cut off the connector, strip the wires and attach them to the indicator using the wiring codes in Table 1.

- Detailed instructions that describe how to prepare the load cell cable for connection to the terminal block inside an indicator.

- The extension cable accessory (P/N: 30429931) cannot be used if the scale will be used in a legal for trade or legally controlled application, because there is no provision to seal the connection between the indicator and platform.

TABLE 1. LOAD CELL CONNECTION


Type1. DxxxRxx Bases
DxxxWxx Bases

FUNCTION	WIRE COLOR
+ Excitation	Green
- Excitation	Black
+ Signal	White
- Signal	Red
+ Sense	Blue
- Sense	Brown
Shield	Yellow

Selecting the Location

Level the base by adjusting the four leveling feet until the bubble in the level indicator is centered.

Note: Ensure that the base is level each time its location is changed.

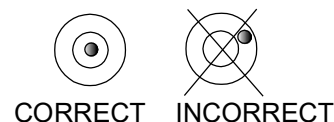MAINTENANCE

Cleaning

The base components should be kept clean and free of excessive material build up.

- A damp cloth with water and a mild detergent may be used to wipe clean the external surfaces – do not use acids, alkalis, strong solvents or abrasive materials.

To maintain surface integrity, do not allow debris to remain on the surface of the painted or stainless steel base frame.

- Remove the pan and use fresh water to clean the base frame surface of any debris. Dry the base with a clean cloth or allow to air dry before replacing the pan. Repeat daily for best care.

Troubleshooting

An Ohaus product service specialist will be available to provide assistance. Outside the USA, please visit our web site, www.ohaus.com to locate the Ohaus office nearest you.

TECHNICAL DATA

Technical Specifications

The technical data is valid under the following ambient conditions:

Ambient temperature: -10°C to +40 °C

Relative humidity: 10% to 90% relative humidity, non-condensing

Altitude: Up to 2,000 m

Operability is assured at ambient temperature -10°C to +40 °C

Note: Not all models may be available in your area. Check with your local Ohaus representative.

TABLE 1 SPECIFICATIONS

MODEL	D6RQDR	D15RQDR	D15RQDL	D30RQDL	D60RQDL	D60RTDL	D150RTDL
Capacity	3 / 6 kg	6 / 15 kg	6 / 15 kg	15 / 30 kg	30 / 60 kg	30 / 60 kg	60 / 150 kg
Approved Resolution	OIML 2 x 3000e						
Safe Overload Capacity	150% of capacity						
Pan Dimensions	305 x 305 mm 12 x 12 in		400 x 400 mm 15.7 x 15.7 in			400 x 500 mm 15.7 x 19.6 in	
Base Construction	Stainless Steel platform with painted steel frame and rubber leveling feet						
Repeatability (std.deviation)	2d						
Linearity	±2d						
Load Cell Capacity	11 kg	22 kg	22 kg	50 kg	100 kg	100 kg	250 kg
Load Cell Cable	2 m L x 6-wire						
Load Cell Type	350 Ohm, aluminum, single point						
Load Cell Excitation	5-15V DC/AC						
Load Cell Rated Output	2mV/V						
Load Cell Protection	IP67						
Net Weight	6 kg		11.5 kg			17 kg	
Shipping Weight	7 kg		13.5 kg			19 kg	

TABLE 2 SPECIFICATIONS

MODEL	D60RQDX	D150RQDX	D60RQDV	D150RQDV	D300RQDV	D150RTDV	D300RTDV	D600RTDV
Capacity	30 / 60 kg	60 / 150 kg	30 / 60 kg	60 / 150 kg	150 / 300 kg	60 / 150 kg	150 / 300 kg	300 / 600 kg
Approved Resolution	OIML 2 x 3000e							
Safe Overload Capacity	150% of capacity							
Pan Dimensions	500 x 500 mm 19.6 x 19.6 in		610 x 610 mm 24 x 24 in			600 x 800 mm 23.6 x 31.5 in		
Base Construction	Stainless Steel platform with painted steel frame and rubber leveling feet							
Repeatability (std.deviation)	2d							
Linearity	±2d							
Load Cell Capacity	100 kg	250 kg	100 kg	250 kg	500 kg	250 kg	500 kg	750 kg
Load Cell Cable	2 m L x 6-wire							
Load Cell Type	350 Ohm, aluminum, single point							
Load Cell Excitation	5-15V DC/AC							
Load Cell Rated Output	2mV/V							
Load Cell Protection	IP67							
Net Weight	19 kg		31 kg			39.2 kg		
Shipping Weight	21 kg		34 kg			42.7 kg		

TABLE 3 SPECIFICATIONS

MODEL	D6WQDR	D15WQDR	D15WQDL	D30WQDL	D60WQDL	D60WTDL	D150WTDL
Capacity	3 / 6 kg	6 / 15 kg	6 / 15 kg	15 / 30 kg	30 / 60 kg	30 / 60 kg	60 / 150 kg
Approved Resolution	OIML 2 x 3000e						
Safe Overload Capacity	150% of capacity						
Pan Dimensions	305 x 305 mm 12 x 12 in		400 x 400 mm 15.7 x 15.7 in			400 x 500 mm 15.7 x 19.6 in	
Base Construction	Stainless Steel platform with stainless steel frame and rubber leveling feet						
Repeatability (std.deviation)	2d						
Linearity	±2d						
Load Cell Capacity	11 kg	22 kg	22 kg	50 kg	100 kg	100 kg	250 kg
Load Cell Cable	2 m L x 6-wire						
Load Cell Type	350 Ohm, aluminum, single point						
Load Cell Excitation	5-15V DC/AC						
Load Cell Rated Output	2mV/V						
Load Cell Protection	IP67						
Net Weight	5 kg		13 kg			15.5 kg	
Shipping Weight	6 kg		14 kg			18.5 kg	

TABLE 4 SPECIFICATIONS

MODEL	D60WQDX	D150WQDX	D60WQDV	D150WQDV	D300WQDV	D150WTDV	D300WTDV	D600WTDV
Capacity	30 / 60 kg	60 / 150 kg	30 / 60 kg	60 / 150 kg	150 / 300 kg	60 / 150 kg	150 / 300 kg	300 / 600 kg
Approved Resolution	OIML 2 x 3000e							
Safe Overload Capacity	150% of capacity							
Pan Dimensions	500 x 500 mm 19.6 x 19.6 in		610 x 610 mm 24 x 24 in			600 x 800 mm 23.6 x 31.5 in		
Base Construction	Stainless Steel platform with stainless steel frame and rubber leveling feet							
Repeatability (std.deviation)	2d							
Linearity	±2d							
Load Cell Capacity	100 kg	250 kg	100 kg	250 kg	500 kg	250 kg	500 kg	750 kg
Load Cell Cable	2 m L x 6-wire							
Load Cell Type	350 Ohm, aluminum, single point							
Load Cell Excitation	5-15V DC/AC							
Load Cell Rated Output	2mV/V							
Load Cell Protection	IP67							
Net Weight	16.5 kg		28 kg			36.5 kg		
Shipping Weight	19.5 kg		32.5 kg			41 kg		

TABLE 5 SPECIFICATIONS

MODEL	D15WQR	D15WQL	D30WQL	D60WQL	D60WQX	D150WQX	D60WQV	D150WQV
Capacity	15 kg	15 kg	30 kg	60 kg	60 kg	150 kg	60 kg	150 kg
Approved Resolution	OIML 3000e							
Safe Overload Capacity	150% of capacity							
Pan Dimensions	305 x 305mm 12 x 12 in		400 x 400 mm 15.7 x 15.7 in		500 x 500 mm 19.6 x 19.6 in		610 x 610 mm 24 x 24 in	
Base Construction	Stainless Steel platform with stainless steel frame and rubber leveling feet							
Repeatability (std.deviation)	2d							
Linearity	±2d							
Load Cell Capacity	30 kg	30 kg	50 kg	100 kg	150 kg	300 kg	150 kg	300 kg
Load Cell Cable	2 m L x 6-wire							
Load Cell Type	350 Ohm, stainless steel, single point							
Load Cell Excitation	5-15V DC/AC							
Load Cell Rated Output	2mV/V							
Load Cell Protection	IP67							
Net Weight	5 kg		13 kg		16.5 kg		28 kg	
Shipping Weight	6 kg		14 kg		19.5 kg		32.5 kg	

COMPLIANCE

	<p>This product complies with the applicable harmonized standards of EU Directive 2011/65/EU (RoHS). The EU Declaration of Conformity is available online at www.ohaus.com/ce.</p>
	<p>Disposal</p> <p>In conformance with the European Directive 2012/19/EU on Waste Electrical and Electronic Equipment (WEEE) this device may not be disposed of in domestic waste. This also applies to countries outside the EU, per their specific requirements.</p> <p>Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment.</p> <p>If you have any questions, please contact the responsible authority or the distributor from which you purchased this device.</p> <p>Should this device be passed on to other parties (for private or professional use), the content of this regulation must also be related.</p> <p>For disposal instructions in Europe, refer to www.ohaus.com/weee.</p> <p>Thank you for your contribution to environmental protection.</p>

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P/N 30428871B © 2018 Ohaus Corporation, all rights reserved.
Ohaus Corporation