



# Defender<sup>®</sup> 5000 Indicators Instruction Manual



TD52XW



TD52P



## TABLE OF CONTENTS

|                                             |           |
|---------------------------------------------|-----------|
| <b>1. INTRODUCTION</b> .....                | <b>5</b>  |
| 1.1 Safety Precautions .....                | 5         |
| 1.1.1 Relay Option Safety Precautions ..... | 5         |
| 1.2 Overview of Parts and Controls .....    | 6         |
| 1.3 Control Functions.....                  | 9         |
| <b>2. INSTALLATION</b> .....                | <b>10</b> |
| 2.1 Unpacking .....                         | 10        |
| 2.2 External Connections.....               | 10        |
| 2.2.1 Scale Base with Connector.....        | 10        |
| 2.2.2 RS232 interface Cable to TD52P .....  | 10        |
| 2.2.3 AC Power to TD52P .....               | 10        |
| 2.2.4 AC Power to TD52XW.....               | 10        |
| 2.2.5 Battery Power.....                    | 10        |
| 2.3 Internal Connections.....               | 11        |
| 2.3.1 Opening the Housing .....             | 11        |
| 2.3.2 Scale Base Without Connector .....    | 11        |
| 2.3.3 RS232 Interface Cable to TD52XW ..... | 13        |
| 2.3.4 MICRO SD Card Installation.....       | 14        |
| 2.4 TD52XW Rear Housing Orientation .....   | 14        |
| 2.5 Mounting Bracket .....                  | 14        |
| <b>3. SETTINGS</b> .....                    | <b>15</b> |
| 3.1 Menu Structure .....                    | 15        |
| 3.2 Menu Navigation.....                    | 18        |
| 3.3 Calibration Menu .....                  | 18        |
| 3.3.1 Zero Calibration .....                | 18        |
| 3.3.2 Span Calibration .....                | 19        |
| 3.3.3 Linearity Calibration .....           | 20        |
| 3.3.4 GEO Adjustment .....                  | 21        |
| 3.4 Setup Menu .....                        | 22        |
| 3.4.1 Capacity Unit.....                    | 22        |
| 3.4.2 Range .....                           | 22        |
| 3.4.3 Capacity .....                        | 22        |
| 3.4.4 Graduation.....                       | 23        |
| 3.4.5 Language .....                        | 23        |
| 3.4.6 Power On Zero .....                   | 23        |
| 3.4.7 Power On Unit .....                   | 23        |
| 3.4.8 Key Beep.....                         | 23        |