# 1. SAFETY INFORMATION

## Definition of Signal Warnings and Symbols

<b>WARNING</b>	For a hazardous situation with medium risk, possibly resulting in injuries or death if not avoided.
<b>CAUTION</b>	For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or injuries if not avoided.
<b>Attention</b>	For important information about the product.
<b>Note</b>	For useful information about the product.

## Warning Symbols

General hazard

Electrical shock

## Safety Precautions

**CAUTION:** Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Verify that the local AC power supply voltage is within the input voltage range printed on the AC adapter's ratings label.
- Only connect the AC adapter to a compatible grounded socket.
- Position the instrument such that the AC adapter can be easily disconnected from the socket.
- Position the power cord so that it does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in the user instructions.
- The equipment is for indoor use only.
- Do not operate the equipment in hazardous or explosive environments.
- Only use the equipment in dry locations.
- Only use approved accessories and peripherals.
- Disconnect the equipment from mains power before cleaning or servicing.
- Service should only be performed by authorized personnel.

## Intended Use

Use the instrument exclusively for weighing as described in the operating instructions. Any other type of use and operation beyond the limits of technical specifications without written consent from OHAUS, is considered as not intended.

This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use.

If the instrument is not used according to these operating instructions, the intended protection of the instrument may be compromised and OHAUS assumes no liability.

Please visit [www.ohaus.com/support](http://www.ohaus.com/support) to download user manual.

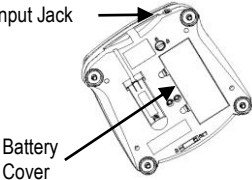
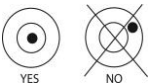
## 2. UNPACKING

Carefully remove your balance and each of its components from the package. Store the packaging material to ensure safe storage and future transport.

Included Components:

- Balance
- Pan
- AC Adapter & Plug
- User Guide
- Warranty Card

## 3. PREPARATION FOR USE

<p><b>Step-1</b></p>	<p><b>Selecting the location</b></p> <p>Place on a firm, steady surface with sufficient space. Avoid excessive air current, vibrations, heat sources, or rapid temperature changes</p>	
<p><b>Step-2</b></p>	<p><b>Transportation Lock</b></p> <p>The Transportation Lock is located under the balance. Rotate the pointer to the unlocked position.</p> <p><b>Note:</b> No shipping lock on Approved models</p>	
<p><b>Step-3</b></p>	<p><b>Connect Power</b></p> <p>Connect the AC adapter to the balance input jack.</p> <p><i>Battery installation</i></p> <p>Install the four "C" batteries (not included) with polarity as shown in the battery compartment.</p>	
<p><b>Step-4</b></p>	<p><b>Leveling the Equipment</b></p> <p>To level the instrument, adjust the feet/leveling wheel so the bubble is centered in the circle of the level indicator. Be sure the equipment is level each time its location is changed.</p>	

**Note:**

After power on, it is recommended to let the balance warm up for at least 5 minutes before using it.

## 4. INITIAL CALIBRATION

This user guide explains how to use the basic functions of the instrument. For information about advanced features, refer to the instruction manual.

When the balance is first installed it should be calibrated to ensure accurate results.

Press and hold Menu until [**MENU**] (Menu) is displayed. When the button is released, the display will display [**C.A.L.**]. Press "Yes" to accept, [**SPAN**] will then be shown. Press "Yes" again to begin the span calibration.

[**--[ ]--**] blinks while zero reading is stored. Next, the display shows the calibration weight value. Place the specified calibration mass on the pan. [**--[ ]--**] blinks while the reading is stored. The balance returns to the previous application mode and is ready for use.

## 5. OPERATION

### 5.1. Controls

Button	Functions
<b>Zero</b> ①	Short Press (when on): Sets display to zero (when off): Turns balance on Long Press (when on): Turns the balance off
<b>Yes</b>	Short Press (in Menu): Selects/accepts displayed setting
<b>Print Units</b> <b>No</b>	Short Press: See Interface Manual for operation description. Long Press: Toggles through active units Short Press (in Menu): Toggles through available settings
<b>Function Mode</b> <b>Back</b>	Short Press: Selects function setting Long Press: Selects active Mode Short Press (in Menu): returns to previous settings
<b>Tare</b> <b>Menu-Cal</b> <b>Exit</b>	Short Press: Enter / clear a Tare value Long Press: Enters User Menu Short Press (in Menu): Quickly exit User Menu

### 5.2. Menu Structure

Sub-Menus:	<i>.C.a.l.</i>	<i>.S.e.t.u.p.*</i>	<i>.M.o.d.e.</i>	<i>.U.n.i.t.*</i>	<i>.E.n.d.</i>
<i>Menu Items:</i>	<i>Span</i> <i>Lin</i>	<i>b.light</i> <i>A-OFF</i> <i>IR</i> <i>Filter</i> <i>AZT</i> <i>Stab</i> <i>Stab.C</i> <i>End</i>	<i>Count</i> <i>Percnt</i> <i>Check</i> <i>End</i>	<i>g, kg, ct, N,</i> <i>oz, ozt,</i> <i>dwt, lb,</i> <i>lb:oz, grn,</i> <i>thk, tsg,</i> <i>ttw, tola,</i> <i>tical</i> <i>End</i>	

\* Note: Available settings vary by models and regions.

### 5.3. Using the Instrument

#### Weighing Application

1. If required, place an empty container on the pan and press **Tare**.
2. Add sample to the pan or container. The display shows the weight of the sample.

## 6. MAINTENANCE

### 6.1. Cleaning

**WARNING:** Electric Shock Hazard. Disconnect the equipment from the power supply before cleaning. Make sure that no liquid enters the interior of the instrument.

**Attention:** Do not use solvents, harsh chemicals, ammonia or abrasive cleaning agents.

The exterior surfaces of the instrument may be cleaned with a cloth dampened with water and a mild detergent.

### 6.2. Troubleshooting

For technical issues contact an Authorized Ohaus Service Agent. Please visit our website [www.ohaus.com](http://www.ohaus.com) to locate the Ohaus office nearest you.

## 7. TECHNICAL DATA

Equipment Ratings:

- Pollution degree 2;
- Installation category II;
- Altitude 2000m;
- Humidity: Maximum 80% for temperatures up to 31°C decreasing linearly to 50% at 40°C; non-condensing;
- Electrical supply: Rated 12VDC 500mA for use with a Certified/Listed power adapter or battery operated;
- Indoor use only;
- Temperature range: 10°C to 40°C; Approved models (M) 0°C to 40°C
- The mains supply voltage fluctuations are not to exceed ± 10 % of the nominal supply voltage.

Model	NV123	NV223	NV323	NV222	NV422	NV622	NV1202	NV2202	NV3202	NV221	NV621	NV1201	NV2201
Capacity (g)	120	220	320	220	420	620	1200	2200	3200	220	620	1200	2200
Readability (g)	0.001	0.001	0.001	0.01	0.01	0.01	0.01	0.01	0.01	0.1	0.1	0.1	0.1
Repeatability (Std. Dev.) (g)	0.002	0.002	0.002	0.01	0.02	0.02	0.02	0.02	0.02	0.1	0.1	0.1	0.1
Linearity (g)	0.003	0.005	0.005	0.02	0.02	0.02	0.03	0.05	0.05	0.2	0.2	0.2	0.2

Model	NVT2201	NVT4201	NVT6201	NVT10201	NVT2200	NVT6200	NVT12000	NVT22000
Capacity (g)	2200	4200	6200	10200	2200	6200	12000	22000
Readability (g)	0.1	0.1	0.1	0.1	1	1	1	1
Repeatability (Std. Dev.) (g)	0.1	0.2	0.2	0.2	1	1	1	1
Linearity (g)	0.2	0.2	0.2	0.2	2	2	2	2

**Approved Models:**

Model	Max	d	e	Approval Class
NVT1601M	1600g	0.5g	0.5g	III
NVT3200M	3200g	1g	1g	III
NVT6400M	6400g	2g	2g	III
NVT16000M	16000g	5g	5g	III

**8. COMPLIANCE**

Compliance to the following standards is indicated by the corresponding mark on the product.

Mark	Standard
	This product complies with the EU Directives 2011/65/EU (RoHS), 2014/30/EU (EMC), 2014/35/EU (LVD) and 2014/31/EU (NAWI). The EU Declaration of Conformity is available online at <a href="http://www.ohaus.com/ce">www.ohaus.com/ce</a> .
	This product complies with the EU Directive 2012/19/EU (WEEE). Please dispose of this product in accordance with local regulations at the collecting point specified for electrical and electronic equipment. For disposal instructions in Europe, refer to <a href="http://www.ohaus.com/weee">www.ohaus.com/weee</a> .
	EN 61326-1

**Verified weighing instruments**

When the instrument is used in trade or a legally controlled application it must be set up, verified and sealed in accordance with local weights and measures regulations. It is the responsibility of the purchaser to ensure that all pertinent legal requirements are met. Weighing Instruments verified at the place of manufacture bear the following supplementary metrology marking on the descriptive plate.

  1259

Weighing Instruments to be verified in two stages have no supplementary metrology marking on the descriptive plate. The second stage of conformity assessment must be carried out by the applicable weights and measures authorities.

If national regulations limit the validity period of the verification, the user of the weighing instrument must strictly observe the re-verification period and inform the weights and measures authorities.

As verification requirements vary by jurisdiction, the purchaser should contact their local weights and measures office if they are not familiar with the requirements.

**INSTRUCTION MANUAL**

For more detailed information about the balance, please refer to the product instruction manual. The instruction manual can be obtained either by scanning the QR code or by going to the product page and from there download it. Go to [www.ohaus.com](http://www.ohaus.com) to register your product's warranty.

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P/N 30467942 E © 2020 Ohaus Corporation, all rights reserved.