



**Defender[®] 6000 Indicators
Instruction Manual**

**Defender[®] 6000 Waagen
Bedienungsanleitung**

**Indicadores Defender[®] 6000
Manual de instrucciones**

**Manuel d'instructions sur les
indicateurs de Defender[®] 6000**

**Indicatori[®] Defender 6000
Manuale di istruzioni**

Table of Contents

| | | |
|-----------|--|-----------|
| 1. | INTRODUCTION | 3 |
| 1.1. | SAFETY PRECAUTIONS | 3 |
| 1.2. | INTENDED USE | 3 |
| 1.3. | OVERVIEW OF PARTS AND CONTROLS | 4 |
| 1.4. | MAINBOARD | 5 |
| 1.5. | CONTROL FUNCTIONS | 6 |
| 2. | INSTALLATION | 9 |
| 2.1 | UNPACKING | 9 |
| 2.2 | EXTERNAL CONNECTIONS..... | 9 |
| 2.2.1 | Scale Base with EasyConnect™ Connector..... | 9 |
| 2.2.2 | Power input to i-DT61PW..... | 9 |
| 2.2.3 | AC Power to i-DT61XWE..... | 9 |
| 2.3 | INTERNAL CONNECTIONS | 9 |
| 2.3.1 | Opening the Housing..... | 9 |
| 2.3.2 | Scale Base without EasyConnect™ Connector..... | 11 |
| 2.3.3 | Communication Interface Cable to i-DT61PW..... | 14 |
| 2.3.4 | RS232 Interface Cable to i-DT61XWE..... | 15 |
| 2.4 | MOUNTING BRACKET..... | 15 |
| 3. | OPERATION | 16 |
| 3.1 | TURNING THE SCALE ON/OFF..... | 16 |
| 3.2 | WEIGHING MODE | 16 |
| 3.2.1 | Enter the Mode and Start Weighing | 16 |
| 3.2.2 | Accumulation and Statistics | 16 |
| 3.2.3 | Check | 17 |
| 3.2.4 | Application Settings | 18 |
| 3.3 | COUNTING MODE | 19 |
| 3.3.1 | Enter the Mode | 19 |
| 3.3.2 | Establish an APW | 19 |
| 3.3.3 | Start Counting | 19 |
| 3.3.4 | Application Settings | 19 |
| 3.4 | PERCENT MODE | 20 |
| 3.4.1 | Enter the Mode | 20 |
| 3.4.2 | Establish a Reference Weight..... | 20 |
| 3.4.3 | Start Percent Weighing | 20 |
| 3.4.4 | Application Settings | 20 |
| 3.5 | DYNAMIC MODE..... | 21 |
| 3.5.1 | Enter the Mode | 21 |
| 3.5.2 | Start Dynamic Weighing | 21 |
| 3.5.3 | Application Settings | 21 |
| 3.6 | FILLING MODE..... | 22 |
| 3.6.1 | Enter the Mode | 22 |
| 3.6.2 | Start Filling | 22 |
| 3.6.3 | Resume and Pause Filling..... | 22 |
| 3.6.4 | Display of the Dot Matrix Screen..... | 22 |
| 3.6.5 | Application Settings | 23 |
| 4. | MENU SETTINGS | 24 |
| 4.1 | MENU NAVIGATION | 24 |
| 4.1.1 | User Menu..... | 24 |
| 4.1.2 | Button Navigation..... | 25 |
| 4.2 | CALIBRATION MENU..... | 26 |
| 4.2.1 | Initial Calibration..... | 26 |
| 4.2.2 | Zero Calibration [ZER]..... | 26 |
| 4.2.3 | Span Calibration [SPAN]..... | 26 |
| 4.2.4 | Linearity Calibration [L #]..... | 27 |
| 4.2.5 | GEO Adjustment [GEO]..... | 27 |
| 4.2.6 | Calibration Test [TEST]..... | 28 |

| | | |
|-----------|---|-----------|
| 4.2.7 | <i>End Cal [End]</i> | 28 |
| 4.3 | SETUP MENU | 29 |
| 4.4 | READOUT MENU | 33 |
| 4.5 | UNIT MENU..... | 35 |
| 4.6 | GLP/GMP MENU | 35 |
| 4.7 | COMMUNICATION | 36 |
| 4.7.1 | <i>RS232 Menu</i> | 36 |
| 4.7.2 | <i>Print Menu</i> | 37 |
| 4.7.3 | <i>RS485 Configuration</i> | 40 |
| 4.7.4 | <i>Ethernet Configuration</i> | 40 |
| 4.7.5 | <i>Analog Configuration</i> | 40 |
| 4.8 | MEMORY | 40 |
| 4.8.1 | <i>Memory menu (for i-DT61PW model)</i> | 40 |
| 4.8.2 | <i>USB memory (for i-DT61XWE model)</i> | 40 |
| 4.8.3 | <i>Alibi Memory (for i-DT61XWE model)</i> | 41 |
| 4.9 | DISCRETE I/O (FOR I-DT61XWE MODEL) | 46 |
| 4.9.1 | <i>I/O Type</i> | 46 |
| 4.9.2 | <i>Input</i> | 47 |
| 4.9.3 | <i>Output</i> | 47 |
| 4.10 | LOCK KEY CONFIGURATION | 47 |
| 4.11 | LIBRARY (FOR I-DT61XWE MODEL)..... | 48 |
| 4.12 | USER (FOR I-DT61XWE MODEL)..... | 50 |
| 4.13 | USB (FOR I-DT61XWE MODEL) | 52 |
| 4.13.1 | <i>USB Flash Drive</i> | 52 |
| 4.13.2 | <i>RFID</i> | 52 |
| 4.13.3 | <i>Barcode</i> | 52 |
| 4.13.4 | <i>Keyboard</i> | 52 |
| 4.13.5 | <i>Wi-Fi/Bluetooth Dongle (Optional)</i> | 53 |
| 5. | LEGAL FOR TRADE | 55 |
| 5.1 | SETTINGS..... | 55 |
| 5.2 | VERIFICATION | 55 |
| 5.3 | SEALING | 55 |
| 6. | MAINTENANCE | 57 |
| 6.1 | CLEANING..... | 57 |
| 6.1.1 | <i>Cleaning for i-DT61PW Model</i> | 57 |
| 6.1.2 | <i>Cleaning for i-DT61XWE Model</i> | 57 |
| 6.2 | TROUBLESHOOTING | 58 |
| 6.3 | SERVICE INFORMATION | 58 |
| 7. | TECHNICAL DATA | 59 |
| 7.1 | SPECIFICATIONS | 59 |
| 7.2 | TABLE OF GEO VALUES..... | 61 |
| 8. | COMPLIANCE | 62 |
| 9. | APPENDICES | 64 |
| 9.1 | APPENDIX A..... | 64 |
| 9.2 | APPENDIX B..... | 66 |

1. INTRODUCTION

This manual contains installation, operation and maintenance instructions for i-DT61PW and i-DT61XWE indicators. Please read it completely before installation and operation.

1.1. Safety Precautions

Definition of Signal Warnings and Symbols

Safety notes are marked with signal words and warning symbols. These show safety issues and warnings. Ignoring the safety notes may lead to personal injury, damage to the instrument, malfunctions and false results.

WARNING For a hazardous situation with medium risk, possibly resulting in severe injuries or death if not avoided.

CAUTION For a hazardous situation with low risk, resulting in damage to the device or the property or in loss of data, or minor or medium injuries if not avoided.

ATTENTION For important information about the product. May lead to equipment damage if not avoided.

NOTE For useful information about the product.

Warning Symbols



General hazard



Explosion hazard



Electrical shock hazard

Safety Precautions



CAUTION: Read all safety warnings before installing, making connections, or servicing this equipment. Failure to comply with these warnings could result in personal injury and/or property damage. Retain all instructions for future reference.

- Before connecting power, verify that the AC adapter's input voltage range and plug type are compatible with the local AC mains power supply.
- Do not position the equipment such that it is difficult to reach the power connection.
- Only connect the power cord to a compatible grounded electrical outlet.
- Make sure that the power cord does not pose a potential obstacle or tripping hazard.
- Operate the equipment only under ambient conditions specified in these instructions.
- The equipment is for indoor use only.
- Do not operate the equipment in hazardous or unstable environments.
- Do not place the equipment upside down on the platform.
- Use only approved accessories and peripherals.
- Disconnect the equipment from the power supply when cleaning.
- Service should only be performed by authorized personnel.



WARNING: Never work in an environment subject to explosion hazards! The housing of the instrument is not gas tight. (Explosion hazard due to spark formation, corrosion caused by the ingress of gases).



WARNING: Electrical shock hazards exist within the housing. The housing should only be opened by authorized and qualified personnel. Remove all power connections to the unit before opening.

1.2. Intended Use

This instrument is intended for use in light industry. It must only be used for measuring the parameters described in these operating instructions. Any other type of use and operation beyond the limits of technical specifications, without written consent from OHAUS, is considered as not intended. This instrument complies with current industry standards and the recognized safety regulations; however, it can constitute a hazard in use. If the instrument is not used according to these operating instructions, the intended protection provided by the instrument may be impaired.

1.3. Overview of Parts and Controls

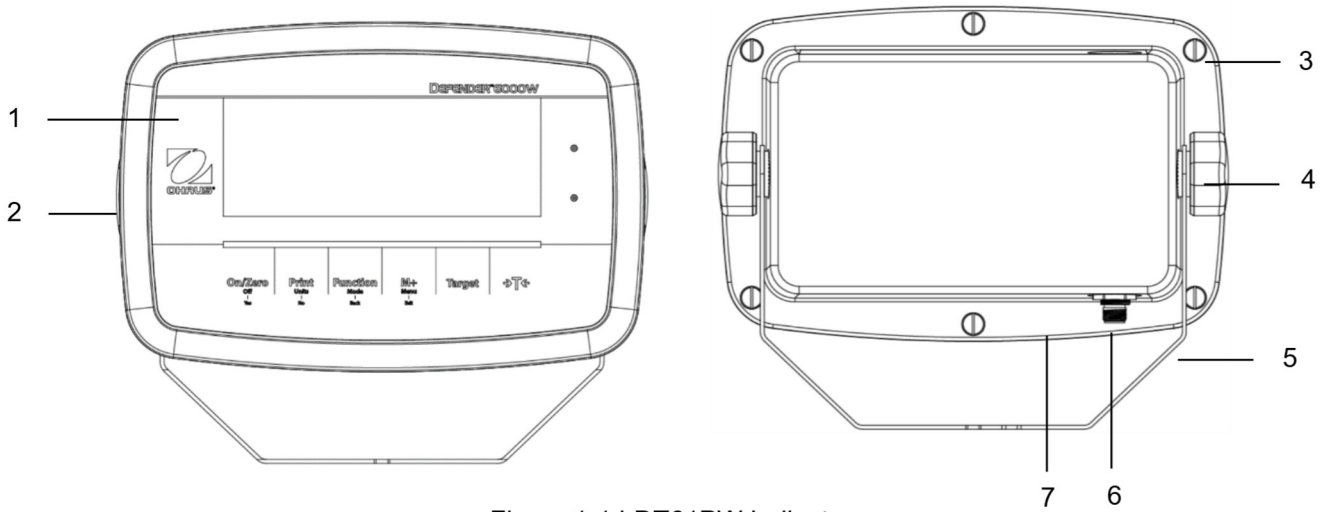


Figure 1-1 i-DT61PW Indicator

| Item | Description |
|------|---------------------|
| 1 | Control Panel |
| 2 | Front Housing |
| 3 | Screws (6) |
| 4 | Adjusting Knobs (2) |
| 5 | Mounting Bracket |
| 6 | Load Cell Connector |
| 7 | Rear Housing |


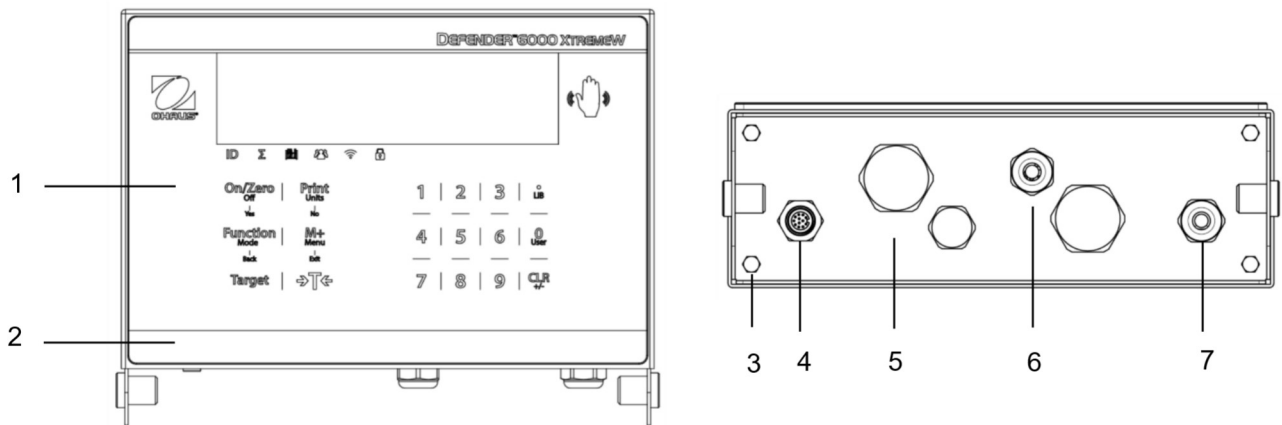


Figure 1-2 i-DT61XWE Indicator

| Item | Description |
|------|--------------------------|
| 1 | Control Panel |
| 2 | Front Housing |
| 3 | Screws (4) |
| 4 | Load Cell Connector |
| 5 | Bottom Housing |
| 6 | Strain Relief for Option |
| 7 | Power cord |

1.4. Mainboard

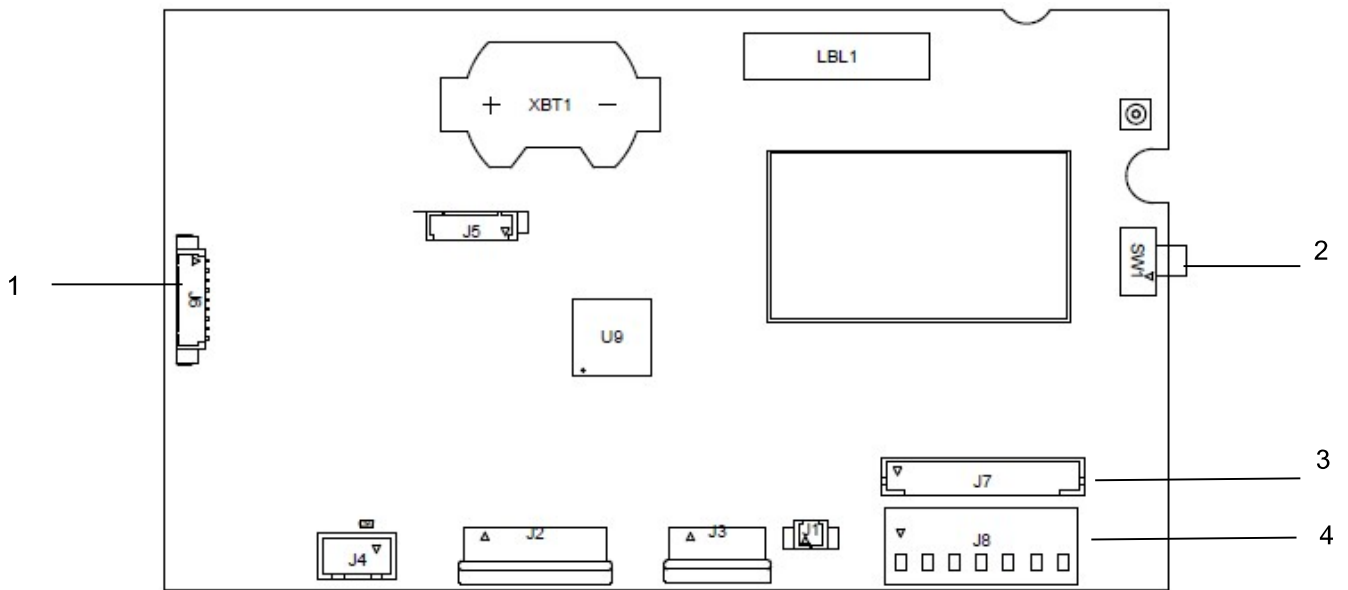

.Figure 1-3 i-DT61PW Mainboard

| Item | Description |
|------|---------------------------------|
| 1 | IR Communication connector (J6) |
| 2 | Security Switch (SW1) |
| 3 | Load Cell connector (J7) |
| 4 | Load Cell Terminal Block (J8) |

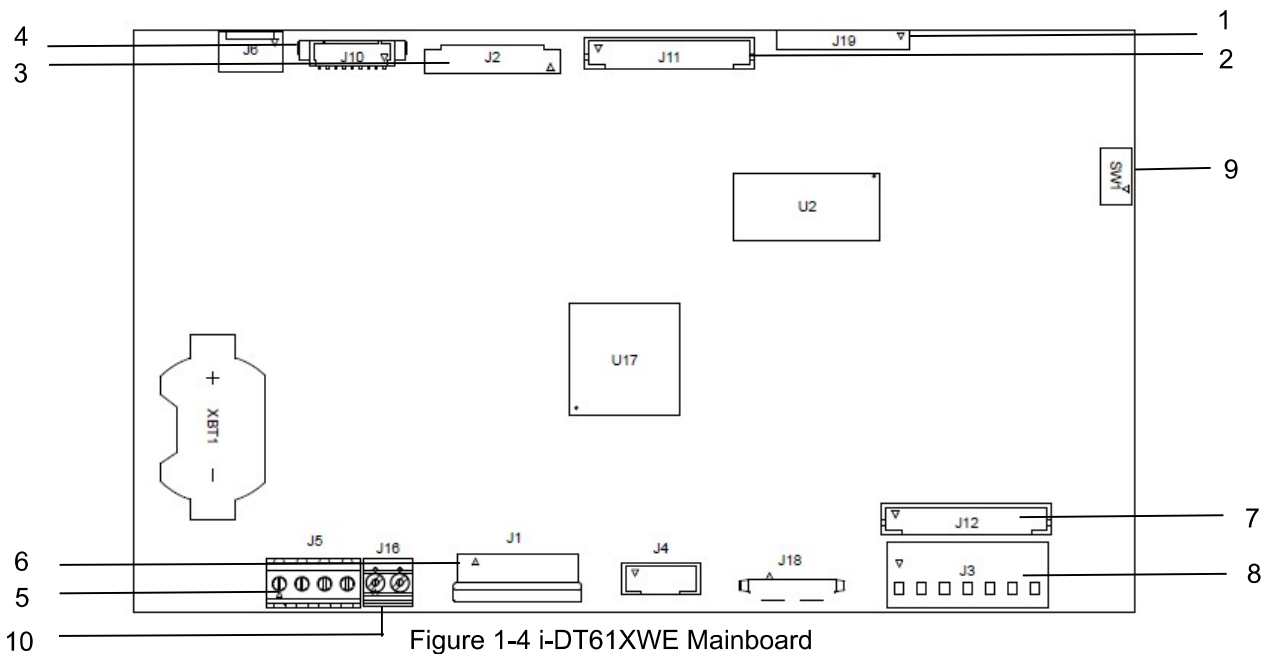


Figure 1-4 i-DT61XWE Mainboard

| Item | Description | Item | Description |
|------|--|------|---------------------------------|
| 1 | Alibi Memory Board connector (J19) | 6 | Keyboard connector (J1) |
| 2 | Discrete I/O/Analog/RS232-RS485-USB Device connector (J11) | 7 | Load Cell connector (J12) |
| 3 | Display Board connector (J2) | 8 | Load Cell Terminal Block (J3) |
| 4 | Ethernet connector (J10) | 9 | Security Switch connector (SW1) |
| 5 | RS232 connector (J5) | 10 | Discrete Input0 (J16) |

1.5. Control Functions



i-DT61PW Control Panel



i-DT61XWE Control Panel

| Button | On/Zero Off Yes | Print Units No | Function Mode Back | M+ Menu Exit | Target | →T← |
|---------------------------------|---|--|--|--|--|---|
| Primary Function (Short Press) | On/Zero If the terminal is Off, press to power on; If the terminal is On, press to set the zero point. | Print Sends the current value to the selected COM ports if AUTOPRINT is disabled. | Function Initiates an application mode. | M+ Accumulates the weight or displays the accumulated information with no load on the pan. | Target Sets under/over limit value for Check. | Tare Enters/clears a tare value; When the accumulation data is displayed, press to clear them. |
| Secondary Function (Long Press) | Off If the terminal is On, press to power off. | Units Changes the weighing unit. | Mode Allows changing the application mode. | Menu Enters the user menu. | Target Shows under/over limit value for Check. | Tare Displays the tare weight. |
| Menu Function (Short Press) | Yes Accepts the current setting on the display. | No Advances to the next menu or menu item. Rejects the current setting on the display and advances to the next available one. | Back Moves back to the previous menu item. | Exit Exits the user menu. Aborts the calibration in progress. | | |

Notes:

- Short Press: press less than 1 second.
- Long Press: press and hold for more than 2 seconds.

Numeric keyboard (i-DT61XWE)

| | | | | |
|---------------------------------|--------------------------------------|---|--|---|
| | | | | |
| Primary Function (Short Press) | 1-9 Enters numeric values. | . Enters decimal point (.). | 0 Enters numeric values 0. | CLR Clears the entered value. Clears an existing APW. When the accumulation data is displayed, press to clear them. |
| Secondary Function (Long Press) | | LIB Searches library items with numeric keys. | User Searches users with numeric keys. | +/- Switches between positive and negative values. |

Note: for i-DT61XWE model, press the and button together for three seconds can lock all buttons. Perform the same procedure again to unlock all buttons. When all buttons are locked, the icon will be lighted.

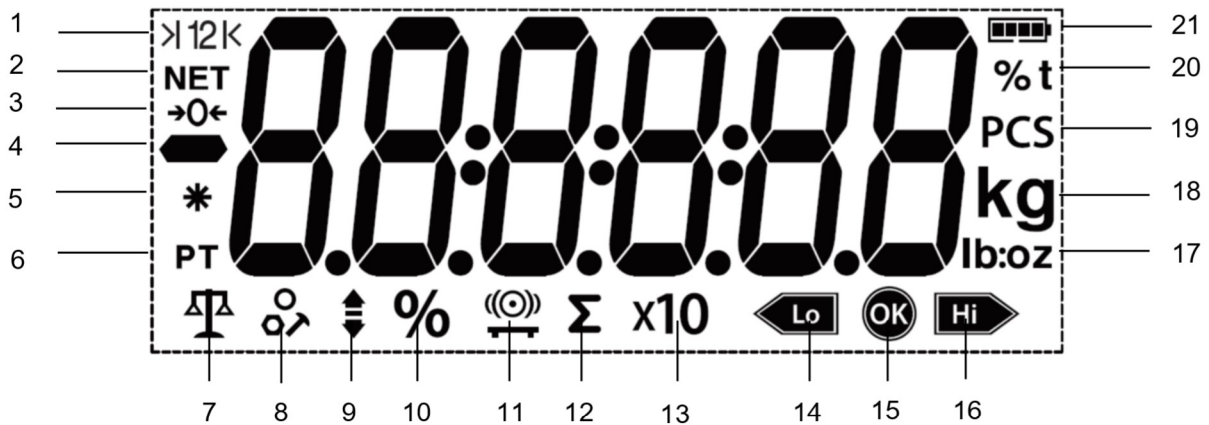


Figure 1-5 i-DT61PW Display

| Item | Description | Item | Description |
|------|---------------------------------|------|---|
| 1 | Range symbol (not used) | 12 | Accumulation symbol |
| 2 | NET symbol | 13 | Resolution extension symbol (not used) |
| 3 | Center of Zero symbol | 14 | Check weighing lower symbol |
| 4 | Negative symbol | 15 | Acceptable symbol |
| 5 | Stable weight symbol | 16 | Check weighing higher symbol |
| 6 | Preset Tare, Tare symbols | 17 | Pound, Ounce, Pound:Ounce symbols |
| 7 | Weighing mode symbol | 18 | Kilogram, gram symbols |
| 8 | Counting mode symbol | 19 | Pieces symbol |
| 9 | Check weighing mode symbol | 20 | Percent symbol, tonne symbol (not used) |
| 10 | Percentage weighing mode symbol | 21 | Battery symbol |
| 11 | Dynamic weighing mode symbol | | |

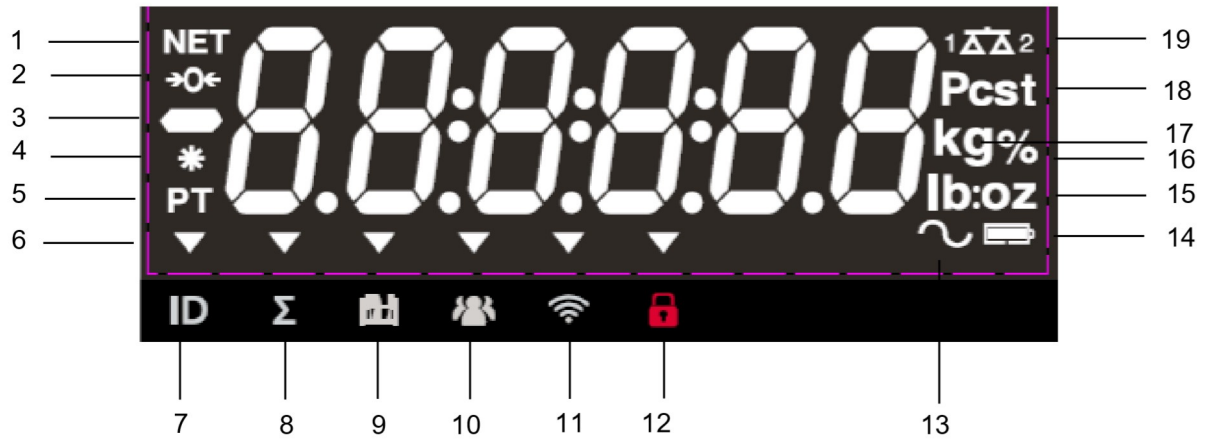


Figure 1-6 i-DT61XWE Display

| Item | Description | Item | Description |
|------|---------------------------|------|--|
| 1 | NET symbol | 11 | Wi-Fi symbol |
| 2 | Center of Zero symbol | 12 | Lock symbol |
| 3 | Negative symbol | 13 | Dynamic (tilde) symbol |
| 4 | Stable weight symbol | 14 | Battery symbol (not use) |
| 5 | Preset Tare, Tare symbols | 15 | Pound, Ounce, Pound:Ounce symbols |
| 6 | Pointer symbols | 16 | Percent symbol |
| 7 | ID symbol | 17 | Kilogram, gram symbols |
| 8 | Accumulation symbol | 18 | Pieces symbol, tonne symbol (not used) |
| 9 | Library symbol | 19 | Scale symbol (not used) |
| 10 | User symbol | | |