VACUO-TEMP

44 000 474

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

SOMMAIRE

INFORMATION GENERALE
SOMMAIRE
LISTE D'EMBALLAGE
ACCESSOIRES
DESCRIPTION DE L'APPAREIL
SPECIFICATIONS TECHNIQUES
INSTALLATION
MODE OPERATOIRE
ENTRETIEN
PIECES DE RECHANGE
GARANTIE
SCHEMA ELECTRIQUE
DECLARATION DE CONFORMITE "CE"

CONTENTS

GENERAL INFORMATION
CONTENTS
PACKING LISTS
ACCESSORIES
EQUIPMENT DESCRIPTION
TECHNICAL FEATURES
INSTALLATION
OPERATION
MAINTENANCE
SPARE PARTS
GUARANTEE
ELECTRICAL DIAGRAM
"EC" CONFORMITY DECLARATION

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PACKING LIST

The standard equipment consist of the following components:

Description	Description	Code / Code	Qté. / Qty.
Dessiccateur "Vacuo-temp"	"Vacuo-temp" vacuum dryer	44 000 474	1
Cloche en verre	Glass bell	44 000 475	1
Guide de l'utilisateur	Instruction manual.	80047	1

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ACCESSORIES

(They must be fitted in the factory)

Denomination	Code
TEMPERED GLASS BELL 15cm high and 23cm useful diameter.	44000475
SILICONE GASKET	44000476
VACUUM PUMP «VACUM» Rotary blade pump with built-in oil return. Motor with thermic protection device and starting up switch. Maximum working temperature: 60°C.	41000003

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EQUIPMENT DESCRIPTION

VACUO-TEMP thermostatic vacuum dryer is designed for drying at adjustable temperatures from 50°C up to 170°C with a temperature fluctuation of $\pm 2\%$. Electronic temperature regulation and digital reading (3 digits) with 1 digit resolution.

It includes Pt100 probe, shielded heating element, vacuum pump rear connection, vacuum valve and tempered glass bell with silicone gasket.

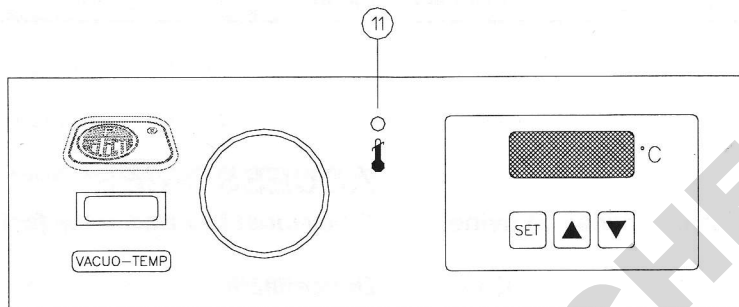
The upper plate is made of special aluminium alloy with surface grinding and sealing gasket. Outer casing is made in AISI 304 stainless steel.

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WORKING OF THE SAFETY THERMOSTAT

If a bad function of the regulation produce an abnormal increase of the temperature, the safety thermostat starts working blocking the heating of the machine, and the indicator lamp (11) of the frontal panel lights up. To reset the thermostat, stop the machine, leave it to cool and start up again. In case that the failure continues, call the authorized technical service.

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TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

Code / Code		44 000 474
Vide maximal / Maximum vacuum		10 ⁻² mm Hg.
Capacité litres / Capacity litres		--
Volume utile litres / Chamber volume litres		3
Dim. intérieures cm. Internal dimensions cm	Ø	23,5
	Prof. / Depth	--
Dimensions extérieures cm. Overall dimensions cm	Haut. / Height	17
	Larg. / Length	28
	Prof. / Width	34
Dispositif pour guides plateau / Shelf runner device		--
Consommation W. / Consumption W		600
Poids Kg. / Weight Kg		9

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INSTALLATION

Put the oven on a flat, horizontal and level surface leaving a space of 10 cm at the back and on each side of the apparatus.

For maximum efficiency, do not place the oven in direct sunlight or other sources of heat.

**ATTENTION ! IMPORTANT POUR VOTRE SECURITE
CAUTION!!! IMPORTANT FOR YOUR SAFETY**

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Do not use the oven for drying or treatment of any products which can give off gases or make explosive or inflammable mixtures.

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Be sure that the voltage supply is the same as the one indicated on the characteristic plate of the apparatus.

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Do not use the apparatus if it is NOT earthed.

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If you have to change the plug bear in mind the following:

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*Blue cable: Neutral.
Brown cable: Phase.
Yellow/green cable: Earth.*

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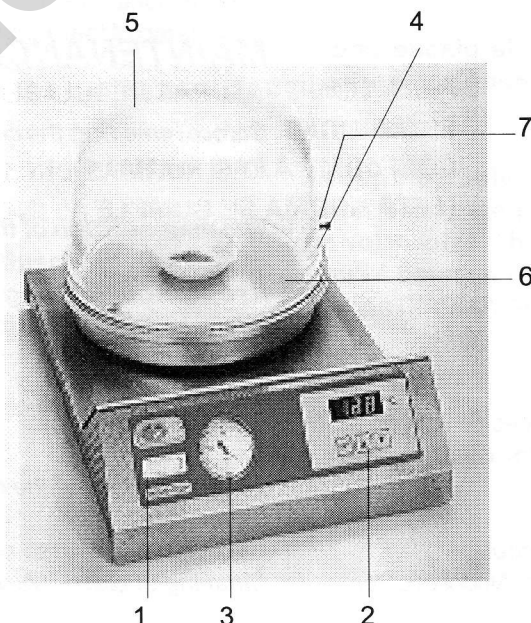
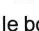
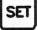
1. Main switch.
2. Temperature control thermostat.
3. Vacuum gauge.
4. Ventilation valve.
5. Tempered glass bell.
6. Hotplate.
7. Nipple.

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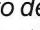
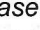
STARTING UP:

1. Press the main switch (1).
2. Select the working temperature by simultaneously pressing buttons  and  to increase or buttons  and  to decrease.
To check the selected temperature, press button  and it will appear on the display.
3. Check that the gasket is adjusted and place the glass bell.
4. Open the valve (4) and connect the vacuum pump in nipple (7).
5. Do the vacuum, when the vacuum gauge shows the desired vacuum, close the valve (4).
6. Disconnect the vacuum pump from the nipple (7).
7. When necessary dry the glass bell, open the valve (4).

Nota: If the material inside the glass bell can produce gases that can damage the membrane of the vacuum pump, it is recommended to place a water jet pump between the vacuum pump and the Vacuo-temp.

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MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

The manipulation of the internal electronic circuits of the oven by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA authorized technical services.

CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol
Cleaning of plastic: Alcohol with cotton duster.

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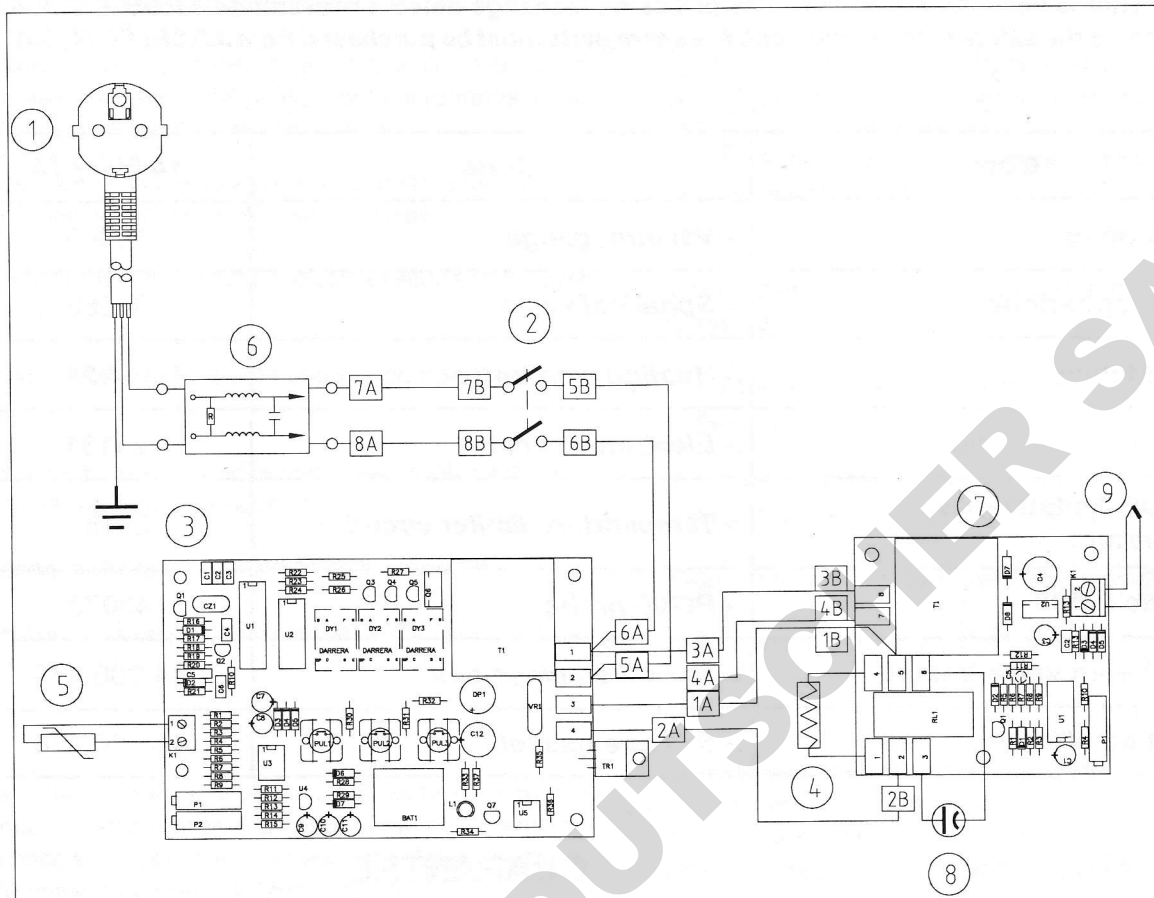
GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use of causes beyond the control of J.P. SELECTA,S.A..

Any manipulation of the apparatus by unauthorized personnel by J.P. SELECTA,S.A. cancels the aguarantee automatically.

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To guarantee the safety of the equipment, the spare parts must be purchased from J.P.SELECTA,S.A.
Agent en France de J.P. SELECTA S.A