PLAQUE CIRCULAIRE	1000718
PLAQUE CIRCULAIRE	1000442
PLAQUE CIRCULAIRE	1000443
SELECTEMP	1000444
PLACFIN	7000497
UNIPLAC	6002312
UNIPLAC	6002314
UNIPLAC	6002316

**PLAQUES CHAUFFANTES**

**HOTPLATES**

**INFORMATION GENERALE**

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**GENERAL INFORMATION**

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

**SOMMAIRE**

INFORMATION GENERALE
SOMMAIRE
LISTE D'EMBALLAGE
ACCESSOIRES
DESCRIPTION DE L'APPAREIL
SPECIFICATIONS TECHNIQUES
INSTALLATION
MODE OPERATOIRE
ENTRETIEN
PIECES DE RECHANGE
SCHEMA ELECTRIQUE
GARANTIE
DECLARATION DE CONFORMITE "CE"

**CONTENTS**

GENERAL INFORMATION
CONTENTS
PACKING LISTS
ACCESSORIES
EQUIPMENT DESCRIPTION
TECHNICAL FEATURES
INSTALLATION
OPERATION
MAINTENANCE
SPARE PARTS
ELECTRICAL DIAGRAM
GUARANTEE
"EC" CONFORMITY DECLARATION

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The standard equipment consist of the following components:

Description	Description	Code / Code	Qté/ Qty.
Plaque circulaire	Circular hotplate	1000718	1
Plaque circulaire	Circular hotplate	1000442	1
Plaque circulaire	Circular hotplate	1000443	1
Plaque circulaire "Selectemp"	Circular hotplate "Selectemp"	1000444	1
Plaque rectangulaire "Placfin"	Rectangular hotplate "Placfin"	7000497	1
Plaque circulaire "Uniplac"	Circular hotplate units "Uniplac"	6002312	1
Tiges verticales Ø12x60	Vertical rods Ø12x60	12103	2
Tige transversale Ø12x45	Horizontal rods Ø12x45	12105	1
Noix Trevin	Bosshead for rods	7008133	2
Clé allen	Head key	9100005	1
Plaque circulaire "Uniplac"	Circular hotplate units "Uniplac"	6002314	1
Tiges verticales Ø12x60	Vertical rods Ø12x60	12103	2
Tige transversale Ø12x77	Horizontal rods Ø12x77	12101	1
Noix Trevin	Bosshead for rods	7008133	2
Clé allen	Head key	9100005	1
Plaque circulaire "Uniplac"	Circular hotplate units "Uniplac"	6002316	1
Tiges verticales Ø12x60	Vertical rods Ø12x60	12103	2
Tige transversale Ø12x105	Horizontal rods Ø12x105	12107	1
Noix Trevin	Bosshead for rods	7008133	2
Clé allen	Head key	9100005	1
Guide de l'utilisateur	Instruction manual	80061	1

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## EQUIPMENT DESCRIPTION

The hotplates are designed to work with a big temperature range, up to 130°C (Placfin), 250°C (Selectemp) up to 400°C (Circular hotplates and Uniplace).

Temperature regulation is done by a hydraulic thermostat (Selectemp and Placfin) or impulse regulators (Uniplace) independent for each plate.

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## TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

		Plaques circulaires Circular hotplates			Plaque circulaire Circular hotplate SELECTEMP	Plaque rectangula- ire Rectangul- ar hotplate PLACFIN	Plaques circulaires Circular hotplate units UNIPLAC		
Code / Code		1 000 718	1 000 442	1 000 443	1000444	7000497	6 002 312	6 002 314	6 002 316
N° de places / N° of plates		1	1	1	1	1	2	4	6
Dim. utiles de la plaque cm. / Plate dimensions cm	Larg. Length	--	--	--	--	23	--	--	--
	Prof. Width	--	--	--	--	22	--	--	--
Ø Plaque en cm. / Ø Plate cm		12	15	19	15	--	12		
Dim. extérieures en cm Overall dimensions cm	Haut. Height	8			6	8	23		
	Larg. Length	20	22	26	30	22	47	77	105
	Prof. Width	20	22	26	22	30	27		
Consom. W. / Consumption W		1000	1500	2000	1500	400	1000	2000	3000
Température max. Maximum temperature		400°C			250°C	130°C	400°C		
Poids Kg / Weight Kg		2.1	3	4.5	3	3.7	8	15	22

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## INSTALLATION

Place the hotplate on a flat horizontal, level surface, leaving a space at the back and sides of the apparatus of about 10cm.

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*For a longer life of the apparatus provided with hotplate, it is recommended that its surface be covered as much as possible by the beaker to be heated. This avoids overheating and deformation of the plate by absorbing more the heat.*

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*Because of that, when using round-bottomed, Soxhlet or Kjeldahl flasks it is necessary to use the adapters in aluminium described in "HOTPLATES" section of our catalog.*

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*Be sure that the voltage supply is the same as the one indicated on the characteristics plate.*

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*Do not use the apparatus if it is not earthed.*

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*If you change the plug bear in mind the following:*

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
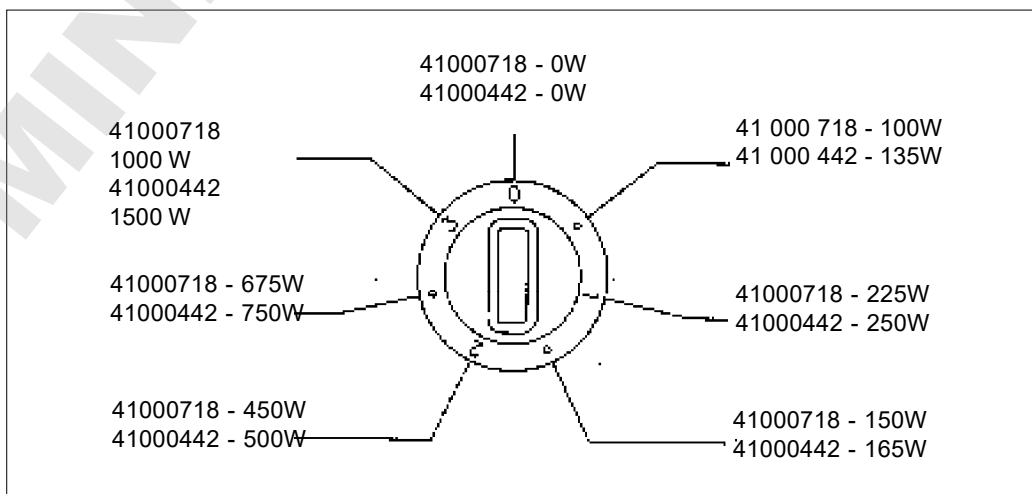
*These heaters include a selector of 6 heating powers settings and stop.*

*In the table below the power in watts is given according to the selector position.*

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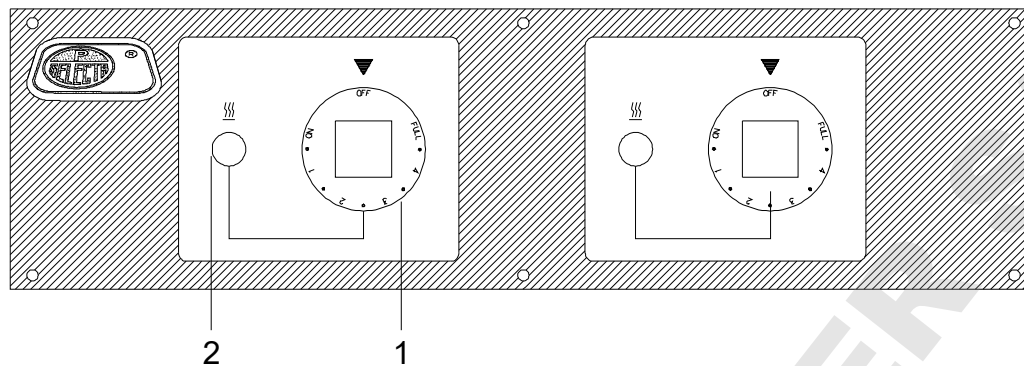## CIRCULAR HOTPLATE UNIPLAC UNITS CONTROL PANEL:

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### CIRCULAR HOTPLATE UNIPLAC UNITS STARTING UP:

1. Move the control knob of temperature (1) from the "OFF" position (disconnected), to the "FULL" position (maximum power).
2. It will be noted that according to the control position, the ON-OFF cycle switch of the indicator lamp for the working of the heater (2) will be longer or shorter, which will permit adjusting to the desired working temperature.
3. To stop, move the control button to "OFF" position.

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### CIRCULAR HOTPLATE SELECTEMP:

Circular hotplate "SELECTEMP" is provided with a temperature regulator thermostat whose sensitive part is in the center of the plate in order to register the temperature of the container to be heated.

Move the thermostat button numbered 1 to 12 to the right and fix it at one point, bearing in mind that the higher the point fixed, the higher the temperature. If after a reasonable time the desired temperature has not been reached, move the button of the thermostat to more or less, as desired.

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### IMPORTANT

Treat the central thermostat of the plate very carefully, do NOT try to scrape it with anything to clean it. Should any liquid spill onto the plate simply clean it with a damp cloth.

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## RECTANGULAR HOTPLATE PLACFIN:

Follow the same instructions as for "SELECTEMP" plate and read the temperature with the thermometer included in the apparatus.

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**WARNING!!! Caution with the hotplate surface, serious burns can be produced, even after switching off the apparatus.**

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## MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

The manipulation of the internal electronic circuits of the apparatus by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA authorized technical services.

### CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol

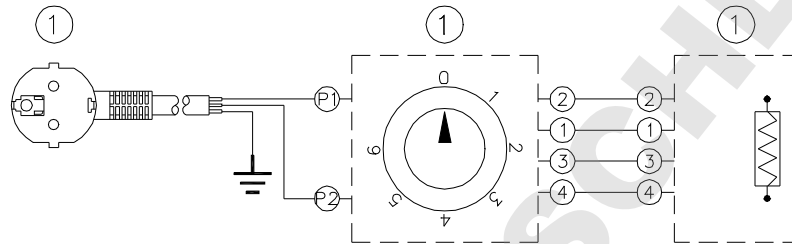
Cleaning of plastic: Alcohol with cotton duster.

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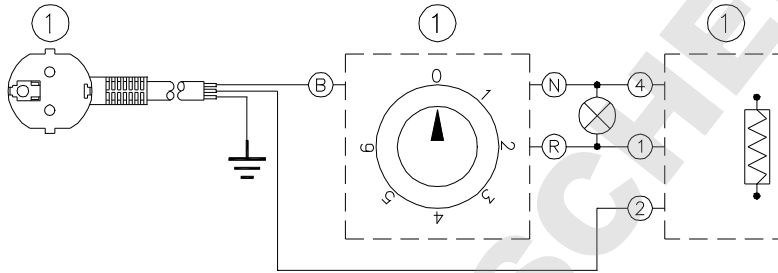
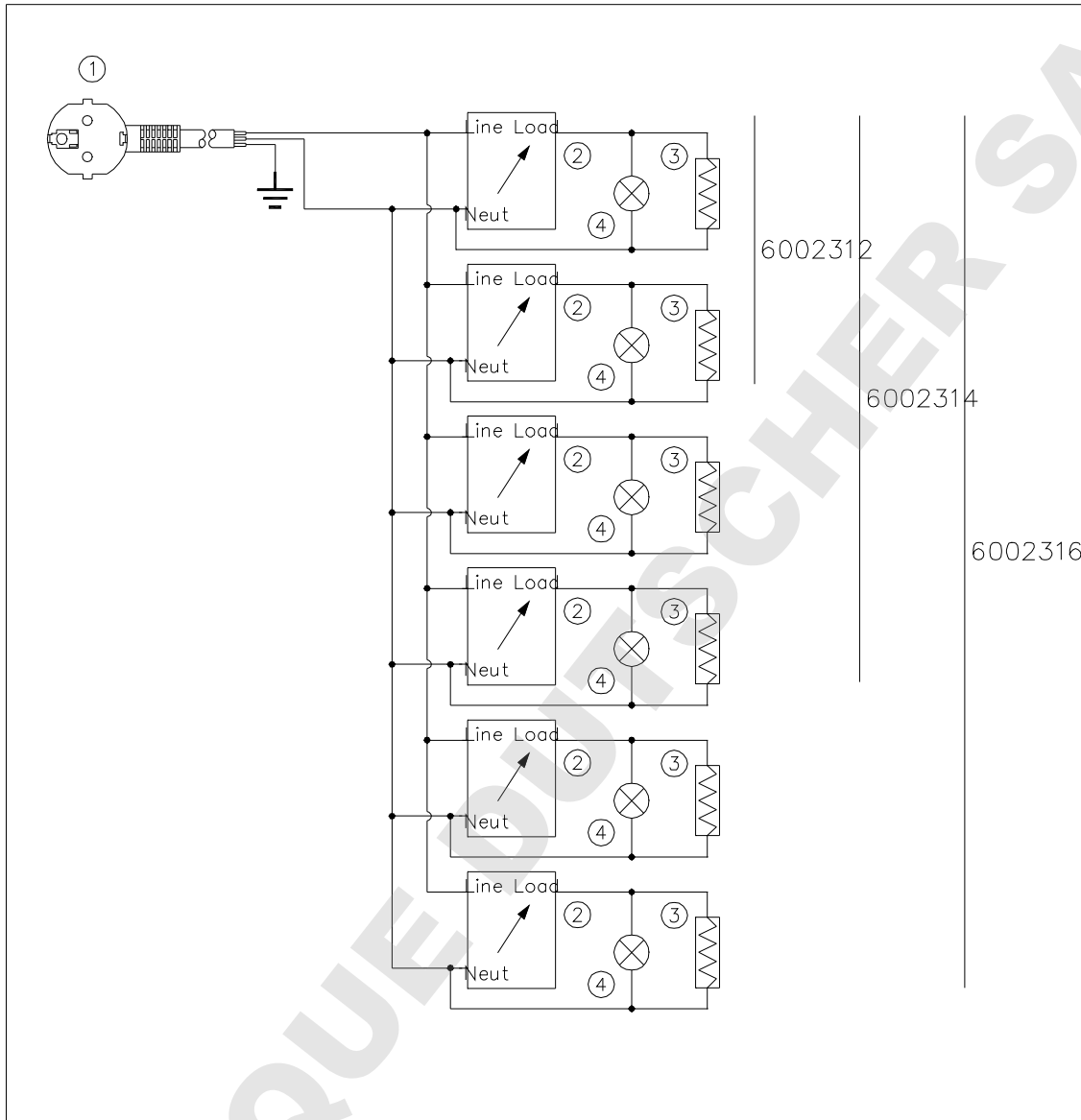
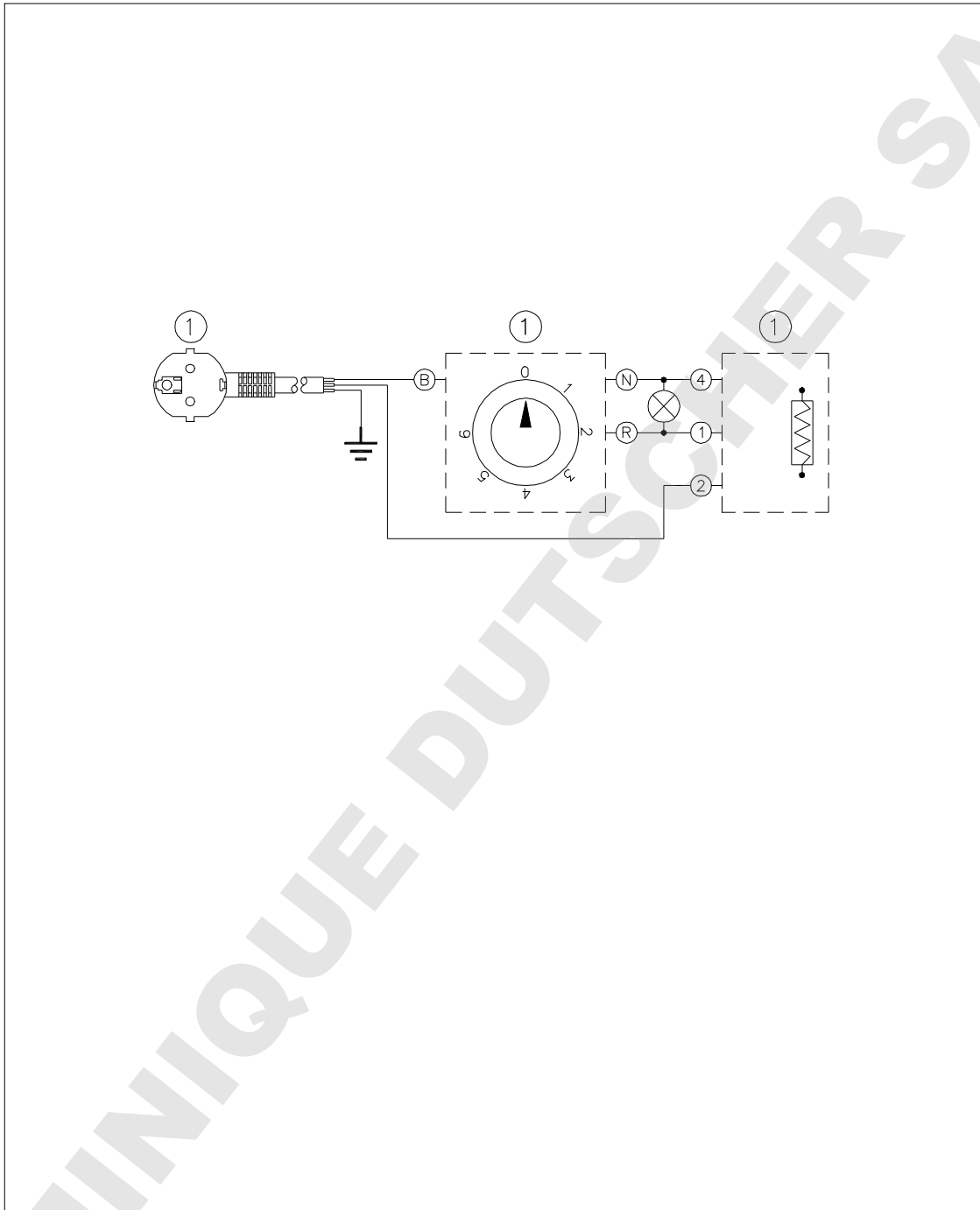
**GUARANTEE**

*This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use of causes beyond the control of J.P. SELECTA, S.A..*

*Any manipulation of the apparatus by unauthorized personnel by J.P. SELECTA, S.A. cancels the guarantee automatically.*

## DECLARATION DE CONFORMITE "CE" "EC" CONFORMITY DECLARATION

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La manipulation de l'appareil par des personnes non agréées par J.P. SELECTA S.A. annule automatiquement la garantie.