IC Bovino Quanti

IC Caprino Quanti

ZE/ICBQ

ZE/ICBQ100

ZE/ICCQ

ZE/ICCQ100

Immunochromatographic test for
quantification of milk species

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www.zeulab.com

SCOPE

IC Bovino Quanti kit is a rapid test to quantify presence of cow's milk in goat or sheep's milk. The test is an immunoassay based on the detection of bovine immunoglobulins (IgG) in milk samples. IRIS reader is required to obtain the quantitative result.

IC Caprino Quanti kit is a rapid test to quantify the presence of goat's milk in sheep's milk. The test is an immunoassay based on the detection of caprine immunoglobulins (IgG) in milk samples. IRIS reader is required to obtain the quantitative result.

Raw milk must be used with these kits.

| KIT COMPONENTS | ZE/ICBQ & ZE/ICCQ | ZE/ICBQ100 & ZE/ICCQ100 |
|----------------------------|------------------------------|------------------------------------|
| Strip tests | 25 | 100 |
| Dilution solution | 1 | 4 |
| Disposable Tips | 50 | 200 |
| Divisible strip of 8 wells | 4 | 13 |
| Well rack | - | 1 |

ADDITIONAL MATERIAL (NOT PROVIDED)

- ✓ Micropipette (20 - 200 μ L)

ASSAY PROCEDURE

1. Take a well from the divisible strip. Place it into the rack.
2. Apply 20 μ L of milk and 200 μ L of dilution solution. Stir gently.
3. Place the strip test into the well following the arrows direction.
4. Let the test progress for 10 minutes at room temperature.
5. Place the strip in the tray. The strip position in the tray is shown in the final scheme.
6. Read the results with IRIS in the following minute after the end of the assay.

IRIS READING

Turn on the IRIS reader.

- App mode:
 - Open the app in the smartphone, connect the devices and follow IRIS instructions.
 - Select “Read strip”.
 - Scan the QR code corresponding to batch of the kit. It appears in the product certificate.
- Key mode:
 - Load the method corresponding to the kit batch into IRIS following IRIS instructions.
 - Select the method “IC Bov Quanti” or “IC Cap Quanti” using the IRIS keyboard and then “Read strip”.

RESULTS

IRIS shows the percentage of milk mixture from the limit of detection (LoD) of the test (see table for each case, Bovino/Caprino) up to 2%.

When the percentage of milk is outside this range, <LoD (see table for each case) or > 2% are displayed.

An assessment of the result is also displayed:

- **Negative:** when the value is lower than the limit of detection.
- **Low positive:** when the sample contains milk mixture in a low percentage, lower than the limit of quantification (LoQ).
- **Positive:** when the result is higher than the limit of quantification (see table for each case).

| Quantification | | Colour | Interpretation |
|---|--|--------|----------------|
| IC Bovino Quanti LoD 0.1% LoQ 0.25% | IC Caprino Quanti LoD 0.25% LoQ 0.5% | | |
| <0.1% | <0.25% | green | Negative |
| 0.1% - 0.25% | 0.25% - 0.5% | orange | Low positive |
| 0.25% - 2% | 0.5% - 2% | red | Positive |
| >2% | >2% | red | Positive |

NOTES

- Use a new disposable tip for each sample.
- Milk samples with a pH below 6.3 have the viscosity altered and can affect the test result.
- The kit is designed to detect traces of milk mixture. As example, 0.125% of mixture implies that a tank of 10 000L is contaminated with 12.5 L of other milk specie.
- To obtain the maximum performance of the test it is necessary to use perfectly calibrated micropipettes.
- Results must be read with IRIS in the following minute after the end of the assay.
- If the blue line does not appear on the white central zone, the test has not been run properly or the reagents have been damaged. The result must be considered INVALID (INV) and repeated. If it happens again, please contact ZEULAB.

STABILITY

Store the test under dry conditions and temperatures between 15 and 25°C. The strips are very sensitive to humidity and must be kept inside the container provided with the desiccant bag. The kit has a shelf life of approximately 12 months. Check the expiration date on the package.

SAFETY

This kit should be used following good laboratory practices. A Safety Data Sheet is available on request.

IC Bovine Quanti and IC Caprine Quanti are in vitro diagnostic tests. In analysis implicating legal processes, the results should be re-evaluated with an official reference method. ZEULAB, S.L. does not assume any legal responsibility.

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