**ZEULAB**

# **PROTEON MILK EXPRESS SW**

**ZE/PR/L10SW**

**Immunocromatographic test for detection of milk in  
working surfaces**

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## SCOPE AND TEST PRINCIPLE

**PROTEON MILK EXPRESS SW** is an immunochromatographic test in strip format, to detect  $\beta$ -lactoglobulin ( $\beta$ -LG) in working surfaces.  $\beta$ -LG together with caseins,  $\alpha$ -lactoalbumin and seroalbumin are the main allergens in milk.

The test is based on the detection of  $\beta$ -LG, which reacts with a specific antibody coated to red particles attached to strip test. The appearance of two lines, red and blue, on the strip test is a positive result.

Limit of detection in food: 0.25 ppm  $\beta$ -LG (2.5 ppm milk proteins or 7.0 ppm powder milk MoniQA)\*

Limit of detection in surfaces: 0.7  $\mu$ g of milk proteins (2.0  $\mu$ g powder milk MoniQA)\*

\***References:** Bobe et al., 2007; Ng-Kwai-Hang, 1986; MoniQA reference material Certificate

## COMPONENTS

	ZE/PR/L10SW
<i><math>\beta</math>-LG STRIP TESTS</i>	10
<i>EXTRACTION BUFFER TUBE (EB) (0.5 ml)</i>	10
<i>SWABS</i>	10
<i>FRAME</i>	10
<i>PRODUCT CERTIFICATE</i>	YES

## SAFETY

This kit should be used following good laboratory practices. Safety clothing should be worn and skin contact with the reagents avoided. Do not ingest. A SAFETY DATA SHEET is available from your local distributor or ZEULAB.

## STABILITY AND STORAGE

Store the test under dry conditions and temperatures between 15 and 25°C. Check the expiration date on the package.

## NOTES

- The strips are very sensitive to moisture, keep them inside the supplied container and fully closed.
- Use a pipette, tube and swab for each sample.
- For manipulating the strip, hold it from the extreme opposite end of the arrows area.
- Swab a representative area of at least 100 cm<sup>2</sup>. The result of the assay is presence or absence of the allergen on the tested surface. It is not possible to extrapolate to ppm allergen values.
- The test limit of detection on surfaces is subjected to the use of the swabs provided by ZEULAB.
- Please, contact ZEULAB for further information.

## TEST PROCEDURE

1. Open an *EXTRACTION BUFFER* tube .
2. Introduce a swab into the tube and soak it in the *EXTRACTION BUFFER*. The swab must be held and manipulated by its opposite end in order to avoid cross contamination. Only ZEULAB swabs are recommended.
3. Rub the swab through all possible directions on the analysed surface. Rotate from time to time the swab's point in contact with the surface.
4. Place the swab in the initial tube with the *EXTRACTION BUFFER* and stir gently into the solution.
5. Cast aside the swab.
6. Introduce the strip test into the tube following the arrows' direction.

7. Wait for 10 min.

8. Read the results:

**NEGATIVE:** Only a blue line is observed in the white central zone of the strip test. This is the control line and it indicates that the test works properly.

**POSITIVE:** Two lines, BLUE and RED, are observed in the white central zone of the strip test.

When no blue line is shown, the result must be considered INVALID. This may be occur when the assay has not been performed properly, when the matrix interferes or when the reagents are damaged. Repeat the assay with a new strip.

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