Biosigma s.r.l.

a Dominique Dutscher Company

Via Valletta, 6 | 30010 Cantarana di Cona (VE), Italy | Tel. ++39 0426 302224 (r.a.) |

Fax ++39 0426 302228 | SMS ++39 348 4077376 |

E-mail info@biosigmaeu.com | http://www.biosigma.com | www.biosigma.it

M521D

20/11/2017

Pag. 1 di 3

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- *Centrifuge tubes made in transparent polypropylene suitable for use in laboratory and molecular biology.*
- *Two volumes 15 ml conical and 50 ml, conical or self standing, are available.*
- *Graduation clearly visible printed in black on the body of tube and imprinted graduation directly into the conical part of the tube.*
- *White writing area to better identify the tubes.*
- *Cap with an interior ring for a perfect seal.*
- *Available in zip bag or in plastic rack packed in zip bag*
- **Sterile SAL 10⁻⁶. DNase, RNase, Pyrogens FREE**
- *The recommended working temperature range for centrifuge tubes is 0 - 40°C. The suitability of these tubes for storage below 0°C depends on both the solution and the storage conditions. It is strongly recommended that a trial run be performed under actual conditions to test the suitability of the tubes for frozen storage*
- *Remove the closures before autoclaving the device. In the absence of this, in an autoclave the differences in pressure that occur can cause the collapse of the container. Make sure that the sterilization method adopted is suitable for all components of the Device (tubes, caps). If not, sterilize them separately each with the most appropriate method. Always keep in mind that sterilization could reduce the mechanical / functional / aesthetic properties of the Device.*

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DESTINATION USE:

Centrifuge tubes with screw caps Sterile ClearLine® are disposable labware intended to be used on laboratory to contain biological liquids for in vitro diagnostic analysis.

FOR IN VITRO USE ONLY

The use and the manipulation of the devices are reserved to the medical and technical staff with the necessary professional qualification.

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CLASSIFICATION: In order to the Directive 98/79/CE concerning medical-diagnostic in vitro devices receipt in Italy with D.L. n. 332 G.U.R.I. n. 269 dated 17th November 2000, Biosigma s.r.l., defines all the components of the above-state family as medical-diagnostic in vitro devices and classifies it as "Other medical-diagnostic in vitro devices", in order to make conformable the attachment III of the directive.

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MATERIALS: The materials utilized for the production of the in vitro labware devices are selected on the base of the properties indicated for the individual destination use. The Centrifuge tubes with screw caps Sterile ClearLine[®] are manufactured by Polypropylene (PP) and Polyethylene (PE).

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IN VITRO DIAGNOSTIC MEDICAL DEVICE CE for professional use only

- Use only for in vitro laboratory analysis,
- Use gloves: contamination risk if contents leaks,
- Keep out of flame or heat sources which might damage the product,
- Do not use after expiry date,
- Do not vary the intended purpose of the product,
- Keep out of reach of children,
- Disposal: use appropriate personal protective equipment and act according to applicable regulations.

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