Biosigma s.r.l.

a Dominique Dutscher Company

Via Valletta, 6 | 30010 Cantarana di Cona (VE), Italy | Tel. ++39 0426 302224 (r.a.) |

Fax ++39 0426 302228 | SMS ++39 348 4077376 |

E-mail info@biosigmaeu.com | http://www.biosigma.com |

M521D

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WATER SAMPLING BOTTLES

Manufactured by LP ITALIAN SPA - Distributed by Biosigma Srl

SQUARE BOTTLES IN VIRGIN PP

Ref	Capacity (ml)	Single pack	Sodium Thiosulfate
391920	500	X	X
391922	500	X	
391932	1000	X	X
391934	1000	X	

Transparency: translucent

Cap colour: natural white

Norms:

.- CE marking: __ N/A

.- food contact: raw material complies with CE regulation 1935/2004

Fields of application:

.- industry

.- microbiology

.- environmental sampling

Dedicated to optimise the sampling of water or food for microbiology analyses, bottles are designed and manufactured to assure functionality, safety and user-friendliness.

Sampling Bottles come sizes: 500 ml 1.000 ml.

Key features, common for all models:

- Manufactured with virgin raw material, in conformity of the environmental norms: PP
- Strong and virtually unbreakable.
- Lightweight, well balanced and easy to handle. Its square cross section and rounded edges assure the best transport and storage optimization, while minimizing their packaging volume.
- Type of opening : "wide neck"
- The volume of the inserted liquid can be easily controlled by means of the clear moulded graduation (see table); accuracy $\pm 2\%$.
- Closure: all bottles are closed by one leakproof screw cap with a special inert inner gasket and with a safety seal.
- Caps: all made in HDPE
- All closures are designed with vertical ribs for easy use and openness, even when wearing gloves
- Traceability : The product label on each box includes the lot number and expiration date. In addition, each bottle bears a pre-pasted label that allows the recording of the main sample identification data :
 - . - Batch number,
 - . - Expiry date
 - . - A unique number, in clear and in form of bar code.

All these allow full traceability of each individual sample.

• Sterilization: based on ISO 11137 - SAL 10^{-6} performed by means of ionizing radiation.

• Shelf life: see product tables.

• The whole manufacturing process, packaging included, is checked according applicable norms and LP's proprietary quality criteria.

WIDE NECK BOTTLES FEATURES

produced in: **PP**

Models with 50 mm nominal opening, facilitate the filling, minimize the risk of contamination during sampling (in accordance with the recommendations of the ISO 19458:2006 - sampling techniques):

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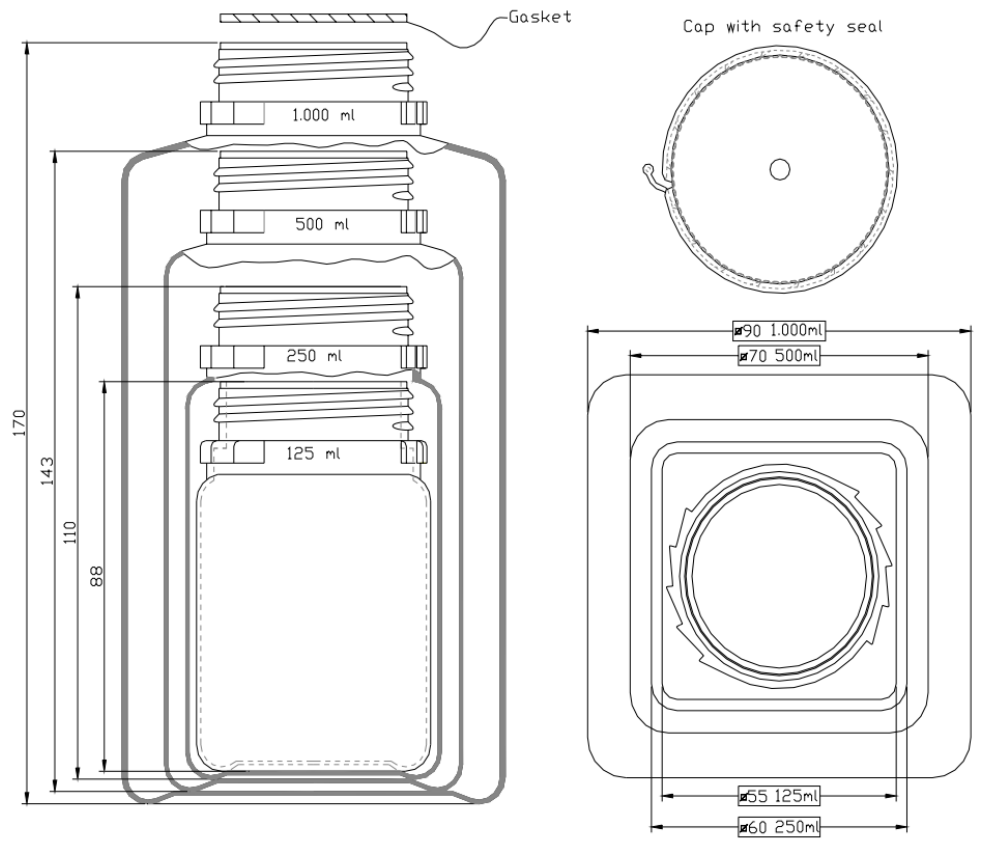
SODIUM THIOSULFATE BUFFERING

Chlorides: with chlorinated water samples, buffering of free or residual chlorine is suggested, to inhibit its bactericide action during transportation and storing. All bottles are available predosed with Sodium Thiosulfate (see "Suggestions") in conformity with:

- ISO 19458:2006 (same as French Standard NFT 90-40): 20 mg/l or

- International Standard ISO 5667-3: 80 mg/l

(Bottles filled to nominal capacity)

In case of highly chlorinated water, by customer request, bottles with higher quantity of Na Thiosulfate are available.

Note: in these cases the standard buffer amount is 100 mg/l, which can be increased up to 120 mg/l (typically used for swimming pool water sampling) with no influence on sample's quality (see "Suggestions").

Sodium Thiosulfate has no influence on the sample, that is why it is possible to use bottles containing Sodium Thiosulfate with non chlorinated samples.

Sterility: Water Sampling Bottles, with and without Na Thiosulfate, are all irradiated to a SAL (sterility assurance level) of 10^{-6} . The sterilization is guaranteed until seal is broken and the cap is opened (see table for shelf life).

Guarantee: screw caps have a tamper evident safety seal. If the cap seal is unbroken closure and

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inner sterility are guaranteed.

SUGGESTIONS FOR BEST USE

1. Microbiological analysis of water samples must always be performed within the shortest possible time after sampling. If the sample has aged too long, many factors can influence its bacterial contents, all of them related to the quality of the water: presence of toxic or nutritive substances for bacterial flora, saltiness, pH and so on.

As a general guideline, suggests to transport and store samples at a temperature of (+4 to +10)°C and to analyse them within 24 hours.

1) Buffering ratio

As a bactericide, sporicide, fungicide and viroicide, a sodium salt (hypochlorite NaClO and/or chlorite NaClO₂) is usually dissolved in water or, more frequently, one mixture of the two. It is not possible to know neither the composition of the mixture, nor the amount of each dissolved salt, therefore it is not possible to supply any indication of how much free chlorine is to be buffered by a predetermined amount of sodium thiosulfate (Na₂S₂O₃).

Furthermore, depending on inactivation dynamics, it is hard to predict which amount of Sodium Thiosulfate is required to inactivate even a known amount of free residual chlorine.

We suggest considering the following indications:

Buffering ratio between thiosulfate and hypochlorite → 1 Mole: 1 Mole

Buffering ratio between thiosulfate and Chlorite → 4 Moles: 1 Mole

In order to know the actual ratio in weight, such relationships must be related to the respective molecular weights, but quantitative indications are not set here because they could be not pertinent or misleading to each single case. We limit ourselves only to give evidence that in one limit case (all chlorite) the amount of Thiosulfate required is approximately quadruple than the one required for the opposite case ("all hypochlorite").

We suggest considering that 18 mg Sodium Thiosulfate will inactivate at least 2 mg/l and up to 5 mg/l of free chlorine residual, which is sufficient for the majority of samples.

DISPOSAL

All components of bottles (bottles, caps and gaskets) are manufactured using virgin raw materials that are completely recyclable and environment friendly; they can be disposed:

- Emptied and cleaned are fully recyclable; or
- Emptied, burned with energy recovery; or
- By incineration, at an appropriate temperature and with an adequate air insufflations. They give residuals of H₂O and CO₂, only; or
- Delivered to a controlled landfill, in accordance with local laws on waste disposal.