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Description: CLEARStab is an innovative system for the freezing and long-term preservation of bacterial and microbial cultures. The system is designed to eliminate the need for freeze-drying or frequent transfers, ensuring optimal preservation of the cultures. The system includes a cryobox containing 80 sterile vial and 1 vial picker. Each vial includes 20 ± 1 porous beads, immersed in 1 ± 0.1 ml of a glycerol-based cryoprotective solution. Sterile SAL 10^{-6} .

Purpose: CLEARStab ensures optimal preservation of bacteria, bacterial spores, and fungi at low temperatures, maintaining their viability and genetic stability. It is ideal for microbiology laboratories, research, and biotechnology production.

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Advantages:

- *Long-term preservation: Maintains cultures for extended periods without loss of viability.*
- *Ease of use: Reduces the need for frequent transfers of cultures.*
- *Versatility: Suitable for a wide range of organisms, including bacteria and fungi.*
- *Space efficiency: Allows for freezing numerous different strains in a compact area.*
- *Product options: Available in versions with cryoprotectant and dry versions without cryoprotectant.*
- *Color variants: CLEARStab offers vials with colored caps and beads in a range of colors—white, red, blue, yellow, and green—to facilitate the use of color coding for different species or strains of bacteria. All vials contain the same cryoprotectant, ensuring consistent and reliable performance. The variety of colors allows for quick and organized identification of different strains, thereby optimizing laboratory work efficiency.*

Materials:

- **Cryobox:** Container designed for low-temperature storage made of polycarbonate.
- **Criotubes:** Tubes made from medical-grade polypropylene with bi-injected polyethylene/rubber caps.
- **Beads:** Porous ceramic
- **Cryoprotectant:** Glycerol-based solution

Features:

- *Storage: Before use, CLEARStab can be stored at room temperature or, preferably, at +4°C for optimal product preservation.*
- *Storage Conditions for Microbiological Cultures: Microorganisms can be stored at different temperatures depending on their specific biological nature. For short periods, storage at -20°C is recommended. For long-term storage, the ideal temperature is -80°C, while for prolonged conservation, the gaseous phase of liquid nitrogen can be used. The selection of the appropriate temperature depends on the biological characteristics of the microorganisms and the operating procedures implemented by the laboratories to ensure maximum viability and integrity of the cultures.*
- *Sterility testing: Incubation for 3 days at 30°C and for 5 days at 25°C showed no growth.*
- *Validity: 24 months.*

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Prodotto Product type	Codice Code	Descrizione Description	Imballaggio Packaging
CLEARStab 1D con codice a barre CLEARStab 1D barcoded	CL1D80/AZ	Tubo 2ml biglie trattate con crioprotettore, Blu <i>2ml tube with treated beads in cryoprotectant, Blue</i>	80
	CL1D80/G	Tubo 2ml biglie trattate con crioprotettore, Giallo <i>2ml tube with treated beads in cryoprotectant, Yellow</i>	80
	CL1D80/RO	Tubo 2ml biglie trattate con crioprotettore, Rosso <i>2ml tube with treated beads in cryoprotectant, Red</i>	80
	CL1D80/V	Tubo 2ml biglie trattate con crioprotettore, Verde <i>2ml tube with treated beads in cryoprotectant, Green</i>	80
	CL1D80/BI	Tubo 2ml biglie trattate con crioprotettore, Bianco <i>2ml tube with treated beads in cryoprotectant, White</i>	80
	CL1D80/MIX	Tubo 2ml biglie trattate con crioprotettore, Assortito <i>2ml tube with treated beads in cryoprotectant, Mixed</i>	80
CLEARStab 2D con codice a barre CLEARStab 2D barcoded	CL2D80/AZ	Tubo 2ml biglie trattate con crioprotettore, Blu <i>2ml tube with treated beads in cryoprotectant, Blue</i>	80
	CL2D80/G	Tubo 2ml biglie trattate con crioprotettore, Giallo <i>2ml tube with treated beads in cryoprotectant, Yellow</i>	80
	CL2D80/RO	Tubo 2ml biglie trattate con crioprotettore, Rosso <i>2ml tube with treated beads in cryoprotectant, Red</i>	80
	CL2D80/V	Tubo 2ml biglie trattate con crioprotettore, Verde <i>2ml tube with treated beads in cryoprotectant, Green</i>	80
	CL2D80/BI	Tubo 2ml biglie trattate con crioprotettore, Bianco <i>2ml tube with treated beads in cryoprotectant, White</i>	80
	CL2D80/MIX	Tubo 2ml biglie trattate con crioprotettore, Assortito <i>2ml tube with treated beads in cryoprotectant, Mixed</i>	80
CLEARStab secco 1D con codice a barre CLEARStab dry 1D barcoded	CLDR1D80/AZ	Tubo 2ml biglie trattate senza crioprotettore, Blu <i>2ml tube with treated beads, No cryoprotectant, Blue</i>	80
	CLDR1D80/G	Tubo 2ml biglie trattate senza crioprotettore, Giallo <i>2ml tube with treated beads, No cryoprotectant, Yellow</i>	80
	CLDR1D80/RO	Tubo 2ml biglie trattate senza crioprotettore, Rosso <i>2ml tube with treated beads, No cryoprotectant, Red</i>	80
	CLDR1D80/V	Tubo 2ml biglie trattate senza crioprotettore, Verde <i>2ml tube with treated beads, No cryoprotectant, Green</i>	80
	CLDR1D80/BI	Tubo 2ml biglie trattate senza crioprotettore, Bianco <i>2ml tube with treated beads, No cryoprotectant, White</i>	80
	CLDR1D80/MIX	Tubo 2ml biglie trattate senza crioprotettore, Assortito <i>2ml tube with treated beads, No cryoprotectant, Mixed</i>	80
CLEARStab secco 2D con codice a barre CLEARStab dry 2D barcoded	CLDR2D80/AZ	Tubo 2ml biglie trattate senza crioprotettore, Blu <i>2ml tube with treated beads, No cryoprotectant, Blue</i>	80
	CLDR2D80/G	Tubo 2ml biglie trattate senza crioprotettore, Giallo <i>2ml tube with treated beads, No cryoprotectant, Yellow</i>	80
	CLDR2D80/RO	Tubo 2ml biglie trattate senza crioprotettore, Rosso <i>2ml tube with treated beads, No cryoprotectant, Red</i>	80
	CLDR2D80/V	Tubo 2ml biglie trattate senza crioprotettore, Verde <i>2ml tube with treated beads, No cryoprotectant, Green</i>	80
	CLDR2D80/BI	Tubo 2ml biglie trattate senza crioprotettore, Bianco <i>2ml tube with treated beads, No cryoprotectant, White</i>	80
	CLDR2D80/MIX	Tubo 2ml biglie trattate senza crioprotettore, Assortito <i>2ml tube with treated beads, No cryoprotectant, Mixed</i>	80
CRYOGEN® Box81	BSM58061B1D/5AZ	Box 1D, 81 posti per provette 1-2 ml, griglia azzurra <i>1D Box, 81 slots for 1-2 ml tubes, blue grid</i>	5
	BSM580612D/5AZ	Box 2D, 81 posti per provette 1-2 ml, griglia azzurra <i>2D Box, 81 slots for 1-2 ml tubes, blue grid</i>	5