

Product Information

AccuBlue™ Broad Range dsDNA Quantitation Solution

Catalog Number: 31009-T, 31009

Unit Size:

31009-T: 200 assays (200 uL microplate assay)

31009: 1000 assays (200 uL microplate assay)

Kit Contents

Component	31009-T	31009
AccuBlue Broad Range Buffer	50 mL 99975	250 mL 99973
AccuBlue Broad Range Dye (100X in DMSO)	1 mL 99974	3 X 1mL 99974
AccuBlue Broad Range Enhancer (100X in water)	1 mL 99943	3 X 1 mL 99943

Storage and Handling

Store kit at 4°C. Protect quantitation solution from light. The kit is stable for at least 6 months from date of receipt when stored as recommended.

Spectral Properties

Ex/Em 350/460nm (in the presence of dsDNA). See Figure 1 for spectra.

Product Description

The AccuBlue Broad Range Quantitation Solution is designed for quantitation of dsDNA in the range of 2 to 2000 ng in a 200 uL assay volume. Unlike absorbance-based measurements, the assay is selective for dsDNA over ssDNA or RNA (Figure 2). The quantitation assay offers the advantages of a wide dynamic range and high sensitivity over other traditional methods of DNA quantitation. The kit contains the AccuBlue Broad Range dsDNA Quantitation Buffer, Dye solution, Enhancer solution, and pre-diluted dsDNA Standards. The assay kit is tolerable to common contaminants such as proteins, salts, organic solvents and detergents. See Table 1 for more information. AccuBlue Broad Range dsDNA Quantitation Solution also is available in a kit with pre-diluted calf thymus dsDNA standards (cat. no. 31007).

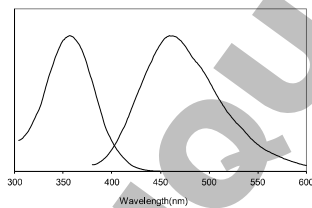


Figure 1: Excitation and emission spectra for AccuBlue Broad Range dsDNA quantitation reagent in the presence of excess dsDNA.

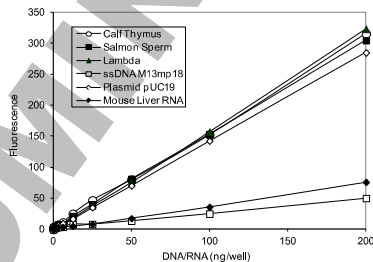


Figure 2. Relative fluorescence intensities of different nucleic acids using the AccuBlue Broad Range dsDNA Quantitation kit.

Assay Protocol

1. Use properly calibrated pipettes and DNase-free pipette tips, tubes and plates for best accuracy. It is recommended to test each DNA standard and each unknown sample in triplicate. If more than one 96 well plate is to be tested in a single assay, it is recommended to include a standard curve on each plate to minimize variability between plates.
2. Prepare DNA standards in 1X TE buffer using the dsDNA of your choice. Prepare a 200 ng/uL stock solution of DNA. Determine the DNA concentration on the basis of absorbance at 260 nm in a cuvette with a 1 cm path length. An A_{260} of 1 corresponds to a 50 ng/uL dsDNA solution. Because most spectrophotometers saturate around an absorbance of 2, dilute the 200 ng/uL stock 1:10 and OD the diluted stock solution. An A_{260} of 0.4 corresponds to a concentration of 20 ng/uL. Factor in the dilution to obtain the concentration of the original stock solution. Perform several 2-fold dilutions of the 20 ng/uL DNA solution to obtain standards ranging between 20 ng/uL and 2 ng/uL. Use 1X TE as the zero DNA standard. For long term storage, we recommend adding sodium azide to a final concentration of 2 mM and storing DNA standards at 4°C.
3. Remove the DNA quantitation kit from storage and warm all components to room temperature before use. AccuBlue Broad Range dye is provided in DMSO, which may freeze during storage at 4°C. You can warm up all kit components in a 37°C water bath; be sure to allow solutions to equilibrate to room temperature before using. Before removing the required volume, mix each component well by shaking or vortexing, and centrifuge vials briefly before opening to minimize reagent loss on the cap.
4. Prepare 200 uL of working solution for each sample to be tested. Working solution should be used within an hour after preparation for best results, but it can be stored and used up to 24 hours later, with only minor loss of accuracy. During storage dye precipitation may occur, but can be resuspended by vortexing. It is recommended to test each DNA standard and each unknown sample in triplicate. If more than one 96 well plate is to be tested in a single assay, it is recommended to include a standard curve on each plate to minimize variability between plates. Combine reagents as shown in Table 2 below and mix well by vortexing or shaking.

Table 1: Preparation of 20 mL AccuBlue Broad Range working solution (scale volumes to prepare only amount required for assay).

Component	Volume
AccuBlue Broad Range Buffer	20 mL
AccuBlue Broad Range Dye (100X)	200 uL
AccuBlue Broad Range Enhancer (100X)	200 uL

5. For each sample, pipette 200 uL of the AccuBlue working solution per well of a black 96-well microplate. To test samples in triplicate, prepare three separate wells for each DNA standard and three separate wells for each unknown DNA sample. Accurate multi-channel pipettes and reagent reservoirs can be used to increase throughput. Black plates are recommended to minimize fluorescence bleed-through between wells. We have found that all black 96-well plates from Greiner Bio-one or Corning give the most consistent signal-to-noise ratio at low DNA concentrations.
6. Add 10 uL of each dsDNA standard and unknown into its own separate well and mix well by pipetting up and down.
7. Incubate the microplate at room temperature for 10 minutes in the dark. The assay plate should be read within 1 hour for best results, but can be stored

Continued next page

and read up to 24 hours after preparation with only minor loss of accuracy.

8. Measure fluorescence using a microplate reader with 350 nm excitation/460 nm emission.
9. Generate a standard curve to determine the unknown DNA concentration (see Figure 3). Average the triplicate values for each sample and subtract the average zero DNA value from each data point. Plot the fluorescence values for the DNA standards on the y-axis and ng/well DNA on the x-axis, and fit a trend line through these points to generate a standard curve with a y-intercept = 0. Use the equation for the standard curve trend line to calculate the amount of unknown DNA in each well ($y = \text{fluorescence}$ and $x = \text{ng DNA per well}$). Note: the standard curve shown in Figure 3 is for reference only. You must generate your own standard curve using your instrument to calculate the amount of DNA in your unknown samples.

Considerations for Data Analysis

Calf thymus DNA can often serve as a reference for most plant and animal DNA because it is double-stranded, highly polymerized and is approximately 58% AT (42% GC). At times it is preferable to use a dsDNA standard similar to the unknown samples (i.e. similar in size, linear vs. circular). We have found that most linear dsDNA yield similar results; however, it is best to compare the concentration of the unknown sample to a more appropriate standard if necessary. If the fluorescence of an unknown sample is higher than the linear range, further dilute the sample and add 10 μL of the diluted sample to perform the assay. For consistency, it is best to use the same volume in all the wells with samples that do not have high levels of contaminating substances.

Fluorescence quantitation by the AccuBlue Broad Range assay is linear from 2 – 2000 ng dsDNA. The dynamic range can be extended to 4000 ng with some loss of linearity (Figure 4).

Due to differences in instruments, check instrument settings to optimize for the best linearity. Some factors that can affect the final linearity and relative fluorescence intensity are: (1) the excitation and emission wavelengths and bandwidths, (2) cut-off filters, (3) sensitivity settings, (4) pipette accuracy, and (5) microplate manufacturers.

The effects of common DNA contaminants such as salts, solvents, detergents and protein on the AccuBlue™ Broad Range assay are listed in Table 2. Please also see our AccuBlue™ High Sensitivity and AccuClear™ Ultra High Sensitivity dsDNA Quantitation Assays (related products), which have different tolerances for certain contaminants compared to AccuBlue™ Broad Range.

Table 2. Effect of common DNA contaminants on AccuBlue Broad Range assay

Compound	Initial concentration in DNA sample	Final concentration in assay (200 μL)	Decrease in Signal
Sodium Chloride	500 mM	25 mM	-5%
Magnesium Chloride	100 mM	5 mM	35%
Sodium Acetate	600 mM	30 mM	50%
Ethanol	20%	1%	4%
Phenol	2%	0.10%	3%
SDS	0.2%	0.01%	32%
SDS	0.02%	0.001%	5%
Triton X-100	0.2%	0.01%	8%
Triton X-100	0.02%	0.001%	-5%
Tween-20	0.1%	0.005%	20%
CTAB*	0.01%	0.0005%	64%
dNTPs	2 mM	100 μM	-1%
BSA**	0.8 mg/mL	0.05 mg/mL	27%

* Average change for 250 ng to 2000 ng samples. Complete loss of signal is seen below 250 ng.

** Average change for 20 ng to 2000 ng samples. Not compatible with quantitation below 20 ng due to increased background at low DNA concentrations

AccuBlue™ Broad Range dsDNA Quantitation Solution
PS006

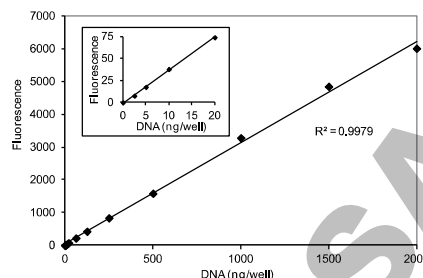


Figure 3: Standard curve of calf thymus DNA assayed using AccuBlue Broad Range Kit and read on a microplate reader (Ex/Em 350/460). Inset shows the lower end of the titration.

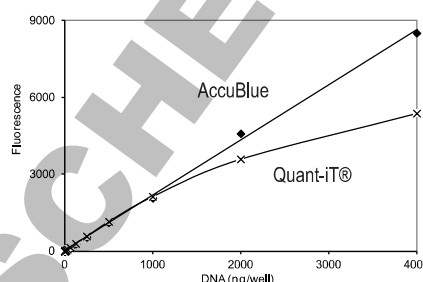


Figure 4: Two-fold dilutions of calf thymus DNA were assayed using AccuBlue or Quant-iT® Broad Range assay kits. AccuBlue has improved linearity and wider dynamic range than the Quant-iT Broad Range.

Related Products

Catalog number	Product
31007C	AccuBlue™ Broad Range dsDNA Standards, Set of 9
31007	AccuBlue™ Broad Range dsDNA Quantitation Kit with 9 DNA Standards (1000 assays)
31060	AccuBlue™ NextGen dsDNA Quantitation Kit (1000 assays)
31028	AccuClear™ Ultra High Sensitivity dsDNA Quantitation Kit with 7 DNA Standards (1000 assays)
31029	AccuClear™ Ultra High Sensitivity dsDNA Quantitation Kit with 1 DNA Standard (2000 assays)
31006	AccuBlue™ High Sensitivity dsDNA Quantitation Kit with 8 DNA Standards (1000 Assays)
31066	AccuGreen™ High Sensitivity dsDNA Quantitation Kit (for Qubit)
41003	GelRed™ Nucleic Acid Gel Stain, 10,000X in water
31041-T	Forget-Me-Not™ qPCR Master Mix, trial size
31043-T	Forget-Me-Not™ Universal Probe Master Mix, trial size

Please visit our website at www.biotium.com for information on our life science research products, including environmentally friendly EvaGreen® qPCR master mixes, fluorescent CF™ dye antibody conjugates and reactive dyes, apoptosis reagents, fluorescent probes, and kits for cell biology research.

Qubit is a registered trademark of Thermo Fisher. AccuBlue, AccuClear, and AccuGreen are trademarks of Biotium, Inc. AccuBlue, AccuClear and AccuGreen are covered by granted and/or pending U.S. and/or international patents. Materials from Biotium are sold for research use only, and are not intended for food, drug, household, or cosmetic use.