Operating Manual

Laboratory autoclave
Steam sterilizer
CertoClav Connect

Content

CE Declaration of Conformity	3
Introduction.....	4
Read Carefully Before Using	4
Proper Use	4
<i>Unwrapping</i>	<i>6</i>
<i>Product contents.....</i>	<i>6</i>
<i>Place of use.....</i>	<i>6</i>
<i>Installation of CertoClav Connect.....</i>	<i>6</i>
<i>Assembly and set up CertoTouch controller</i>	<i>7</i>
<i>Turn on CertoTouch</i>	<i>7</i>
<i>Turn on CertoClav Connect pot.....</i>	<i>7</i>
<i>Choosing the language.....</i>	<i>7</i>
<i>Checking for updates.....</i>	<i>7</i>
<i>Connecting CertoTouch with the pot.....</i>	<i>8</i>
<i>Registering new users.....</i>	<i>8</i>
<i>Logging in without using the CertoCloud.....</i>	<i>9</i>
<i>Logging in using the CertoCloud.....</i>	<i>9</i>
Test run safety instructions	10
Description of the working steps.....	11
<i>Turning on the CertoTouch.....</i>	<i>11</i>
<i>Turning on the CertoClav Connect pot.....</i>	<i>11</i>
<i>Connecting the CertoTouch with the the pot.....</i>	<i>11</i>
<i>Opening lid</i>	<i>12</i>
<i>Pouring in water.....</i>	<i>12</i>
<i>Loading substance</i>	<i>13</i>
<i>Closing lid.....</i>	<i>13</i>
<i>Select and run a program.....</i>	<i>15</i>
<i>Unloading.....</i>	<i>16</i>
<i>Interrupting a cycle</i>	<i>17</i>
<i>If power supply is cut off during operation.....</i>	<i>17</i>
Description of a program.....	18
<i>Process status screen</i>	<i>18</i>
<i>Operation of Cycles.....</i>	<i>18</i>
<i>Sterilization Cycle.....</i>	<i>18</i>
<i>End of the program.....</i>	<i>19</i>
Description of the options	19
<i>Flexible media temperature sensor.....</i>	<i>19</i>
Maintaining of the autoclave	20
<i>Condensate container</i>	<i>20</i>
<i>Cleaning the device housing</i>	<i>20</i>
<i>Real-time-clock battery.....</i>	<i>20</i>
<i>Gasket</i>	<i>20</i>
Troubleshooting	21
Technical data.....	22
<i>Certificates.....</i>	<i>23</i>
<i>Technical data CertoTouch</i>	<i>23</i>
Indication labels	24
Warranty	24

CE Declaration of Conformity

PRODUCT: CERTOCLAV AUTOCLAV

This is to certify, that the **CERTOCLAV** Product Range Connect, Essential, MultiControl and MultiControl 2 complies with the following regulations of the European Community:

2014/35/EU Low Voltage Directive

Associated Standards:

EN 61010-2-040:2015 Safety requirements for electrical equipment for measurements, control and laboratory Part 2-040: Particular requirements for sterilizers and washer-disinfectors used to treat medical materials.

EN 61010-1:2010 Safety requirements for electrical equipment for measurements, control and laboratory use – Part 1: General requirements

The CertoClav pressure equipment is constructed according Low Voltage Directive.

2014/30/EU Electromagnetic compatibility

Associated Standards:

EN 61326-1:2013 Electrical Equipment for measurements, control and laboratory use - EMC Requirements-

Part 1: General requirements (provided replacement: EDIN EN 61326-1:2018)

EN 61000-3-2:2014 Electromagnetic Compatibility (EMC) — Part 3- 2: Limits —

Limits for harmonic current emissions (equipment input current equal to or less than 16 A per phase)

EN 61000-3-3:2013 Electromagnetic compatibility — Part 3- 3: Limits — Limits for voltage fluctuation and flicker on the public low-voltage system (equipment having an input current equal to or less 16 A per phase) and not subject to conditional connection.

2011/65/EU Restriction of the use of certain hazardous substances (RoHS) Directive

Associated Standards:

EN 50581:2012 Technical documentation for the assessment of electrical and electronic products with respect to the restriction of hazardous substances

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Quality of Production:

CertoClav produces according modern production methods and with high quality materials. Experience for many years and a competent team ensure high quality of products and professional service. All units are 100 % tested on function and quality. CertoClav manufactures in Austria only.

Manufacturer:

CERTOCLAV Sterilizer GmbH

Georg Grinninger-Str. 37

A-4050 Traun / Austria

Traun, 27/09/2018

Michael Dirix
CEO

Introduction

This document has been carefully compiled. If you have any questions or required information is not covered in the manual, please contact:

CertoClav Sterilizer GmbH

Georg Grinninger Str. 37

A-4050 Traun

Austria

Tel.: +43 (0) 7229 689

Email: support@certoclav.com

Read Carefully Before Using

Please read this manual carefully before using the device to ensure safe and proper use.

Please consider that the following instructions insure your own safety and proper use of the device.

Proper Use

The CertoClav Connect small-steam-sterilizer is used to sterilize materials which can withstand temperatures up to 140°C and 2.7 bar of relative pressure. This autoclave is mainly used to sterilize microbiological materials in the laboratory:

- Glass
- Instruments
- Liquid substances and serums
- Media
- Rubber
- Filters
- Synthetic materials

Because of the freely programmable nature of this autoclave it can also be used for material tests and research purposes.

- ⚠ Use exclusively the CertoTouch-controller to operate your autoclave.
- ⚠ The lid, chamber, gasket and panel are extremely hot immediately after the completion of operation. Do not touch the equipment or you may get burned.
- ⚠ Keep this manual in a visible, easily accessible location.
- ⚠ The autoclave should not be installed in explosive areas
- ⚠ Never connect the power cord to a power supply other than one of the rated voltage. Connection to such a power supply can cause fire or electric shock.
- ⚠ Only for laboratory use by trained personnel. Do not use this autoclave for medical purposes.
- ⚠ This autoclave can be serviced only by our authorised dealers. Only original parts may be installed.
- ⚠ Remove the power plug before removing the electronic housing or the bottom.
- ⚠ Do not use the autoclave without water. The amount of water needed for sterilization: minimally 0.7 Litre, maximally 2.5 Litre of demineralized soft water.
- ⚠ Please do not sterilize tightly closed containers, especially sealed glass bottles without pressure equalisation.
- ⚠ Please, transport only with emptied chamber.
- ⚠ Do not open the lid under pressure with force.
- ⚠ Under no circumstance let the steam in the direction of another person. Risk of a burn. Please use either the provided waste steam container or connect an external waste steam chamber under own responsibility with a hose (inside diameter 10mm) to let out the hot steam.
- ⚠ If filled with glass bottles, open the lid only after the temperature of the liquid has dropped under 80°C. Danger of over boil of the liquids or of an explosion of the closed bottles.
- ⚠ Under no circumstance put explosive or flammable materials in the autoclave.
- ⚠ If the electrical supply is damaged immediately remove the device from the power supply. It can be used again only after a qualified person has repaired the autoclave.
- ⚠ Never put objects in the ventilation slot nor pour liquids in the ventilation slot. Danger of electric shocks and short circuits.
- ⚠ Manipulating or rebuilding the device or the programs can lead to malfunction.
- ⚠ To avoid corrosion in the chamber and the conduit system do not sterilize heavy saline substances.

Unwrapping

- Take the device out of the carton by holding the handles on the lid and lifting it out. Furthermore, remove the CertoTouch controller, power cord and equipment.
- Place the device on stable ground.
- Check if the information on the label coincides with the delivery note.

Product contents

The following items were delivered with the autoclave:

- CertoClav Connect autoclave 18L
- Stainless steel base plate with water level indicator
- Power cord
- CertoTouch control terminal with stand bracket, 4 screws and a power cord
- Operating manual CertoClav Connect
- Test and guarantee certificate

Should one of those parts be missing, please contact your distributor or us directly at:

support@certoclav.com or per telephone +43 (0) 7229 689

Should you need additional equipment or a warranty packet, you can order it per telephone, per email or via your distributor or in the online shop at:

<https://www.certoclav.com/shop/>

Place of use

- ⚠ Avoid temperatures below 5 ° and beyond 35 °C and air moisture beyond 80 %.
- ⚠ The operating terminal is splash-proof. Please guard the autoclave from any splash water.
- ⚠ Avoid places where convulsions and vibrations appear and do not use the device outdoors.
- ⚠ Take care that the device is not used near steam sensitive laboratory devices (for example microscopes).
- ⚠ Avoid places where the ventilation slot might be blocked.

Installation of CertoClav Connect

The voltage values required for this device are to be observed.

If you do not know what values the power supply supplies, ask your electrician or at the local electrical utility.

- Turn on the device by connecting the power cord to the power supply and turn the rocker switch on the left side to |.
- If the device was locked, it will now unlock automatically.
- Push and turn the red lever on the right lid handle clockwise.
- Turn the lid anti-clockwise to open the autoclave.
- Take out the baskets and other enclosed equipment.
- Assemble the CertoTouch controller next to the autoclave as described in its operating manual

Assembly and set up CertoTouch controller

- Please mount the stand bracket to CertoTouch-controller with 4 screws. The screws are included in the packaging with the stand bracket.
- If provided, connect the extension cable directly to the CertoTouch controller. There is only one suitable connector.
- Plug the enclosed power supply unit either to the extension cable or directly to the CertoTouch.
- Place the CertoTouch controller only on a levelled, stable surface. A suitable viewing angle can be adjusted by means of the standing bracket.

Turn on CertoTouch

Please press the -button at the front right of CertoTouch for at least 3 seconds.

Turn on CertoClav Connect pot

Turn on by turning the rocker switch to | . If the switch lights green, then the device is turned on. This is necessary to be able to connect the CertoTouch with the CertoClav Connect pot via Bluetooth.

Choosing the language

- 1) Go to the settings menu by clicking the -button at the upper right corner of the screen.
- 2) In the Menu – settings open the language settings by clicking on the “Language” button.
- 3) In the detail view at the upper right corner of the screen click the current language settings.
- 4) A scrolling list of supported languages can be seen now. Please choose here by clicking the desired one.

Checking for updates

- 1) Please make sure CertoTouch is connected to the Internet.
- 2) Go to the Settings menu by clicking the -button at the top right corner of the screen.
- 3) Please enter the “Device” section by selecting the entry “Device” in the left list.
- 4) In the detailed view on the right, please scroll to “Check for updates” and select this entry.
- 5) A progress bar with the download progress will pop up. Once all updates have been downloaded, they will be installed. If a new software version is available, an installation wizard will guide you through the installation process.

Connecting CertoTouch with the pot

- 1) Please navigate to the Connection Manager. For this purpose, please log out and press the


-button at the login window.

- 2) Please, press the  -button in order to determine all reachable CertoClav controllers. This may take a few seconds.
- 3) In this list, you can see the serial numbers of all reachable CertoClav controllers.
- 4) Please choose the controller you want to connect with by pressing that serial number in the list. The serial number of each controller can be found on the left or right side of the controller box.
- 5) CertoTouch will connect itself to the selected CertoClav autoclave. The connection is successfully established, if a “connected” status message appears in the selected list sentry. Only after a connection has been successfully established, it is possible to return to the login-window. Please log in there with your email and password in order to use the CertoTouch controller.

Registering new users

- 1) Navigate to the login window. For this, a logged in user must log out first.
- 2) Click on the  - button at the top right of the screen.
- 3) Please confirm the popup notification in order to create a new user.

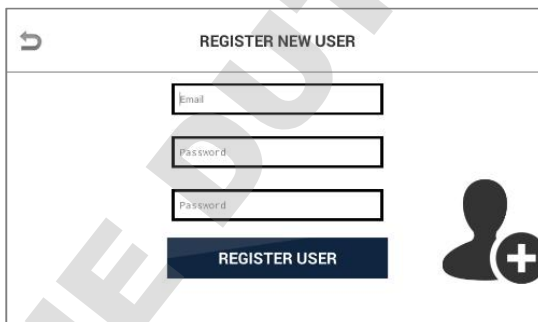- 4) A new window for the user registration will appear.
- 5) Enter the user's email address. Note: The “@” symbol can be reached in the keyboard by pressing the “SYM” (symbols) button.
- 6) Please choose a password with at least 4 characters and retype it.
- 7) Confirm the user registration, by clicking the “Register user” button.

Logging in without using the CertoCloud

- 1) Make sure that the online mode is deactivated. The online mode can be enabled or disabled in the settings menu  .
- 2) Please select the email of the user from the drop-down menu.
- 3) Please type the user password.
- 4) Click the “Login” button.

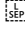
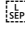
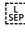
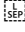
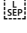
Logging in using the CertoCloud

- 1) In order to take advantage of the CertoCloud, please create a CertoCloud account first. You can create a CertoCloud account at www.certocloud.com.
- 2) Please connect the CertoTouch with the internet.
- 3) On the CertoTouch, the online mode must be enabled. The online mode can be enabled or disabled in the settings menu  .
- 4) Navigate to the login window. For this, a logged in user must log out first.
- 5) Please select the email of the user from the drop-down list. The email must be the same email as the one of the CertoCloud account(www.certocloud.com). If this email address is not available in the dropdown list, refer to the chapter “Register a new user”.
- 6) Please enter the CertoCloud password of the user.
- 7) Click the “Login” button.
- 8) Your account is now connected with the CertoCloud. On www.certocloud.com you can access you autoclave, define custom programs and access your protocols.

Test run safety instructions

Please follow the following safety instructions:

Please take into consideration all of the instructions and warnings noted on the device.

- ⚠ Do not put your face or hands close to the chamber when lifting the lid after a program has finished; steam will gush out of the chamber. 
- ⚠ Never put your fingers above the steam chamber. The outflowing steam may cause burns.
- ⚠ Open the lid using the handle just a few centimetres first, to let the steam out. Keep safety distance from the device. Afterwards, carefully open the lid fully.
- ⚠ Please, keep foreign materials (metals, liquids) from the ventilation, because a current voltage can be induced inside the autoclave, which can cause short circuits or electric shocks. Do not forcibly bend, twist, tie or extend the power cord. Do not place heavy objects on the cord. A damaged cord or exposed wire can cause fire or electric shock. 
- ⚠ Close the lid after confirming that no foreign matter is adhering to the section contacting the lid gasket. Foreign matter in this section can cause vapor leaks. When using a waste processing bag or other kind of bag and disinfecting, place the bag in the metal mesh holder and then insert it into the chamber. Using the bag “as is” can cause excessive temperatures, pressures, lack-of-water, etc. 
- ⚠ Please take care not to pinch your hands while closing the lid.
- ⚠ The lid, chamber, gasket and panel are extremely hot immediately after completing a program. Do not touch the equipment or you may get burned. Put on heat insulating gloves before removing a substance from the chamber. Do not put hands into the chamber until the steam has been vented. 
- ⚠ Some time is required for liquids to cool. Be sure to check that the temperature has dropped sufficiently before unloading a liquid from the chamber or burns can result. 
- ⚠ Do not remove the exhaust bottle before water in the bottle has sufficiently cooled. 
- ⚠ If any abnormality occurs (e.g. abnormal sounds, smells, smoke), immediately shut off the power. After checking to see that the abnormal condition does not continue, call our authorized distributor in your region. 
- ⚠ When transporting the device keep the red lever on the handle on “Close” at all times. It prevents the lid from opening on its own.
- ⚠ Do not pour anything except demineralized water for sterilizing into the autoclave.

The first test run serves as a functionality control. It should discover the damage that may have occurred during the transport. Furthermore, it demonstrates how to properly handle the device.

Description of the working steps

Turning on the CertoTouch

Please press the -button at the front right of the CertoTouch for at least 3 seconds.

Turning on the CertoClav Connect pot

Turn on the device by turning the switch rocker on | . If the rocker lights green, the device is on. If the temperature in the chamber is under 80°C, the device unlocks itself automatically.

Connecting the CertoTouch with the the pot

- 1) Please navigate to the connection manager. For this purpose, please log out and press the

-button at the login window.

- 2) Please press the -button in order to determine all reachable CertoClav controllers. This may take a few seconds.
- 3) In this list, you can see the serial numbers of all reachable CertoClav controllers. Please choose the controller you want to connect with, by pressing that serial number in the list. The serial number of each controller can be found at the left or right side of the controller box.
- 4) CertoTouch will connect to the selected CertoClav controller. The connection is successfully established, if a “connected” status message appears on the selected list entry. Only when a connection is successfully established, it’s possible to return to the login-window. Please log in there with your email and password in order to use the CertoClav controller.

Opening lid

Ensure that the autoclave is powered on. As soon as the program was finished or the temperature and pressure in the chamber are low enough, the electrical locker unlocks the autoclave automatically. Rotate the lid anti-clockwise to take it off.

- ⚠ SAFETY HINT: If loading contains glass bottles wait until chamber temperature is below 80°C, for liquids wait until chamber temperature has reached 70°C or until liquid temperature has reached 80°C. Danger of boiling over of fluids.

Pouring in water

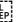
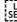
- ⚠ Do not pour anything except demineralized water. Please use soft water since the properties of the water affect the life time of the device.

Please follow the following instructions:

Pour in demineralized soft water in the sterilization chamber until you can see the water through the openings of the base plate. On the base plate is a mark for the minimum amount of water. Needed amount of water: minimally 0.7 Litre, maximally 2.5 Litre.


Loading substance

Please follow the following instructions:

- ⚠ Close the lid after confirming that no foreign matter is adhering to the section contacting the lid gasket. Foreign matter in this section can cause vapor leaks. When using a waste processing bag or other kind of bag and disinfecting, place the bag in the metal mesh holder and then insert it into the chamber. Using the bag "as is" can cause excessive temperatures, pressures, lack-of-water, etc.
- ⚠ Be careful not to pinch hands when closing the lid. Always use the base plate.
- ⚠ Be sure that the temperature has dropped to at least 50°C in the chamber before the beginning of the next cycle.
- ⚠ If a waste disposal bag is used in sterilization, open the bag enough that the bag is not in contact with the inside surface of the chamber. Insufficient sterilization may be caused if the bag is closed during sterilization. When the bag is opened excessively, steam is prevented from circulating in the chamber. This may also result in insufficient sterilization. 
- ⚠ Do not pile specimen on top of one another. When the chamber is overly packed, steam fails to penetrate all points, resulting in an incomplete sterilization. 
- ⚠ In sterilizing liquids such as chemicals and media pay attention to the quantity of the liquid in relation to its container. For an Erlenmeyer flask, the amount of chemical should be approx. 3/4 of the capacity of the container; for a test tube, the appropriate quantity of chemical is approx. half of the capacity of the container. Too much chemical may result in overflow from the container during the temperature rising or cooling process. 
- ⚠ Use container caps that are loose fitting and allow the passage of air. Containers may break if venting is not possible. 
- ⚠ In the case of dissolution of agar medium, its quantity should be 2 liters or less per container. Two liters or more of agar medium may not be completely dissolved. 

Closing lid

Please check that the lid seal is correctly inserted. The signature "Place This Side Facing Down" must be readable. See figure below. Then close the lid in exactly the opposite way, as it was previously opened.

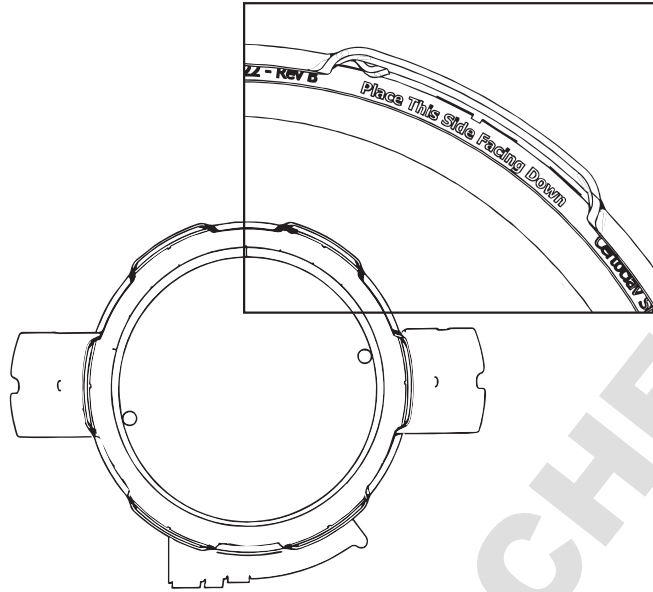DOMINIQUE DUTSCHER SAS

Select and run a program

- 1) Log into your account first. After a user has logged in, the main menu appears.
- 2) Visit the STERILISATION-window by navigation to: MAIN-MENU->STERILIZATION.
- 3) Find the correct program by scrolling the list of available programs.
- 4) By tapping on the start button of the specific program, the user can start the program.
- 5) If activated, please follow the STEP-BY-STEP guide until the program starts. Note: The STEP-BY-STEP guide can be disabled or enabled in the Settings->Sterilisation->STEP-BY-STEP-GUIDE section.

For further instructions on how to use the CertoTouch please refer to the provided operating manual.

Unloading

In order to open the device and take out the sterilized goods must the temperature in the chamber drop to under 80°C. This serves as a safety measure to prevent burns.

As soon as the cycle ends, a message "Please open the lid" appears on the screen. The program counts as successfully finished only after opening of the lid.

- ⚠ Keep the face and hands away from the chamber when lifting the lid after operations are complete; steam will gush out of the opening of the chamber.
- ⚠ After operation is finished, the lid, chamber, packing and panel are hot. To protect yourself from burns, do not touch with bare hands. Put on heat insulation gloves before taking a substance out of the chamber. Do not put your hands into the chamber until the steam has been vented.
- ⚠ It takes a lot of time for the liquid to cool. Be sure to check that the temperature has dropped sufficiently before unloading the liquid from the chamber, or you may get burned.

To open the lid please follow the instructions:

Ensure that the autoclave is powered on. As soon as the program was finished or the temperature and pressure in the chamber are low enough, the electrical locker unlocks the autoclave automatically. Rotate the lid anti-clockwise to take it off.

- ⚠ SAFETY HINT: If loading contains glass bottles wait until chamber temperature is below 80°C, for liquids wait until chamber temperature has reached 70°C or until liquid temperature has reached 80°C. Danger of boiling over of fluids.

Interrupting a cycle

- ⚠ If the cycle is interrupted during the sterilization of liquids, the lid is electrically unlocked after the temperature in the chamber has dropped to under 80°C.
- ⚠ In case of power loss, the lid locks itself.


If power supply is cut off during operation

If the power supply is cut off due to power failure or the like, operation is interrupted. When power supply is restored, the equipment will be back in the standby state (state before operation). In this case, repeat operations from the beginning. The autoclave can not be opened when turned off.

Description of a program

Process status screen

If a program is running, the MONITOR window is visible. In the monitor window, all important information about the current cycle is displayed. The following information can be found (additionally marked in the screenshot below).

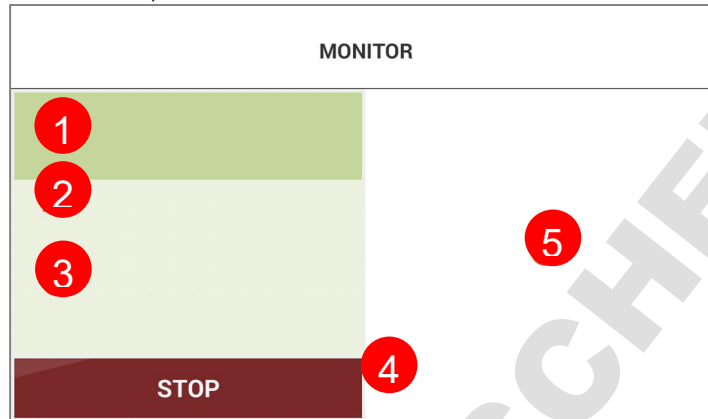- 1 Name of the running program.
- 2 Status of the running program. This can be "Running", "Finished", "Prepare to start", "Not running".
- 3 Description of the running program.
- 4 Current temperature in [°C] and pressure in [kPa], measured in the chamber.
- 5 Graphical overview.

Operation of Cycles

If activated in the program, the autoclave will first be aired.

Any air remaining in the chamber makes the temperature distribution in the chamber uneven. This hinders temperature rise (sterilization). To cope with this, a microcomputer-controlled automatic exhaust valve vents out virtually 100 % of the air in 10 minutes at 115°C. In the chamber remains water steam. At the end the valve closes.

Sterilization Cycle

A constant temperature (pressure) is maintained during the set sterilization time period. If the temperature in the chamber drops 1°C or more or rises 2°C or more from the set value due to any trouble, the program is interrupted.

-  In sterilization of petri dishes or empty containers, the air remaining in the container expands and may increase the pressure extraordinarily within the chamber. If the pressure in the chamber exceeds the saturated steam pressure, the automatic exhaust valve opens and discharges the air in the chamber into the exhaust bottle.

End of the program

After all the cycles have been finished, a message "Please, open the lid" comes forth. When the lid is opened, the program is finished.

- ⚠ If the touch screen is not touched for 30 minutes, the display becomes black. By touching the touch screen will the display turn on again.

Description of the options

Flexible media temperature sensor

In general

The flexible media temperature sensor measures the temperature in the media that need to be sterilized. Because of the delay of boiling, for example by sterilization of liquids, it is reasonable to use a media temperature sensor. Additionally, you can control the exact temperatures, which the media was exposed to, after the sterilization cycle.

- ⚠ Do not forcibly bend, twist, tie or extend the cable of the temperature sensor. Do not place heavy objects on the cord. A damaged cord or exposed wire can cause malfunction.
- ⚠ When sterilizing plastic petri dishes, please, take care that the media temperature sensor does not contact any part of the petri dishes. Because the plastic sclerotizes when the temperature drops, it becomes impossible to separate those parts from the temperature sensor.

Use

After loading the chamber, set the media sensor in the middle of the media you want to sterilize. The media temperature during the program always has precedence over the chamber temperature. For example, the sterilization duration begins only after the media has reached the sterilization temperature.

Deactivating of the media sensor

In the program settings (see CertoTouch operating manual) it can be specified whether the media sensor should be ignored or used.

Maintaining of the autoclave

Condensate container

During every sterilizations process the water gathers in the condensate container. Empty the condensate container after each cycle. Check before each new cycle, if the condensate container has been emptied.

- ⚠ Do not empty the condensate container before it has cooled sufficiently.
- ⚠ The steam exhaust valve closes itself if the condensate container was not installed.

Cleaning the device housing

- ⚠ Please, do not use any aggressive detergents to clean the device housing.

Real-time-clock battery

The devices intern battery supplies a real-time-clock with power. As soon as the battery dies, there will be a warning showed on the screen that must be confirmed. The average life expectancy of the battery is 5 years. The autoclave can also be operated with a dead battery. Without the real-time-clock it may come to inaccuracies up to a couple of minutes at longer user-defined programs. To change the battery please contact us. When your autoclave is not connected to the network or the internet there is a possibility of fluctuations in time and date.

Gasket

A brittle gasket could lead to leakage of steam. The average life expectancy of a gasket is 4 years. When the gasket needs to be replaced or you wish to buy a new one, please contact us.

Troubleshooting

Should any malfunction occur in the autoclave, the error detection circuit will be triggered to assure safety. Once the circuit is activated, an error number appears on the digital display. The alarm will stop only after the error has been confirmed.

Error code	Error description	Next step
Error 1	Faulty ventilator	Replace the ventilator
Error 2	The lid is not closed	Please contact the service-team. Please, close the lid prior to starting a program. Follow the step-by-step video.
Error 3	Too much water in the chamber	Please, let some water out of the chamber. There should be maximally 2,5 Litre water in the chamber during the cycle. If the error does not disappear, please contact the service-team.
Error 4	Too little water in the chamber	Please, fill the chamber with minimally 0,7 Litre water in the chamber before the start of a program. If the error does not disappear, please contact the service-team.
Error 5	The lid cannot be locked	Please, turn off the device and try once more. If the error does not disappear, please contact the service-team.
Error 6	Sterilization heater trouble alarm	Temperature in the chamber cannot be achieved. Either the autoclave is empty of water or the heater is faulty. Please, contact the service-team.
Error 7	Over – pressure alarm	Please, contact the service-team.
Error 8	Over – temperature alarm	Please, contact the service-team. Possible defects: temperature sensor is faulty or the heater malfunctions.
Error 9	Pressure out of permissible range	Unexpected pressure fluctuations in the chamber. Please, do not put any tightly closed containers in the chamber.
Error 10	Temperature in the chamber out of permissible range	Please, contact the service-team. Temperature fluctuations during the cycle. Possible causes: heater or temperature sensor broken.
Error 11	Temperature of the substance out of permissible range	Please, contact the service-team. Temperature fluctuations during the cycle. Possible causes: heater defect, temperature sensor defect.
Error 13	Temperature band not adhered to	Please, contact the service-team. Possible combined error of Error 10 and Error 11
Error 14	Cycle aborted because of a fatal error.	If a fatal error during the sterilization occurs, and the program must be aborted, then will an additional error be showed.
Error 15	Cycle aborted because of user's interference.	No action needed. The program was manually cancelled. Please, wait until the temperature in the chamber has dropped to under 80°C.
Error 16	Tablet-PC and autoclave connection terminated	Please set the Touch controller closer to the autoclave and renew the connection. You should not move the Touch controller further that 6m from the autoclave.

Technical data

Measures and weight			
Material	Stainless steel 304L	Certificates	CE
Dimensions			
Chamber Volume	18 Litre	Device height	60 cm
Chamber diameter	26 cm	Device width	44 cm
Chamber height	35 cm	Device depth	40 cm
Chamber usable height	30 cm	Device weight	15 kg
Operating data			
Max. operating temperature	140 °C	Power supply	230V, 50/60 Hz
Max. operating pressure	2,7 Bar (39 psi)	Power	1300 W
Test pressure	4,1 Bar (59 psi)		

Heating duration

Amount of load	Heating from 25 °C to 121 °C
Little	15 – 20 Minutes
A lot	25 – 40 Minutes

Cooling duration

Amount of load	Cooling duration at faster steam exhaust from 121 °C to 80°C	Cooling duration at slower steam exhaust from 121 °C to 80°C
Little	20 – 35 minutes	40 – 60 minutes
A lot	40 – 60 minutes	80 – 120 minutes

Correlations table

T in °C	p in bar (kgf/cm ²)	T in °C	p in bar (kgf/cm ²)
100	0,03	118	0,90
101	0,07	119	0,96
102	0,11	120	1,02
103	0,15	121	1,09
104	0,19	122	1,15
105	0,23	123	1,22
106	0,28	124	1,29
107	0,32	125	1,37
108	0,37	126	1,44
109	0,41	127	1,52
110	0,46	128	1,59
111	0,51	129	1,67
112	0,56	130	1,75
113	0,61	131	1,84
114	0,67	132	1,92
115	0,72	133	2,01
116	0,78	134	2,10
117	0,84	135	2,19

Certificates

CE certified, please see chapter EC declaration of conformity.

Technical data CertoTouch

Display screen	7" LED background light
Resolution	480 pixels
Touch panel	capacitive
Brightness	450cd/m ²
Aspect ratio	16:9
CPU	Freescale iMX.53 x 800MHz / 1.0GHz
Core	ARM Cortex-A8
Memory card connectors	Micro-SD card und USB-Stick optional
LAN (cable)	Ethernet 10/100M (RJ45) optional
W-LAN (without cable)	IEEE 802.11
Bluetooth	Bluetooth 2.1 Host
Other Interfaces	CertoClav barcode scanner port CertoClav label printer port
Audio	Earphone slot Speaker
Speaker	1W
Power consumption	9W
Working temperature	-10°C to 65°C
Storage temperature	-25°C to 80°C
Power supply	DC 9V to 36V
Dimensions (LWD)	220x132x36.5mm
Weight	740g
Degree of protection	All-round splash water protection

Indication labels

	Directive 2012/19/EU (WEEE): the product should not be disposed of with your other household waste. It is your responsibility to dispose of your waste equipment separately from the municipal waste stream.
	WARNING: Read the operating manual before usage
	WARNING: Hot area

Warranty

This autoclave was quality, safety and functionality checked by our company. We provide a 2-year warranty on all parts, that are not consumables, from the day of the delivery.

Please keep the separate warranty card for future reference.

An optional 5-year extension to the warranty can be purchased.

No warranty in case of inappropriate use.

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German Version Page 1

Content

1	Validity of the operating instructions	3
2	General safety instructions.....	3
3	Introduction	3
3.1	Field of application	3
3.2	Technical data.....	4
3.3	Requirements for transport and storage:.....	4
4	Scope of delivery	4
5	Installation of the CertoTouch.....	5
6	Basic operation	6
6.1	Switching on the CertoTouch	6
6.2	Switching off the touch screen	6
6.3	Stand-by.....	6
7	Connection to the autoclave	7
7.1	Connecting to an MultiControl 2 Touch	7
7.2	Establish connection with a CertoClav Connect.....	7
7.3	Establish connection with a CertoClav VacuumPro.....	7
8	User registration and login	8
8.1	Sterilization program selection on screen.....	10
8.1.1	Status screen.....	10
8.2	Support menu	11
8.2.1	Tutorial videos	11
8.2.2	Send e-mail to CertoSupport.....	11
8.2.3	TeamViewer remote support and screen sharing	12
8.3	Logging menu	13
8.3.1	View logs.....	13
8.3.2	Print out protocols.....	13
8.3.3	Example protocol print out.....	14
9	Settings	15
9.1	User administration	15
9.2	Network settings.....	15
9.2.1	Connection to a WLAN network.....	15
9.2.2	Connection to a LAN network.....	16
9.2.3	Network Communication	16
9.2.4	Connection of the autoclave to the local CertoServer (for Connect and VacuumPro)	16
9.3	Sterilization settings	17

9.3.1	Export of all protocols to USB flash drive	17
9.3.2	Export of all protocols to SD card	18
9.3.3	Export of all protocols via network.....	18
9.3.4	Label Designer (only available for VacuumPro and Connect)	18
9.4	System settings.....	19
9.4.1	Install updates via the internet.....	19
9.4.2	Install updates via USB flash drive.....	19
9.4.3	Install updates via SD card.....	20
9.4.4	Restore factory settings.....	20
9.4.5	Change language.....	21
10	Error messages	21
11	Transport and storage	21
11.1	Storage conditions.....	21
12	Warranty.....	21

13 Validity of the operating instructions

	These operating instructions are only valid for the CertoTouch Tablet PC.
	Please read this manual before using the device for the first time
	Strictly adhere to all safety instructions in this operating manual.
	Keep these operating instructions within easy reach.
	If there are questions or problems with the device, contact your dealer or the manufacturer.

14 General safety instructions

-  Only use the equipment in accordance with these instructions. Store these operating instructions in a place that is easily accessible to all users.
-  Use a power outlet with a protective ground conductor and make sure that the protective ground conductor is well grounded.
-  Operate the device exclusively with the supply voltage declared on the type plate.
-  Do not touch the power plug or socket with wet hands.
-  Do not load, bend or twist the cables. Do not place heavy objects on the cables.
-  If you notice any unusual odors or noises during operation, immediately disconnect the power supply and contact your local dealer or the manufacturer via e-mail.
-  If you are not going to operate the device for a long time, please disconnect the power supply.
-  Only connect original accessories.
-  Do NOT modify the device.
-  The device may only be used in laboratories. It is not suitable for domestic use.

15 Introduction

The CertoTouch controller can be connected to the CertoClav autoclaves of the MultiControl, Connect and VacuumPro series via cable connection. CertoTouch offers you extended possibilities for operation, programming, monitoring and logging of your autoclave.

These include the editing of programs, the storage of cycle protocols locally on the touch PC or in CertoCloud, the graphic view of time, temperature and pressure progression on the display.

15.1 Field of application

This touch PC is exclusively intended for the use with a CertoClav laboratory autoclave in a laboratory setting. For example, to control the products CertoClav Connect, CertoClav MultiControl 2 and CertoClav Vacuum Pro.

15.2 Technical data

Display screen	7" LED backlight
Resolution	800×480 pixels
Touch screen	Capacitive
Brightness	450cd/m ²
Picture format	16:9
CPU	Freescale iMX.53x 800MHz / 1.0GHz
Core	ARM Cortex-A8
Memory	RAM: 512MB DDR3 (Optional 1G) ROM: 4GB NAND FLASH (Optional 8G)
Memory card connections	Micro-SD card and USB stick
LAN	Ethernet 10/100M (RJ45)
Bluetooth	Bluetooth 2.1 Host and Slave
Further interfaces	CertoClav Barcode Scanner Slot
Audio	External Audio Jack loudspeaker in front
Speaker	1W
max. Consumption	9W
Acceptable working temperature	-10 – 65 °C
Acceptable storage temperature	-25 – 8 °C
Power supply	DC 9-36V
Dimensions (LBT)	220 × 132 × 36.5 mm
Weight	740g

15.3 Requirements for transport and storage:

- Ambient temperature: 5~40 °C
- Relative air humidity: ≤80%.
- No corrosive gases in the environment

16 Scope of delivery

After you have opened the packaging, please check the product for transport damage. If you find any damage, please contact your dealer and do not operate the device for safety reasons. Please also check if the full scope of delivery is included.

Description	Quantity
Operating instructions	1
Mobile touch screen	1
Power cable	1
Base for touch screen with four fastening screws	1
Interface box for touch screen	1
RS232 cable	1
Touch Pen	1

17 Installation of the CertoTouch

Mount the base to the touch screen using the four screws provided.

Now, the touch screen with the base should look like the illustration on the right.


Connect the touch screen to the interface box with the supplied RS232 cable.

Connect the power cable of the touch screen to the interface box.


Connect the interface box to the corresponding autoclave as described in the autoclave manual.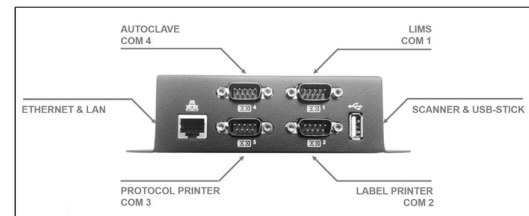


Only place the touch screen on flat, stable surfaces. You can also place the waterproof touch screen on top of or next to the horizontal autoclave, or next to the vertical models. A suitable viewing angle can be set using the base.

18 Basic operation

Click and wipe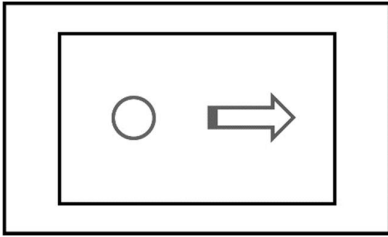


Long click and drag

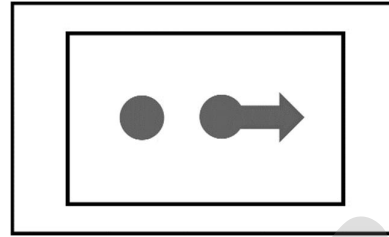18.1 Switching on the CertoTouch

To turn on the device, press the -button on the front of the device for at least 3 seconds. If a login window appears on the screen after a few seconds, the touch screen is correctly connected to the autoclave.

If a message about a connection termination with the autoclave appears on the screen, the touch screen and autoclave are not connected correctly. Check whether the autoclave is switched on and connected to the COM4 port of the interface box

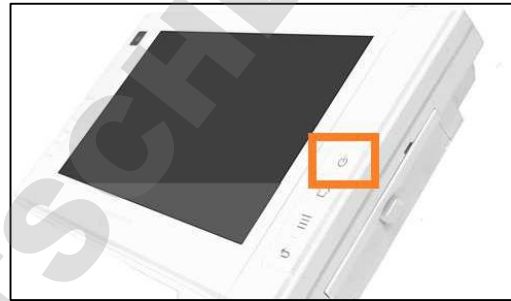First switch on the autoclave at the green power switch before switching on the CertoTouch to ensure that the connection can be established correctly.

18.2 Switching off the touch screen

Press the -button on the front panel (to the right of the screen) for at least 3 seconds. A notification will then appear on the screen. Select "Shutdown" to turn off the screen.

18.3 Stand-by

Briefly press the -button next to the screen. The backlight of the screen will turn off. Briefly press the -button again to reactivate the touch screen.

19 Connection to the autoclave

19.1 Connecting to an MultiControl 2 Touch

If you use a CertoTouch in combination with a MultiControl2 controller, connect it as follows:

7. The CertoTouch is connected to the interface box as described in chapter 5.
8. The MC2 controller is connected to the interface box with the RS232 cable. For the control, the cable goes into the printer port. For the interface box, the COM 4 port is used.
9. Switch on the autoclave first.
10. Then switch on the MC2 control unit.
11. Finally, switch on the CertoTouch and wait until the connection with the controller is established. This is done automatically but may take a moment.
12. Then create a user as described in chapter 8.

19.2 Establish connection with a CertoClav Connect

If you use a CertoTouch in combination with a CertoClav Connect, connect it as follows:


5. The CertoTouch is connected to the interface box as described in chapter 5.
6. The pot is connected to the interface box with the RS232 cable. With the interface box, the COM 4 port is used.
7. First switch on the autoclave, then the CertoTouch and wait until the connection is established. This is done automatically but may take a moment.
8. Then create a user as described in chapter 8.

19.3 Establish connection with a CertoClav VacuumPro


If you use a CertoTouch in combination with a CertoClav VacuumPro, connect it as follows:

5. The CertoTouch is connected to the interface box as described in chapter 5.
6. The VacPro is connected to the interface box with the RS232 cable. With the interface box, the COM 4 port is used.
7. First switch on the autoclave, then the CertoTouch and wait until the connection is established. This is done automatically but may take a moment.
8. Then create a user as described in chapter 8.


20 User registration and login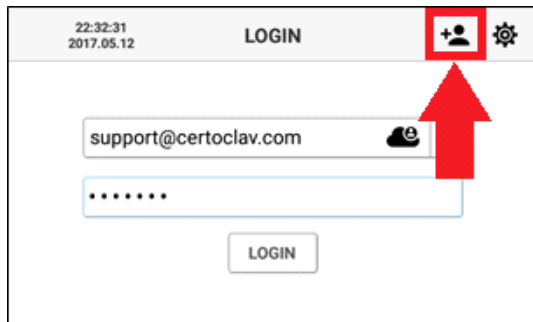

When the touch screen is turned on, the login window appears. To create a new user, click on the plus symbol at the upper right edge of the screen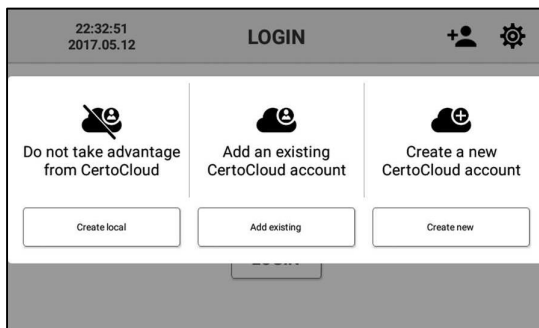

To create a local account without server connection, click the "Local" button on the left side of the selection window. To add an existing CertoCloud user account with server connection click the middle button. To create a new CertoCloud account with server connection, click the "Create New" button on the right side of the selection window.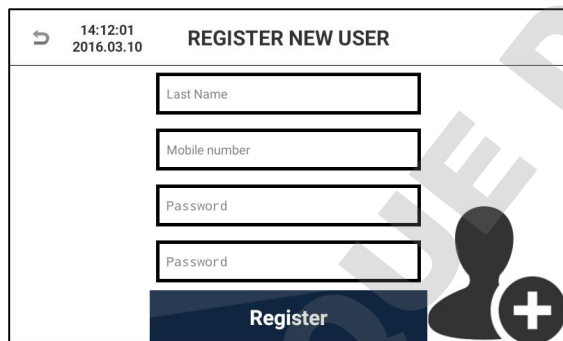

To create an account, enter your user data. Text fields can be selected and edited by touch. Enter your phone number including country code. This is required if you want to receive status notifications from your autoclave via text message. Enter a password of your choice with at least 4 characters twice. To cancel the registration, click on the "Return" button in the upper left corner of the screen. To complete the registration, click on the "Register" button.

To add an existing CertoCloud account, enter your email address and password. Then click "Add Account". To cancel the registration, press the "Return" button in the upper left corner of the screen.


To create a CertoCloud account, enter your user data into the provided text fields. Please enter your mobile phone number including country code. To confirm your email address and phone number, you will receive an e-mail and a text message. Enter the password of your choice twice and then click the "Register" button.

To sign in to a recently created account, first select your e-mail address from the drop-down menu and then enter your password. If you want to use a server-connected account, make sure that online mode is enabled. To enable online mode, go to Settings → Device → Online Mode.

When you log into your CertoCloud account for the first time, you will be asked for an activation key. You received the activation key via text message and e-mail after you created your CertoCloud account. Enter this key in the text box and click "Activate Account". Now your CertoCloud account is activated and you can login to your CertoCloud account.


20.1 Sterilization program selection on screen

- After logging in, navigate to the sterilization menu.
- In the middle 4 programs are visible. Scroll down the list to see all available programs.
- Click a program to select it. Confirm that you want to start the program.
- If enabled in the settings, step-by-step videos are shown before the program starts. This is not available with all devices.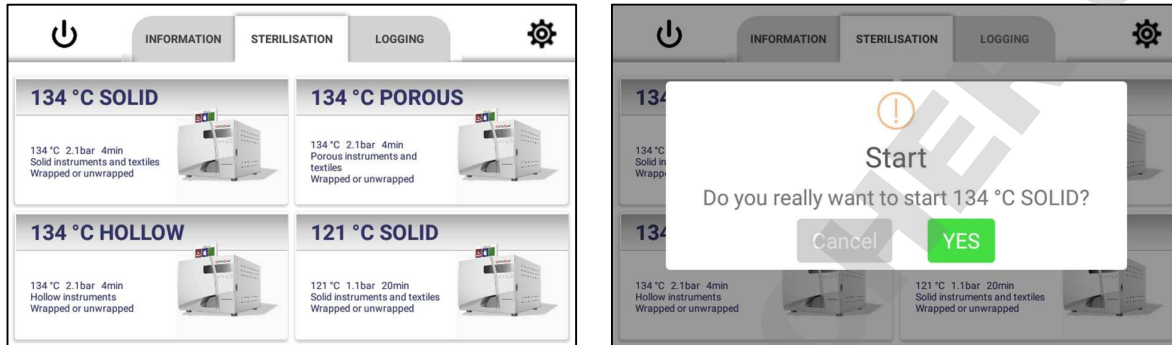

20.1.1 Status screen

During the current cycle, the status screen is visible. On the left side, the information about the current cycle is listed. On the right side of the screen, the current steam temperature and pressure are displayed. If the liquid program has been started, the liquid temperature is also displayed. To access the graphical view of the temperature and pressure graph, hold and drag the autoclave icon to the left. The graphical record will then be displayed.

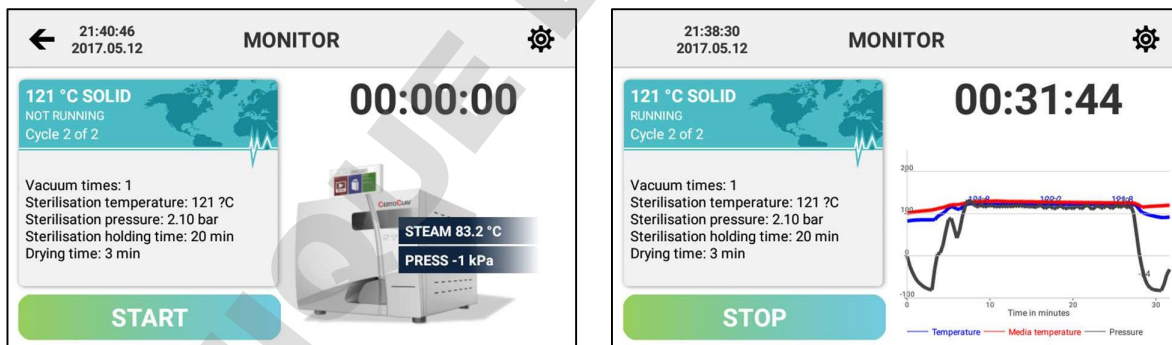
20.2 Support menu

Navigate to the Information tab to display the Information menu. Here you can watch training videos and contact CertoSupport directly from the touch screen.


20.2.1 Tutorial videos

To scroll through the list of all available videos, click the video section of the information tab. Select the desired video to view it in full screen mode.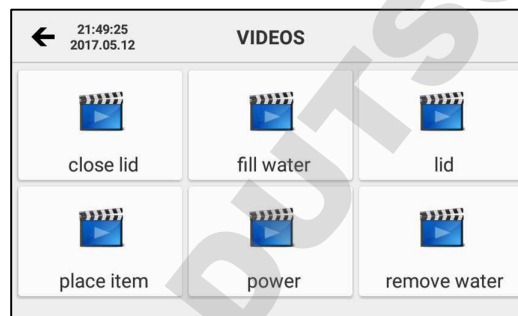

20.2.2 Send e-mail to CertoSupport


To send an e-mail to CertoSupport directly from the device, the device must be connected to the internet. To connect the device to the Internet, you can plug a network cable into the Ethernet port or establish a wireless connection to your wireless router. Click the "SUPPORT" button to open the email window.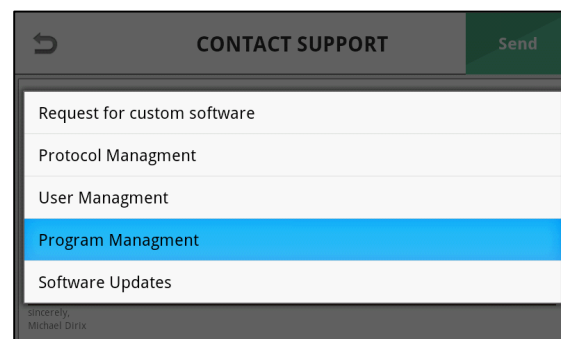


Select the subject of your message to CertoSupport from the drop-down list. Then enter your message to CertoSupport in the message text field. To send the e-mail, click the "Send" button.

20.2.3 TeamViewer remote support and screen sharing


To share the screen with CertoSupport using TeamViewer, navigate to the information tab and click the “Support” button. Remote support is only available after making an appointment via e-mail or telephone.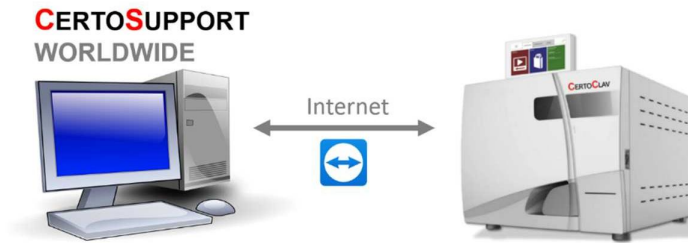

As soon as you start a TeamViewer session, a window appears which gives you the opportunity to roughly describe the problem once more. Then click “Start”.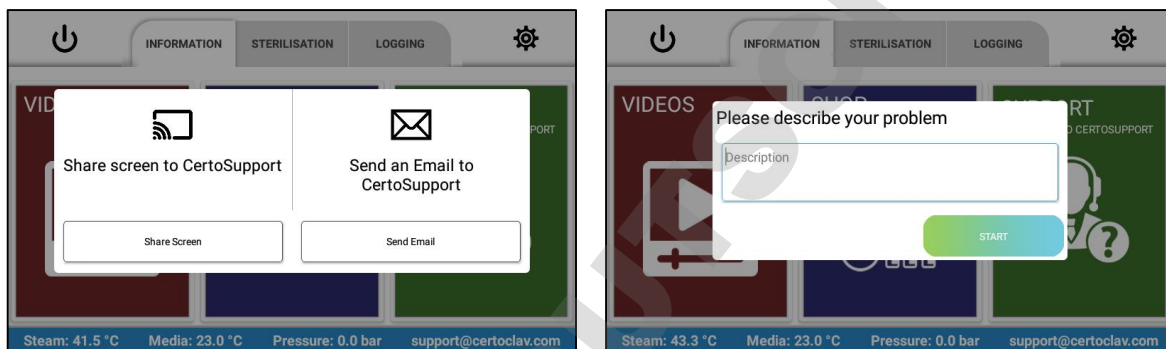


After you have clicked “Start”, wait until CertoSupport joins the TeamViewer session. Once CertoSupport has joined the session, you will be asked for your permission to share the screen. Click “Allow”. Now the technician can access the menu of your device. To cancel the screen sharing, drag the TeamViewer icon to the X icon at the bottom of the screen.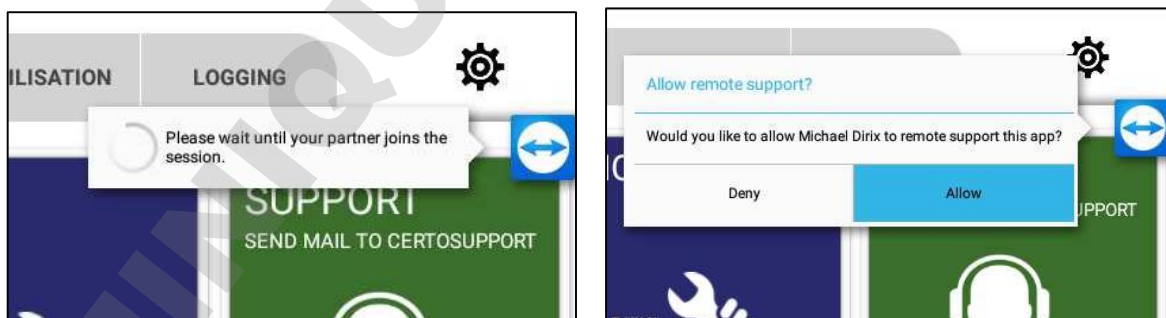

20.3 Logging menu


20.3.1 View logs

To view sterilization protocols, navigate to the “Logging” tab. A list of all recorded protocols is displayed on the left side of the screen. The cloud icon in the protocol list indicates whether the protocols have been backed up to the CertoCloud or to your local CertoServer. Press the drop down menu above the protocol list to sort the protocols by specific criteria. To view the details and graphical view of a protocol, select the protocol by touching the list item. The graphical history of the protocol is now displayed on the right side of the window. To display the protocol as a list, click on the list icon above the graph.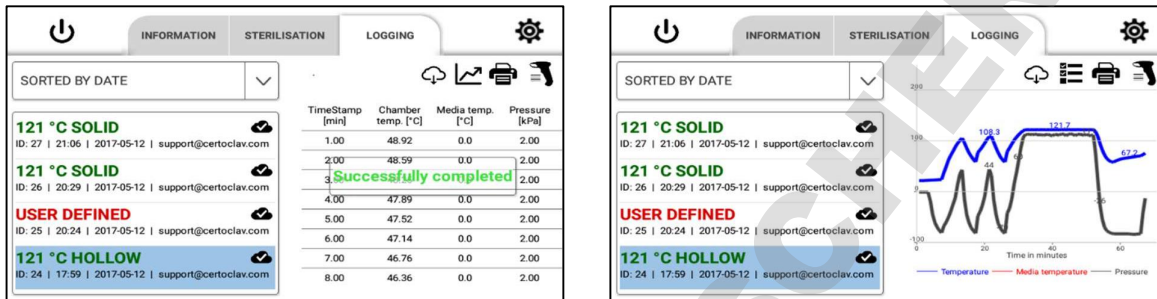

20.3.2 Print out protocols

- 6) Make sure a log printer is connected to the device.
- 7) Navigate to the “Logging” tab
- 8) On the left side of the screen, select the protocol you want to print.
- 9) Click the print icon in the upper right corner of the screen.
- 10) Select the “Print protocol” option

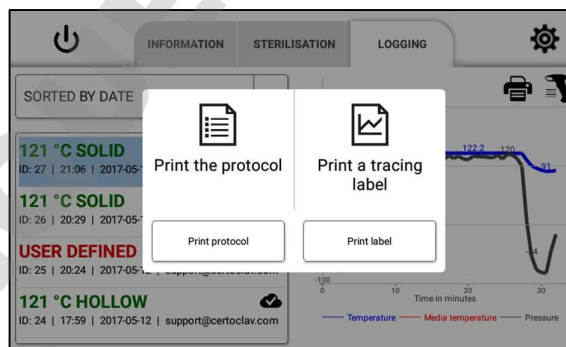20.3.3 Example protocol print out

```

Header line 1
Header line 2
Header line 3

Protocol      CertoClav Vacuum Pro Series
Model:       VAC-PRO-22L
S/N:        22ABCDEFGHJ
Program:     134°C HOLLOW
Program description:
  3 x venting
  134 °C   2.1bar   30min
  Hollow instruments
  wrapped or unwrapped
Media:       NO
Cycle no:   971
Start time: 12.01.15 17:42:00
End time:   12.01.15 18:16:00
Result:     PASSED
hh:mm:ss    Temp °C    Press [bar]
17:42:00    27.5      1.00
17:42:20    27.5      0.50
17:42:40    27.6      0.18
17:43:00    80.4      0.80
[...]
18:15:00    65.9      0.18
18:15:20    65.4      0.18
18:15:40    68.0      0.50
18:16:00    70.2      1.00
Program finished successfully
END

Signed by _____
Confirmed by _____

```

Optional

21 Settings

Press the  icon at the top right of the screen to enter the settings menu.

The following settings can be made here:

21.1 User administration


All registered users are now displayed on the right side of the window. Here you can change the user data and passwords with the “Edit” button and remove a user account from the Touch PC with the “Delete” button.

21.2 Network settings


21.2.1 Connection to a WLAN network

4. Click on the “Network” entry.
5. If the Wi-Fi checkbox is not set, click the empty checkbox to enable Wi-Fi.
6. Then click “Manage WLAN connections” to display a list of available WLAN networks. Select the relevant WLAN network and enter access keys and other parameters if necessary.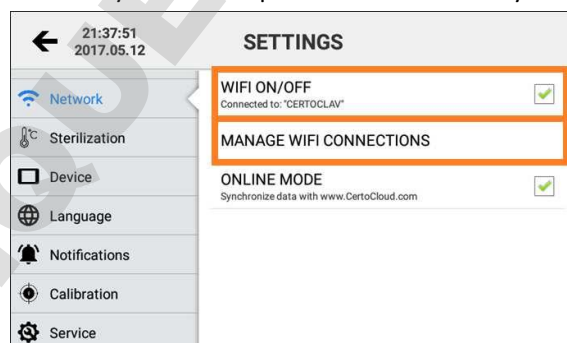

21.2.2 Connection to a LAN network


The autoclave can also be connected to a LAN network via ethernet cable. Connect the network cable to the ethernet port of the interface box.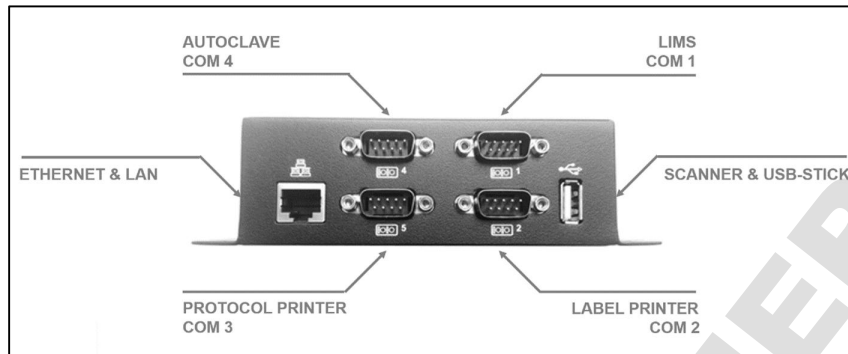

21.2.3 Network Communication



If the autoclave data is to be synchronized with the CertoCloud (www.certocloud.com) or the local CertoServer (<http://certocloud>), the online mode must be activated.

3. Select the "Network" entry on the left side of the settings window.
4. Click the "Online Mode" entry to enable or disable data synchronization with CertoCloud. **Note:** If the online mode is disabled, your protocols are only stored locally on the autoclave. Once you reactivate online mode, all protocols, stored locally up to that point, will be uploaded to CertoCloud and are available online in your backup storage.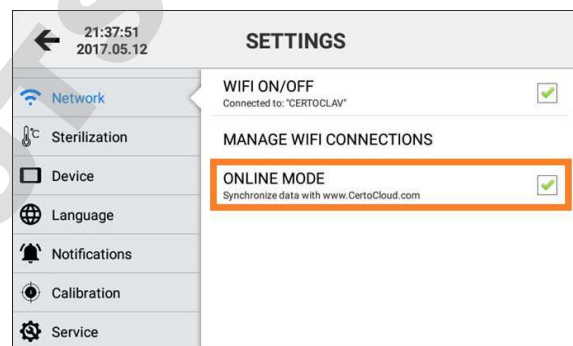


21.2.4 Connection of the autoclave to the local CertoServer (for Connect and VacuumPro)

7. Connect the CertoServer to the Interface-Box or the in-house network.
8. Connect the interface box to the same network as the CertoServer.
9. Navigate to the button: Settings → Network → Select server
10. Select the option "Add server from local network".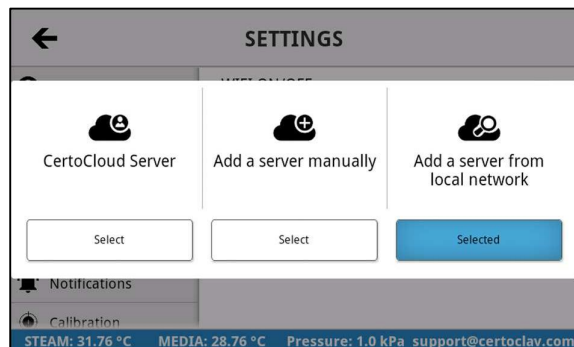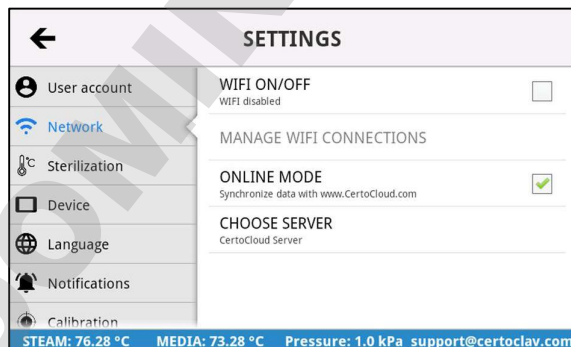

11. The autoclave searches for the CertoServer in the local network.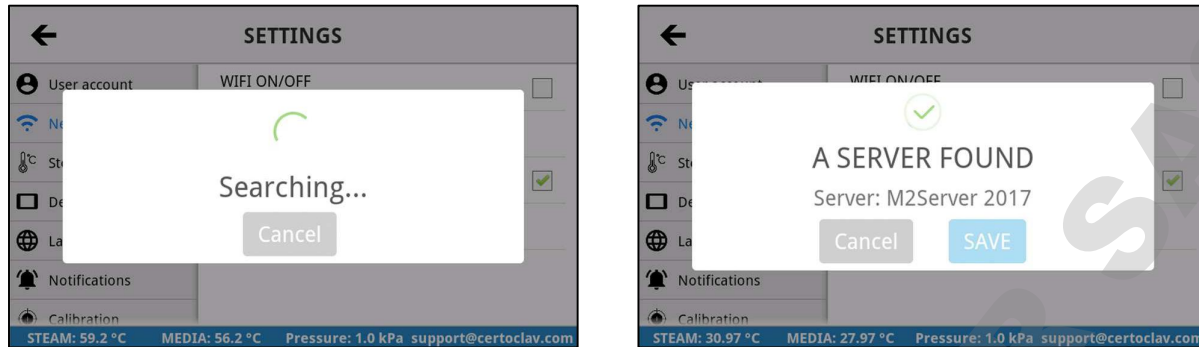


12. After the CertoServer has been found, confirm with the “Save” button.

21.3 Sterilization settings

21.3.1 Export of all protocols to USB flash drive




The touch PC can store up to 1000 protocols in its internal memory. To export all protocols to USB flash drive, insert a USB flash drive at the interface box. Navigate to Settings → Sterilization → Export protocols to USB flash drive.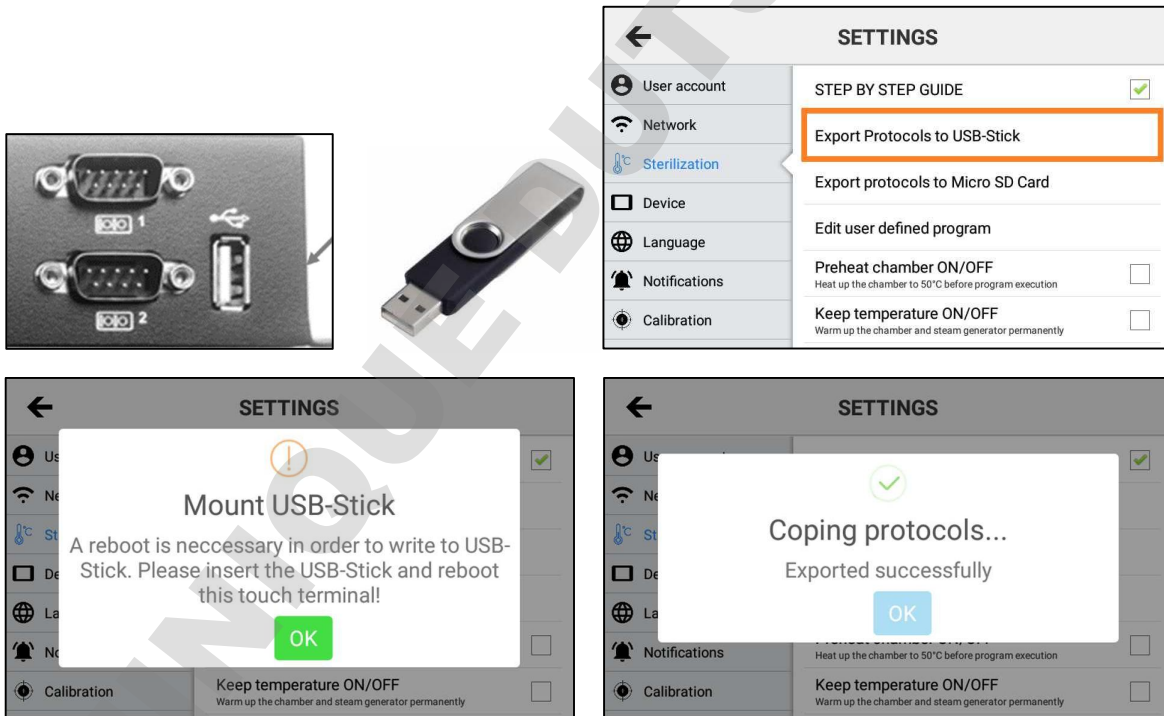

If the protocol export fails, your USB flash drive was not recognized immediately. In this case, leave the USB flash drive inserted in the device and restart the touch PC. Now try to export the protocols again.


21.3.2 Export of all protocols to SD card

4. Insert a micro SD card into the hinged slot on the right side of the touch screen.
5. Press the list entry "Export all protocols to Micro-SD card"
6. If the protocol export fails, restart the touch PC with the SD card inserted. Now try to export the protocols again.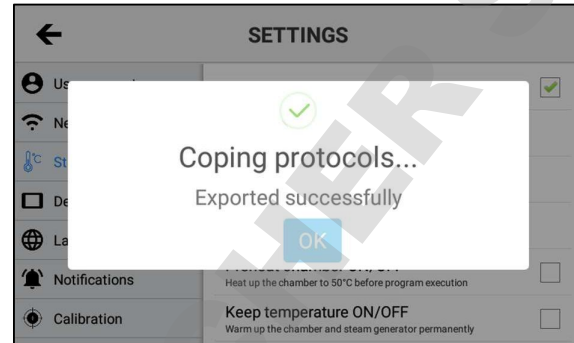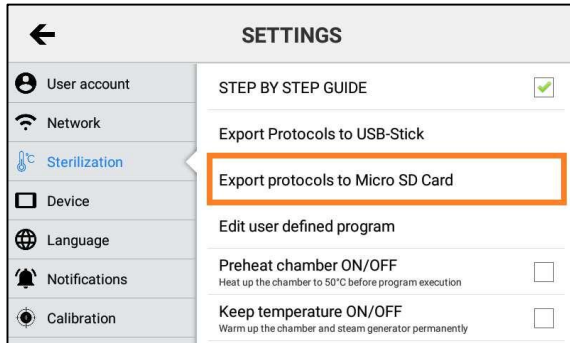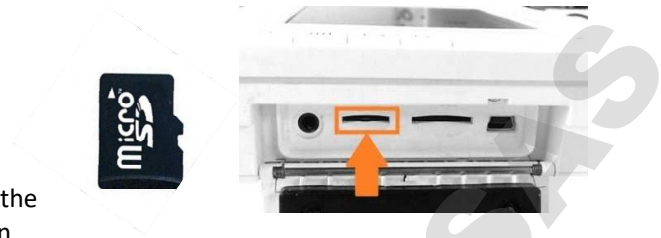



21.3.3 Export of all protocols via network

2. Log into the autoclave with your CertoCloud or CertoServer account. Make sure online mode is activated. If the login is successful, all protocols will be uploaded to CertoCloud. The export was successful, if the Cloud icon is displayed with a check mark.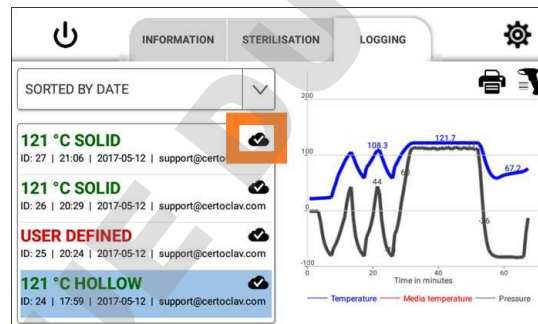

21.3.4 Label Designer (only available for VacuumPro and Connect)


To mark articles with barcodes, you can use the integrated label designer. The label designer can be opened under Settings → Sterilization → Label Designer.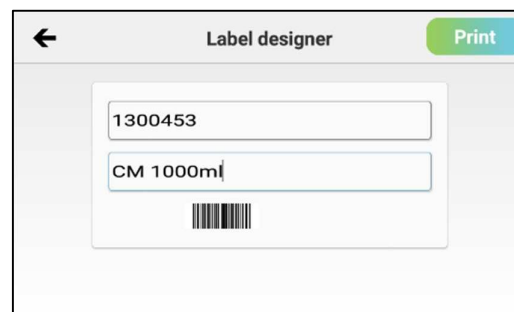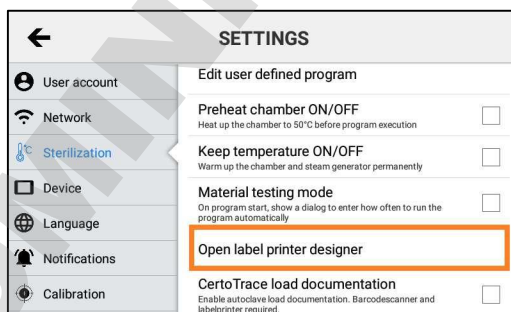

21.4 System settings


21.4.1 Install updates via the internet

5. Make sure the autoclave is connected to the Internet.
6. Click the "Device" entry.
7. In the detail view on the right side, scroll "Check for updates" and click this entry.
8. A progress bar appears with the download. Once all updates have been downloaded, they will be installed automatically. If a new software version is available, an installation wizard will guide you through the installation process.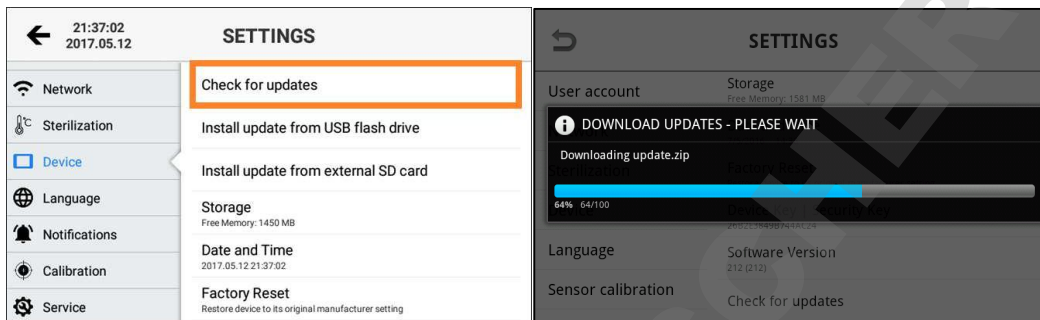


21.4.2 Install updates via USB flash drive

6. Download the latest update from www.certoclav.com or ask for the latest update for your device via e-mail to support@certoclav.com. Please inform the support team about the serial number of the autoclave. The serial number can be found on the type plate of the autoclave or under Settings → Device → Serial number. Copy the "update.zip" to the main directory of a USB flash drive and insert this USB flash drive into the interface box.
7. Click the "Device" entry.
8. In the detailed view on the right-hand side, scroll down to the entry "Installation of updates via USB flash drive" and click this entry.
9. If the USB flash drive is not recognized, restart the touch PC and try again.
10. Follow the instructions on the screen to install the appropriate updates.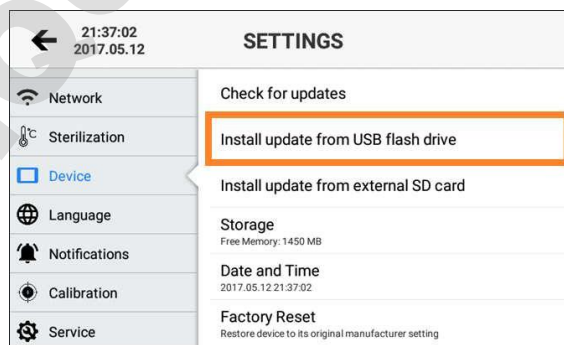

21.4.3 Install updates via SD card

- Download the latest update from www.certoclav.com or ask for the latest update for your device by e-mail to support@certoclav.com. Please inform the support team about the serial number of the autoclave. The serial number can be found on the autoclave's type plate or under Settings → Device → Serial number. Copy the "update.zip" to the root directory of a micro SD card and insert this SD card into the slot on the right side of the touch PC.
- Click the "Device" entry.
- In the detailed view on the right-hand side, scroll "Installing updates via SD card" and click this entry.
- If the SD card is not recognized, restart the touch PC and try again.
- Follow the instructions on the screen to install the appropriate updates.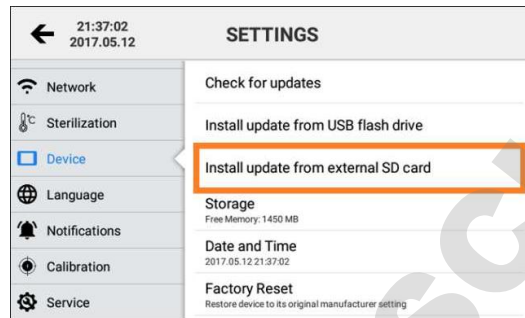

21.4.4 Restore factory settings


When you reset the touch PC to factory defaults, all protocols and user accounts that have not been previously exported and/ or synchronized with CertoCloud are permanently deleted.

- Click the "Device" entry.
- In the detail view on the right side, scroll to "Restore Factory Defaults" and click this entry.
- The touch PC has been reset. All stored user accounts and protocols have been permanently deleted.

21.4.5 Change language

4. In the settings menu, open the language menu by clicking on the “Language” button.
5. Click “current language” in the detail view on the right side of the screen
6. You see a scrollable list of supported languages. Select a language by clicking it.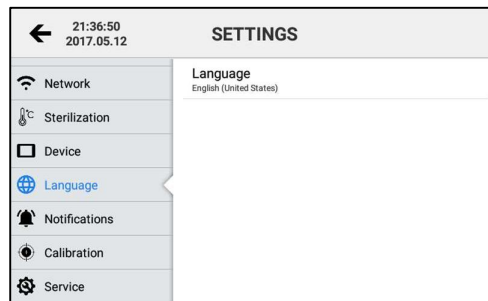

Finally, click “Save” to save the changes.


22 Error messages

If an error occurs during a program cycle, the autoclave emits an acoustic signal and displays an error code on the screen. For detailed error descriptions please reference the main operating instructions of the respective device.

23 Transport and storage

23.1 Storage conditions

The CertoTouch should be stored at a room temperature between 5 °C and a maximum of 40 °C. The relative humidity must not exceed 80%. Make sure that the storage area is well ventilated and does not contain any corrosive gases.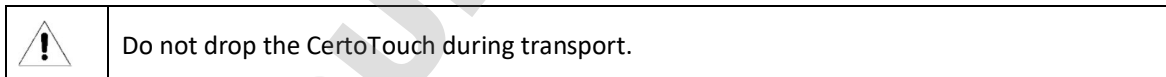

24 Warranty

Warranty: 2 years. The right to free service is forfeited if any of the following conditions apply:

- The damage was caused by incorrect installation and/or operation.
- The damage was caused by negligent behavior.
- The damage was caused by force majeure such as inadmissible supply voltage or fire.