
Cole-Parmer®

BH-250 Series

Analog Block Heaters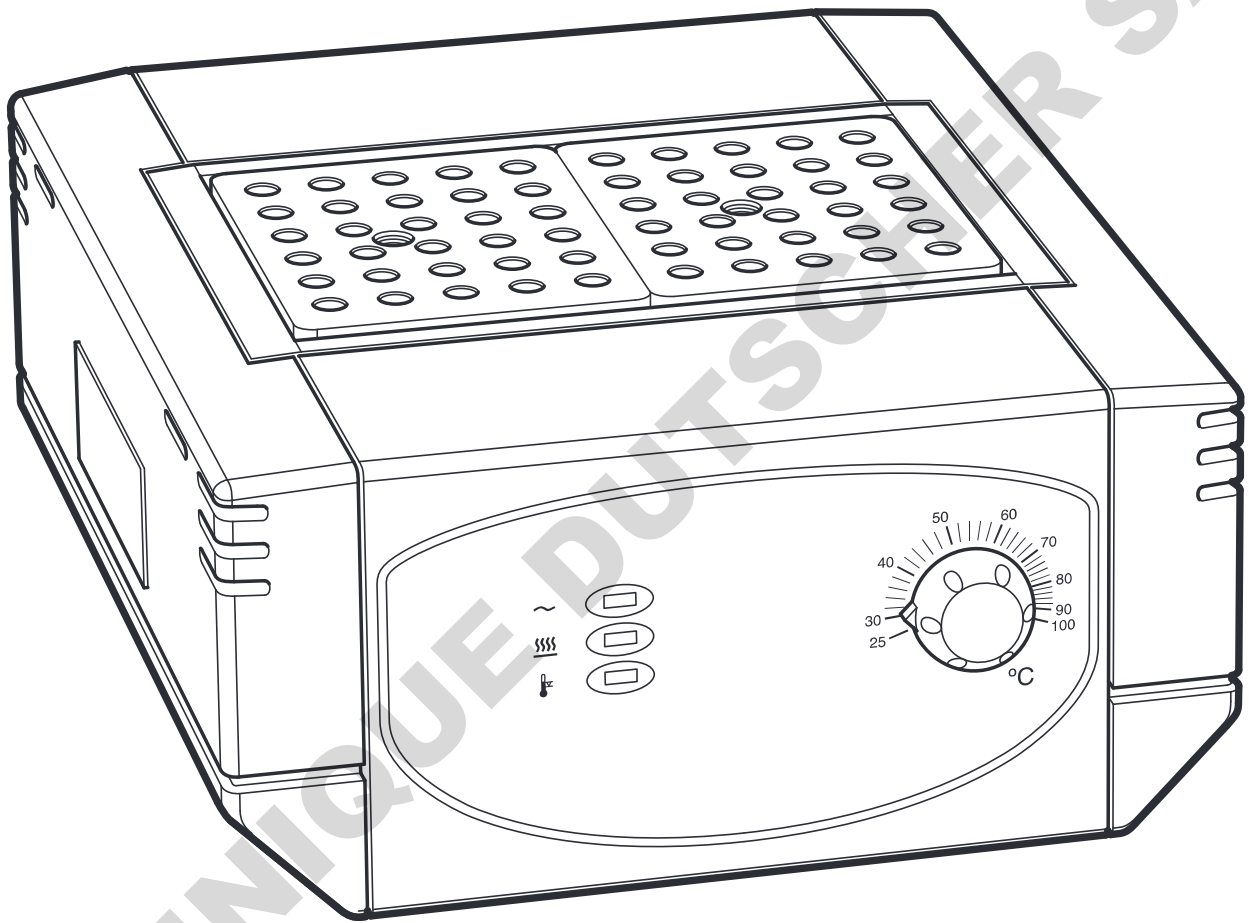

Instruction Manual

Version 1.3

Cole-Parmer®
essentials

CONTENTS**PAGE**

INTRODUCTION	3
BEFORE USE	4
SAFETY and INSTALLATION	5
English	5
Français	7
Deutsch	9
Español	11
Italia	13
SPECIFICATIONS	15
OPERATION	16
Installation	16
The front panel controls	17
Setting the operating temperature	18
After use	18
ADDITIONAL INFORMATION	19
Fault finding	19
Fuses	19
Safety over-temperature cut-out	19
User maintenance	19
Technical support and servicing	20
BLOCK INSERTS	21
REPLACEMENT PARTS	22
DECLARATION OF CONFORMITY	23

INTRODUCTION

Please read all the information in this booklet before using the unit.

Antylia Scientific Block Heaters provide a safe, dry, constant temperature source in the laboratory. The units can be used for incubation, boiling, inactivation, wet ashing, sample concentration, enzyme analysis and many other clinical and industrial purposes.

The units cover the temperature range from ambient +25°C to 100°C or 200°C (depending on model) using machined aluminium alloy blocks as the heat transfer medium. Independent indicator lights indicate mains power, heater action and over-temperature cut-out. An ON/OFF switch is mounted on the rear panel, together with the mains connector and a double fuse holder.

The units are constructed of polyphenylene sulphide (PPS) plastic, which is strong yet lightweight and can resist sustained temperatures as high as 260°C. Even though the units heat up rapidly, venting and highly efficient insulation ensure that the cases remain cool enough to handle even at maximum operating temperatures.

Each unit is supplied with an extraction tool for easy removal of the insert blocks.

The operating instructions in this Operator's Manual cover the following range of digital Blockheaters:

Model number	Description	Voltage, frequency
BH-250-2	Block Heater ambient to 100°C, accepts 2 insert blocks	230V, 50-60Hz
BH-250-2-115	Block Heater ambient to 100°C, accepts 2 insert blocks	115V, 50-60Hz
BH-250-3-HT	Block Heater ambient to 200°C, accepts 3 insert blocks	230V, 50-60Hz
BH-250-3-HT-115	Block Heater ambient to 200°C, accepts 3 insert blocks	115V, 50-60Hz

The features of individual models are outlined in the Specifications section of this manual.

Note: This manual covers the analog Block Heater range. The digital Block Heaters BH-250D-2, BH-250D-3, BH-250D-4, BH-250DC-2, BH-250D-2-HT and BH-250D-3-HT are covered in a separate user manual.

BEFORE USE

Before using the Block Heater please make sure you have read this manual carefully. If there is any doubt relating to the proper use of this equipment, the staff at Antylia Scientific Ltd. or your supplier will be happy to assist you.

UNPACKING

When unpacking the unit please ensure that the following have been removed from the packaging:


Block Heater

6101308 Extraction tool


Mains cables (UK and EU or US)

Operator's manual

The user is advised to keep the original packaging in case the instrument ever needs to be returned for service or repair. Antylia Scientific Ltd. accepts no responsibility for damage incurred unless the unit is correctly packed and transported in its original packaging.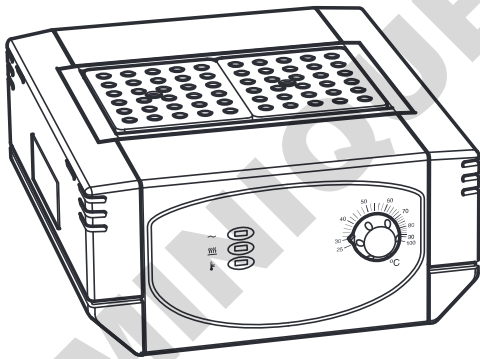

BH-250-2 (230V)
BH-250-2-115 (115V)

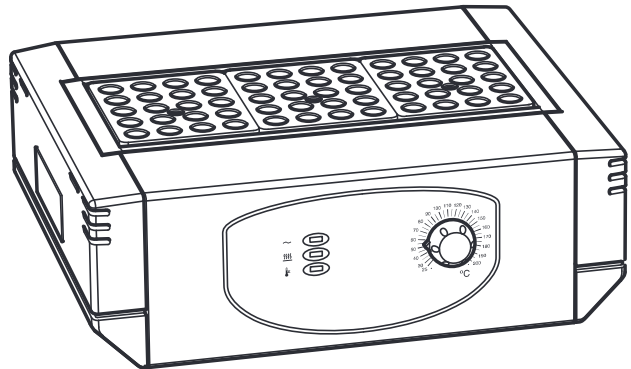BH-250-3-HT (230V)
BH-250-3-HT-115 (115V)

The Block Heater range

SAFETY AND INSTALLATION

Please read all the information in this manual before using the Block Heater.

WARNING

HIGH TEMPERATURES ARE DANGEROUS: they can cause serious burns to operators and ignite combustible material.

Antylia Scientific have taken great care in the design of these units to protect operators from hazards, but operators should pay attention to the following points:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS.
- DO NOT use combustible substances near hot objects.
- DO NOT operate the instrument in the vicinity of inflammable liquids or gases.
- DO NOT place any liquid directly into the instrument.

OPERATOR SAFETY

All operators of Antylia Scientific equipment must have available the relevant literature needed to ensure their safety. It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures. If the equipment is used in a manner not specified by Antylia Scientific Ltd. the protection provided by the equipment to the operator may be impaired.

All Antylia Scientific instruments are designed to conform to international safety requirements and are fitted with an over-temperature cut-out.

If a safety problem should be encountered, switch off the unit at the mains socket and remove the plug from the electricity supply.

INSTALLATION

The instrument should be carried using both hands. Never move or carry the instrument when in use or connected to the mains electricity supply.

1. All Antylia Scientific instruments are supplied with a power cable; this may be integral or plug-in.

2. Before connecting the instrument to the mains electricity supply, check the voltage against the rating plate (located on the back of the unit). **Please note that the unit must be earthed to ensure proper electrical safety.** Connect the mains cable to a suitable plug according to the table below.

<i>Connections</i>	<i>230V, 50/60Hz</i>	<i>115V, 50/60Hz</i>
Live	Brown	Black
Neutral	Blue	White
Earth	Green/yellow	Green

Note that units marked 230V on the rating plate work at 207V; units marked 115V work at 104V. In both cases, however, the heating rate will degrade by approximately 8%.

3. Place the unit on a suitable flat bench or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction.
4. Plug the mains cable into the socket on the back of the instrument.
5. Switch on the instrument. Symbols on or near the power switch of the unit have the following meanings:
 - I Mains Switch On
 - O Mains Switch Off

Replacement cable

Should the mains lead need replacement, a cable of 1mm² of harmonized code H05VV-F connected to an IEC320 plug should be used. **IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN.**

ENVIRONMENTAL CONDITIONS

The Block Heater is designed operate under the following conditions:

- Indoor use
- Ambient temperature range +5°C to +40°C
- Altitude to 2000m
- Relative humidity up to 80%
- Mains supply fluctuations not exceeding 10%
- Over voltage category II IEC 60364-4-443
- Pollution degree 2
- Use with a minimum distance all-around of 200mm from walls or other items

Note: The control specifications are quoted at an ambient temperature of 20°C. The specification may deteriorate outside an ambient temperature of between 10°C and 30°C.

The instrument has been tested for radio frequency interference and is certified under EN61326.

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
SPECIFICATIONS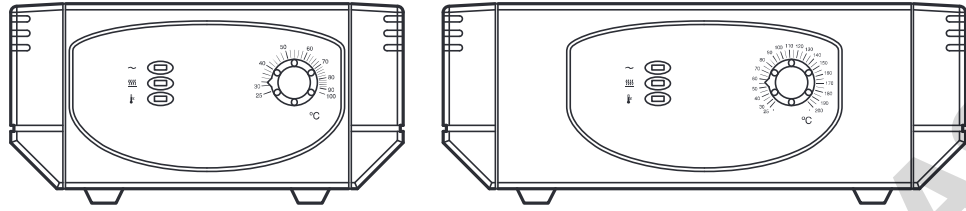

Specification

	BH-250-2	BH-250-3-HT
Blocks	2	3
Temperature Setting	Rotary dial	Rotary dial
Working range	Ambient +5°C to 100°C	Ambient +5°C to 200°C
Temperature Scale	2°C graduation	2°C graduation
Set point accuracy	±2°C	±2°C
Stability @ 40°C	±0.05°C	±0.1°C
Stability @ 100°C	±0.1°C	±0.15°C
Max variation between blocks	0.2°C at 40°C	0.2°C at 40°C
Heat up 30-37°C	8 minutes	8 minutes
Heat up 30-56°C	12 minutes	30 minutes
Dimensions (WDH)	202 x 260 x 105mm	279 x 260 x 105mm
Electrical supply	230V or 115V 50-60Hz	230V or 115V 50-60Hz
Power	300W	450W
Shipping weight	4kg	6kg
Warranty	3 years	3 years

Note: The control specifications are quoted for an ambient temperature of 20°C for units fitted with 2, 3 or 4 insert blocks type F3506. The specification will be closely held if the ambient temperature is in the range 10°C to 30°C. Outside this range the quoted figures may deteriorate but the unit will still work safely.


Note: Different block inserts may take different amounts of time to equilibrate to the same temperature.

OPERATION

INSTALLATION

See also the sections on Safety and Installation.


1. Place the unit on a suitable flat bench or in a fume cupboard if required, ensuring that the air inlet vents on the underside are free from obstruction.
2. Select the insert aluminium alloy block(s) appropriate for the application. A list of available blocks is given in the section 'Accessories'. Ensure that both the underside of the block(s) and the top of the hotplate are clean; efficient heat conduction between these two surfaces is essential.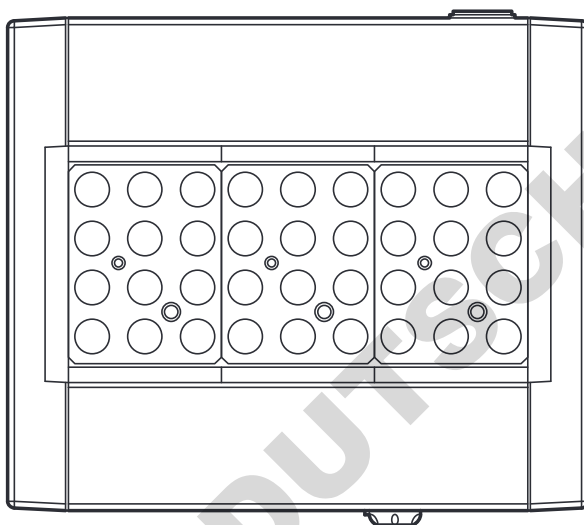

Top view showing three insert blocks in place

3. Place the insert blocks onto the hotplate in the well of the unit and place the tubes containing the sample liquid in the blocks. **Note:** Aluminium block inserts are designed to allow for thermal expansion up to 200°C. For this reason there will be a small air space around the blocks at ambient temperature. The heat source is the base of the Block Heater. For best results push all the block inserts to the right hand side so they are touching.
4. The heater design, temperature sensor and control circuit give good temperature control and uniformity, but it is important that there is a close fit of the tubes in the block to allow efficient heat transfer.
5. If less than the maximum number of insert blocks are fitted it is recommended that plastic half blocks (part code F4466) are fitted to the space either side of the aluminium block. The half blocks have the effect of reducing heat loss, thus improving temperature stability.

6. Plug the mains cable into the socket in the rear of the unit. Connect to the mains electricity supply with the plug provided or one wired correctly for your supply. Switch the power ON from the switch located next to the mains input lead. The front display will then light up.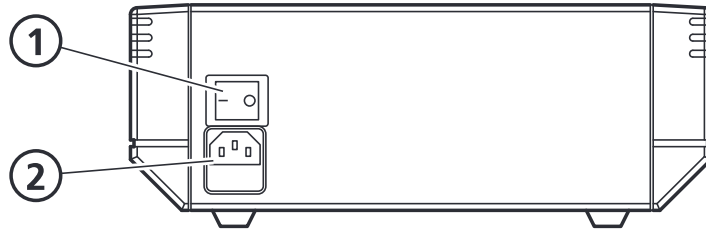


Rear view

1. Power ON/OFF switch.
2. Mains cable socket.

THE FRONT PANEL CONTROLS

The front panel controls consist of three indicators and a rotary dial for controlling the temperature.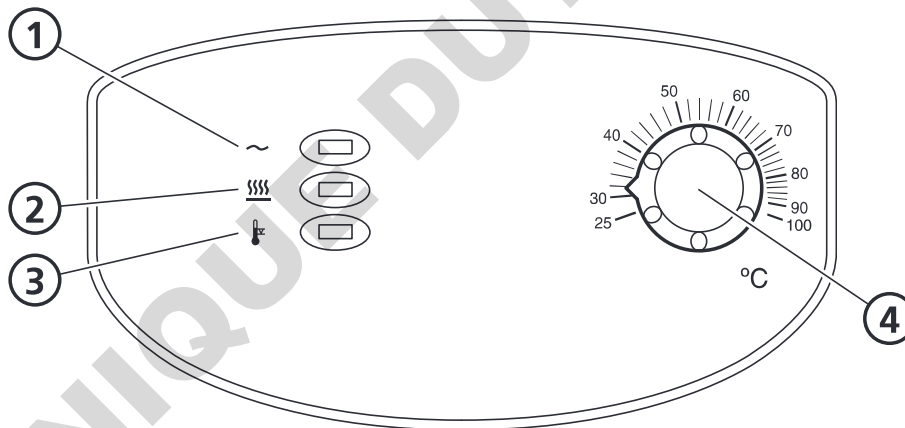

The front panel controls

1. Power indicator. When lit, this shows that there is power to the unit.
2. Heater indicator. When this is flashing, it indicates that the heater is heating the unit or controlling the temperature. If the Block is above the set temperature then the indicator will be off.
3. Over-temperature indicator. If this is lit, it indicates that the unit has developed a fault and the over-temperature cut-out has been triggered. This will cause the heater to be switched off and the unit will begin to cool even if the heater light is on. The unit should be returned to Antylia Scientific or its authorised agent for investigation.
4. Temperature rotary dial. Turn the dial to the desired temperature. BH-250-2 displays to 100°C, BH-250-3-HT displays to 200°C

SETTING THE OPERATING TEMPERATURE

1. Turn the temperature rotary dial to the required temperature. Note that the actual temperature of the block critically depends on the type of block used and the quality of the thermal contact between the base of the block and the unit.
2. The heater (and heater indicator) comes on if the set temperature is higher than the current block temperature.
3. When the set temperature is approached, the heater indicator will begin to flash and when the temperature is stabilised the indicator will flash slowly.
4. Due to variations in heat losses with different designs of insert block, the actual temperature may vary. The units are calibrated using model F3506. Running a unit with only one insert block and without the plastic half blocks will reduce the maximum achievable temperature by a few degrees.


If you need to control the temperature to a greater accuracy than the instrument's set point, place a thermometer in the special hole in one of the blocks. For greater accuracy still, you can place the thermometer in the sample liquid in one of the test tubes. In either of these cases it may be necessary to re-adjust the temperature control knob or preset adjustment to achieve the precise temperature required. Allow the temperature of the unit to stabilise after each adjustment.

5. There will be a time lag between the heater platten and the insert block achieving the set temperature due to thermal contact between them. Guide times are shown in the specification tables.

AFTER USE

When you have finished heating samples, remember that parts of the unit – the tubes, blocks and associated accessories – may be very hot. Take the precautions listed in the section Safety and Installation. We recommend that the blocks should be allowed to cool to 70°C or below before being removed from the heating unit.

To remove an insert block, screw the extractor tool into the threaded hole and lift the block out vertically. Never leave the extractor tool in the block while it is being used in the Block Heater.

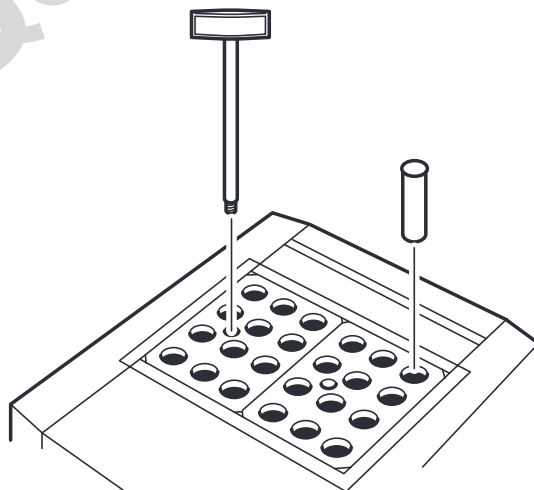Removing an insert block with the extractor tool

ADDITIONAL INFORMATION

Note that this equipment should only be dismantled by properly trained personnel. Removing the outer cover exposes potentially lethal mains voltages. There are no user serviceable parts within this equipment.

FAULT FINDING

Should you have any problems with your Block Heater which cannot be easily remedied, you should contact your supplier and return the unit if necessary. Please include details of the fault observed and remember to return the unit in its original packing. Antylia Scientific Ltd. accepts no responsibility for damage to units which are not properly packed for shipping: if in doubt, contact your supplier, giving the full serial number of the unit.

FUSES

If the display on the front panel is not lit, one of the two fuses may have blown. Check that there is no external cause, such as a faulty plug or lead. Check both fuses and replace the faulty fuse with a new one of the correct value (fuse values are given on the label next to the power inlet). Note that fuses should only be replaced by a qualified electrician.

The holder for the two fuses is built into the mains input socket. First remove the power cable and then gently prise the fuse drawer open with a flat-bladed screwdriver or similar tool. Each fuse can be removed by using the screwdriver as a lever.

Exchange the faulty fuse in the fuse holder for a working fuse of the correct value. Finally, replace the fuse drawer in the fuse compartment and push the drawer shut. Fuses which blow repeatedly indicate a serious fault and you should return the unit to your supplier for repair.

Model	Fuses for 230V	Fuses for 115V
BH-250-2	2 x F2 Amp	2 x F4 Amp
BH-250-3-HT	2 x F4 Amp	2 x F6.30 Amp

SAFETY OVER-TEMPERATURE CUT-OUT

In the event of a fault occurring which could cause the heater to over-heat, a thermal fuse will cut in and remove power from the heater. This will occur when the temperature reaches approximately 110°C in the BH-250-2 model and 210°C in the BH-250-3-HT model.

USER MAINTENANCE

Before cleaning your unit, disconnect it from the power supply and allow to cool to below 50°C. The outer case of the Block Heater may be cleaned with a cloth dipped in water or ethanol (methanol can also be used). No part of the case or cover should be immersed in the solvents. Do not use aggressive solvents such as acetone or abrasive cleaners. Do not use bleach or strong acids.

Before using any cleaning or decontamination method except those recommended here, the responsible person should check with Antylia Scientific that the proposed method will not damage the equipment.

TECHNICAL SUPPORT AND SERVICING

If you require further technical or application assistance please contact Antylia Scientific Ltd at:

E-mail: cptechsupport@antylia.com

Phone: +44 (0)1785 810433

For servicing information please contact:

Antylia Scientific Ltd.

Service Department,

Beacon Road,

Stone,

Staffordshire,

ST15 0SA,

E-mail: cpervice@antylia.com.

Phone: +44 (0)1785 810475

We are continually striving to improve our products and software. If you have any comments and suggestions on how we can do things better please send them to us at: cptechsupport@antylia.com

DOMINIQUE DUTSCHER SAS

HEATER BLOCK INSERTS

Insert blocks are made of aluminium alloy and must be ordered separately from the heater units.

Part code	Nominal tube size	Number of holes	Hole diameter (mm)	Hole depth (mm)	Base profile	Block dimensions (mm)
F3501	Plain block (single)	None	-	-	-	95 x 75 x 50
F3502	6mm	30	6.5	38	Conical	95 x 75 x 50
F3503	10mm	20	10.5	48	Conical	95 x 75 x 50
F3504	12mm	20	12.75	48	Conical	95 x 75 x 50
F3505	13mm	20	13.8	48	Conical	95 x 75 x 50
F3506	15mm	12	15.75	48	Conical	95 x 75 x 50
F3507	16mm and 15ml Falcon tubes	12	16.75	48	Conical	95 x 75 x 50
F3508	19mm	8	19.75	48	Conical	95 x 75 x 50
F3509	25mm	6	25.75	48	Conical	95 x 75 x 50
F3510	10mm Cuvettes	2 channels	13 x 95	38	Flat	95 x 75 x 50
F4460	Plain block (single)	Thermometer only	-	-	-	95 x 75 x 50
F4461	7mm and 9mm	20 + 10	7.0 and 9.5	38	Conical	95 x 75 x 50
F4462	24mm	6	24.75	48	Conical	95 x 75 x 50
F4463	26mm	6	26.75	48	Conical	95 x 75 x 50
F4464	1.5ml Microfuge tubes	20	10.9	35	Tapered	95 x 75 x 50
F4465	0.5ml Microfuge tubes	30	7.9	27	Tapered	95 x 75 x 50
F4466	Half block plastic spacers	-	-	-	-	98.8 x 40.2 x 18.5
F4470	2.0ml Microfuge tubes	20	10.8	35	Conical	95 x 75 x 50
F4471	0.2ml Microfuge tubes	72	4.4	13.8	Tapered	95 x 75 x 50

REPLACEMENT PARTS

The following replacement parts can be obtained from Antylia Scientific Ltd. or your Antylia Scientific dealer:

Product code	Description
6101308	Extraction tool
HH179(S)	UK mains lead with plug, 230V
HH180(S)	European mains lead with plug, 230V
FCABLEUS	US mains lead with plug, 115V
F4466	Plastic half block
6101325	Screw driver for presets

DOMINIQUE DUTSCHER SAS

CE UK CA This product meets the applicable CE Directives and UKCA Legislation for radio frequency interference and may be expected not to interfere with, or be affected by, other equipment with similar qualifications. We cannot be sure that other equipment used in its vicinity will meet these standards and so we cannot guarantee

that interference will not occur in practise. Where there is a possibility that injury, damage or loss might occur if equipment malfunctions due to radio frequency interference, or for general advise before use, contact the manufacturer.

Declaration of Conformity is available to view online at www.coleparmer.com

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T: +1 847 549 7600

Cole-Parmer®
essentials

Antylia Scientific Ltd.
Beacon Road,
Stone,
Staffordshire,
ST15 0SA,
United Kingdom