Temperature

Controlled Shaking

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

Operating Manual

Page 2

Overhead Shaker Hei-MIX Reax 2

Operating manual must be read before initial start-up.

Please follow the safety instructions provided.

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About this Document	4
Version and variants	4
About this manual	4
Symbols and keywords.....	5
Safety Instructions	6
General safety instructions.....	6
EU Declaration of Conformity.....	6
Intended use.....	6
Installation / electrical safety	7
Qualifications of personnel	7
Obligations of the operator	7
Installation site	7
Modifications to the device.....	8
Safety of personnel.....	8
Safety during use	9
Disposal	9
Device Description	10
Device overview	10
Overall view.....	10
Control panel	10
Start-Up	11
Setting up device	11
Switch device on and off.....	11
Connect/disconnect power cord	11
Inserting Adaptor	12
Operation	13
Installing vessels.....	13
Installing vessels: Universal adaptor	13
Installing vessels: Adaptor 1	13
Starting rotation.....	14
Cleaning and Maintenance	16
Cleaning.....	16
Maintenance	16
Troubleshooting	17
Possible errors and how to resolve	17

Assembly	18
Electrical connection	18
Disassembly and Storage	19
Dismantling, transportation and storage	19
Dismantling	19
Transportation and storage	19
Accessories and Spare Parts	20
Scope of delivery	20
Accessories	20
Attachments	20
Technical data	20
Service	21
Contact / Technical Service	21
Warranty	22
Confirmation of condition	23

DOMINIQUE DOTSCHER SAS

Copyright

This operating manual is protected by copyright and is exclusively intended for the internal use of the purchaser.

Handing over this operating manual to third parties, reproducing it in any type or form – in part or in full – and utilizing and/or communicating its contents is prohibited without prior written approval from the manufacturer.

Any violations will result in an obligation to provide compensation for damages. Other claims remain unaffected.

➤ Version and variants

Version

This manual describes the function, operation and maintenance of the Hei-MIX Reax 2 overhead shaker.

Version	Alteration date
1.2	08/2018

Variants

The device is available in different variants. Certain features or functions are only available in specific product variants. These variants are described in this operating manual.

➤ About this manual

This operating manual is an integral part of the device described here.

- Carefully read through this operating manual and observe all of the safety instructions and warnings notices.
- Ensure that every user has carefully read the operating manual before using the device for the first time.
- Store this operating manual so that it is accessible to all users at all times.
- Pass the operating manual on to any subsequent owner.

The current version of this operating manual can also be found at:

[https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers \(Reax 2\)](https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers (Reax 2))

In addition, also observe the local regulations at the installation site.

➤ Symbols and keywords

Standardized symbols and signal words are used in this operating manual to warn against any dangers and provide important instructions. These instructions must be strictly observed to avoid accidents and damage.

The following terms and basic symbols are used:

Symbol	Additional signal words / explanation
Warning signs 	The yellow triangle indicates hazardous situations. It is used in combination with the following signal words: DANGER: Indicates a hazardous situation which, if not avoided, will result in serious injury or death. WARNING: Indicates a hazardous situation which, if not avoided, may result in serious injury or death. CAUTION: Indicates a hazardous situation which, if not avoided, may result in property damage and minor or moderate injury.
Prohibitory signs 	Prohibited: The red circle indicates a situation that should be avoided under all circumstances and which, if not avoided, may result in serious injury or death.
Mandatory signs 	Must be observed: The blue circle indicates important information. Please observe this information to avoid any property damage.

Other symbols used:

Symbol	Description
→	Handling instruction, action required
✓	Result of action
▪	List of information
▪	
▪	
a.	List of variants
b.	
c.	

➤ General safety instructions

EU Declaration of Conformity

This device complies with the following EC-Directives*:

- 2006/42/EC Machinery Directive
- 2014/30/EU Electromagnetic Compatibility Directive

*Also see the attachment "EU Declaration of Conformity".

The device has been manufactured according to state-of-the-art technology and in compliance with recognized safety regulations. However, risks may still arise during installation, operation and maintenance.

→ Please ensure that the operating manual is available at all times.

The device may only be used under the following conditions:

- Only operate the device if it is in a technically perfect condition.
- Only operate the device if you are properly aware of the risks and required safety measures.
- Only operate the device in accordance with the instructions given in this operating manual.
- If any information is ambiguous or missing, ask your superior or contact the manufacturer.
- Do not operate anything on the device without authorization.
- Only use the device in accordance with its intended use.

Intended use

The device may only be operated by authorized personnel.

The device is suitable for the following uses:

- Shaking
- Mixing
- Emulsifying
- Suspending
- Separating

The device may be operated in research laboratories, other laboratories and production facilities in the following sectors:

- Chemistry
- Pharmacy
- Biology
- Environmental analytics
- Basic research
- Similar research laboratories

Reasonable foreseeable misuse

Any use which deviates from the device's intended use is considered to be improper. The manufacturer is not liable for damage that occurs as a result. The risk is borne by the operator alone.

Installation / electrical safety

- The device may only be connected up if the available power supply voltage corresponds to the information stated on the rating plate for the device.
- The power supply connection must be easily accessible at all times.
- An RCD circuit breaker (residual current circuit breaker) must be fitted to the electrical system in the building so that the device is disconnected from the power circuit in the event of a fault.
- Repairs may only be performed by electricians authorised by Heidolph Instruments.
- Never operate the device with a damaged power cord.
- Always turn the device OFF and disconnect the power cord before carrying out any maintenance or repair work.

Qualifications of personnel

- The device may only be operated by trained personnel.
- The device may only be operated by persons who have been instructed and supervised in its proper use by trained specialist personnel.
- The device may only be operated by specialist personnel who are above the legal minimum age.
- Other persons may only work on the device under the constant supervision of experienced and trained specialist personnel.
- This operating manual must be read and understood by all persons working with the device.
- Personnel must receive safety training that ensures responsible and safe working practices.

➤ Obligations of the operator

Installation site

- The device must be installed in a suitable location.
- The device must be installed on a firm stable and temperature resistant surface.
- Ensure that the device and all of its components are easily accessible at all times.
- Maintain a sufficient safety clearance to, in particular, any moving and/or hot equipment components.
- It is not permitted to place or store any objects such as accessories, tools or chemicals within this clearance area during operation.
- All screw connections must be securely tightened.
- It is not permitted to operate the device near to highly flammable or explosive substances.
- Operate the device in conjunction with an extractor hood, if working with potential harmful media (see DIN EN 14175 and DIN 12924).

Safety Instructions

- The device is designed for indoor use only and under the following ambient conditions:

Ambient temperature	5 - 31 °C at 80 % relative humidity
	32 - 40 °C decreasing linearly to a maximum 50 % relative humidity
Installation altitude	0 - 2,000 m above sea level
Contamination level	2
Overvoltage category	II
Permissible supply deviations	± 10 %

- If the device is operated in corrosive atmospheres, the service life of the device will decrease based on the concentration, duration and frequency of the exposure to the corrosive atmosphere e.g. concentrated hydrochloric acid (HCl).

Modifications to the device

- It is not permitted to make any unauthorised modifications or changes to the device.
- Do not attach or install any parts that have not been approved by the manufacturer.
- Unauthorized modifications or changes will void the EC Declaration of Conformity for the device and operation of the device will no longer be permitted.
- The manufacturer is not liable for any damage, dangers or injuries that result from unauthorized modifications and changes or due to the non-observance of the instructions in this manual.

Safety of personnel

- Ensure that the device is only operated by qualified specialist personnel and trained employees.
- Observe the following instructions to avoid any personal injuries and property damage:
 - Laboratory regulations
 - Accident prevention regulations
 - Hazardous Substance Act
 - Other generally accepted occupational health and safety regulations
 - Local regulations

Safety Instructions

➤ Safety during use

- Wear the appropriate clothing when working on the device (protective clothing, protective glasses and, if necessary, safety gloves).
- Do not use the device in potentially explosive areas. The device is not protected against explosion. There is no explosion or ATEX protection available.
- Do not carry out work with naked flames in the vicinity of the device (risk of explosion).
- Do not operate or assemble devices in the vicinity which are emission or radiation sources (electromagnetic waves) for the frequency range ($3 \cdot 10^{11}$ Hz to $3 \cdot 10^{15}$ Hz).
- Do not operate or assemble appliances in the vicinity of the device which constitute emission or radiation sources for ionizing radiation or in the ultrasonic range.
- Do not operate the device where adiabatic compression or shock waves might occur (shock wave combustion).
- Do not use substances, where the energy input due to mixing might pose a hazard.
- Do not spill liquids over the device or any parts of it.
- Remove any accidentally spilled liquids immediately.
- Eliminate errors immediately.
- Always switch the device OFF after use.
- Do not use abrasive material to clean the surface. Only wipe with damp cloth.

➤ Disposal

- Check the device components for hazardous substances and solvents.
- Clean all components before disposal.
- Dispose of the device in accordance with the relevant national regulations.
- Dispose of the packaging material in accordance with the appropriate national regulations.

➤ Device overview

Overall view

Control panel

➤ Setting up device

Warning: **Risk of slipping device!**

As a result of vibrations generated during operation the device might slide off the table top.

The surface must be smooth and clean, should be properly leveled and has to feature an anti-skid surface coat.

- Locate the shaker on a stable, horizontal surface.
- Clean surface and feet with a damp cloth and ethanol regularly.

➤ Switch device on and off

The ON/OFF switch is situated on the front panel.

Switch device on

- Press the ON/OFF switch on the left side.
- ✓ The device is switched on.

ON/OFF switch: device switched on

Switch device off

- Press the ON/OFF switch on the right side.
- ✓ The device is switched off.

ON/OFF switch: device switched off

➤ Connect/disconnect power cord

The three-wire socket is firmly connected to the device.

The connection is at the rear side of the device.

- ✓ The toggle switch ON/OFF is in the "Off" position.
- Connect the power cord with a grounded socket.
- Disconnect the plug if the device is not in use.

➤ Inserting Adaptor

There are two adaptors available to meet with a wide range of applications:

- „Universal adaptor“ for any kind of vessels or racks
- „Adaptor 1“ for test tubes

A square bolt is situated on one side of the adaptor.

On the opposite you will find a rounded pin.

Universal adaptor

Square bolt

Rounded pin

Adaptor 1

Square bolt

Rounded pin

→ Insert the square bolt into the carrier at the right side of the housing.

→ Draw the knurled grip.

✓ The rounded pin is placed in the receptacle.

→ Release knurled grip.

Knurled grip

Receptacle

➤ Installing vessels

Installing vessels: Universal adaptor

The universal adaptor allows for fixing vessels with ground-in stoppers or screwed caps of different sizes as well as complete test tube racks. Simultaneously fixed vessel must be of the same height.

Dimensions of vessels suitable for usage:

- height min./max. 45 mm/160 mm
- width max. 230 mm

→ Loosen the two quick-clamping nuts by turning them once or twice.

✓ When they are in an inclined position, the quick-clamping nuts can be moved along the threaded bolt.

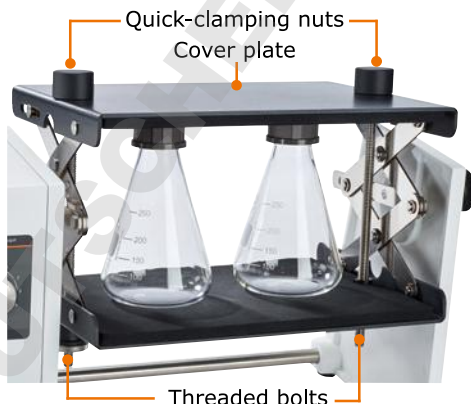
→ Lift the cover plate level so not to jam it.

✓ The universal adaptor can be opened to a width from 45-160 mm, depending on the case of application.

→ Place vessels or the test tube rack (see underneath) into the adaptor, as shown in the illustration.

→ Align the cover plate parallel to the ground plate, press it down and hold pressed.

→ Move the quick-clamping nuts along the threaded bolt and tighten the cover plate by turning them once or twice.

Installing vessels: Adaptor 1

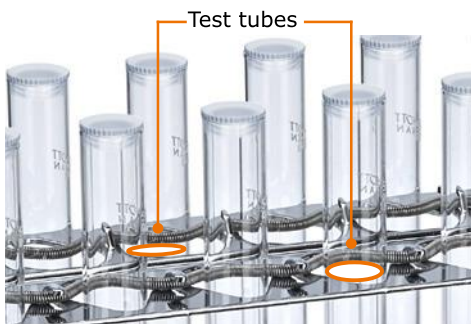
Adaptor 1 is a test tube rack, which is to be fixed into the universal adaptor. It can hold up to 20 test tubes with diameters from 10 -18 mm.

A tension ring runs above the gap of the test tubes in two levels.

→ Pull the tension ring on the 1st level above the gap towards the middle of the adaptor.

→ Move the test tube in front of the tension spring through the gap, so that the tension ring pushes the test tube against the narrow side of the gap to the outside.

→ Move the test tube further through the 2nd level and repeat to tensioning the spring.

➤ Starting rotation

**Caution:**
Risk of breakage and spillage

If the vessels are not attached securely they may fall out and spill the sample.

Prior to switching the device on ensure that all vessels are securely fastened.

**Warning:**
Risk of poisoning

Open vessels will result in samples splashing.

Close all vessels tight especially if sample is dangerous or toxic.

Adjust rotation speed step by step until you have reached the required rpm settings.

Wear safety glasses and especially with dangerous and toxic samples appropriate protective clothing.

**Warning:**
Risk of injury!

High speeds in combination with high loads and/or a tall assembly may lead to an unstable device. As result the device may start moving and slide off the table top.

When working with with high loads and/or a tall assembly, increase the speed with care and watch for the devices stability. Ensure the device is installed in such a way as it can not collide with the surrounding environment or slip off its work surface.

**Warning:**
Risk of injury! Risk of entanglement!

Rotating parts could cause injuries to the user if their hair, clothing or jewellery comes into contact with or become entangled in these parts.

- Never touch any rotating parts.
- Do not wear any loose clothing, e.g. belts or ties, which could become entangled in the rotating parts.
- Do not wear long necklaces or bracelets.
- If you have long hair, wear a suitable head covering that keeps the hair in place.

A single flask should be arranged in the middle of the shaker plate.
Several flasks should be distributed equally on the plate.

A power cut will stop the device.
With power return the device must be switched on manually.

The rotation of the adaptor is steplessly adjustable in the range from 20-100 rpm through the control knob. Rotation starts as soon as device is switched on.

Operation

An electronic overspeed monitor provides for constant rotation.

- Turn device on with the ON/OFF switch.
- Set rotation speed with analog knob.
- When the ON/OFF switch is switched off the adaptor can be turned to the desired position by hand.

When using shakers inside conditioning cabinets, make reference to ambient conditions as stipulated in the "Technical data" section.

DOMINIQUE DUTSCHER SAS

➤ Cleaning

All surfaces of the device allow for cleaning with a damp cloth and if required a mild soap lotion.

Warning:
Risk of electric shock!

If the device is still connected to power during cleaning, you might get in contact with live parts due to accidentally entering liquids.

Prior to cleaning:

- ensure that the device is switched OFF
- ensure that the plug is pulled from the electric socket.

Caution:
Damage to the device

Accidentally entering liquids will cause the device to fail .
Only wipe with a damp cloth. Never use a soaked cloth.

Attention:
Risk of damage to the surfaces

All of the surfaces may be damaged by improper cleaning.

Never use:

- Chlorine bleach or any cleaning agent containing chlorine
- Solvent-bearing agents (e.g. acetone)
- Ammonia
- Abrasive cleaning agents such as scouring agents, cleaning wool or agents with metallic particles

As a result of vibrations generated during operation the device might slide off the table top.

To ensure a stable position clean surface and feet with a damp cloth and ethanol regularly.

➤ Maintenance

The device is maintenance-free. Any necessary repair must be performed by an authorized Heidolph distributor.

Please contact Heidolph Instruments or your local Heidolph Instruments distributor.

➤ Possible errors and how to resolve

Malfunctions	Possible reason	Troubleshooting
On-/Off switch does not enlighten	No power	<ul style="list-style-type: none"> ▪ Check size and shape of plug and compatibility to the mains ▪ Check circuit breakers
Device stops rotating	Thermal motor circuit breaker triggered by motor overload.	Wait 20 minutes, decrease overall load
	<ul style="list-style-type: none"> ▪ Mechanical parts broken ▪ Electronic failure 	Contact your local Heidolph Instruments distributor

If you experience a malfunction which can not be resolved, please contact your authorized Heidolph Instruments distributor immediately.

Electrical connection

Repairs to the device must **only be carried out by a qualified electrician approved by Heidolph Instruments**. Improper repairs could result in serious dangers. Please contact your specialist dealer if the device needs to be repaired.

Warning: Risk of electric shock

The device must be connected to an earthed power socket.

The rated voltage for the device and the supply voltage must be the same. The rated voltage for the device can be found on the rating plate on the rear of the device.

The device is earthed via the supplied power cord when delivered. If the original plug is replaced, it is essential that the new plug is connected to a protective earth!

Rating plate

Power cord

Colour coding for the electrical connection:

Europe		North America	
GREEN/ YELLOW	PE: Protective earth (earth)	GREEN	PE: Protective earth (earth)
BLUE	N: Neutral conductor	WHITE	N: Neutral conductor
BROWN	P: Phase	BLACK	P: Phase

The device is connected to the power supply with the supplied power cord. The device socket is located on the rear of the device.

If you want to use the device in a country with a different type of plug:

- The supplied plug may only be replaced by a qualified electrician.
- If you want to use an adapter, it must be earthed and approved in accordance with local regulations.
- The cable, adapter and plug must correspond at least to the power stated on the rating plate.

➤ Dismantling, transportation and storage

Dismantling

**Warning:
Risk of injury!**

While the device is plugged in you might be at harm if the device is accidentally switched to ON mode.

Prior to dismantling device ensure that the device is switched off and the plug is pulled from the electric socket.

- ✓ The device is switched off.
- Disconnect plug from power socket.
- Disconnect plug from socket on the device.
- Remove all vessels.
- Disassemble attachments.

Transportation and storage

- Store and transport the device and its components only if they were emptied, cleaned and dismantled.
- To carry the device wear safety gloves and lift the device from below.
- Store and transport the device and parts of device in the original packing material or alternatively in an adequate container to prevent damage.
- Seal the package carefully against unauthorized or accidental opening.
- Store the device in a dry and frost-free place.

Improper storage and transportation may result in damages to the system and the mechanical components.

Avoid any kind of shocks during transportation.

➤ Scope of delivery

Component	Quantity	Product number
Hei-MIX Reax 2*	1	541-21001-00
Universal adaptor	1	549-30000-00
Operating Manual English / German	1	01-005-001-49
Warranty registration / Confirmation of condition	1	01-006-002-78
EC Declaration of Conformity	1	01-001-025-06
Power cord	1	depending on country

* The product number is valid for 230 V devices in the european comdevicey. Please contact Heidolph Instruments for further product numbers on country specific varieties.

➤ Accessories

Component	Quantity	Product number
Adaptor 1 for 20 test tubes	1	509-08100-00

Additional accessories are shown in the general catalog or on our website at www.heidolph-instruments.com

➤ Technical data

Hei-MIX Reax 2

Standard supply voltage	230-240 V/50 Hz or 115 V/60 Hz
Supply power (W)	33
Protective class (IEC 61140)	1
Protection class (IEC 60529)	IP 21
Sound pressure level (dB(A)) (based on IEC 61010)	< 45
Drive	shaded pole motor
Overload	overheat protection, self reset
Speed control	electronic
Motion	overhead rotation
Speed range (rpm)	20 - 100
Rotation speed setting	analog
Dimensions (mm) (w x h x d)	510 x 235 x 180
Effective area	230 x 160
Weight (kg)	5.2
Load capacity (kg)	1

➤ Contact / Technical Service

Questions / Repair work

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please get in touch with the following address.

For repairs please call Heidolph Instruments or your local authorized Heidolph Instruments distributor.

Warning! **Danger of poisoning!**

Contaminated devices can lead to severe injury or death of our employees!

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact substance
- take adequate protective action towards our parts receiving and service personnel
- mark the pack in accordance with Ordinance on Hazardous Substances

A "Confirmation of Condition" form can be found at the end of this operating manual.

→ Prior to shipping the device for repair, complete a copy of this form and submit it in advance.

Contact details:

Heidolph Instruments Germany

Heidolph Instruments GmbH & Co. KG
Technical Service
Walpersdorfer Str. 12
D-91126 Schwabach / Germany

Tel.: +49 – 9122 – 9920-74

Fax: +49 – 9122 – 9920-84

E-Mail: service@heidolph.de
www.heidolph.com

Heidolph Instruments North America

Phone: 1-866-650-9604
E-mail: service@heidolph.com
www.heidolphNA.com

Heidolph Instruments United Kingdom

Phone: 01799 – 5133-20
E-mail: service@radleys.co.uk
www.heidolph-instruments.co.uk

All other countries

You will find contact details of your local Heidolph distributor at
www.heidolph.com

➤ Warranty

Heidolph Instruments provides a 3 year warranty for the products described here (excluding glass and wearing parts) if you register using the warranty card enclosed or online (www.heidolph.com). The warranty period begins with the date of registration. The serial number is also valid without registration.

This warranty covers material and manufacturing defects.

In the event of a material or manufacturing defect, the device shall be repaired or replaced free of charge under the terms of the warranty.

Heidolph Instruments shall not assume liability for any damage incurred as a result of improper handling or transport.

Warranty claim?

- ➔ Please notify Heidolph Instruments or your local Heidolph distributor if you wish to make a warranty claim.

➤ Confirmation of condition

→ In the case of repair, copy and complete the Confirmation of condition form and send it to Heidolph Instruments.

1. Details about the device

Product number _____

Serial number _____

Reason for repair _____

2. Has the device been cleaned, decontaminated/sterilized?

Yes _____ No _____

3. Is the device in a condition which does not represent any health threats for the staff of our service department?

Yes _____ No _____

If not, which substances has the device come into contact with?

4. Legally binding declaration

The customer is aware of being legally liable to Heidolph Instruments for any damages arising from incomplete and incorrect information.

Date _____ Signature _____

Company stamp

Please note

The shipper is responsible for the return of the goods in well packed condition, suitable for the mode of transport.

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This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
© Heidolph Instruments GmbH & Co. KG