Temperature

Controlled Shaking

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

Operating Manual

Page 2

Platform Shakers Hei-MIX Unimax 1010/Promax 1020/
Duomax 1030/Polymax 1040

Operating manual must be read before initial start-up.
Please follow the safety instructions provided.

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About this Document	4
Version and variants	4
About this manual	4
Symbols and keywords.....	5
Safety Instructions	6
General safety instructions.....	6
EU Declaration of Conformity.....	6
Intended use.....	6
Installation / electrical safety	7
Qualifications of personnel	7
Obligations of the operator	7
Installation site	7
Modifications to the device.....	8
Safety of personnel.....	8
Safety during use	9
Disposal	9
Device Description	10
Device overview	10
Hei-MIX Duomax 1030 overall view.....	10
Hei-MIX Polymax 1040 overall view.....	10
Hei-MIX Duomax 1030 / Polymax 1040 control Panel.....	10
Hei-MIX Unimax 1010 overall view.....	11
Hei-MIX Promax 1020 overall view.....	11
Hei-MIX Unimax 1010 / Promax 1020 control panel.....	11
Start-Up	12
Setting up device	12
Connect/disconnect power cord	12
Switch unit on and off	13
Operation	14
Mounting vessels.....	14
Secure vessels on shaker plate	14
Shaking mode.....	14
Hei-MIX Duomax 1030 / Polymax 1040.....	15
Hei-MIX Unimax 1010 / Promax 1020	16

Cleaning and Maintenance	18
Cleaning.....	18
Maintenance	18
Troubleshooting	19
Possible errors and how to resolve	19
Assembly	20
Electrical connections.....	20
Mounting attachments.....	21
Erlenmeyer attachment	21
Attachment separatory funnels (Promax 1020 preferred)	22
Frame with tension rollers.....	23
Disassembly and Storage	24
Dismantling, transportation and storage	24
Dismantling	24
Transportation and storage	24
Accessories and Spare Parts	25
Scope of delivery.....	25
Accessories.....	25
Attachments	26
Technical data	26
Range / admissible load.....	27
Service	28
Contact / Technical Service	28
Warranty	29
Confirmation of condition.....	30

Copyright

This operating manual is protected by copyright and is exclusively intended for the internal use of the purchaser.

Handing over this operating manual to third parties, reproducing it in any type or form – in part or in full – and utilizing and/or communicating its contents is prohibited without prior written approval from the manufacturer.

Any violations will result in an obligation to provide compensation for damages. Other claims remain unaffected.

➤ Version and variants

Version

This manual describes the function, operation and maintenance of the Hei-MIX Unimax 1010/Promax 1020/Duomax 1030/Polymax 1040 platform shakers.

Version	Alteration date
1.0	07/2019

Variants

The device is available in different variants. Certain features or functions are only available in specific product variants. These variants are described in this operating manual.

➤ About this manual

This operating manual is an integral part of the device described here.

- Carefully read through this operating manual and observe all of the safety instructions and warnings notices.
- Ensure that every user has carefully read the operating manual before using the device for the first time.
- Store this operating manual so that it is accessible to all users at all times.
- Pass the operating manual on to any subsequent owner.

The current version of this operating manual can also be found at:

[https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers \(Uni- Pro- Duo- Polymax 1000\)](https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers (Uni- Pro- Duo- Polymax 1000))

In addition, also observe the local regulations at the installation site.

➤ Symbols and keywords

Standardized symbols and signal words are used in this operating manual to warn against any dangers and provide important instructions. These instructions must be strictly observed to avoid accidents and damage.

The following terms and basic symbols are used:

Symbol	Additional signal words / explanation
Warning signs 	The yellow triangle indicates hazardous situations. It is used in combination with the following signal words: DANGER: Indicates a hazardous situation which, if not avoided, will result in serious injury or death. WARNING: Indicates a hazardous situation which, if not avoided, may result in serious injury or death. CAUTION: Indicates a hazardous situation which, if not avoided, may result in property damage and minor or moderate injury.
Prohibitory signs 	Prohibited: The red circle indicates a situation that should be avoided under all circumstances and which, if not avoided, may result in serious injury or death.
Mandatory signs 	Must be observed: The blue circle indicates important information. Please observe this information to avoid any property damage.

Other symbols used:

Symbol	Description
→	Handling instruction, action required
✓	Result of action
▪ ▪ ▪	List of information
a. b. c.	List of variants

➤ General safety instructions

EU Declaration of Conformity

This device complies with the following EC-Directives*:

- 2006/42/EC Machinery Directive
- 2014/30/EU Electromagnetic Compatibility Directive

*Also see the attachment "EU Declaration of Conformity".

The device has been manufactured according to state-of-the-art technology and in compliance with recognized safety regulations. However, risks may still arise during installation, operation and maintenance.

→ Please ensure that the operating manual is available at all times.

The device may only be used under the following conditions:

- Only operate the device if it is in a technically perfect condition.
- Only operate the device if you are properly aware of the risks and required safety measures.
- Only operate the device in accordance with the instructions given in this operating manual.
- If any information is ambiguous or missing, ask your superior or contact the manufacturer.
- Do not operate anything on the device without authorization.
- Only use the device in accordance with its intended use.

Intended use

The device may only be operated by authorized personnel.

The device is suitable for the following uses:

- Shaking
- Mixing
- Emulsifying
- Suspending
- Separating
- Dissolving
- Staining

The device may be operated in research laboratories, other laboratories and production facilities in the following sectors:

- Chemistry
- Pharmacy
- Biology
- Environmental analytics
- Basic research
- Similar research laboratories

Unintended use/Reasonable foreseeable misuse

Any use which deviates from the device's intended use is considered to be improper. The manufacturer is not liable for damage that occurs as a result. The risk is borne by the operator alone.

When used in medical and food sectors the operator alone is responsible for compliance with the applicable regulations and standards.

Installation / electrical safety

- The device may only be connected up if the available power supply voltage corresponds to the information stated on the rating plate for the device.
- The power supply connection must be easily accessible at all times.
- An RCD circuit breaker (residual current circuit breaker) must be fitted to the electrical system in the building so that the device is disconnected from the power circuit in the event of a fault.
- Repairs may only be performed by electricians authorised by Heidolph Instruments.
- Never operate the device with a damaged power cord.
- Always turn the device OFF and disconnect the power cord before carrying out any maintenance or repair work.

Qualifications of employees

- The device may only be operated by trained personnel.
- The device may only be operated by persons who have been instructed and supervised in its proper use by trained specialist personnel.
- The device may only be operated by specialist personnel who are above the legal minimum age.
- Other persons may only work on the device under the constant supervision of experienced and trained specialist personnel.
- This operating manual must be read and understood by all persons working with the device.
- Personnel must receive safety training that ensures responsible and safe working practices.

➤ Obligations of the operator

Installation site

- The device must be installed in a suitable location.
- The device must be installed on a firm and stable surface.
- Ensure that the device and all of its components are easily accessible at all times.
- Maintain a sufficient safety clearance to, in particular, any moving and/or hot equipment components.
- It is not permitted to place or store any objects such as accessories, tools or chemicals within this clearance area during operation.
- All screw connections must be securely tightened.
- It is not permitted to operate the device near to highly flammable or explosive substances.
- Operate the device in conjunction with an extractor hood, if working with potential harmful media (see DIN EN 14175 and DIN 12924).

Safety Instructions

- The device is designed for indoor use only and under the following ambient conditions:

Ambient temperature	5 - 31 °C at 80 % relative humidity
	32 - 40 °C decreasing linearly to a maximum 50 % relative humidity
Installation altitude	0 - 2,000 m above sea level
Contamination level	2
Overvoltage category	II
Permissible supply deviations	± 10 %
- If the device is operated in corrosive atmospheres, the service life of the device will decrease based on the concentration, duration and frequency of the exposure to the corrosive atmosphere e.g. concentrated hydrochloric acid (HCl).

Modifications to the device

- It is not permitted to make any unauthorised modifications or changes to the device.
- Do not attach or install any parts that have not been approved by the manufacturer.
- Unauthorized modifications or changes will void the EC Declaration of Conformity for the device and operation of the device will no longer be permitted.
- The manufacturer is not liable for any damage, dangers or injuries that result from unauthorized modifications and changes or due to the non-observance of the instructions in this manual.

Employee safety

- Ensure that the device is only operated by qualified specialist personnel and trained employees.
- Observe the following instructions to avoid any personal injuries and property damage:
 - Laboratory regulations
 - Accident prevention regulations
 - Hazardous Substance Act
 - Other generally accepted occupational health and safety regulations
 - Local regulations

➤ Safety during use

- Wear the appropriate protective clothing when working on the device (protective glasses and, if necessary, safety gloves).
- Do not use the device in potentially explosive areas. The device is not protected against explosion. There is no explosion or ATEX protection available.
- Do not carry out work with naked flames in the vicinity of the device (risk of explosion).
- Do not operate or assemble devices in the vicinity which are emission or radiation sources (electromagnetic waves) for the frequency range ($3 \cdot 10^{11}$ Hz to $3 \cdot 10^{15}$ Hz).
- Do not operate or assemble appliances in the vicinity of the device which constitute emission or radiation sources for ionizing radiation or in the ultrasonic range.
- Do not operate the device where adiabatic compression or shock waves might occur (shock wave combustion).
- Do not use substances, where the energy input due to mixing might pose a hazard.
- Avoid putting pressure on the display when you are not operating the device.
- Do not spill liquids over the device or any parts of it.
- Remove any accidentally spilled liquids immediately.
- Eliminate errors immediately.
- Do not use abrasive material to clean the surface. Only wipe with damp cloth.
- Always switch the device OFF after use.

➤ Disposal

- Check the device components for hazardous substances and solvents.
- Clean all components before disposal.
- Dispose of the device in accordance with the relevant national regulations.
- Dispose of the packaging material in accordance with the appropriate national regulations.

Device Description

➤ Device overview

Hei-MIX Duomax 1030 overall view

Example:
culture bottles

Hei-MIX Polymax 1040 overall view

Example:
petri dishes

Hei-MIX Duomax 1030 / Polymax 1040 control Panel

Analog knob:
set speed

Analog knob:
set time

Device Description


Hei-MIX Unimax 1010 overall view

Hei-MIX Promax 1020 overall view

Hei-MIX Unimax 1010 / Promax 1020 control panel

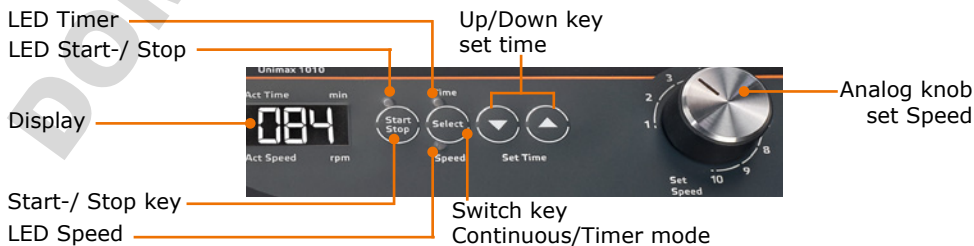➤ Setting up device

Warning:
Risk of slipping device!

As a result of vibrations generated during operation the device might slide off the table top.

The surface must be smooth and clean and should be properly leveled.

- Locate the shaker on a stable, horizontal surface.
- Clean surface and feet with a damp cloth and ethanol regularly.

➤ Connect/disconnect power cord

Connect power cord

The power cord comes with a three-wire plug and a recess on the bottom side.

Unit plug

Power cord

The socket is located on the back side of the unit.

- ✓ The device is switched OFF.
- Connect plug to the socket with the recess facing down.
- Push plug up against socket until securely attached.
- Connect plug to power socket.

Unit socket

Disconnect power cord

- Disconnect plug from power socket first.
- Then disconnect plug from socket on unit.

➤ **Switch unit on and off**

The ON/OFF switch is situated on the front panel.

Switch unit on


- ➔ Press the On/Off switch on the lower side.
- ✓ The green control light switches on.
- ✓ The unit is switched on.

Switch unit off

- ➔ Press the On/Off switch on the upper side.
- ✓ The green control light switches off.
- ✓ The unit is switched off.

Duomax 1030 / Polymax 1040

ON/OFF switch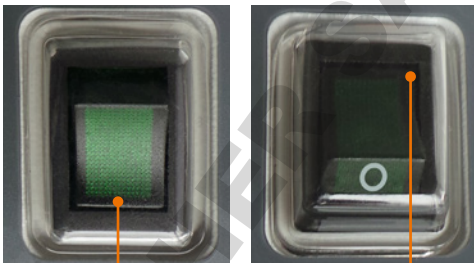

Unit switched on

Unit switched off

The ON/OFF switch is situated at the lower left side of the device.

Switch unit on


- ➔ Press the On/Off switch at the front side.
- ✓ The green control light switches on.
- ✓ The unit is switched on.

Switch unit off

- ➔ Press the On/Off switch at the back side.
- ✓ The green control light switches off.
- ✓ The unit is switched off.

Unimax 1010 / Promax 1020

ON/OFF switch

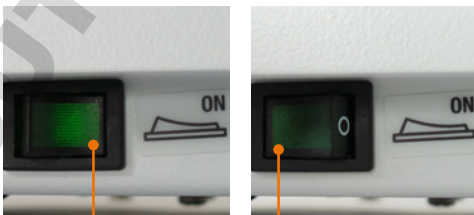Unit switched on

Unit switched off

DOMINIQUE DUTSCH

➤ Mounting vessels

Secure vessels on shaker plate

Adaptors for Erlenmeyer flasks, separating funnels and an adaptor with tension rollers for universal use are available.

- To secure vessels on shaker plate, use the optional adaptors (see chapter „Spare parts and accessories“ and chapter „Assembly“).
- At low shaking speeds, low-profile vessels like petri dishes may be placed directly on the non-skid rubber plate.

A single flask should be arranged in the middle of the shaker plate.
Several flasks should be distributed equally on the plate.

➤ Shaking mode

Caution: Risk of breakage and spillage

If the vessels are not attached securely they may fall out and spill the sample.

Prior to switching the unit on ensure that all vessels are securely fastened.

Warning: Risk of poisoning

Open vessels and too high vibration frequencies may result in samples splashing.

- Close all vessels tight especially if sample is dangerous or toxic.
- Adjust vibration frequency step by step until you have reached the required rpm settings.
- Wear safety glasses and especially with dangerous and toxic samples appropriate protective clothing.

Warning: Risk of injury!

High speeds in combination with high loads and/or a tall assembly may lead to an unstable device. As result the device may start moving and slide off the table top.

When working with with high loads and/or a tall assembly, increase the speed with care and watch for the devices stability. Ensure the device is installed in such a way as it can not collide with the surrounding environment or slip off its work surface.

At high loads (load bearing capacity of shaker plate) and high frequency of shaking, always make reference to load graph as applicable for your item (refer to „Technical data“).

When using shakers inside conditioning cabinets, make reference to ambient conditions as stipulated in the „Technical data“ section.

Hei-MIX Duomax 1030 / Polymax 1040

(with mechanical timer)

You can choose from two shaking options:

- Continuous mode: The unit starts and stops shaking as soon as you switch it on or off.
- Timer mode: The time for shaking can be set between 0 and 120 minutes. The unit stops shaking when the time has elapsed.

Warning: Risk of injury!

Returning power after a power cut off will restart shaking mode. There is a risk of injury in the vicinity of moving parts of device.

Before working with the device or in the vicinity ensure all controls are switched to "0" or "OFF" mode, when device is not moving.

Continuous mode

- ✓ Select mode switch is positioned to continuous mode „0”.
- Turn item on with ON/OFF switch.
- Set speed with analog knob „Speed”.
- ✓ Actual speed appears in the 3-digit digital display.
- Turn speed off.
- Turn item off with ON/OFF switch.

Timer mode

- Select timer mode „**Timer**” with select mode switch.
- Set time with analog knob „Time”.
- Set speed with analog knob „Speed”.
- Turn item on with ON/OFF switch.
- ✓ When time has elapsed an acoustic signal will be audible and shaking action will stop.
- Turn speed off.
- Turn item off with ON/OFF switch.

A power cut will stop shaking mode. The timer will continue running even in case of power supply failure.

Hei-MIX Unimax 1010 / Promax 1020 (with digital timer and digital speed display)

You can choose from two shaking options:


- Continuous mode: The unit starts and stops shaking as soon as you switch it on or off.
- Timer mode: The time for shaking can be set between 0 and 999 minutes. The unit stops shaking when the time has elapsed.

A power cut will stop the unit.
With power return the device is switched on automatically. Shaking must be started manually.

Continuous mode

- Turn unit on with ON/OFF switch.
- Set time to „000” by pressing keys ▲ and ▼ together.
- Start continuous operation by hitting the Start/Stop key.
- ✓ Start and speed LEDs enlighten.
- ✓ Shaking frequency is displayed (4-digit display).
- Set shaking frequency with knob „Speed”.
- Hit Start/Stop key to stop shaking.

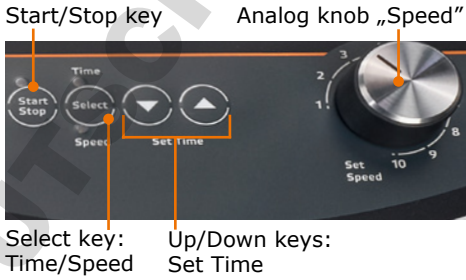DOMINIQUE DIETHELM

Timer mode

- Set time between 0 and 999 minutes with Set Time keys ▲ or ▼.
- ✓ Short hits on the keys change time setting slowly.
- ✓ Depressing keys is going to change time rapidly.
- Set speed with analog knob.
- Start shaking with Start/Stop key.
- ✓ Start and speed LEDs enlighten.
- ✓ Time lapse appears in a 3-digit display.
- Discontinue current shaking action with the Start/Stop key.
- ✓ Actual time won't change.
- Hit Start/Stop key to continue shaking at the moment it had been discontinued.
- ✓ In case of a power supply failure or interruption (item turned off), this feature is lost.
- Hit Select key during timer-controlled operation to change display from time to shaking frequency and vice versa.
- ✓ Speed LED turns on.
- ✓ Timer LED flashes to indicate timer operation mode.
- ✓ Timer LED also flashes, when timer-controlled operation was discontinued with the Start/Stop key.

Repetition

Repeating last timer controlled operation does not require new time setting, as long as you did not touch the master switch.

Edit time during shaking operation

Time setting may be changed during operation, while timer is running and even if timer controlled operation is discontinued.

The pre-programmed time remains in memory.

➤ Cleaning

All surfaces of the unit allow for cleaning with a damp cloth and if required a mild soap lotion.

Warning:
Risk of electric shock!

If the device is still connected to power during cleaning, you might get in contact with live parts due to accidentally entering liquids.

Prior to cleaning:

- ensure that the device is switched OFF
- ensure that the plug is pulled from the electric socket.

Caution:
Damage to the device

Accidentally entering liquids will cause the device to fail .
Only wipe with a damp cloth. Never use a soaked cloth.

Attention:
Risk of damage to the surfaces

All of the surfaces may be damaged by improper cleaning.

Never use:

- Chlorine bleach or any cleaning agent containing chlorine
- Solvent-bearing agents (e.g. acetone)
- Ammonia
- Abrasive cleaning agents such as scouring agents, cleaning wool or agents with metallic particles

As a result of vibrations generated during operation the device might slide off the table top.

To ensure a stable position clean surface and feet with a damp cloth and ethanol regularly.

➤ Maintenance

The unit is maintenance-free. Any necessary repair must be performed by an authorized Heidolph distributor.

Please contact Heidolph Instruments or your local Heidolph Instruments distributor.

Troubleshooting

➤ Possible errors and how to resolve

Malfunctions	Possible reason	Troubleshooting
On-/Off switch does not enlighten	No power	<ul style="list-style-type: none"> ▪ Check size and shape of plug and compatibility to your electrical socket ▪ Check circuit breakers ▪ Check device fuse (see below)
	Light diode defective	Contact your local Heidolph Instruments distributor
Shaking movement stops	Pre-set timer elapsed	Restart timer, if necessary
	Thermal motor circuit breaker triggered by motor overload.	Wait 20 minutes, decrease overall load
	<ul style="list-style-type: none"> ▪ Motor defective ▪ Control electronics defective 	Contact your local Heidolph Instruments distributor

If you experience a malfunction which can not be resolved, please contact your authorized Heidolph Instruments distributor immediately.

Electrical connections

- Fuses must **only** be replaced by a **professional electrician**.
- **Repairs** to the device must **only be carried out by a qualified electrician approved by Heidolph Instruments**. Improper repairs could result in serious dangers. Please contact your specialist dealer if the device needs to be repaired.

Warning: Risk of electric shock

The device must be connected to an earthed power socket.

The rated voltage for the device and the supply voltage must be the same. The rated voltage for the device can be found on the rating plate on the rear of the device.

The device is supplied with a grounded power cord. In case the original plug is replaced, the new plug must have a protective conductor!

If the device is still connected to power while changing the fuse, you might get in contact with live parts.

Prior to replacing the fuse ensure that the unit is switched off and pull the plug from the electric socket. Only use original spare fuses from Heidolph Instruments.

After changing a fuse check the device for safe operating condition according to IEC 61010-1.

Rating plate

Unit socket

Fuse

Color code for electrical connections:

Europe

GREEN/YEL-LOW	PE: Protective conductor (Earth)
BLUE	N: Neutral conductor
BROWN	P: Phase

North America

GREEN	PE: Protective conductor (Earth)
WHITE	N: Neutral conductor
BLACK	P: Phase

The device is connected to the mains with the power cord supplied. The power connection is located on the rear of the housing.

For countries where a different plug than the standard plug supplied is used:

- The plug supplied may only be changed by a professional electrician.
- If you wish to use an adapter, ensure that it is grounded and approved in accordance with the local regulations.
- Wiring, adapter and plug must at least meet the electrical power stated on the rating plate.


➤ Mounting attachments

Erlenmeyer attachment

Components Erlenmeyer attachment:
(Example)

- 1 attachment
- 4 knurled bolts

- Locate attachment on rubber plate.
- Secure attachment at the four corners of the shaker plate with 4 knurled bolts.
- Load Erlenmeyer flasks in cut-outs.
- ✓ For proper seating, the cut-outs feature a preformed packing.

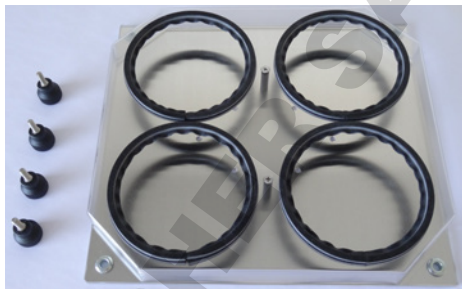4 x Knurled bolts

Lightly tilt Erlenmeyer flasks to fit them under the clamp.

Attachment separatory funnels (Promax 1020 preferred)

Caution

Risk of breakage! Risk of injury!

Separatory funnels are very delicate especially at the tip and may break when inserted without care.

Handle separatory funnels carefully when clamping.

Components Erlenmeyer attachment:

- 1 attachment
- 4 knurled bolts
- Locate attachment on rubber plate.
- Secure attachment at the four corners of shaker plate with 4 knurled bolts.
- Load separatory funnel outlet pipe first in place.
- Unlock pressure pad by pulling knurled knob.
- Introduce opposite end of funnel in rubber grommet and slowly release pressure pad until funnel is locked in place.
- ✓ With this sort of clamping, funnel's ground plug is held in place firmly.

Outlet pipe

Knurled knob


Frame with tension rollers


Frame with tension rollers allow to safely attach whatever sort of vessels and racks on the shaker platform.

Components frame with tension roller:

- 4 bolts
- 1 box-end spanner
- 2 holding bars
- 2 tension rollers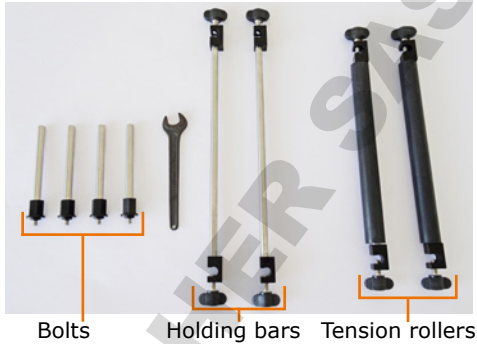

- Thread 4 ea. bolts in the respective holes of platform and pull tight with box-end spanner from hardware bag.
- Slip holding bar over bolts and secure with wing bolts.
- Slip tension rollers over holding bars, push against the vessel to be clamped, and secure with wing bolts.
- Attaching one more row of vessels requires installation of extra tension rollers.
- Lock tension rollers safely in place and pull all screws of the universal mount tight as specified.

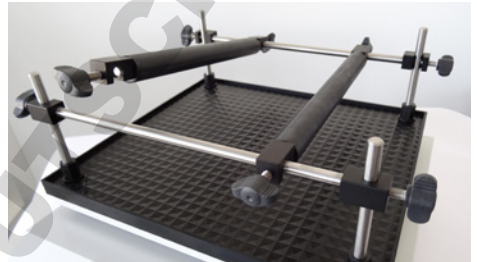➤ Dismantling, transportation and storage

Dismantling

Warning:
Risk of injury!

While the device is plugged in you might be at harm if the device is accidentally switched to ON mode.

Prior to dismantling device ensure that the device is switched off and the plug is pulled from the electric socket.

- ✓ The device is switched off.
- Disconnect plug from power socket.
- Disconnect plug from socket on the device.
- Remove all vessels.
- Disassemble attachments.

Transportation and storage

- Store and transport the device and its components only if they were emptied, cleaned and dismantled.
- To carry the device wear safety gloves and lift the unit from below.
- Store and transport the device and parts of device in the original packing material or alternatively in an adequate container to prevent damage.
- Seal the package carefully against unauthorized or accidental opening.
- Store the device in a dry and frost-free place.

Improper storage and transportation may result in damages to the system and the mechanical components.

Avoid any kind of shocks during transportation.

➤ Scope of delivery

Component	Variant	Quantity	Product number
Hei-MIX	Unimax 1010*	1	543-12310-00
	Promax 1020*	1	543-22332-00
	Duomax 1030 tilt angle 5° *	1	543-32205-00
	Duomax 1030 tilt angle 10° *	1	543-32210-00
	Polymax 1040 tilt angle 5° *	1	543-42205-00
	Polymax 1040 tilt angle 10° *	1	543-42210-00
Operating Manual English / German		1	01-005-005-67
Warranty registration / Confirmation of condition		1	01-006-002-78
EC Declaration of Conformity		1	01-001-025-06
Power cord		1	depending on country

* The product numbers are valid for 230 V devices in the european community. Please contact Heidolph Instruments for further product numbers on country specific varieties.

➤ Accessories

Component	Quantity	Product number
Fuse for 230 V (device needs 2 fuses)	1	14-002-015-23
Fuse for 230 V (device needs 2 fuses)	1	14-002-015-45
Attachment with 2 tension rollers	1	549-70000-00
Tension rollers (additional)	1	549-71000-00
Attachment for up to 22 Erlenmeyer flasks, 25 ml	1	549-72000-00
Attachment for up to 14 Erlenmeyer flasks, 50 ml	1	549-73000-00
Attachment for up to 9 Erlenmeyer flasks, 100 ml	1	549-74000-00
Attachment for up to 5 Erlenmeyer flasks, 250 ml	1	549-75000-00
Attachment for up to 4 Erlenmeyer flasks, 500 ml	1	549-76000-00
Attachment for up to 2 Erlenmeyer flasks, 1000 ml	1	549-77000-00
Attachment for separatory funnels 4 x 50 ml / 4 x 100 ml	1	549-78000-00

Additional accessories are shown in the general catalog or on our website at www.heidolph-instruments.com.


➤ Technical data

Protective class (IEC 61140)	1 Ⓢ			
Protection class (IEC 60529)	IP 40			
Device fuses	M 1,25 A (230 V) oder T 2,0 A (115 V)			
Sound pressure level (dB(A)) (based on IEC 61010)	< 50			
Operating mode	timer or continuous			
Platform size	290 x 258			
Weight (kg)	8.0			
Load capacity (kg)	5			
Model Hei-MIX	Unimax 1010	Promax 1020	Duomax 1030	Polymax 1040
Speed range (rpm)	30 - 500	30 - 250	2 - 50	2 - 50
Motion	orbital	reciprocating	rocking	wave
Orbit (mm)	10	32	-	
Tilt angle (°)	-		5 / 10	
Dimensions (w x h x d) (mm)	320 x 125 x 375		320 x 185 x 375	320 x 195 x 375
Standard supply voltage	100 - 240 V (50/60 Hz)		250 V/50 Hz or 115 V/60 Hz	
Drive	EC-motor		asynchronous motor	
Supply power (W)	50		115	
Overheat protection	-		self-resetting	
Speed setting	digital		analog	


Range / admissable load


Unimax 1010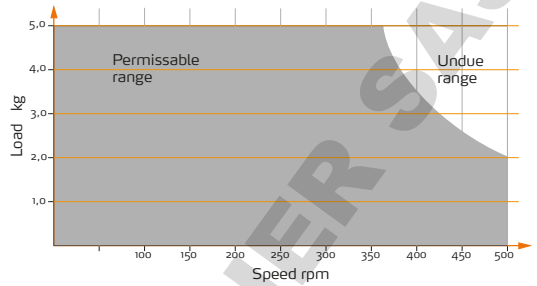

Promax 1020

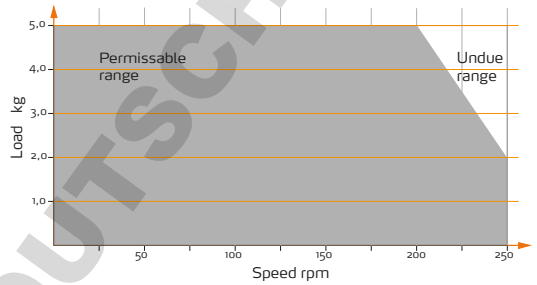Duomax 1030,
Polymax 1040

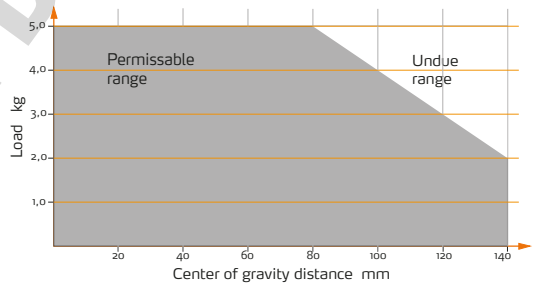DOMINIQUE DUTSCHER SAS

➤ Contact / Technical Service

Questions / Repair work

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please get in touch with the following address.

For repairs please call Heidolph Instruments or your local authorized Heidolph Instruments distributor.

Warning! **Danger of poisoning!**

Contaminated devices can lead to severe injury or death of our employees!

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact substance
- take adequate protective action towards our parts receiving and service personnel
- mark the pack in accordance with Ordinance on Hazardous Substances

A "Confirmation of Condition" form can be found at the end of this operating manual.

→ Prior to shipping the device for repair, complete a copy of this form and submit it in advance.

Contact details:

Heidolph Instruments Germany

Heidolph Instruments GmbH & Co. KG
Technical Service
Walpersdorfer Str. 12
D-91126 Schwabach / Germany

Tel.: +49 – 9122 - 9920-74

Fax: +49 – 9122 - 9920-84

E-Mail: service@heidolph.de

www.heidolph.com

Heidolph Instruments North America

Phone: 1-866-650-9604

E-mail: service@heidolph.com

www.heidolphNA.com

Heidolph Instruments United Kingdom

Phone: 01799 - 5133-20

E-mail: service@radleys.co.uk

www.heidolph-instruments.co.uk

All other countries

You will find contact details of your local Heidolph distributor at www.heidolph.com

➤ Warranty

Heidolph Instruments provides a 3 year warranty for the products described here (excluding glass and wearing parts) if you register using the warranty card enclosed or online (www.heidolph.com). The warranty period begins with the date of purchase. The serial number is also valid without registration.

This warranty covers material and manufacturing defects.

In the event of a material or manufacturing defect, the device shall be repaired or replaced free of charge under the terms of the warranty.

Heidolph Instruments shall not assume liability for any damage incurred as a result of improper handling or transport.

Warranty claim?

- ➔ Please notify Heidolph Instruments or your local Heidolph distributor if you wish to make a warranty claim.

DOMINIQUE DUTSCHER MAS

➤ Confirmation of condition

→ In the case of repair, copy and complete the Confirmation of condition form and send it to Heidolph Instruments.

1. Details about the device

Product number _____

Serial number _____

Reason for repair _____

2. Has the device been cleaned, decontaminated/sterilized?

Yes _____ No _____

3. Is the device in a condition which does not represent any health threats for the staff of our service department?

Yes _____ No _____

If not, which substances has the device come into contact with?

4. Legally binding declaration

The customer is aware of being legally liable to Heidolph Instruments for any damages arising from incomplete and incorrect information.

Date

Signature

Company stamp

Please note

The shipper is responsible for the return of the goods in well packed condition, suitable for the mode of transport.

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This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
© Heidolph Instruments GmbH & Co. KG