Temperature Controlled Shaking

Leading Safety Standards

Superior Ease of Use

Reduced Cost of Ownership

Operating Manual

Plattform Shaker Hei-MIX Titramax 100/101/1000,
Vibramax 100/110 and Rotamax 120.

Operating manual must be read before initial start-up.

Please follow the safety instructions provided.

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Seite 28

About this Document	4
Version and variants	4
About this manual	4
Symbols and keywords.....	5
Safety Instructions	6
General safety instructions.....	6
EU Declaration of Conformity.....	6
Intended use.....	6
Installation / electrical safety	7
Qualifications of employees.....	7
Obligations of the operator	7
Installation site	7
Modifications to the device.....	8
Employee safety	8
Safety during use	9
Disposal	9
Device Description	10
Device overview	10
Vibramax 100 overall view	10
Vibramax 110 overall view	10
Titramax 100/101/1000 overall view	11
Rotamax 120 overall view	11
Control panel	11
Start-Up	12
Setting up device	12
Connect/disconnect power cord	12
Switch unit on and off	13
Operation	14
Mounting vessels.....	14
Secure vessels on shaker plate.....	14
Shaking mode.....	16
Continuous mode.....	17
Timer mode	17

Cleaning and Maintenance	18
Cleaning.....	18
Maintenance	18
Troubleshooting	19
Possible errors and how to resolve	19
Assembly	20
Electrical connections	20
Disassembly and Storage	21
Dismantling, transportation and storage	21
Dismantling	21
Transportation and storage	21
Accessories and Spare Parts	22
Scope of delivery.....	22
Accessories.....	22
Attachments	23
Technical data	23
Performance range / admissible load	24
Service	25
Contact / Technical Service	25
Warranty	26
Confirmation of condition.....	27

Copyright

This operating manual is protected by copyright and is exclusively intended for the internal use of the purchaser.

Handing over this operating manual to third parties, reproducing it in any type or form –in part or in full – and utilizing and/or communicating its contents is prohibited without prior written approval from the manufacturer.

Any violations will result in an obligation to provide compensation for damages. Other claims remain unaffected.

➤ Version and variants

Version

This manual describes the function, operation and maintenance of the Hei-MIX Titramax 100/101/1000, Vibramax 100/101 and Rotamax 120 platform shakers.

Version	Alteration date
1.2	04/2019

Variants

The device is available in different variants. Certain features or functions are only available in specific product variants. These variants are described in this operating manual.

➤ About this manual

This operating manual is an integral part of the device described here.

- Carefully read through this operating manual and observe all of the safety instructions and warnings notices.
- Ensure that every user has carefully read the operating manual before using the device for the first time.
- Store this operating manual so that it is accessible to all users at all times.
- Pass the operating manual on to any subsequent owner.

The current version of this operating manual can also be found at:

[https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers \(Hei-MIX Titra- Vibra- Rota- 100\)](https://heidolph-instruments.com/en/service/downloads/operation-manuals#Shakers & Mixers (Hei-MIX Titra- Vibra- Rota- 100))

In addition, also observe the local regulations at the installation site.

➤ Symbols and keywords

Standardized symbols and signal words are used in this operating manual to warn against any dangers and provide important instructions. These instructions must be strictly observed to avoid accidents and damage.

The following terms and basic symbols are used:

Symbol	Additional signal words / explanation
<p>Warning signs</p> 	<p>The yellow triangle indicates hazardous situations. It is used in combination with the following signal words:</p> <p>DANGER: Indicates a hazardous situation which, if not avoided, will result in serious injury or death.</p> <p>WARNING: Indicates a hazardous situation which, if not avoided, may result in serious injury or death.</p> <p>CAUTION: Indicates a hazardous situation which, if not avoided, may result in property damage and minor or moderate injury.</p>
<p>Prohibitory signs</p> 	<p>Prohibited:</p> <p>The red circle indicates a situation that should be avoided under all circumstances and which, if not avoided, may result in serious injury or death.</p>
<p>Mandatory signs</p> 	<p>Must be observed:</p> <p>The blue circle indicates important information. Please observe this information to avoid any property damage.</p>

Other symbols used:

Symbol	Description
→	Handling instruction, action required
✓	Result of action
▪	List of information
▪	
▪	
a.	List of variants
b.	
c.	

➤ General safety instructions

EU Declaration of Conformity

This device complies with the following EC-Directives*:

- 2006/42/EC Machinery Directive
- 2014/30/EU Electromagnetic Compatibility Directive

*Also see the attachment "EU Declaration of Conformity".

The device has been manufactured according to state-of-the-art technology and in compliance with recognized safety regulations. However, risks may still arise during installation, operation and maintenance.

→ Please ensure that the operating manual is available at all times.

The device may only be used under the following conditions:

- Only operate the device if it is in a technically perfect condition.
- Only operate the device if you are properly aware of the risks and required safety measures.
- Only operate the device in accordance with the instructions given in this operating manual.
- If any information is ambiguous or missing, ask your superior or contact the manufacturer.
- Do not operate anything on the device without authorization.
- Only use the device in accordance with its intended use.

Intended use

The device may only be operated by authorized personnel.

The device is suitable for the following uses:

- Shaking
- Mixing
- Emulsifying
- Suspending
- Separating
- Dissolving
- Staining

The device may be operated in research laboratories, other laboratories and production facilities in the following sectors:

- Chemistry
- Pharmacy
- Biology
- Environmental analytics
- Basic research
- Similar research laboratories

Unintended use/Reasonable foreseeable misuse

Any use which deviates from the device's intended use is considered to be improper. The manufacturer is not liable for damage that occurs as a result. The risk is borne by the operator alone.

When used in medical and food sectors the operator alone is responsible for compliance with the applicable regulations and standards.

Installation / electrical safety

- The device may only be connected up if the available power supply voltage corresponds to the information stated on the rating plate for the device.
- The power supply connection must be easily accessible at all times.
- An RCD circuit breaker (residual current circuit breaker) must be fitted to the electrical system in the building so that the device is disconnected from the power circuit in the event of a fault.
- Repairs may only be performed by electricians authorised by Heidolph Instruments.
- Never operate the device with a damaged power cord.
- Always turn the device OFF and disconnect the power cord before carrying out any maintenance or repair work.

Qualifications of employees

- The device may only be operated by trained personnel.
- The device may only be operated by persons who have been instructed and supervised in its proper use by trained specialist personnel.
- The device may only be operated by specialist personnel who are above the legal minimum age.
- Other persons may only work on the device under the constant supervision of experienced and trained specialist personnel.
- This operating manual must be read and understood by all persons working with the device.
- Personnel must receive safety training that ensures responsible and safe working practices.

➤ Obligations of the operator

Installation site

- The device must be installed in a suitable location.
- The device must be installed on a firm and stable surface.
- Ensure that the device and all of its components are easily accessible at all times.
- Maintain a sufficient safety clearance to, in particular, any moving and/or hot equipment components.
- It is not permitted to place or store any objects such as accessories, tools or chemicals within this clearance area during operation.
- All screw connections must be securely tightened.
- It is not permitted to operate the device near to highly flammable or explosive substances.
- Operate the device in conjunction with an extractor hood, if working with potential harmful media (see DIN EN 14175 and DIN 12924).

Safety Instructions

- The device is designed for indoor use only and under the following ambient conditions:

Ambient temperature	5 - 31 °C at 80 % relative humidity
	32 - 40 °C decreasing linearly to a maximum 50 % relative humidity
Installation altitude	0 - 2,000 m above sea level
Contamination level	2
Overvoltage category	II
Permissible supply deviations	± 10 %

- If the device is operated in corrosive atmospheres, the service life of the device will decrease based on the concentration, duration and frequency of the exposure to the corrosive atmosphere e.g. concentrated hydrochloric acid (HCl).

Modifications to the device

- It is not permitted to make any unauthorised modifications or changes to the device.
- Do not attach or install any parts that have not been approved by the manufacturer.
- Unauthorized modifications or changes will void the EC Declaration of Conformity for the device and operation of the device will no longer be permitted.
- The manufacturer is not liable for any damage, dangers or injuries that result from unauthorized modifications and changes or due to the non-observance of the instructions in this manual.

Employee safety

- Ensure that the device is only operated by qualified specialist personnel and trained employees.
- Observe the following instructions to avoid any personal injuries and property damage:
 - Laboratory regulations
 - Accident prevention regulations
 - Hazardous Substance Act
 - Other generally accepted occupational health and safety regulations
 - Local regulations

Safety Instructions

➤ Safety during use

- Wear the appropriate clothing when working on the device (protective clothing, protective glasses and, if necessary, safety gloves).
- Do not use the device in potentially explosive areas. The device is not protected against explosion. There is no explosion or ATEX protection available.
- Do not carry out work with naked flames in the vicinity of the device (risk of explosion).
- Do not operate or assemble devices in the vicinity which are emission or radiation sources (electromagnetic waves) for the frequency range ($3 \cdot 10^{11}$ Hz to $3 \cdot 10^{15}$ Hz).
- Do not operate or assemble appliances in the vicinity of the device which constitute emission or radiation sources for ionizing radiation or in the ultrasonic range.
- Do not operate the device where adiabatic compression or shock waves might occur (shock wave combustion).
- Do not use substances, where the energy input due to mixing might pose a hazard.
- Do not spill liquids over the device or any parts of it.
- Remove any accidentally spilled liquids immediately.
- Eliminate errors immediately.
- Always switch the device OFF after use.
- Avoid putting pressure on the display when you are not operating the device.
- Do not use abrasive material to clean the surface. Only wipe with damp cloth.

➤ Disposal

- Check the device components for hazardous substances and solvents.
- Clean all components before disposal.
- Dispose of the device in accordance with the relevant national regulations.
- Dispose of the packaging material in accordance with the appropriate national regulations.

➤ Device overview

Vibramax 100 overall view

Vibramax 110 overall view

Titramax 100/101/1000 overall view

Rotamax 120 overall view

Control panel

➤ Setting up device

Warning:
Risk of slipping device!

As a result of vibrations generated during operation the device might slide off the table top.

The surface must be smooth and clean and should be properly leveled.

- Locate the shaker on a stable, horizontal surface.
- Clean surface and feet with a damp cloth and ethanol regularly.

➤ Connect/disconnect power cord

Connect power cord

The power cord comes with a three-wire plug and a recess on the bottom side.

Unit plug

Power cord

The socket is located on the back side of the unit.

- ✓ The device is switched OFF.
- Connect plug to the socket with the recess facing down.
- Push plug up against socket until securely attached.
- Connect plug to power socket.

Unit socket

Disconnect power cord

- Disconnect plug from power socket first.
- Then disconnect plug from socket on unit.

➤ Switch unit on and off

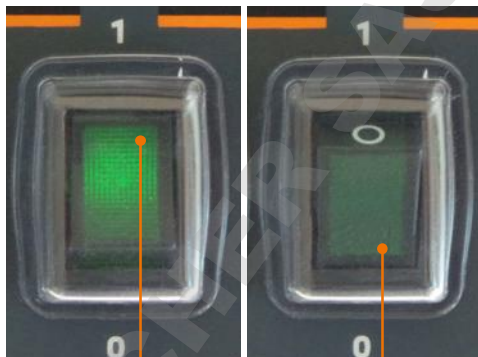
The ON/OFF switch is situated on the front panel.

Switch unit on

- Press the On/Off switch on the upper side.
- ✓ The green control light switches on.
- ✓ The unit is switched on.

Switch unit off

- Press the On/Off switch on the lower side.
- ✓ The green control light switches off.
- ✓ The unit is switched off.

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➤ Mounting vessels

Secure vessels on shaker plate

Use optional attachment to secure vessels on shaker plate:

- To secure vessels on shaker plate, use optional adaptors (see chapter „Spare parts and accessories“ and chapter „Assembly“).
- At low shaking speed, low-profile vessels like petri dishes may be placed directly on the non-skid rubber plate.

A single flask should be arranged in the middle of the shaker plate.
Several flasks should be distributed equally on the plate.

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- Place microtiter plates directly in indentations in the rubber mat.
- ✓ Microtiter plates are positively locked in these indentations.
- ✓ The rubber mat can be removed easily for cleaning or replacement.

VIBRAMAX 100 / ROTAMAX 120 / attachment with clamping rollers

Vessels of different sizes can be secured by an optional attachment with clamping rollers. Vessels clamped at the same time by two clamping rollers must have the same diameter.

- Secure attachment on the shaker plate with the wing bolts supplied.
- Place vessels on the anti-skid rubber mat.
- Clamp vessels on both sides with the clamping rollers.

Operation

VIBRAMAX 110 / attachment for test tubes

- Place a single test tube on the rubber mat by hand and softly push tube downwards.

Caution:
Possible risk of exertion

There is a possible risk of exertion due to physical force required to holding test tube in place.

Do not use the device any longer than recommended in the directive 2002-44 EC on vibration strain.

Several test tubes may be shaken at the same time, using the optional test tube attachments.

- Secure attachment on the shaker table with the wing bolts supplied.
- Plug test tubes into "collets".
- ✓ The test tubes will be secured by the collets.

➤ Shaking mode

Warning:
Risk of injury!

Returning power after a power cut off will restart shaking mode. There is a risk of injury in the vicinity of moving parts of device.

Before working with the device or in the vicinity ensure all controls are switched to "0" or "OFF" mode, when device is not moving.

Caution:
Risk of breakage and spillage

If the vessels are not attached securely they may fall out and spill the sample.

Prior to switching the unit on ensure that all vessels are securely fastened.

Warning:
Risk of poisoning

Open vessels and too high vibration frequencies may result in samples splashing.

- Close all vessels tight especially if sample is dangerous or toxic.
- Adjust vibration frequency step by step until you have reached the required rpm settings..
- Wear safety glasses and especially with dangerous and toxic samples appropriate protective clothing.

Warning:
Risk of injury!

High speeds in combination with high loads and/or a tall assembly may lead to an unstable device. As result the device may start moving and slide off the table top.

When working with with high loads and/or a tall assembly, increase the speed with care and watch for the devices stability. Ensure the device is installed in such a way as it can not collide with the surrounding environment or slip off its work surface.

At high loads (load bearing capacity of shaker plate) and high frequency of shaking, always make reference to load graph as applicable for your item (see chapter "Technical data", section "Performance range, permissible load").

When using shakers inside conditioning cabinets, make reference to ambient conditions (see chapter "Safety, Operating company's obligations", section "Installation site").

Operation

When using shakers inside conditioning cabinets, make reference to ambient conditions (see chapter "Safety, Operating company's obligations", section "Installation site").

You can choose from two shaking options:

- Continuous mode: The unit starts and stops shaking as soon as you switch it on or off.
- Timer mode: The time for shaking can be set between 0 and 120 minutes. The unit stops shaking when the time has elapsed.

Continuous mode

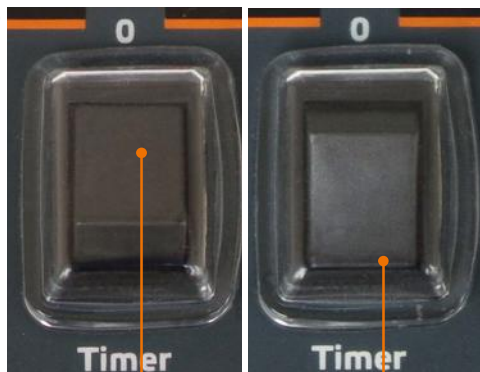
- ✓ Select mode switch is positioned to continuous mode „0”.
- Turn item on with ON/OFF switch.
- Set speed with rotating knob „Speed”.
- ✓ Actual speed appears in the 3-digit digital display.
- Reduce speed with rotating knob.
- Turn item off with ON/OFF switch.

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Timer mode

- Select mode „Timer” with select mode switch.
- Set time with rotating knob „Timer”.
- Set speed with rotating knob „Speed”.
- Turn item on with ON/OFF switch.
- ✓ When time has elapsed an acoustic signal will be audible and shaking action will stop.
- Reduce speed with rotating knob.
- Turn item off with ON/OFF switch.

Select mode switch: "continuous"

Select mode switch: "timer"

› Cleaning

All surfaces of the unit allow for cleaning with a damp cloth and if required a mild soap lotion.

Warning:
Risk of electric shock!

If the device is still connected to power during cleaning, you might get in contact with live parts due to accidentally entering liquids.

Prior to cleaning:

- ensure that the device is switched OFF
- ensure that the plug is pulled from the electric socket.

Caution:
Damage to the device

Accidentally entering liquids will cause the device to fail .
Only wipe with a damp cloth. Never use a soaked cloth.

Attention:
Risk of damage to the surfaces

All of the surfaces may be damaged by improper cleaning.

Never use:

- Chlorine bleach or any cleaning agent containing chlorine
- Solvent-bearing agents (e.g. acetone)
- Ammonia
- Abrasive cleaning agents such as scouring agents, cleaning wool or agents with metallic particles

As a result of vibrations generated during operation the device might slide off the table top.

To ensure a stable position clean surface and feet with a damp cloth and ethanol regularly.

› Maintenance

The unit is maintenance-free. Any necessary repair must be performed by an authorized Heidolph distributor.

Please contact Heidolph Instruments or your local Heidolph Instruments distributor.

➤ Possible errors and how to resolve

Malfunctions	Possible reason	Troubleshooting
On-/Off switch does not enlighten	No power	<ul style="list-style-type: none"> ▪ Check size and shape of plug and compatibility to your electrical socket ▪ Check circuit breakers
	Light diode defective	Contact your local Heidolph Instruments distributor
Unit stops shaking	Pre-set timer elapsed	Restart timer, if necessary
	Thermal motor circuit breaker triggered by motor overload.	Wait 20 minutes, decrease overall load
	<ul style="list-style-type: none"> ▪ Mechanical parts broken ▪ Electronic failure 	Contact your local Heidolph Instruments distributor

If you experience a malfunction which can not be resolved, please contact your authorized Heidolph Instruments distributor immediately.

Electrical connections

- Fuses must **only** be replaced by a **professional electrician**.
- **Repairs** to the device must **only be carried out by a qualified electrician approved by Heidolph Instruments**. Improper repairs could result in serious dangers. Please contact your specialist dealer if the device needs to be repaired.

Warning: Risk of electric shock

The device must be connected to an earthed power socket.

The rated voltage for the device and the supply voltage must be the same. The rated voltage for the device can be found on the rating plate on the rear of the device.

The device is supplied with a grounded power cord. In case the original plug is replaced, the new plug must have a protective conductor!

If the device is still connected to power while changing the fuse, you might get in contact with live parts.

Prior to replacing the fuse ensure that the unit is switched off and pull the plug from the electric socket. Only use original spare fuses from Heidolph Instruments.

After changing a fuse check the device for safe operating condition according to IEC 61010-1.

Rating plate

Fuse

Color code for electrical connections:

Europe

GREEN/YEL-LOW	PE: Protective conductor (Earth)
BLUE	N: Neutral conductor
BROWN	P: Phase

North America

GREEN	PE: Protective conductor (Earth)
WHITE	N: Neutral conductor
BLACK	P: Phase

The device is connected to the mains with the power cord supplied. The power connection is located on the rear of the housing.

For countries where a different plug than the standard plug supplied is used:

- The plug supplied may only be changed by a professional electrician.
- If you wish to use an adapter, ensure that it is grounded and approved in accordance with the local regulations.
- Wiring, adapter and plug must at least meet the electrical power stated on the rating plate.

➤ Dismantling, transportation and storage

Dismantling

**Warning:
Risk of injury!**

While the device is plugged in you might be at harm if the device is accidentally switched to ON mode.

Prior to dismantling device ensure that the device is switched off and the plug is pulled from the electric socket.

- ✓ The device is switched off.
- Disconnect plug from power socket.
- Disconnect plug from socket on the device.
- Remove all vessels.
- Disassemble attachments.

Transportation and storage

- Store and transport the device and its components only if they were emptied, cleaned and dismantled.
- To carry the device wear safety gloves and lift the unit from below.
- Store and transport the device and parts of device in the original packing material or alternatively in an adequate container to prevent damage.
- Seal the package carefully against unauthorized or accidental opening.
- Store the device in a dry and frost-free place.

Improper storage and transportation may result in damages to the system and the mechanical components.

Avoid any kind of shocks during transportation.

➤ Scope of delivery

Component	Variant	Quantity	Product number
Hei-MIX (one of these variants)	Titramax 100*	1	544-11200-00
	Titramax 101*	1	544-11300-00
	Titramax 1000*	1	544-12200-00
	Vibramax 100*	1	544-21200-00
	Vibramax 110*	1	544-31200-00
	Rotamax 120*	1	544-41200-04
Operating Manual English / German		1	01-005-002-34
Warranty registration / Confirmation of condition		1	01-006-002-78
EC Declaration of Conformity		1	01-001-025-06
Power cord		1	depending on country

* The product numbers are valid for 230 V devices in the european community. Please contact Heidolph Instruments for further product numbers on country specific varieties.

➤ Accessories

Component		Quantity	Product number
Fuse for 230 V (device needs 2 fuses)		1	14-002-015-23
Fuse for 115 V (device needs 2 fuses)		1	14-002-015-45
Vibramax 100 Rotamax 120	Attachment with 2 clamping rollers	1	549-81000-00
	Clamping roller (additional)	1	11-008-007-08
Vibramax 110	Attachment for 49 test tubes, 12 mm Ø	1	549-82000-00
	Attachment for 36 test tubes, 16 mm Ø	1	549-83000-00

Additional accessories are shown in the general catalog or on our website at www.heidolph-instruments.com

➤ Technical data

Plarform shakers Hei-MIX

Standard supply voltage	230-240 V 50/60 Hz oder 115 V/60 Hz)
Device fuses	M 1,25 A (230 V) oder T 2,0 A (115 V)
Protective class (IEC 61140)	1 ⊕
Protection class (IEC 60529)	IP30
Sound pressure level (dB(A)) (based on IEC 61010)	< 50
Overheat protection	self-resetting
Speed control	analog
Operation mode	timer / continuous

Modell Hei-MIX


	Titramax 100/101	Titramax 1000	Vibramax 100	Vibramax 110	Rotamax 120
Dimensions (wxdxh) (mm)	245x310x125	320x375x125	245x310x125	245x310x146	245x310x125
Load capacity (kg)	2	5	2	2	2
Drive	shaded pole motor			asyn- chronus motor	shaded pole motor
Motion	circular vibration			circular vibration	rotation
Orbit (mm)	1,5 / 3	1,5	3	1,5	20
Rotation speed (rpm)	150 - 1.350			150 - 2.500	20 - 300
Supply power (W)	31			46	33
Weight (kg)	5,5	6,5	5,5	5,0	5,5


Performance range / admissible load

Hei-MIX Titramax 100/101/1000

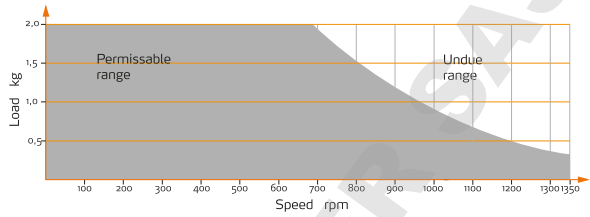
Hei-MIX Vibramax 100

Hei-MIX Vibramax 110

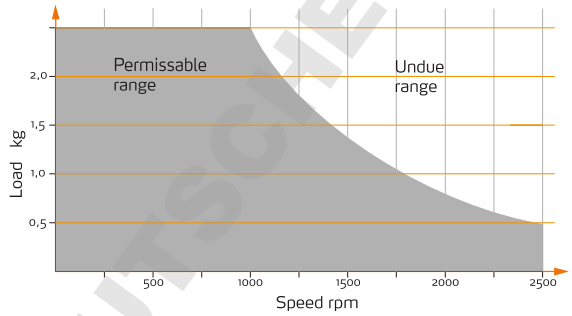Hei-MIX Rotamax 120

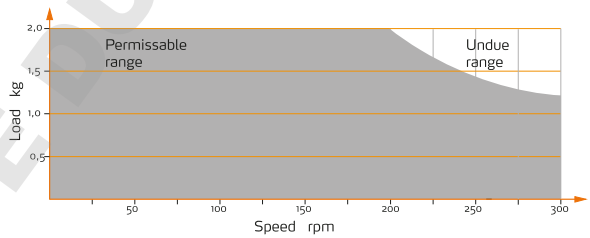➤ Contact / Technical Service

Questions / Repair work

If any aspect of installation, operation or maintenance remains unanswered in the present manual, please get in touch with the following address.

For repairs please call Heidolph Instruments or your local authorized Heidolph Instruments distributor.

Warning! **Danger of poisoning!**

Contaminated devices can lead to severe injury or death of our employees!

When shipping items for repair that may have been contaminated by hazardous substances, please:

- advise exact substance
- take adequate protective action towards our parts receiving and service personnel
- mark the pack in accordance with Ordinance on Hazardous Substances

A "Confirmation of Condition" form can be found at the end of this operating manual.

→ Prior to shipping the device for repair, complete a copy of this form and submit it in advance.

Contact details:

Heidolph Instruments Germany

Heidolph Instruments GmbH & Co. KG
Technical Service
Walpersdorfer Str. 12
D-91126 Schwabach / Germany

Tel.: +49 – 9122 - 9920-74
Fax: +49 – 9122 - 9920-84

E-Mail: service@heidolph.de
www.heidolph.com

Heidolph Instruments North America

Phone: 1-866-650-9604
E-mail: service@heidolph.com
www.heidolphNA.com

Heidolph Instruments United Kingdom

Phone: 01799 - 5133-20
E-mail: service@radleys.co.uk
www.heidolph-instruments.co.uk

All other countries

You will find contact details of your local Heidolph distributor at www.heidolph.com

➤ Warranty

Heidolph Instruments provides a 3 year warranty for the products described here (excluding glass and wearing parts) if you register using the warranty card enclosed or online (www.heidolph.com). The warranty period begins with the date of purchase. The serial number is also valid without registration.

This warranty covers material and manufacturing defects.

In the event of a material or manufacturing defect, the device shall be repaired or replaced free of charge under the terms of the warranty.

Heidolph Instruments shall not assume liability for any damage incurred as a result of improper handling or transport.

Warranty claim?

- Please notify Heidolph Instruments or your local Heidolph distributor if you wish to make a warranty claim.

➤ Confirmation of condition

→ In the case of repair, copy and complete the Confirmation of condition form and send it to Heidolph Instruments.

1. Details about the device

Product number _____

Serial number _____

Reason for repair _____

2. Has the device been cleaned, decontaminated/sterilized?

Yes _____ No _____

3. Is the device in a condition which does not represent any health threats for the staff of our service department?

Yes _____ No _____

If not, which substances has the device come into contact with?

4. Legally binding declaration

The customer is aware of being legally liable to Heidolph Instruments for any damages arising from incomplete and incorrect information.

Date

Signature

Company stamp

Please note

The shipper is responsible for the return of the goods in well packed condition, suitable for the mode of transport.

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This is a translation of the original operating manual. We reserve the right to make any technical modifications without prior notification. This documentation is not subject to revision; the latest version of this documentation can be found on our homepage. The copyright for all texts and images is held by Heidolph Instruments GmbH & Co. KG.
DOMINIQUE DUTSCHER SAS
© Heidolph Instruments GmbH & Co. KG