


### Short explanation to close the lid at the A-BOX

- Put all Microplates inside the A-Box.
- Put the lid on, there is a short and a long side of the lid.
- Position the lid onto the square box, make sure the clamps are position correctly.
- Turn the knobs as per drawing. Turn A=B a few times, then C and Da few times then A+B again etc. until the lid is fixed upon the jar.
- Verify the lid is positioned correctly and give small turns to all knobs for the last time again 1 or 2 times.

Connect the Snap Shut coupling to the device BioMax jar tube or use a GasPak to achieve anaerobic conditions.

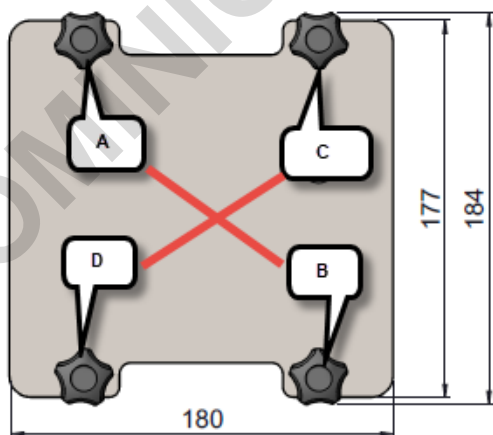

 - BOX

