

LUNA-FX™ Calibration Beads Kit

F73101

Product Description

Appearance	Colorless beads in clear liquid
Bead content	PMMA
Excitation/emission	460 nm / 520 nm, 555 nm / 605 nm

Storage

4°C in the dark

✓ F73102
LUNA-FX™
Calibration GF Beads

✓ F73103
LUNA-FX™
Calibration RF Beads

The LUNA-FX™ Calibration Beads are specialized to calibrate the LUNA-FX7™ Automated Cell Counter.

Directions for Use

1. Each time mix the LUNA-FX™ Calibration Beads thoroughly to ensure a homogeneous suspension before loading into a slide.
2. Load 10 µL of the LUNA-FX™ GF Calibration Beads into the chamber A, and 10 µL of the LUNA-FX™ RF Calibration Beads into the chamber B of a LUNA PhotonSlide™.
3. Wait for 30 seconds for the beads to settle down.
4. Insert the slide face up and chamber A first into the slide port. Follow the calibration directions of the LUNA-FX7™ Automated Cell Counter.



HEADQUARTERS

FL 3
28 Simindaero 327beon-gil, Dongan-gu
Anyang-si, Gyeonggi-do 14055
South Korea

Tel: +82 (31) 478-4185

USA

7700 Little River Turnpike STE 207
Annandale, VA 22003
USA

Tel: +1 (703) 622-4660, +1 (703) 942-8867

EUROPE

1 allée Lavoisier 59650 Villeneuve d'Ascq
France

Tel: +33 (0)3 74 09 44 35

www.logosbio.com

Disclaimer

This product is for research use only.
Please consult the material safety data sheet for information regarding hazards and safe handling practices.

Additional information is available on our website www.logosbio.com.
© Logos Biosystems 2020. All rights reserved.

LUNA-FX™ Calibration Beads Kit

F73101

Product Description

Appearance	Colorless beads in clear liquid
Bead content	PMMA
Excitation/emission	460 nm / 520 nm, 555 nm / 605 nm

Storage

4°C in the dark

✓ F73102
LUNA-FX™
Calibration GF Beads

✓ F73103
LUNA-FX™
Calibration RF Beads

The LUNA-FX™ Calibration Beads are specialized to calibrate the LUNA-FX7™ Automated Cell Counter.

Directions for Use

- Each time mix the LUNA-FX™ Calibration Beads thoroughly to ensure a homogeneous suspension before loading into a slide.
- Load 10 µL of the LUNA-FX™ GF Calibration Beads into the chamber A, and 10 µL of the LUNA-FX™ RF Calibration Beads into the chamber B of a LUNA PhotonSlide™.
- Wait for 30 seconds for the beads to settle down.
- Insert the slide face up and chamber A first into the slide port. Follow the calibration directions of the LUNA-FX7™ Automated Cell Counter.



HEADQUARTERS

FL 3
28 Simindaero 327beon-gil, Dongan-gu
Anyang-si, Gyeonggi-do 14055
South Korea

Tel: +82 (31) 478-4185

USA

7700 Little River Turnpike STE 207
Annandale, VA 22003
USA

Tel: +1 (703) 622-4660, +1 (703) 942-8867

EUROPE

1 allée Lavoisier 59650 Villeneuve d'Ascq
France

Tel: +33 (0)3 74 09 44 35

www.logosbio.com

Disclaimer

This product is for research use only.
Please consult the material safety data sheet for information regarding hazards and safe handling practices.

Additional information is available on our website www.logosbio.com.
© Logos Biosystems 2020. All rights reserved.