
Gebrauchsanleitung | Operating manual |  
Mode d'emploi | Instrucciones de manejo |  
Istruzione | Instruções | 操作手册

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Titrierapparate | Burettes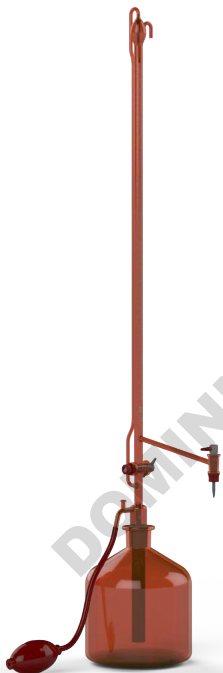

# Impressum

BRAND GMBH + CO KG

Otto-Schott-Str. 25

97877 Wertheim (Germany)

T +49 9342 808 0

F +49 9342 808 98000

info@brand.de

www.brand.de

## Languages

|                               |     |
|-------------------------------|-----|
| Gebrauchsanleitung.....       | 2   |
| Operating manual .....        | 19  |
| Mode d'emploi .....           | 35  |
| Instrucciones de manejo ..... | 51  |
| Istruzione.....               | 68  |
| Instruções.....               | 86  |
| 操作手册 .....                    | 103 |

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DOMINIQUE DUTSCHER SAS

# Table of contents

|  |           |
|--|-----------|
| <b>1 Introduction</b> .....                          | <b>20</b> |
| 1.1 Scope of delivery .....                          | 20        |
| 1.2 Terms of use .....                               | 20        |
| <b>2 Safety instructions</b> .....                   | <b>22</b> |
| <b>3 Functional and operational components</b> ..... | <b>23</b> |
| <b>4 Operation</b> .....                             | <b>24</b> |
| 4.1 Preparing the burette .....                      | 24        |
| 4.2 Filling the burette.....                         | 24        |
| 4.3 Deaerating the burette .....                     | 24        |
| 4.4 Titration .....                                  | 25        |
| 4.5 Post-titration .....                             | 26        |
| <b>5 Cleaning</b> .....                              | <b>27</b> |
| <b>6 Recommendation</b> .....                        | <b>27</b> |
| <b>7 Ordering Information</b> .....                  | <b>28</b> |
| 7.1 Accessories .....                                | 28        |
| 7.2 Spare parts .....                                | 29        |
| <b>8 Warranty</b> .....                              | <b>31</b> |
| <b>9 Repairs</b> .....                               | <b>32</b> |
| 9.1 Sending for repair .....                         | 32        |
| <b>10 Disposal</b> .....                             | <b>34</b> |

# 1 Introduction

## 1.1 Scope of delivery

BLAUBRAND® or SILBERBRAND model Pellet-type burette with PTFE valve stopcock or PTFE key; models with or without intermediate stopcock (PTFE key); supply bottle is packaged separately, if applicable; BLAUBRAND® models with quality certificate.

## 1.2 Terms of use

- Please carefully read the operating manual before using the device for the first time.
- The operating manual is part of the device and must be kept in an easily accessible place.
- Be sure to include the operating manual if you transfer possession of this device to a third party.
- You can find up-to-date versions of the operating manual on our website: [www.brand.de](http://www.brand.de).

### 1.2.1 Hazard levels

The following signal words identify possible hazards:

| Signal word | Meaning                                 |
|-------------|---|
| DANGER      | Will lead to serious injury or death.   |
| WARNING     | May lead to serious injury or death.    |
| CAUTION     | May lead to minor or moderate injuries. |
| NOTICE      | May lead to property damage.            |

### 1.2.2 Symbols

| Symbol  | Meaning        |
|---|----------------|
|  | Hazardous area |

## 1.2.3 Format

| Layout         | Meaning                                   | Layout | Meaning                |
|----------------|---|--------|------------------------|
| <b>1. Task</b> | Indicates a task.                         | >      | Indicates a condition. |
| a., b., c.     | Indicates the individual steps of a task. | ⇒      | Indicates a result.    |

DOMINIQUE DUTSCHER SAS

## 2 Safety instructions

### **Please read carefully!**

The can be used in combination with hazardous materials, work processes and equipment. However, the operating manual cannot cover all of the safety issues that may occur in doing so. It is the user's responsibility to ensure compliance with the safety and health regulations and to specify the corresponding restrictions before use.

1. Follow the general instructions for hazard prevention and the safety regulations. E.g., wear protective clothing, eye protection and protective gloves.
2. Follow the instructions given by the reagent manufacturer.
3. Every user must read this operating manual before using the instrument.
4. Always perform work in a way that does not endanger the user or any other person.
5. Use the instrument only for titrating and only within the limits of its material resistance.
6. Never use force.
7. To unstick stuck stopcocks or ground joint connections, follow the cleaning instructions (see Cleaning, p. 27).
8. Use only original accessories and original replacement parts.

## 3 Functional and operational components

**1** Automatic zero point setting; located beside it is the overflow.

**3** Supply bottle

**5** NS 29/32 ground joint connection, olive and vents

**2** Burette stopcock, pictured here is a PTFE key. Other models have a valve stopcock.

**4** Rubber blowball\*

**6** Intermediate stopcock; located behind it is the filling tube.

\* This item is not included in the scope of delivery.

## 4 Operation

### 4.1 Preparing the burette

- a. Check whether the stopcocks are easy to turn.
- b. Lightly grease the ground joint connection to the supply bottle with stopcock grease, or use a PTFE ground joint sleeve.

### 4.2 Filling the burette

#### WARNING

#### Media leakage

In general, do not fill while the intermediate stopcock is open. If overfilled, media will escape through the overflow at the top of the burette.

- a. Place the burette onto the filled supply bottle.
- b. Connect the rubber blowball to the side tubing connector.
- c. Close the intermediate stopcock (model with intermediate stopcock).
- d. Hold the vent closed with a finger. Squeeze the blowball to fill the burette.
- e. When the liquid level is approx. 5 mm above the zero mark, lift your finger off the vent.
  - ⇒ This ends the process.
  - ⇒ The burette will automatically set to zero.

### 4.3 Deaerating the burette

- a. To remove trapped air, briefly open and close the burette stopcock.

- b. After the first filling operation, a small air bubble may remain on the burette stopcock.
- c. To remove this bubble, hold the burette at an angle and lightly tap a finger on the spot where the bubble is located.

## 4.4 Titration

### Preventing volume errors

- a. Begin titration from the zero point.  
⇒ This prevents volume errors.
- b. Turn the valve stopcock gently when closing. Excessive force may lead to deformation of the valve spindle, resulting in an unwanted rise of the liquid column in the burette tube (reading error).

### Observing wait times

#### NOTICE

If titration is carried out, the wait time can be omitted since the time required to reach the titration end point is normally longer than the specified wait time.

| Class | Wait time in accordance with DIN EN ISO 385 |
|-------|---|
| AS    | 30 s  |
| B     | Not applicable                              |

When calibrating Class A burettes, the wait time according to DIN EN ISO 385 must be adhered to.

## 4.5 Post-titration

On burettes with an intermediate stopcock, the content of the burette can be allowed to flow back into the supply bottle. It should be noted that liquid residue on the wall may partially evaporate.

### **Preventing volume errors**

To prevent volume errors, fill and empty the burette one to two times before the next titration.

### **Storing the burette**

Clean the burette if you will not be using it for several days. If it is not cleaned, stopcocks and ground joint connections may get stuck.

## 5 Cleaning

### CAUTION

#### **Risk of injury from broken glass!**

Never use excessive force to unstick stuck stopcocks or ground joint connections.

### Cleaning

If the burette does not discharge without dripping or if it sets the zero point too low, the burette must be cleaned. To clean, fill the burette with a mild alkaline cleaner and rinse it out with water several times after the prescribed soaking time.

### Unsticking stuck parts

- a. Completely immerse stuck stopcocks or ground joint connections in a cleaning solution and allow to soak overnight.
- b. Carefully unstick stuck stopcocks or ground joint connections while wearing suitable safety gloves.
- c. If unsuccessful, send the instrument in for repair.

## 6 Recommendation

We recommend using a PTFE ground joint sleeve for the ground joint connection to the supply bottle, as these do not require grease for sealing.

## 7 Ordering Information

### 7.1 Accessories

#### Rubber blowball

| Description     | Order No. |
|-----------------|-----------|
| Rubber blowball | 234 00    |

#### Ground joint sleeve

| Description              | Packaging unit | Order No. |
|--------------------------|----------------|-----------|
| PTFE ground joint sleeve | 10 pc          | 514 22    |

#### Ground joint socket

| Description                       | Order No. |
|-----------------------------------|-----------|
| PTFE ground joint socket NS 29/32 | 514 67    |

#### Joint clip NS 29/32

| Description     | Order No. |
|-----------------|-----------|
| Stainless steel | 556 18    |
| POM (red)       | 556 46    |

#### Supply bottle 2000 ml, NS 29/32

| Description            | Order No. |
|------------------------|-----------|
| Boro 3.3, clear        | 233 10    |
| Soda-lime glass, clear | 126 965   |
| Boro 3.3, amber        | 233 20    |
| Soda-lime glass, amber | 127 065   |

## 7.2 Spare parts

### 7.2.1 Clear glass models

#### Burette stopcock, PTFE (10, 25, 50 ml)

| Description                            | Order No. |
|--|-----------|
| Burette stopcock, PTFE (10, 25, 50 ml) | 821 70    |

#### Burette stopcock, PTFE with housing and tip

| Description  | Order No. |
|--|-----------|
| Burette stopcock, PTFE with housing and tip 10 ml, Boro 3.3, clear | 821 20    |
| Burette stopcock, PTFE with housing and tip 25 ml, Boro 3.3, clear | 821 21    |
| Burette stopcock, PTFE with housing and tip 50 ml, Boro 3.3, clear | 821 22    |

#### Intermediate stopcock, PTFE key 4NS/19

| Description                            | Order No. |
|--|-----------|
| Intermediate stopcock, PTFE key 4NS/19 | 812 65    |

## 7.2.2 Amber glass models

### Burette stopcock, PTFE key 3NS/12 incl. PTFE-tip (25, 50 ml)

| Description  | Order No. |
|--|-----------|
| Burette stopcock, PTFE key 3NS/12 incl. PTFE-tip (25, 50 ml) | 811 67    |

### Burette stopcock, PTFE key 3NS/12 incl. PTFE-tip and housing

| Description  | Order No. |
|--|-----------|
| Burette stopcock, PTFE key 3NS/12 incl. PTFE-tip and housing (25, 50 ml, Boro 3.3 amber) | 819 27    |

### Intermediate stopcock, PTFE key 4NS/14

| Description                     | Order No. |
|---------------------------------|-----------|
| Intermediate stopcock, PTFE key | 812 70    |

## 8 Warranty

We shall not be liable for the consequences of improper handling, use, servicing, operating or unauthorized repairs of the device or for the consequences of normal wear and tear, especially of wearing parts such as pistons, seals, valves and the breakage of glass. The same applies for failure to follow the instructions of the operating manual. We are not liable for damage resulting from disassembly beyond that described in the operating manual or if non-original spare parts or components have been installed.

### **USA and Canada:**

Find more warranty information on [www.brandtech.com](http://www.brandtech.com).

## 9 Repairs

### 9.1 Sending for repair

#### NOTICE

Transporting of hazardous materials without a permit is a violation of federal law.

---

#### **Clean the instrument thoroughly and decontaminate!**

- When returning products, please enclose a general description of the type of malfunction and the media used. If information regarding media used is missing, the instrument cannot be repaired.
- Shipment is at the risk and the cost of the sender.

#### **Outside USA and Canada**

Complete the “Declaration on Absence of Health Hazards” and send the instrument to the manufacturer or supplier. Ask your supplier or manufacturer for the form. The form can also be downloaded from [www.brand.de](http://www.brand.de).

#### **Within USA and Canada**

Please clarify the requirements for the return delivery with BrandTech Scientific, Inc **before** sending the instrument in for service.

Return only cleaned and decontaminated instruments to the address provided with the Return Authorization Number. Place the Return Authorization number so that it is clearly visible on the outside of the package.

## Contact addresses

**Germany:**

BRAND GMBH + CO KG  
Otto-Schott-Str. 25  
97877 Wertheim (Germany)  
T +49 9342 808 0  
F +49 9342 808 98000  
info@brand.de  
www.brand.de

**USA and Canada:**

BrandTech® Scientific, Inc.  
11 Bokum Road  
Essex, CT 06426-1506 (USA)  
T +1-860-767 2562  
F +1-860-767 2563  
info@brandtech.com  
www.brandtech.com

**India:**

BRAND Scientific Equipment Pvt. Ltd.  
303, 3rd Floor, 'C' Wing, Delphi  
Hiranandani Business Park,  
Powai  
Mumbai-400 076 (India)  
T +91 22 42957790  
F +91 22 42957791  
info@brand.co.in  
www.brand.co.in

**China:**

BRAND (Shanghai) Trading Co., Ltd.  
Guangqi Culture Plaza  
Room 506, Building B  
No. 2899, Xietu Road  
Shanghai 200030 (P.R. China)  
T +86 21 6422 2318  
F +86 21 6422 2268  
info@brand.com.cn  
www.brand.cn.com

## 10 Disposal

Before disposal, observe the relevant national disposal regulations and ensure that the product is disposed of properly.

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BRAND Scientific Equipment Pvt.Ltd.