


TFD 4

FOAMING DEGREASING DETERGENT

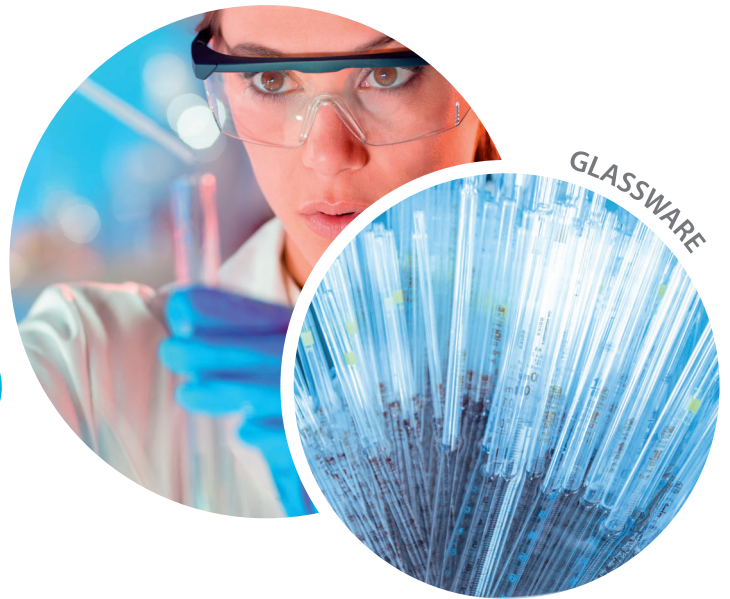

■ PRODUCT BENEFITS

- A free alkalinity and a wetting ability allow a powerful surfactant and sequestering activity
- Perfect rinsability even in presence of hard water
- Action through a saponification activity instead of a simple emulsification
- Perfect dissolution of paraffin
- Stronger detergent action for ultrasound vats.



FOAMING DEGREASING DETERGENT

■ INSTRUCTIONS FOR USE

- Wear gloves, goggles, apron.

■ Classic washing:

- 1- Prepare a TFD 4 solution (3 to 5%) in water at room temperature
 - 2- Immerse material for 15 minutes
 - 3- Scrub if necessary
 - 4- Thoroughly rinse with water
- Renew the action if necessary*

■ Reinforced cleaning:

- 1- Prepare a TFD 4 solution (5 to 10%) in hot water (about 40°C)
 - 2- Immerse material for 15 minutes
 - 3- Thoroughly rinse with water
- Renew the action if necessary*

■ PRECAUTIONS OF USE

See Safety Data Sheet (SDS 010104)

■ FEATURES

- Aspect: **Liquid**
- Color: **Whitish**
- pH at 3%: **12.1**
- pH at 5%: **12.2**
- pH at 10%: **12.5**
- Pure pH: **13.3**
- Expiration date: **4 years**
- Conservation after opening: **6 months**

■ COMPOSITION

- Non-ionic and anionic surfactants
- Sequestrants

■ PACKAGING

- 1 Box of 4 x 5L.....ref. **01010420**
- 2 1 x 20L.....ref. **01010421**



Hazardous product - Follow precautions of use. Always read label and product information before use.