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Fournisseur: Merck

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Anaerocult® A for microbiology

Reagent for the generation of an anaerobic medium in an anaerobic jar.

Anaerocult® C system

The Anaerocult® C system is used specifically for the incubation of microaerophiles and capnophiles. The special formulation of the reagent mixture is designed so that the cultivation jar reaches gas concentrations of 8-10% vol. CO<sub>2</sub> and 5-6% vol. O<sub>2</sub>. At an incubation temperature of 42°C, the temperature in the jar does not increase further because, unlike systems using a catalyst, the amount of energy released is low, so the microorganisms are not harmed by the heat.

Anaerocult® C is used to generate oxygen-depleted and CO<sub>2</sub>-enriched atmosphere in a 2.5 L anaerobe jar for culturing *Campylobacter* species and other microorganisms with fastidious requirements (e.g. *Neisseria* sp., *Capnocytophaga* sp., *Eikenella corrodens*, *Haemophilus* sp.).

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Gas generator system for the incubation of one to four petri dishes in an anaerobic atmosphere.

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For microbiology gas generator system for incubation of one to two Petri dishes in a oxygen-depleted and CO<sub>2</sub>-enriched atmosphere.

Anaerocult® P for microbiology

Reagent for the generation of an anaerobic medium for one petri dish.

Informations sur les accessoires: Anaerotest® strips are designed as an indicator strip to indicate an anaerobic atmosphere. Anaerobic atmosphere favors growth of anaerobic bacteria and this strip can indicate the absence/presence of oxygen in the environment. The methylene blue dye in the absence of oxygen undergoes a chemical reduction reaction and turns colorless thereby indicating a total anaerobic atmosphere. The reaction zone on the test strip returns to normal (blue color) upon re-introduction of oxygen in the environment.

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