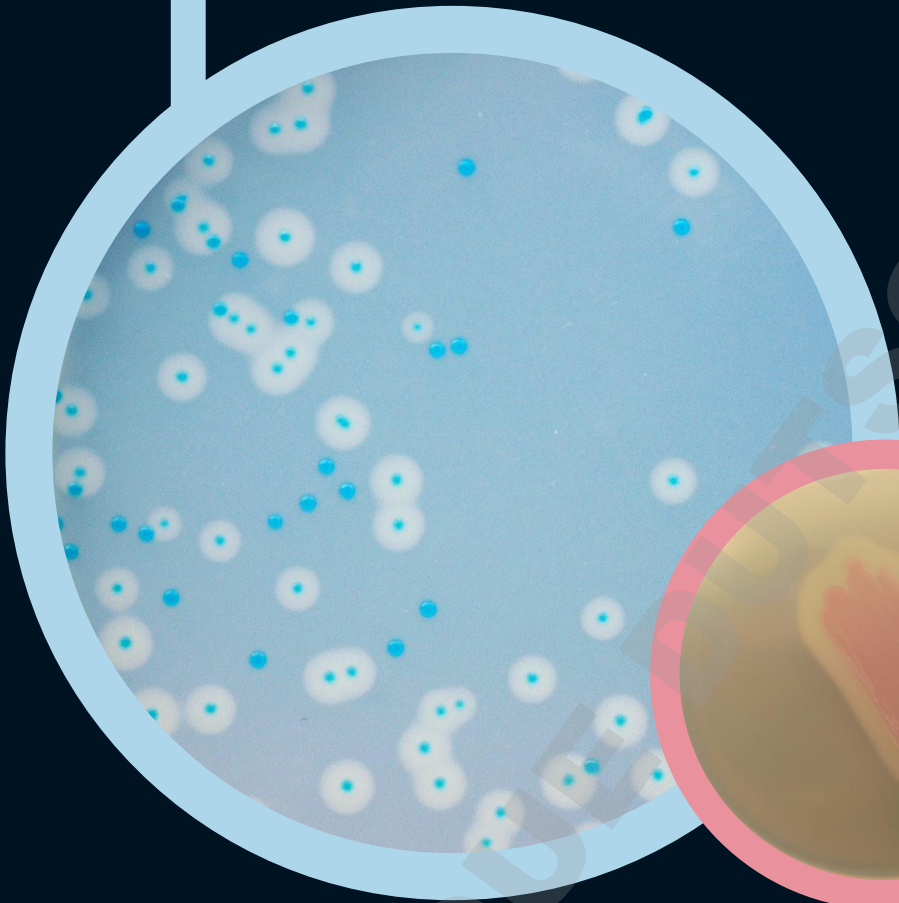
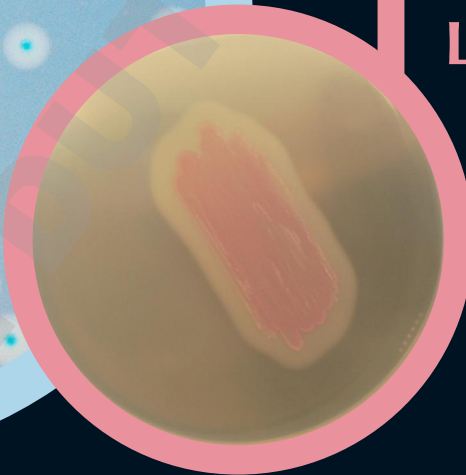
# ● CHROMagar™ Listeria

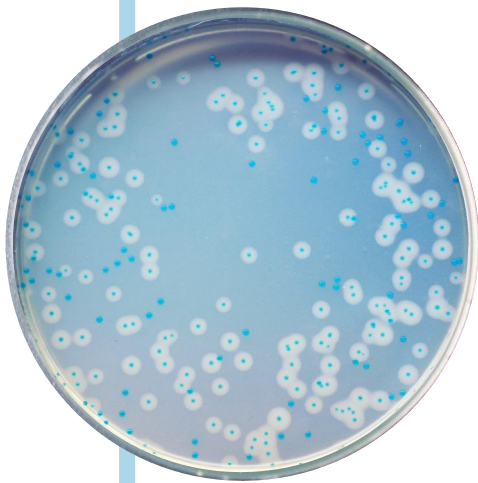
# ● CHROMagar™ Identification Listeria

For detection, enumeration and confirmation  
of *Listeria monocytogenes*

# ● CHROMagar™ Listeria Method

www.CHROMagar.com


## Plate Reading

- *L.monocytogenes*  
→ blue, diameter less than 3mm, regular and white halo

## Quality Control Strains

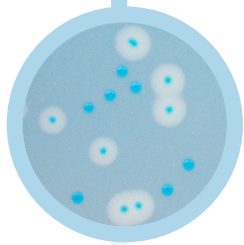- L. monocytogenes* WDCM 00021 = ATCC® 13932  
→ blue with halo
- L. innocua* CIP 8012 = ATCC® 33091  
→ blue without halo
- E. faecalis* ATCC® 29212  
→ inhibited
- E. coli* ATCC® 25922  
→ inhibited

## Plate Reading

- *L.monocytogenes*  
→ rose surrounded by a white halo

## Quality Control Strains

- L. monocytogenes* WDCM 00021 = ATCC® 13932  
→ mauve with halo
- L. innocua* CIP 8012 = ATCC® 33091  
→ mauve without halo
- L. ivanovii* ATCC® 19119  
→ colourless with halo
- B. cereus* CIP 5832 = ATCC® 14893  
→ colourless with big halo

## CHROMagar™ Listeria Method

*Listeria monocytogenes* is a widespread bacteria, present in the soil, sewage or faecal matter. Its ability to form listerial biofilms on contact surfaces makes it difficult to eliminate. This pathogen can cause serious food poisoning and is therefore frequently a microbial Q.C. target in food processing facilities to avoid food contamination. Contamination can occur at all steps of the food manufacturing chain from raw materials to place of consumption.

CHROMagar has developed a **rapid method for the detection of *L.monocytogenes*** in food stuff, comprising the following simple steps (1) enrichment in half frazer broth for 24h, (2) isolation in CHROMagar™ Listeria, (3) confirmation of *Listeria monocytogenes* species in CHROMagar™ Identification Listeria.

CHROMagar™ Listeria Method vs ISO 11290 Method:

|                              | CHROMagar Listeria Method | ISO 11290 Method (classical) |
|------------------------------|---------------------------|------------------------------|
| Number of Enrichments        | <b>1 for 24h</b>          | 2 = total 72h                |
| Number of plates Incubation  | <b>1 plate for 24h</b>    | 4 plates for 24h & 48h       |
| Number of confirmatory tests | <b>1 test</b>             | 6 tests                      |
| <b>Time to result:</b>       |                           |                              |
| Negative results             | <b>After 2 days</b>       | After 7 days                 |
| Positive results             | <b>After 3 days</b>       | After 11 days                |

## CHROMagar™ Listeria

For detection and numeration of *Listeria monocytogenes*

### Medium Performance

#### CLEAR READING

Since *L. monocytogenes* and *L. innocua* have similar biochemical properties, they cannot be differentiated on traditional media (Palcam, Oxford). CHROMagar™ Listeria helps to easily differentiate *Listeria monocytogenes* from other *Listeria* directly at the isolation step: the colonies are blue and surrounded by a white halo due to a specific phospholipase activity.

## CHROMagar™ Identification Listeria

For confirmation of *L.monocytogenes* species from suspect colonies on CHROMagar Listeria

### Medium Performance

#### SIMPLICITY/QUICK RESULTS

Classical confirmatory tests for the *L.monocytogenes* species include many tedious and time consuming steps (purification + catalase + haemolysis + dextrose + rhamnose + xylose...) CHROMagar™ Identification Listeria simplifies the species confirmatory step and reduces the workload, hence improves the efficiency of the laboratory.

A single spot of a suspect colony out of **CHROMagar™ Listeria** directly put onto **CHROMagar™ Identification Listeria**, will provide confirmation of *L. monocytogenes* species within 24 hours.

ATCC® is a registered trademark of the American Type Culture Collection

## Order References

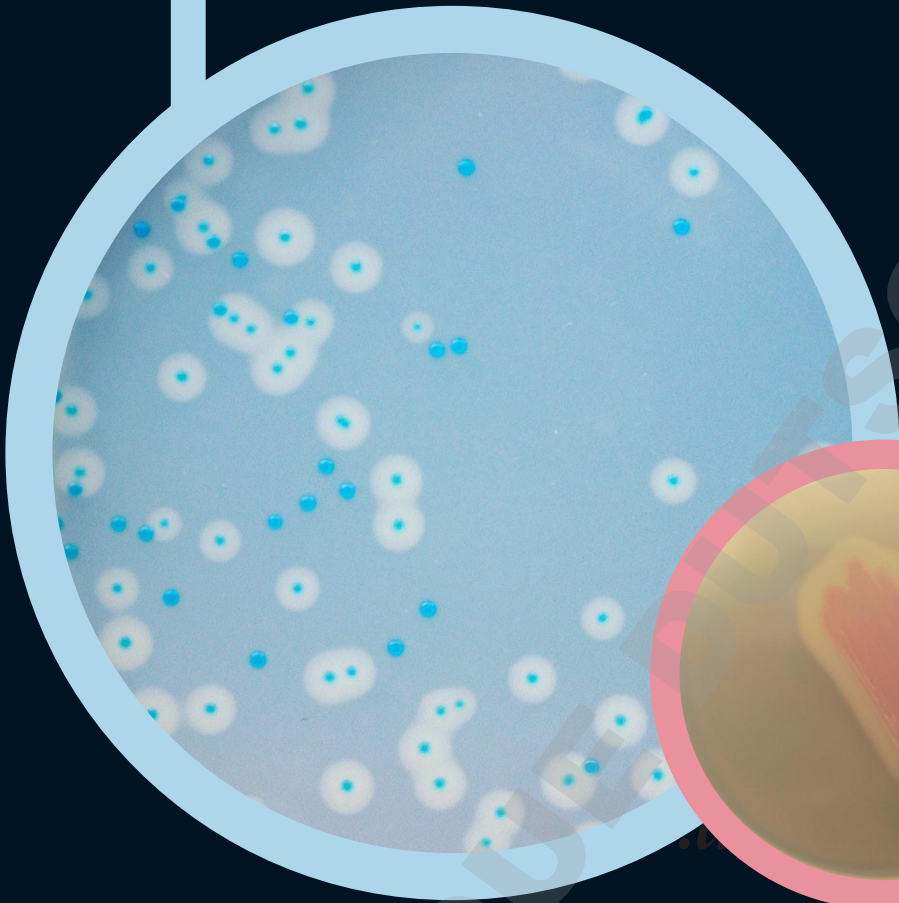
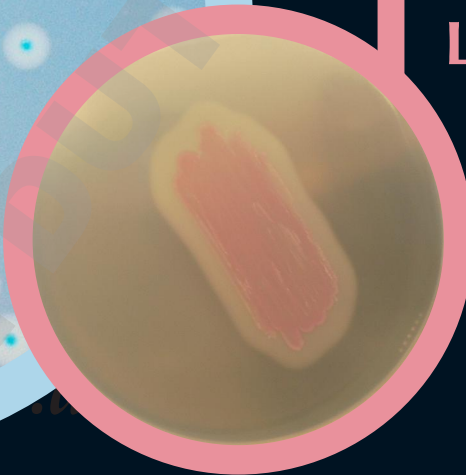
Please use these product references when contacting your local distributor:

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