52 31

Garantia

E

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GB pH electrode

Introduction

The **52 31** is a puncture electrode with stainless steel body. The unique electrode that can be used in any position.

Description / Instructions

- 1 Storage protector with some drops of Crisolylt G (KCl + glycerine). Unscrew to remove it.
- 2 316 stainless steel body. Do not buckle it, the inside body is glass made.
- 3 Diaphragms, 1 ceramic + 1 open.
- 4 Membrane sensitive. Do not rub it. Safety electrolyte, in case of breakage, the affected area can be located easily.

Specifications

| | |
|---------------------|----------------------|
| Measuring range | 2 ... 14 pH |
| Operating temp. | 0 ... 60 °C |
| Reference | encapsulated Ag/AgCl |
| Diaphragms | 1 ceramic + 1 open |
| Electrolyte | solid |
| Body material | glass / 316 s. steel |
| Min. puncture depth | 15 mm |

Applications: Cheeses, meats, fish, fruits, bread doughs, etc.

Limits: pH values < 2. Temperatures > 60 °C.

Warning

- **Do not introduce the electrode directly** in thick samples or samples containing solid materials, bones, etc. It is necessary to open a path using a stainless steel perforator.
- **Cleaning:** between samples clean the electrode with cotton soaked in alcohol. At the end, clean with lukewarm soapy water. Never use cleaning solutions with pH < 2.
- **Store** the electrode in the protector with Crisolylt G.

Certificate of quality

The attached electrode **52 31** has passed the quality test and follows the next specifications:

- Asymmetry potential < ± 15 mV.
- Sensitivity, pH 4...7 (at 25 °C) > 98%.
- Response time, pH 4...7 < 20 s.