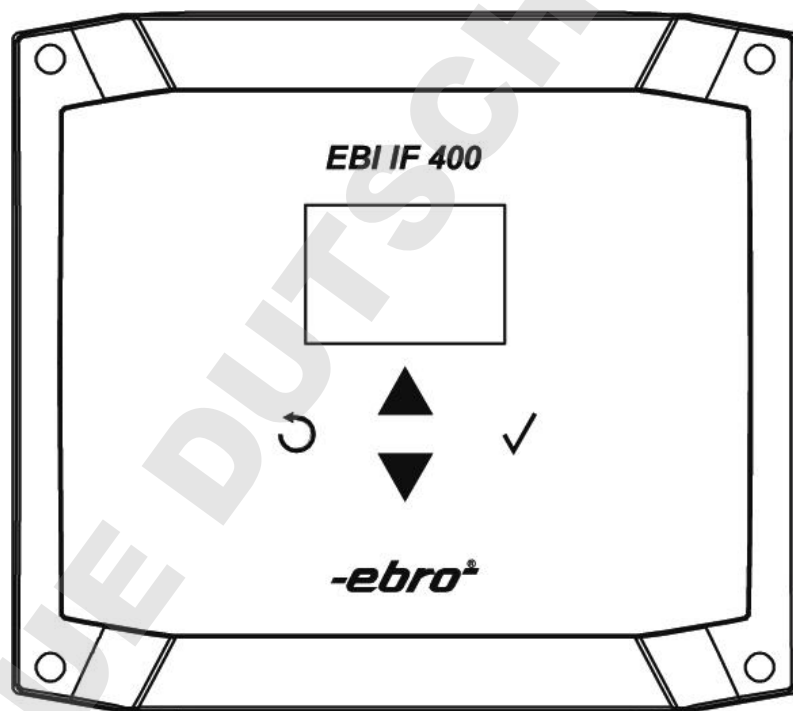
**-ebro-**  
a xylem brand

### **EBI 25-System**

Basisstation  
Interface  
Interface

EBI IF-400

## Inhalt

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## Contents

Overview .....	5
Safety instructions .....	7
Unpacking / scope of delivery .....	9
Operating the IF 400 .....	11
Connect antenna.....	11
Connecting power supply unit.....	13
PC connection.....	13
System Menu .....	15
Display .....	17
What to do, if.....	19
Cleaning and maintenance .....	25
Disposal .....	25
Technical data .....	27
Approvals.....	33
Operation.....	33

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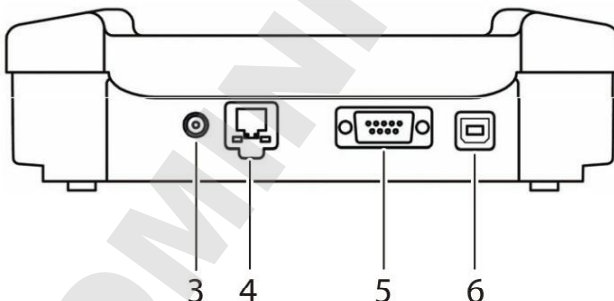
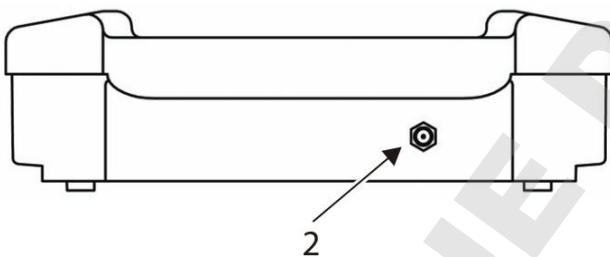
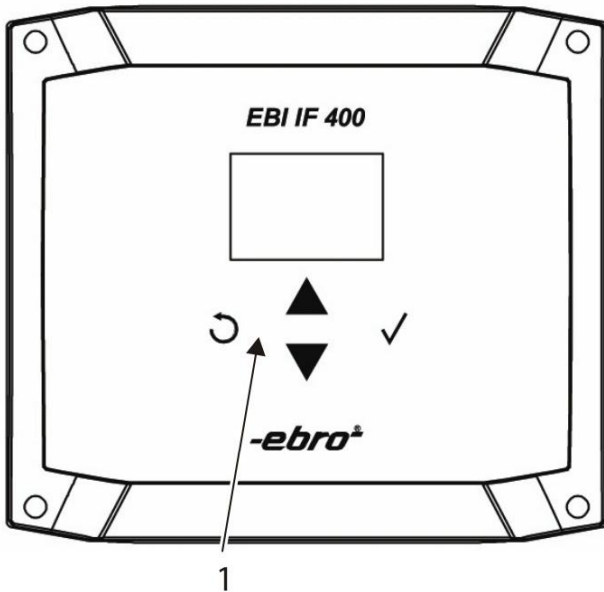
## Overview

The EBI IF-400 interface is used as receiving unit for data loggers of the EBI 25 series. The interface receives and stores measuring values from the EBI 25 loggers; the measuring values will be transmitted to a PC via USB or Ethernet port.

Up to 50 EBI 25 loggers can be registered on one interface.

Several base stations can be used within one EBI 25 wireless logger system in order to be able to use more EBI 25 loggers or to provide for a better radio coverage of large buildings.

In order to be able to work with the EBI 25 system, you will need the Winlog.wave or Winlog.web software

Structure of the interface.

- 1 Display and keyboard
- 2 Connector for antenna
- 3 Connector for power supply
- 4 Ethernet connection
- 5 Service port
- 6 USB port

For more information about how to program the data logger, and how to extract the data, please refer to the software manual of Winlog.wave or Winlog.web.

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## Safety instructions

Do not use the device in areas in which there is a danger of explosions. Danger of death!

During installation, please remember to keep the power supply easily accessible (in case you need to disconnect it quickly)!

Do not use the device in an environment hotter than 60 °C!

Please avoid direct contact with water.

The device emits wireless signals in a frequency band in the range between 779 and 928 MHz. This may cause interference to other technical devices! For this reason, only use the data loggers in areas where the wireless emission does not represent a risk!

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## Unpacking / scope of delivery

Check the device packaging and contents to ensure it is complete and undamaged. You must also check that the contents of the packaging match your order.

The delivery includes these parts:

- This user manual
- Base station EBI IF-400
- USB cable
- Power supply
- Rod antenna (AL 117)

The delivery may also include various accessories.

If you have any reason for complaint, do not hesitate to contact us. You will find our contact data on the back of this manual.

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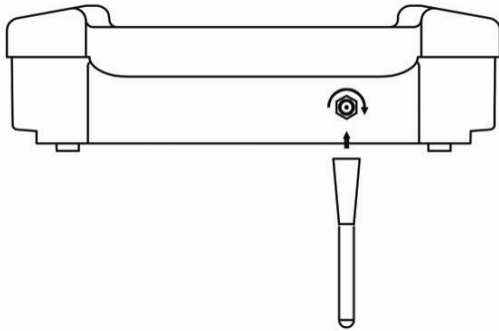
## Operating the IF 400

### Connect antenna

- Mount the antenna (AL 117) on the antenna connector on the top side of the IF 400.

Please do not use any force when mounting the antenna.

In order to maintain operational safety, please do exclusively mount antennas delivered by ebro.

The interface is delivered with a rod antenna by default.

For difficult radio environments – e.g. long distances or radio transmission through walls, you can use the Ground Plane antenna AL 116 which is optionally available.

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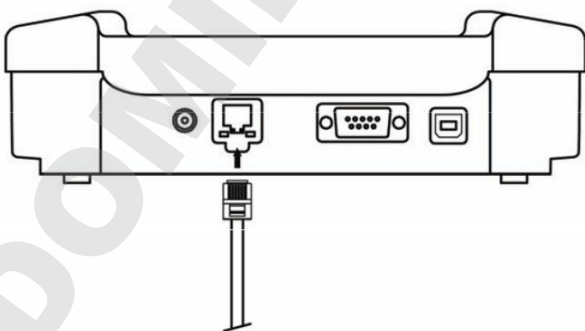
## Connecting power supply unit

Connect the power supply to the interface.

### NOTE:

In order to maintain operational safety, please do exclusively connect power supplies delivered by ebro.

As soon as the interface is supplied with power, the display will show date and time.

The internal clock of the interface is synchronized with the current time by the PC software.

## PC connection

### USB connection

A USB cable is delivered with the IF 400. Connect the plug with the rectangular cross section (Type A) with the PC. The plug with the square cross section (Type B) matches with the interface.

### Ethernet connection

Connect an RJ45 network cable to the interface and connect the interface to your network.

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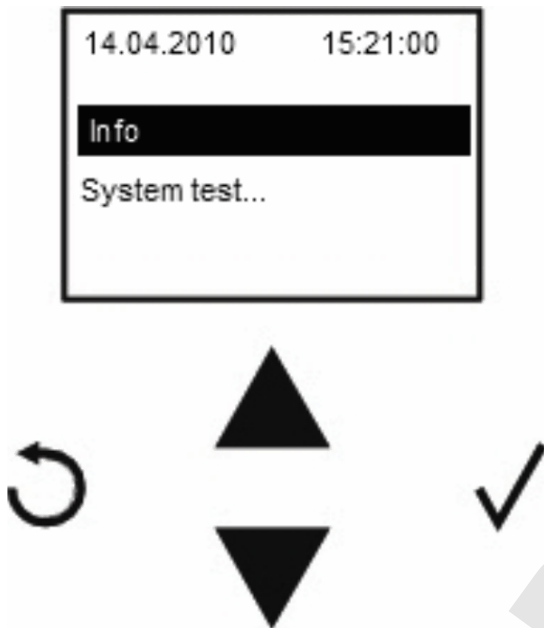
-

## System Menu

The configuration of the interface is done with the software Winlog.wave or Winlog.web.

However, you can read back status information directly from the device and perform a hardware check.

- Push ✓ in order to start the system menu.
- Use the buttons ▲ and ▼ in order to select the menu items.
- Push ✓ in order to execute the selected menu item.
- Push ↻ to exit the menu.

You have 2 menu items to choose from:

- Info  
Shows the firmware and hardware version of the device.
- Hardware check  
Runs a test of the device.
  - If the test is finished successfully, OK will be shown on the display.
  - If the test fails, the corresponding error code will be shown on the display.
- RF Settings
  - Furthermore, the area code and the area in which the base station should be operated, but also the frequency are displayed.

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## Display

During operation, the display of the interface shows the current date and time as well as the serial number and the time of the data logger that last reported to the interface. The current transmitted temperature or temperature and humidity measured value is also displayed.

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## What to do, if...

### ... the display remain dark:

In this case, the interface is not supplied with power.

- Please check if the power supply is connected properly to power socket and interface.

### ... no data logger is identified:

- Register the data logger to the base station via software Winlog.wave or Winlog.web.
- Check the base station by running the hardware check.
- Check if the antenna is connected properly.
- Check if the data loggers are still in the reception range of the base station. If necessary, please reduce the distance between logger and base station or use the Ground Plane Antenna AL 116.

### ... the hardware check fails:

- Re-start the interface: disconnect and then re-connect the power supply.
- Run the hardware check again. If the hardware check fails again, please make a note of the error code and get in contact with the ebro support.

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**... the software cannot access via USB:**

- Try to connect the interface to another USB port of your PC.  
The interface is registered on the PC as so-called USB-HID Device. You do not need to install any drivers.
- Please start the Windows Device Manager in order to check if the interface was identified by the PC.
- Click on the + symbol in the section Enter devices.
- Double-click on the first entry with USB-HID (Human Interface Device) and click on Details.
- If the device entity identification starts with USB\VID\_19BA, this is the entry for the IF 400.
- If necessary, please check all further USB-HID entries.
- If there is no adequate entry, disconnect and then re-connect the USB cable.

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**... the software cannot access via Ethernet:**

- Please check if the network cable is connected properly to the IF 400.
- There are two state LED at the network connector:
  - The right (yellow) LED shows the link state. If this LED flashes, there is a connection to the network. If the link LED does not flash, check the network connection.
  - The left (green) LED shows the data communication of the network connection. When connected properly, this LED flashes depending on the data communication.
- Check the network settings (IP address, gateway, Subnet Mask). These settings are conducted through the software Winlog.web. Default IP address of the IF 400 is 192.168.25.25, and default Subnet Mask is 255.255.255.0.

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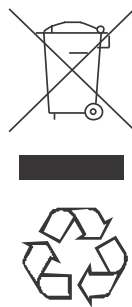
## Cleaning and maintenance

Clean the device with a slightly dampened cloth.

Never use solvents (such as acetone) for cleaning because these may attack the plastic.

## Disposal

If the device becomes no longer fit for purpose, it must be disposed of in a suitable, environmentally-friendly manner.

Do not, under any circumstances, simply dispose of it in domestic garbage. You must return it to the manufacturer.

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Dry ambient conditions we set for trouble-free operation expected.

During operation of the interface at negative temperatures the function of the display and keypad can be disturbed.

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## Approvals

### Wireless:

Please refer the operating range of EBI 25 logger to the relevant area code.

#### Area Code: 999 (also without labeling)

This device may only be used in European Union, Switzerland and Norway, where EN 300 220 is valid.

used frequency: 868.45 MHz

#### Area Code: 052

This device may only be used in Turkey.

used frequency: 868.45 MHz

#### Area Code: 075

This device may only be used in Russian Federation.

77-16/1109/3756

used frequency: 868.45 MHz

#### Area Code: 664

This device may only be used in India.

registration number: NR-ETA-/3671

No.J-22022/182016-RLO(NR)/0233

used frequency: 866.00 MHz

#### Area Code: 701

This device may only be used in Malaysia.

SIRIM / MCMC: E 046578

used frequency: 922.40 MHz

#### Area Code: 706

This device may only be used in Singapore.

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DOMINIQUE DUTSCHER SAS

**Area Code: 720**

This device may only be used in People's republic of China.  
used frequency: 868.45 MHz

**Area Code: 740**

This device may only be used in Hong Kong.  
certification Number: HK0021600072  
used frequency: 821.10 MHz

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## Approvals

The conformity certificate confirms that this product complies to CE guidelines 2014/30/EU.

## Operation

Operation is subject to the following two conditions:

- (1) May only be used within the permitted area. see. Chapter Approvals
- (2) This device may not cause harmful interference, and
- (3) This device must accept any interference received, including interference that may cause undesired operation.

This device has been designed to operate with an internal antenna. Other Antennas are strictly prohibited for use with this device.

Changes or modifications of the device could void the user's authority to operate the equipment. (Exception: Replacing the battery as described in the user manual.)

User manuals can be downloaded from the Internet at [www.ebro.com](http://www.ebro.com)

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Xylem Analytics Germany Sales GmbH & Co. KG
Xylem Analytics Germany GmbH