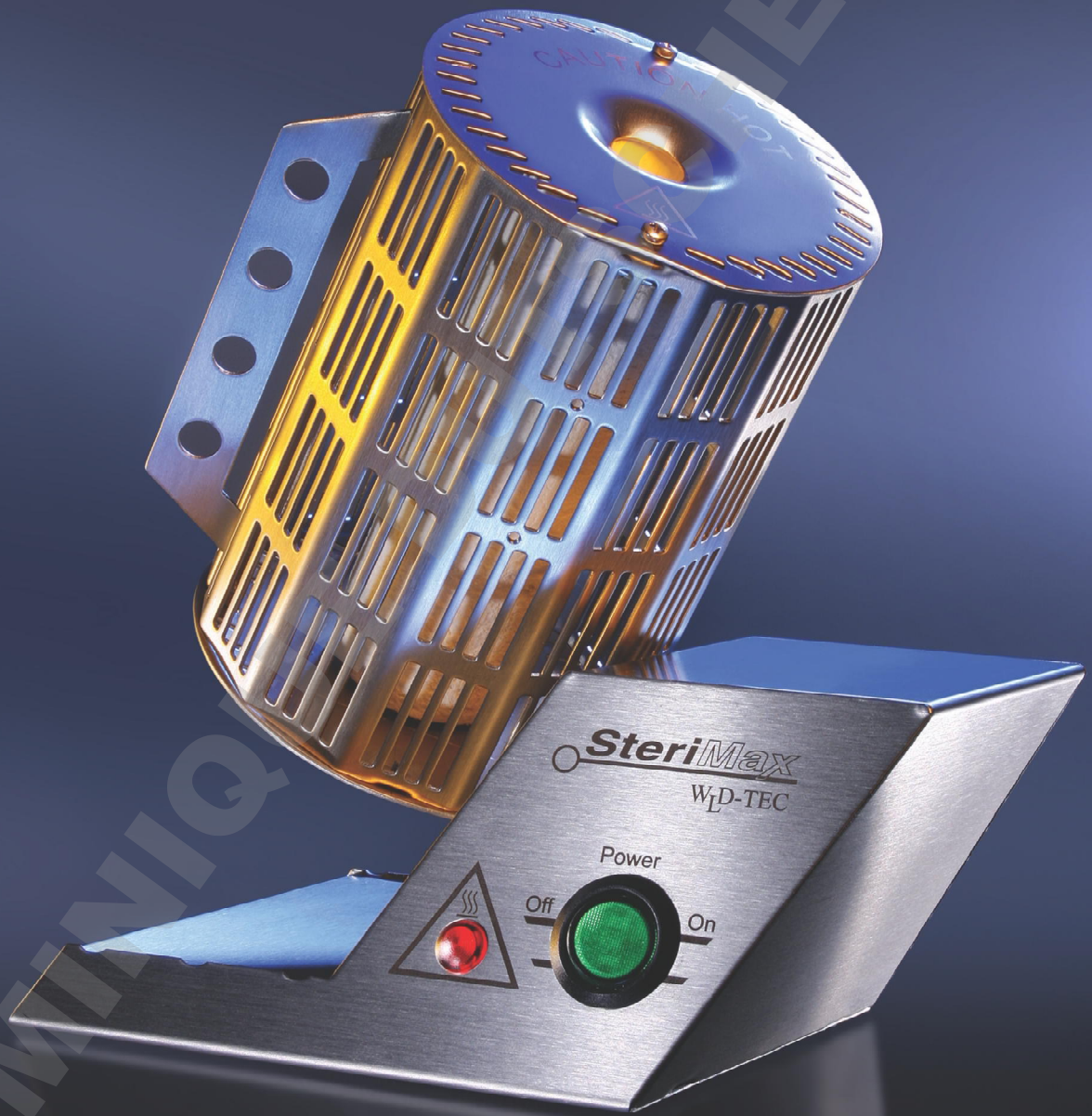
SteriMax

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Annealing without Flame

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Double Tube System
Ceramic / Quartz glass

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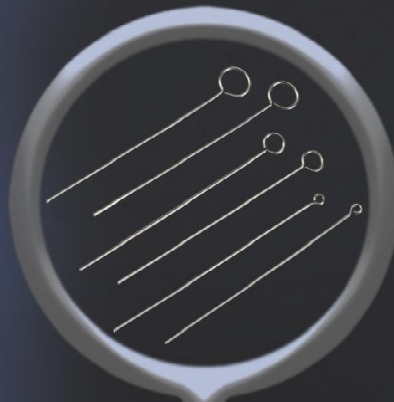
**Holding device
for WLD-TEC
inoculation
loop holders only
suitable for all angles**

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**Universal
support for all
inoculation
loop holders
for angular pos. 1 - 3**

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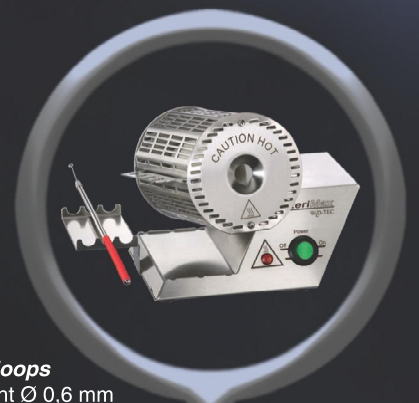
Tray
for left side, made
of stainless steel
Art.-Nr./Art.-No.
6.000.341

Annealing without Flame!

The electrical sterilization system SteriMax is most suitable for the sterilization of inoculation loops, needles and instruments. It is ideal for laboratories and safety cabinets where the use of gases and open flames is not permitted.

Robust

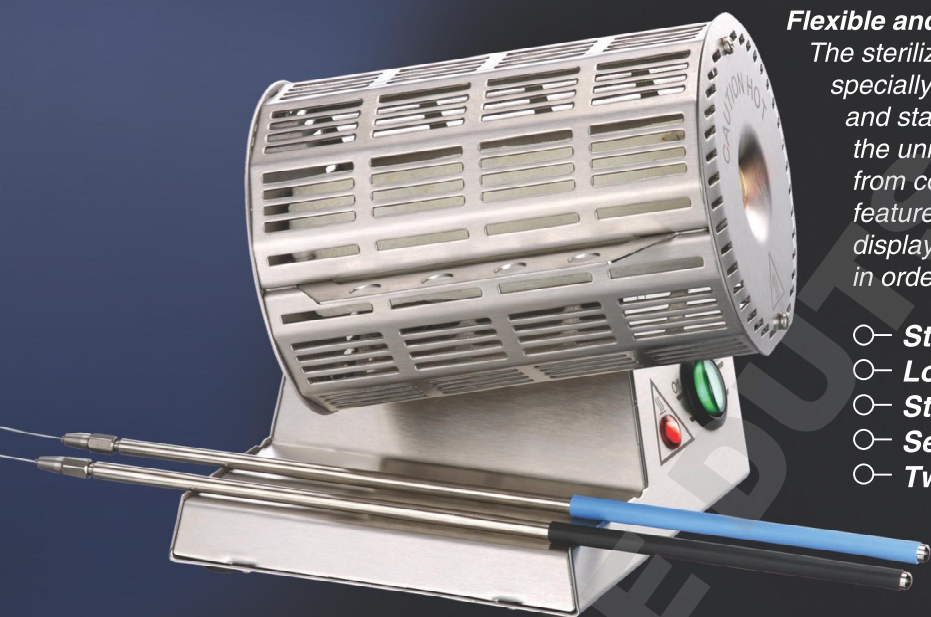
The unique double tube system, which is made of robust quartz glass and wear-resistant ceramic, reaches its optimal temperature of 900°C (1650°F) after only a few minutes. To sterilize the inoculation loop, simply insert it into the ceramic annealing tube. After 5 - 7 seconds, the inoculation loop is sterilized¹.

A back side spray protector is integrated for more safety when handling with pathogenic material. Even clogging is not an issue. The ceramic annealing tube can be removed easily for in-depth cleaning. A 100% stainless steel design makes the SteriMax resistant to even the most extreme laboratory conditions.


Flexible and safe

The sterilization tube can lock itself into place, using the specially designed angular adjusting device. The low and stable housing facilitates ergonomic operation; the unique design protects the working surface from contamination. Exceptional passive safety feature: After prolonged use, the residual heat display signals a hot surface of the sterilization tube in order to protect the user from burns.


- **Stainless steel design**
- **Long - lasting double tube system**
- **Sterilization period 5-7 seconds¹**
- **Self - locking angular adjustment**
- **Two sizes available**

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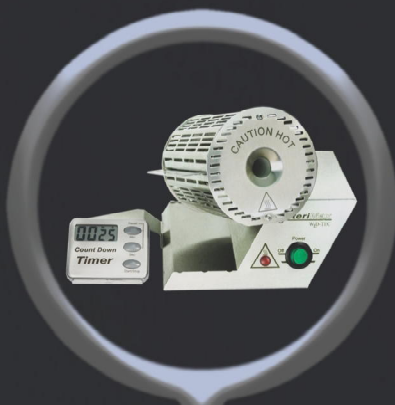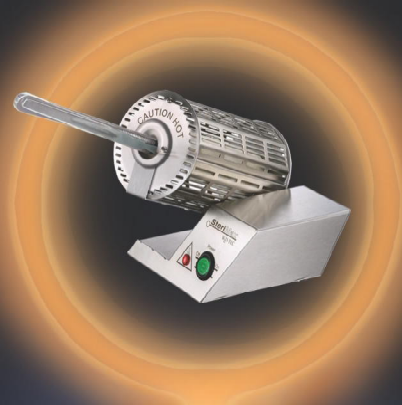**Quartz glass heating element
 wear-persistent**

The range:

with 7 angular adjustments
 Residual heat display
 1 inoculation loop holder and inoculation loop
 Tray for 3 instruments, screwdriver
 2-year warranty

Double Tube System:

Diameter Ø 15 mm, length 118 mm:

for 230 V mains voltage **Art.-Nr. 5.001.000**
 for 120 V mains voltage **Art.-Nr. 5.002.000**
 for 100 V mains voltage **Art.-Nr. 5.003.000**

Double Tube System:

Diameter Ø 20 mm, length 161 mm:

for 230 V mains voltage **Art.-Nr. 5.101.000**
 for 120 V mains voltage **Art.-Nr. 5.102.000**
 for 100 V mains voltage **Art.-Nr. 5.103.000**

Technical Data:

Double Tube System
 Ø 15 mm: length 118 mm
 Ceramic annealing tube Ø 15 mm
 Quartz glass heating element Ø 21 mm
 Ø 20 mm: length 161 mm
 Ceramic annealing tube Ø 20 mm
 Quartz glass heating element Ø 26.6 mm
 7
 Angular adjustments:
 Sterilization temperature: 900 °C - 950 °C
 (1650 °F - 1742 °F)
 Power supply Ø 15 mm: 160 W
 Power supply Ø 20 mm: 225 W
 Heating up time: 10 - 16 minutes
 Sterilization time: 5 - 7 seconds¹
 Heat transfer: Average warming of a stainless
 steel surface, located above in
 motionless air and at an ambient
 temperature of 20 °C (131 °F)
 Distance: 20 cm: 55 °C (131 °F)
 30 cm: 38 °C (100 °F)
 40 cm: 32 °C (90 °F)
 Safety features: residual heat display
 Casing: stainless steel, UV and
 solvent resistant
 Weight Ø 15 mm: 960 g
 Measurements Ø 15 mm: 175 x 129 x 139 mm (w x h x d)
 Weight Ø 20 mm: 1060 g
 Measurements Ø 20 mm: 175 x 129 x 169 mm (w x h x d)
 Licenses: CE: EN 61326-1, EN 61010-1,
 EN 61010-2-010
 EU guidelines: 2014/30/EU,
 2014/35/EU, 2011/65/EU

Sales department:

Spandauer Weg 1
 D-37085 Göttingen
 Telefon: +49 (0)551 / 793789
 Telefax: +49 (0)551 / 793707

Production & Service:

Halle-Kasseler-Straße 49
 D-37318 Arenshausen
 Telefon: +49 (0)36081 / 68940
 Telefax: +49 (0)36081 / 68942