


## VF12 Glass Filtration Set

### Unpacking

Carefully check if the shipping carton is damaged before unpacking. Please check all supplied items according to the standard package list showed as the right photo. If you have any questions, please contact your local dealer and provide with serial number & shipping carton for help.

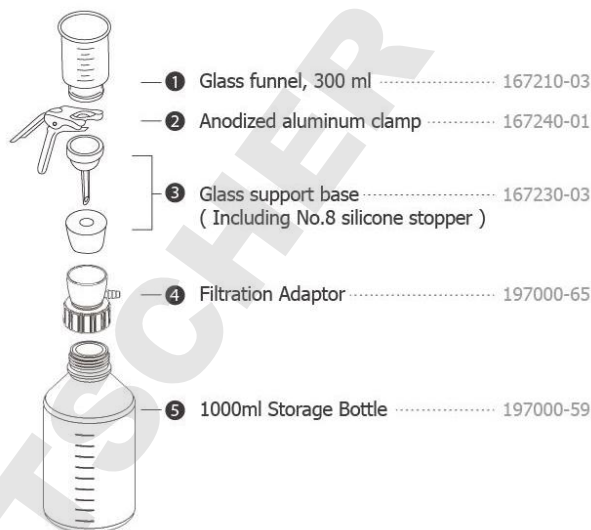

### Important Notice

- The product is suitable for GL45 storage bottle and fit to ID  $\phi$  4~8 mm tube.
- Made of PP material filtration adaptor resistant to high temperature up to 130°C and autoclavable.
- Do not sterilize the product with flame.

### Usage Guide

1. Please ensure that 『 PTFE Gasket 』 has been installed properly in the filtration adaptor.
2. Screw the adaptor onto storage bottle tightly and place No. 8 stopper on the upper receiving port.
3. Connect glass support base to adaptor with No.8 stopper.
4. Put 47mm membrane on the glass support base.
5. Put the glass funnel on the glass support base then clamp them together.
6. Connect a matched tube between its hose barb and vacuum source to start filtration.

### Ordering Information

|            |                            |
|------------|----------------------------|
| 167200-12: | VF12 Glass Filtration Set  |
| 197000-65: | Filtration adaptor         |
| 167210-03: | 300ml Glass Funnel (47mm)  |
| 167240-01: | Anodized Aluminum Clamp    |
| 167230-03: | 47mm Glass Support Base    |
| 197000-59: | 1000ml GL45 storage bottle |