


VF5 Glass Filtration Set

Unpacking

Carefully check if the shipping carton is damaged before unpacking. Please check all supplied items according to the standard package list showed as the right photo. If you have any questions, please contact your local dealer and provide with serial number & shipping carton for help.

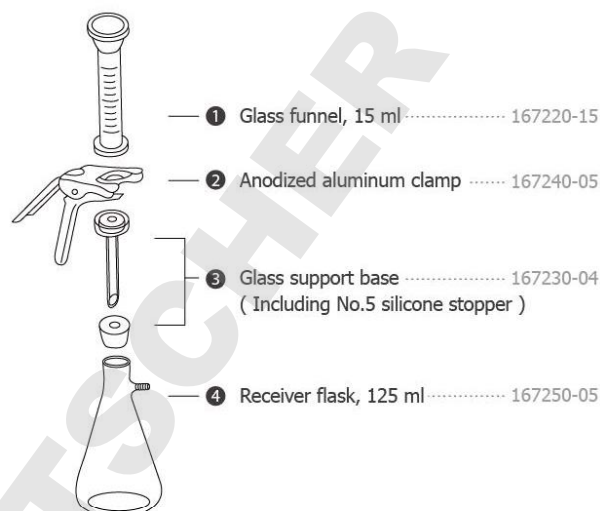

Usage Guide

1. Put the glass support base and silicone stopper together.
2. Connect the glass support base to the flask with silicone stopper.
3. Put a 25mm membrane on the glass support base.
4. Put the glass funnel on the glass support base then clamp them together.

***** Please confirm all steps above are completed before doing filtration test.**

Ordering Information

167200-05:	VF5 Glass filtration set
167220-15:	15ml Glass funnel (25mm)
167240-05:	Anodized aluminum clamp
167230-04:	25mm Glass support base
167250-05:	125ml Receiver flask