


VF7 Glass Filtration Set

Unpacking

Carefully check if the shipping carton is damaged before unpacking. Please check all supplied items according to the standard package list showed as the right photo. If you have any questions, please contact your local dealer and provide with serial number & shipping carton for help.

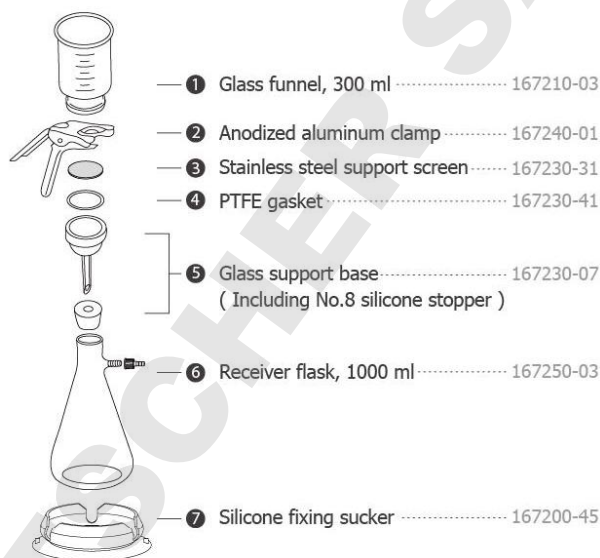

Usage Guide

1. Put the fixing sucker on a flat table.
2. Press down the fixing sucker from the middle to squeeze the air outward, and make sure the "holding base" is connected to the table tightly.
(It is easier to remove the fixing sucker from table by taking the "Handle")
3. Put the glass flask on the "Holding base" and let the "Holding side" cover the glass flask.
4. Put the glass support base and silicone stopper together, and connect to the flask.
5. Screw the screwthread connector onto the hose barb.
6. Put the PTFE gasket and the SS support screen into the support base then put 47mm membrane filter on the SS support screen.
7. Put the glass funnel on the glass support base then clamp them together.

*** Please confirm all steps above are completed before doing filtration test.

Ordering Information

167200-07:	VF7 Glass filtration set
167210-03:	300ml Glass funnel (47mm)
167240-01:	Anodized aluminum clamp
167230-31:	47mm SS support screen
167230-41:	47mm PTFE gasket
167230-07:	47mm Glass support base
167250-03:	1000ml Receiver flask
167200-45:	Silicone fixing sucker (for 1000ml flask)