


VF8 Glass Filtration Set

Unpacking

Carefully check if the shipping carton is damaged before unpacking. Please check all supplied items according to the standard package list showed as the right photo. If you have any questions, please contact your local dealer and provide with serial number & shipping carton for help.

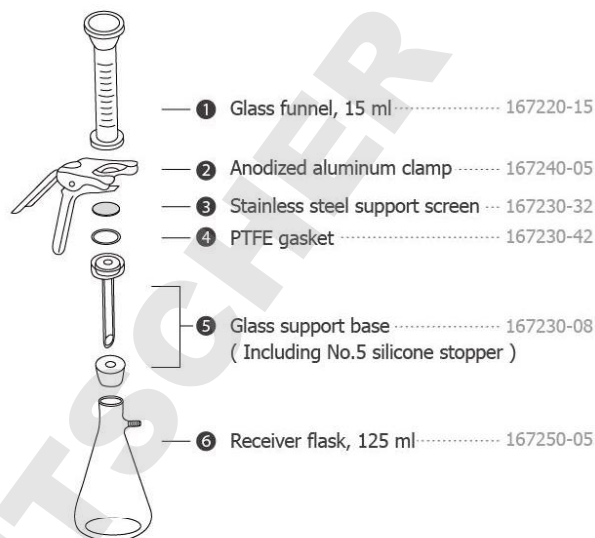

Usage Guide

1. Put the glass support base and silicone stopper together, and connect to the flask.
2. Put the PTFE gasket and the SS support screen into the support base.
3. Put 25mm membrane filter on the SS support screen.
4. Put the glass funnel on the glass support base then clamp them together.

***** Please confirm all steps above are completed before doing filtration test.**

Ordering Information

167200-08:	VF8 Glass filtration set
167220-15:	15ml Glass funnel (25mm)
167240-05:	Anodized aluminum clamp
167230-32:	25mm SS support screen
167230-42:	25mm PTFE gasket
167230-08:	25mm Glass support base
167250-05:	125ml Receiver flask