

eppendorf

General Lab Products EU



Endless Possibilities

Eppendorf Microcentrifuges



Meet the Next Generation in Microcentrifugation

We introduced our very first centrifuge in 1964 as an integral part of the Eppendorf Microliter System, and revolutionized laboratory practices worldwide. To this day, the Eppendorf name remains synonymous with creative design, innovative technology, and reliable performance that stand the test of time.

The Eppendorf approach to product development is about listening to our customer's needs, understanding daily laboratory routines, and anticipating future trends. The result is a comprehensive line of microcentrifuges that serve the multiple applications you encounter in your lab – whether you just have to spin a few tubes at a time or need the extra versatility for handling plates and larger volume tubes directly on your lab bench.

Our goal is to provide you with a product solution that not only leaves you thoroughly satisfied but that is also fun to work with. That is why Eppendorf microcentrifuges are designed to go beyond speed and capacity to offer you unparalleled ergonomic operation and superior temperature management. You will not only experience enhanced performance, but also exceptional joy in their use.

With this sound fundamental experience Eppendorf is proud to serve you as your expert partner in microcentrifugation.

The Benchmark in Ergonomic Operation

To make your daily routines faster and easier we strongly focus on the ergonomic aspects of our products. Today Eppendorf microcentrifuges set the standards in terms of design and user-friendliness. These key success factors have also been awarded the »excellent! Universal Design« certificate of the TÜV Nord organization.



Virtually silent operation

Eppendorf centrifuges are designed for whisper quiet operation to benefit your work environment – you can work right next to them while they are running.



Soft-touch lid closure

All our centrifuges are equipped with a soft-touch one-finger lid closure. Just lower the lid gently and it will lock automatically. No more pressing or slamming down the lid.



Intuitive operation

Large, full-digit displays, and intuitively labeled keys make operation super simple.




Eppendorf QuickLock® rotors

Our aerosol-tight Eppendorf QuickLock rotors close with only ¼ of a turn. This not only saves you time but also takes repetitive stress off your wrist.

Superior Temperature Management

Eppendorf's commitment is to provide you centrifuges that offer the cooling performance you need for successful experiments. Our refrigerated microcentrifuges are best in class when it comes to energy efficiency and sample protection. But that's not all: Non-refrigerated Centrifuge 5424 features a unique cooling concept. Instead of ventilating air directly through the rotor bowl it employs fans to guide air around the outside of the rotor chamber where specially designed cooling fins ensure an optimal heat exchange to minimize sample warming. The additional benefits: The lid of Centrifuge 5424 seals the rotor chamber which leads to a significant noise reduction and prevents particles from being blown into the lab environment.

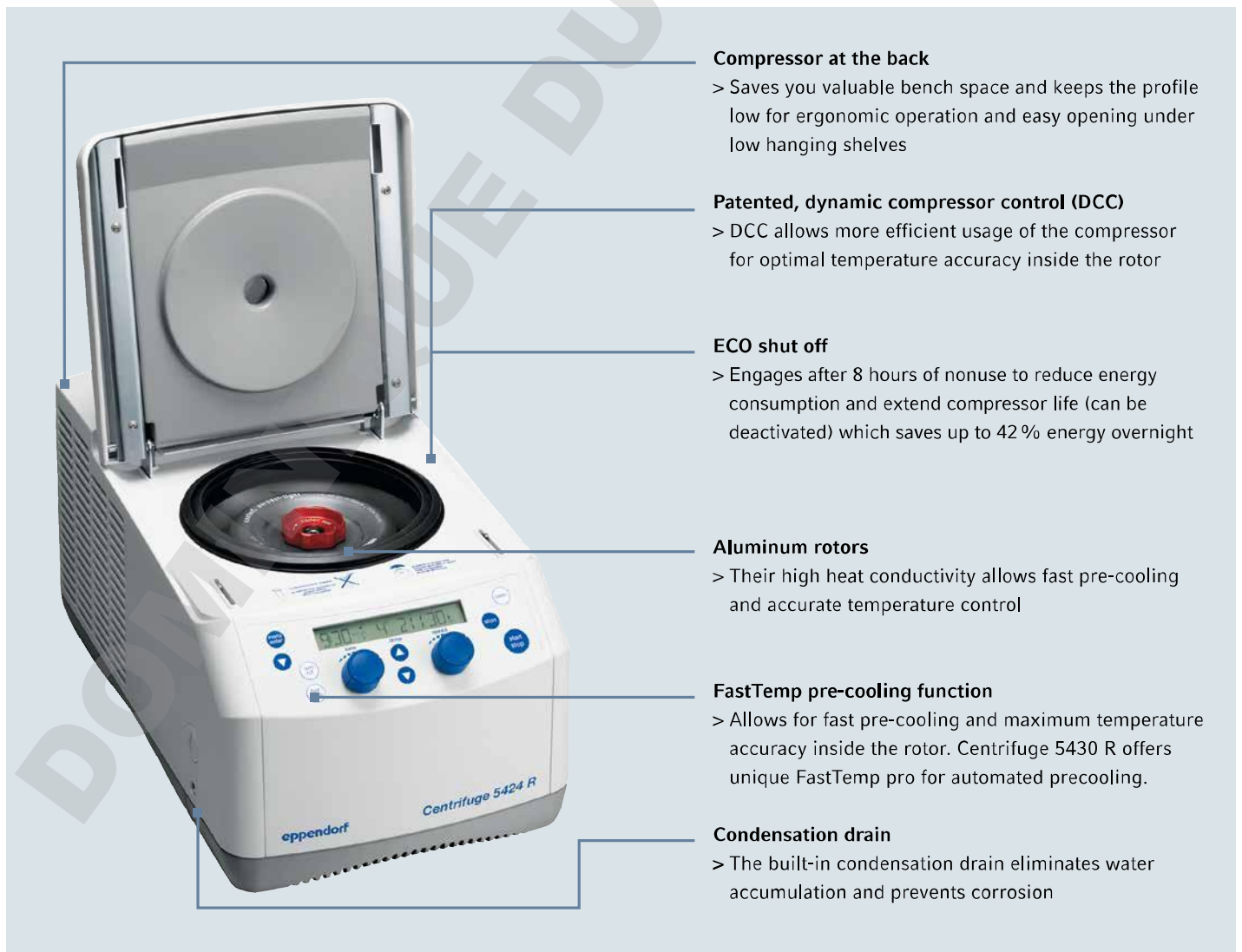


»Our Eppendorf microcentrifuges provide exactly what we expect. Not only speed and cooling performance, but also ease of use and a quality that is second to none! Thus they significantly increase our efficiency and give us more time to focus on.«

Dr. Per Hoffmann
Institute for Human Genetics
Biomedical Center,
Bonn, Germany



Energy Saving Right from the Start

Save energy with every new generation of Eppendorf centrifuges

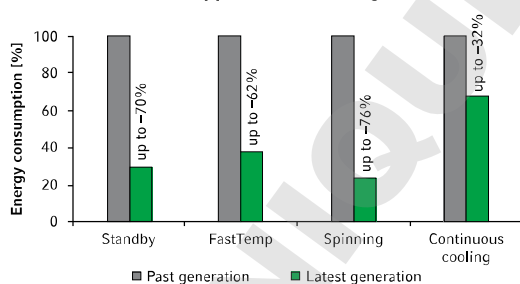


As a commitment to the environment and future generations we constantly look for ways to improve the eco-friendliness of our products. This already started many years ago when we switched all our refrigerated models to CFC-free refrigerants with zero ozone depletion potential.

We continue to further reduce energy consumption of our centrifuges by optimizing motor systems, building lighter weight rotors, and using highly efficient compressor systems. All our centrifuges follow our philosophy by providing several eco-friendly product features – for example:

- > Use of highly efficient, brushless induction motors
- > Unique ECO shut-off deactivates the compressor after 1 to 8 hours of nonuse for up to 88 % energy savings
- > New standby module (for Centrifuge 5427 R and Centrifuge 5430 R) additionally reduces baseline energy consumption to a minimum
- > FastTemp pro® (Centrifuge 5430 R) allows for automated pre-cooling based on preprogrammable time and date which saves up to 90 % of energy (see page 13 for more details)
- > Rotors made of 100 % recyclable aluminum
- > Recyclable packaging material

Energy savings overnight with ECO shut-off and FastTemp pro

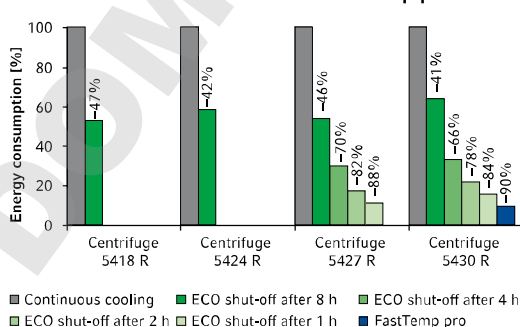



MiniSpin®/MiniSpin® plus

Entry-level centrifuges

Powerful and easy-to-use MiniSpin and MiniSpin plus microcentrifuges are small enough so that each workstation can be equipped with a »personal« centrifuge.

Ordering information

Description	Order no.
MiniSpin® , with 12 × 1.5/2.0 mL rotor	5452 000.010
MiniSpin® plus , with 12 × 1.5/2.0 mL rotor	5453 000.015
Rotor 2 × PCR-strip , with lid (F-55-16-5-PCR)	5452 727.007
Adapter for all 1.5/2.0 mL microcentrifuge rotors	
Adapter for 0.2 mL PCR tubes (set of 6)	5425 715.005
Adapter for 0.4 mL tubes (set of 6)	5425 717.008
Adapter for 0.5 mL tubes and 0.6 mL Microtainer® (set of 6)	5425 716.001

All listed models are 230 V/50–60 Hz.

Product features

- > Capacity: 12 × 1.5/2.0 mL, 2 × PCR strip
- > Speed up to 12,100/14,100 × g (13,400/14,500 rpm)
- > Extremely compact footprint
- > Digital display for reproducible setting of time and speed
- > Rotor, inner centrifuge lid and lid latch are made of metal for maximum operational safety
- > Soft-touch lid closure for ergonomic lid locking
- > Lid opens automatically at the end of the run to prevent sample warming and to allow easy access to samples

Additional features of MiniSpin plus

- > Automatic rpm/rcf conversion
- > Extended time to 99 min or continuous centrifugation



Centrifuge 5418 R

Medium capacity

The Centrifuge 5418 R with its 18-place capacity rotors are specifically designed for low to medium capacity in 1.5/2.0 mL tubes and Microtainer®.

Ordering information

Description	Order no.
Centrifuge 5418 R (refrigerated)	
with 18 × 1.5/2.0 mL aerosol-tight QuickLock rotor	5401 000.013
Adapter for all 1.5/2.0 mL microcentrifuge rotors	
Adapter for 0.2 mL PCR tubes (set of 6)	5425 715.005
Adapter for 0.4 mL tubes (set of 6)	5425 717.008
Adapter for 0.5 mL tubes and 0.6 mL Microtainer® (set of 6)	5425 716.001

All listed models are 230 V/50–60 Hz.

Product features

- > Capacity: 18 × 1.5/2.0 mL
- > Speed up to 16,873 × *g* (14,000 rpm)
- > Compact footprint to fit on crowded lab benches
- > OptiBowl® design allows for whisper quiet operation without rotor lid to benefit your work environment
- > Soft-touch one-finger lid closure for ergonomic operation
- > Aerosol-tight rotor for safe centrifugation of hazardous samples
- > Eppendorf QuickLock® rotor allows fast and ergonomic lid locking
- > Temperature range from 0 °C to 40 °C
- > FastTemp function for fast pre-cooling and maximum temperature accuracy inside the rotor
- > ECO shut-off engages after 8 hours of nonuse to reduce energy consumption (47 % overnight) and extends compressor life (can be deactivated)



Centrifuge 5424 R

The laboratory standard

The Centrifuge 5424 R is the laboratory standard. With its 24-place capacity and speed up to $21,130 \times g$ it is perfectly equipped for all applications in Eppendorf and PCR tubes. When working with Miniprep spin column kits – the Eppendorf Kit rotor® with raised rim solves the issue of open tube caps shearing off during the elution step.

Centrifuge 5424 R features state of the art cooling technology to ensure energy efficiency and high temperature accuracy for maximum sample protection. For ease-of-operation and time savings, Centrifuge 5424 R cools down to 4°C in only 8 minutes.

Product features

- > Capacity: $24 \times 1.5/2.0 \text{ mL}$, $4 \times \text{PCR strips}$
- > Speed up to $21,130 \times g$ (15,000 rpm)
- > OptiBowl design allows for whisper quiet operation without rotor lid
- > Soft-touch one-finger closure for ergonomic operation
- > Temperature range from -10°C to 40°C
- > ECO shut-off engages after 8 hours of nonuse to reduce energy consumption (42% overnight) and extend compressor life (can be deactivated in menu)
- > Built-in condensation drain to eliminate water accumulation



Aerosol-tight rotor for 24 × 1.5/2.0 mL tubes provides speed up to 21,130 × *g*/15,000 rpm. (FA-45-24-11; Order no. 5424 702.007).



Eppendorf Kit rotor for 18 × spin columns and 1.5/2.0 mL tubes for speed up to 18,111 × *g*/15,000 rpm. This rotor is designed with an extended rim that protects tube caps from shearing off (FA-45-18-11-Kit; Order no. 5424 706.002).



Aerosol-tight rotor for 24 × 1.5/2 mL tubes with PTFE-coating for high chemical resistance. Provides speed up to 21,130 × *g*/15,000 rpm (FA-45-24-11-Special; Order no. 5424 700.004).



4 × PCR-strip rotor for centrifugation up to 18,615 × *g*/15,000 rpm (F-45-32-5-PCR; Order no. 5424 704.000).

Ordering information

Description	Order no.	Order no.
	Knob	Keypad
Centrifuge 5424 R (refrigerated)		
with 24 × 1.5/2.0 mL aerosol-tight rotor	5404 000.413	5404 000.014
without rotor	5404 000.618	5404 000.219
Adapter for all 1.5/2.0 mL microcentrifuge rotors	Order no.	
Adapter for 0.2 mL PCR tubes (set of 6)	5425 715.005	
Adapter for 0.4 mL tubes (set of 6)	5425 717.008	
Adapter for 0.5 mL tubes and 0.6 mL Microtainer® (set of 6)	5425 716.001	

All listed models are 230 V/50–60 Hz.



Centrifuge 5427 R

High-throughput refrigerated microcentrifuge

Centrifuge 5427 R was designed with high-throughput applications in mind. With its aerosol-tight 48-place rotor it handles even the toughest challenges in a safe and reliable manner to provide you the efficiency boost you need. All aerosol-tight rotors, including the new rotors for Eppendorf Tubes® 5.0 mL as well as the 24 × 1.5/2.0 mL swing-bucket rotor, are equipped with Eppendorf QuickLock lids for easy rotor lid opening and closing.

Choose the Centrifuge 5427 R if a 24-place microcentrifuge (e.g., Centrifuge 5424 R) doesn't meet your needs regarding tube capacity and if you don't need the extra versatility our cross-over Centrifuge 5430 R can offer you.

Product features

- > Capacity: 48 × 1.5/2.0 mL, 12 × 5.0 mL
- > Speed up to 25,000 × g (16,220 rpm)
- > High-capacity 48-place rotor
- > Aerosol-tight Eppendorf QuickLock rotor lids allow for fast and ergonomic lid locking.
- > Whisper quiet operation due to optimized sound design
- > Soft-touch one-finger closure for ergonomic operation
- > Temperature range from -10°C to 40°C
- > FastTemp program optimized for fast 11 min pre-cooling
- > ECO shut-off can be programmed to engage after 1, 2, 4 or 8 hours of nonuse for energy savings up to 88 % and extended compressor life (see page 5 for more details)



> Watch video clips of Centrifuge 5427 R on our dedicated website for mobile devices or www.eppendorf.com/centrifugation



Rotor for 48 × 1.5/2.0 mL tubes
provides speed up to 18,210 × *g*/12,700 rpm
(FA-45-48-11*; Order no. 5409 710.001).
Also available with basic plastic lid
(F-45-48-11; Order no. 5409 712.004).



Rotor for 30 × 1.5/2.0 mL tubes
provides speed up to 20,800 × *g*/14,000 rpm
(FA-45-30-11*; Order no. 5409 706.004).
Also available with basic plastic lid
(F-45-30-11; Order no. 5409 708.007).



Rotor for 24 × 1.5/2.0 mL tubes
provides speed up to 25,000 × *g*/16,220 rpm
(FA-45-24-11*; Order no. 5409 702.009).



Swing-bucket rotor for 24 × 1.5/2.0 mL tubes
spins up to 16,050 × *g*/12,700 rpm.
Swing-bucket design is ideal for phase
separation (S-24-11-AT*;
Order no. 5409 715.003).



**Eppendorf Kit rotor for 24 × spin columns
and 1.5/2.0 mL tubes.** Designed with an
extra high rim to support open tube lids
during centrifugation. Spins at max.
19,100 × *g*/13,200 rpm (FA-45-24-11-Kit*;
Order no. 5409 704.001).



6 × PCR strip rotor with speed up to
11,710 × *g*/10,500 rpm (F-45-48-5-PCR;
Order no. 5409 714.007).



Rotor for 12 × 5.0 mL Eppendorf Tubes
provides speed up to 20,600 × *g*/14,000 rpm
(FA-45-12-17*; Order no. 5409 700.006).
Adapters for smaller tubes available:
> 1.5/2.0 mL, pk/4 (Order no. 5820 768.002)
> Cryogenic tubes, pk/4
(Order no. 5820 769.009)

Ordering information

Description	Order no.
Centrifuge 5427 R	
with 48 × 1.5/2.0 mL aerosol-tight QuickLock rotor	5409 000.210
with 30 × 1.5/2.0 mL aerosol-tight QuickLock rotor	5409 000.530
with 12 × 5.0 mL aerosol-tight QuickLock rotor	5409 000.630
without rotor	5409 000.012
Adapter for all 1.5/2.0 mL microcentrifuge rotors	
Adapter for 0.2 mL PCR tubes (set of 6)	5425 715.005
Adapter for 0.4 mL tubes (set of 6)	5425 717.008
Adapter for 0.5 mL tubes and 0.6 mL Microtainer® (set of 6)	5425 716.001

All listed models are 230 V/50–60 Hz.

* Aerosol-tight and Eppendorf QuickLock-sealing



Centrifuge 5430/5430 R

Microcentrifuge with multipurpose capabilities

Unique cross-over Centrifuges 5430 and 5430 R combine the best features of a microcentrifuge (small footprint) and multipurpose centrifuge (versatility) in one instrument. These centrifuges spin rotors for Eppendorf Tubes and PCR strips as you would expect from any microcentrifuge. But that's not all. In a compact size, Centrifuges 5430 and 5430 R also accommodate a fixed-angle rotor for 15/50 mL conical tubes, Vacutainer®, 10 to 50 mL OakRidge tubes, cryo vials as well as a swingbucket rotor for MTP and PCR plates. Until now, this has only been possible with large multipurpose centrifuges.

Product features


- > Capacity: 48 × 1.5/2.0 mL, 6 × 50 mL, 2 × MTP
- > Speed up to 30,130 × g (17,500 rpm)
- > Soft-touch closure for ergonomic operation
- > Menu-driven, multi lingual operation menu (English, German, French, Spanish) with large backlit display
- > Saves up to 50 user-defined programs – 5 program keys for easy access to routine programs

Additional features of refrigerated Model 5430 R

- > Temperature range from –11 °C to 40 °C
- > FastTemp program for fast pre-cooling and maximum temperature accuracy inside the rotor
- > ECO shut-off can be programmed to engage after 1, 2, 4 or 8 hours of nonuse for up to 84 % energy savings and extended compressor life

Easy Pre-cooling with FastTemp pro®

In addition to the standard FastTemp pre-cooling program Eppendorf Centrifuge 5430 R features a unique software option called FastTemp pro. This program allows for automated pre-cooling based on pre-programmable time and date. FastTemp pro can be set to a specific date or as a repetitive event during several days every week. Simply turn the centrifuge into standby mode when you leave your lab and let FastTemp pro take care of pre-cooling in the morning. This not only makes operation super easy, it also saves you up to 90 % of energy over night.

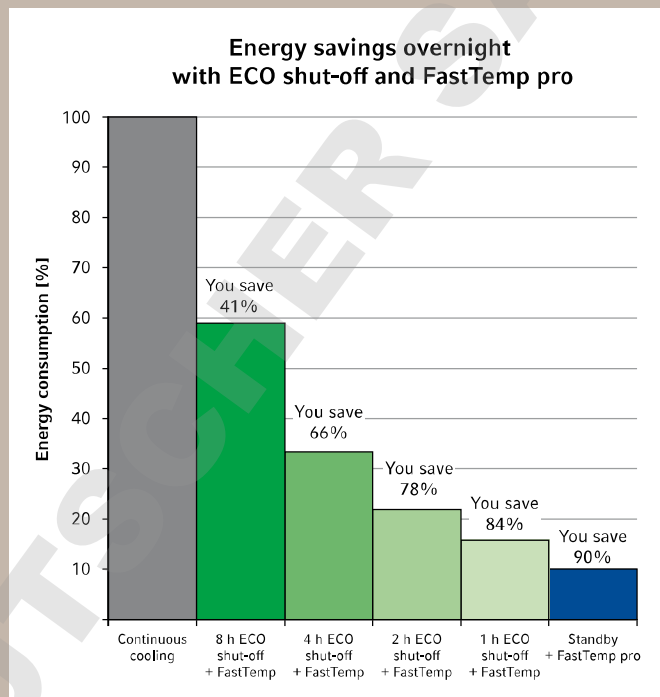

Up to 90 % energy savings with FastTemp pro
Energy savings of ECO shut-off (with 8, 4, 2, and 1 hour setting) and FastTemp pro compared to continuous overnight cooling (17 hours).

Ordering information

Description	Order no.	Order no.
	Knob	Keypad
Centrifuge 5430		
with 30 × 1.5/2.0 mL aerosol-tight QuickLock rotor	5427 000.415	5427 000.210
without rotor	5427 000.610	5427 000.015
Centrifuge 5430 R (refrigerated)		
with 30 × 1.5/2.0 mL aerosol-tight QuickLock rotor	5428 000.410	5428 000.010
without rotor	5428 000.610	5428 000.205
Adapter for all 1.5/2.0 mL microcentrifuge rotors	Order no.	
Adapter for 0.2 mL PCR tubes (set of 6)	5425 715.005	
Adapter for 0.4 mL tubes (set of 6)	5425 717.008	
Adapter for 0.5 mL tubes and 0.6 mL Microtainer® (set of 6)	5425 716.001	

All listed models are 230 V/50–60 Hz.



Versatility for Your Needs

Rotors for Centrifuge 5430/5430 R



Rotor for 48 × 1.5/2.0 mL tubes spins up to 18,210 × g/12,700 rpm (FA-45-48-11*/***; Order no. 54277 540.081). Also available with basic plastic lid (F-45-48-11***; Order no. 54277 550.041).



Rotor for 30 × 1.5/2.0 mL tubes spins up to 20,817 × g/14,000 rpm (FA-45-30-11*/***; Order no. 5427 753.001). Also available with basic plastic lid (F-45-30-11***; Order no. 5427 712.003).



High-speed rotor for 24 × 1.5/2.0 mL tubes spins up to 30,130 × g/17,500 rpm for fast separation results. (FA-45-24-11-HS**/**; Order no. 5427 710.000).



Swing-bucket rotor for 24 × 1.5/2.0 mL tubes spins up to 16,050 × *g*/12,700 rpm. Ideal for phase separation (S-24-11-AT*/****; Order no. 54277 570.071).



Eppendorf Kit rotor for 24 × spin columns and 1.5/2.0 mL tubes. Designed with an extra high rim to support open tube lids during centrifugation. Spins at max. 19,090 × *g*/13,200 rpm (FA-45-24-11-Kit*; Order no. 5427 752.005).



Rotor for 16 × 5.0 mL Eppendorf Tubes for speed up to 21,190 × *g*/14,200 rpm (FA-45-16-17*/****; Order no. 54277 500.021). Adapters for smaller tubes available: > 1.5/2.0 mL, pk/4 (Order no. 5820 768.002) > Cryogenic tubes, pk/4 (Order no. 5820 769.009)



8 × PCR strip rotor with removal adapters that can be used as a pipetting rack; maximum speed is 13,543 × *g*/11,800 rpm (F-45-64-5-PCR; Order no. 5427 714.006, includes adapters).



18-place rotor for Cryovial® and HPLC tubes. Speed up to 8,324 × *g*/8,900 rpm (F-45-18-17-Cryo; Order no. 5427 705.007 incl. cryo adapters).



2-place swing-bucket rotor for MTP, PCR and deepwell plates of up to 29 mm height; speed up to 2,204 × *g*/4,680 rpm. (A-2-MTP; Order no. 5427 700.005).



6 × 15/50 mL conical tube rotor accommodates 12 × Vacutainer, 6 × 15/50 mL conical tubes. Max. speed of 7,815 × *g*/7,830 rpm (F-35-6-30; Order no. 5427 716.009, includes adapters for 15 mL and 50 mL conical tubes).

Adapter options 6 × 15/50 mL conical tube rotor

Adapter	Rotor capacity ¹	Tube Dimensions Ø × L (min-max)	Max. RCF	For bore hole	Order no. (set of 2)
13 mm round	12	13 × 65–89 mm	6,443 × <i>g</i>	Small	5427 740.007
			6,100 × <i>g</i>	Large	5427 742.000
13 mm round	12	13 × 90–110 mm	7,471 × <i>g</i>	Small	5427 741.003
			7,129 × <i>g</i>	Large	5427 743.006
16 mm round	12	16 × 75–105 mm	7,005 × <i>g</i>	Small	5427 732.004
			7,005 × <i>g</i>	Large	5427 734.007
17 mm round	12	17 × 90–125 mm	7,745 × <i>g</i>	Small	5427 735.003
			7,403 × <i>g</i>	Large	5427 738.002
Eppendorf Tubes® 5.0 mL	12	17.5 × 80 mm	6,237 × <i>g</i>	Small	5427 746.005
			7,129 × <i>g</i>	Large	5427 747.001
15 mL conical	6	17.5 × 100–125 mm	7,540 × <i>g</i>	Small	5427 726.004
30 mL round	6	26 × 92–112 mm	7,087 × <i>g</i>	Large	5427 736.000
50 mL round	6	29 × 95–125 mm	7,581 × <i>g</i>	Large	5427 737.006
50 mL conical ^{2,3}	6	29.5 × 100–125 mm	7,197 × <i>g</i>	Large	5427 727.000 ²
Centriprep®	6	29.5 × 100–125 mm	7,567 × <i>g</i>	Large	5427 723.005

* Aerosol-tight and Eppendorf QuickLock-sealing

**Aerosol-tight

***PTFE coated for chemical resistance

****Please contact your local Eppendorf organization or dealer about required software update for your existing device(s).

¹ Please perform a manual test to determine the actual loading capacity.

² Three sets of 2 are included within each 6 × 15/50 mL rotor package.

³ Adapter accommodates skirted 50 mL conical tubes.

A System Striving for Perfection

Eppendorf has always listened to customers and worked together with them on improving existing workflows. Since 1964 Eppendorf is offering system solutions that make everyday lab routines faster and easier. Whether it is spinning down samples, transferring liquids,

storing samples at -80°C , or incubating them at up to 100°C , the perfect combination of Eppendorf instruments, Eppendorf Tubes®, plates and pipette tips will help you to excel your results and give you the best working experience possible.



High-speed centrifuge rotors and high centrifugation resistance of our tubes allow you to get your results even quicker.



Adapters for most common swing-bucket and fixed-angle rotors make it easy to integrate the Eppendorf Tubes into existing centrifuges.



Optimized thermoblocks guarantee accurate temperature control and efficient mixing in Eppendorf ThermoMixer® and MixMate®.



For sample preparation and storage, we offer the ideal racks and boxes for our Eppendorf Tubes.



Broad range of pipettes from $0.1\ \mu\text{L}$ to $10\ \text{mL}$. The complementary set of epT.I.P.S.® allows for stress-free and contamination-free work.



A broad range of adapters make for easy integration into our automated pipetting system epMotion®.



Eppendorf Tubes® 0.5/1.5/2.0 and 5.0 mL

The perfect addition

Eppendorf Tubes combine all the experience from 50 years of constant improvement and development. The original Eppendorf Tube gives you outstanding stability during centrifugation, perfect sealing and evaporation protection.

> The Eppendorf 5.0 mL System at www.eppendorf.com/5ml

Product features

- > Hinged lid of Eppendorf Safe-Lock tubes prevent accidental lid opening during incubation
- > High centrifugation resistance (up to 30,000 × g) prevents tube breakage
- > Precise lid sealing for lowest evaporation rates during storage
- > Manufactured free from slip agents, plasticizers and biocides – substances that have been shown to leach from plastic consumables into the sample and negatively affect assay results
- > Available in Protein and DNA/RNA LoBind quality

Ordering information

Eppendorf Tubes®		0.5 mL, Safe-Lock	1.5 mL, Safe-Lock	2.0 mL, Safe-Lock	5.0 mL ³
Eppendorf Quality	clear	0030 121.023	0030 120.086	0030 120.094	0030 119.401
		500 pcs. (1 × pk/500)	1,000 pcs. (1 × pk/1,000)	1,000 pcs. (1 × pk/1,000)	200 pcs. (2 × pk/100)
PCR clean¹	clear	0030 123.301	0030 123.328	0030 123.344	0030 119.460
		500 pcs. (1 × pk/500)	1,000 pcs. (1 × pk/1,000)	1,000 pcs. (1 × pk/1,000)	200 pcs. (2 × pk/100)
Biopur^{®2}	clear	0030 121.570	0030 121.589	0030 121.597	0030 119.479
		50 pcs. (50 × pk/1)	100 pcs. (100 × pk/1)	100 pcs. (100 × pk/1)	50 pcs. (50 × pk/1)
Protein LoBind (PCR clean)¹	clear	0030 108.094	0030 108.116	0030 108.132	0030 108.302
		100 pcs. (2 × pk/50)	100 pcs. (2 × pk/50)	100 pcs. (2 × pk/50)	100 pcs. (2 × pk/50)
DNA LoBind (PCR clean)¹	clear	0030 108.035	0030 108.051	0030 108.078	0030 108.310
		250 pcs. (5 × pk/50)	250 pcs. (5 × pk/50)	250 pcs. (5 × pk/50)	200 pcs. (4 × pk/50)

¹ PCR clean: Batch-certified free from DNA, DNase, RNase, and PCR inhibitors.

² Biopur[®]: Batch-certified sterile and free from DNA, DNase, RNase, PCR inhibitors, ATP, and pyrogens/endotoxins. Individually packaged.

³ Also available in Sterile (Batch-certified sterile and free from pyrogens/endotoxins): 0030 119.487, 200 pcs. (10 × pk/20)

Eppendorf Microcentrifuges



Model	MiniSpin®/MiniSpin® plus	5418 R
Positioning	Entry-level	Medium capacity
Max. capacity	12 × 1.5/2.0 mL/12 × 1.5/2.0 mL	18 × 1.5/2.0 mL
Max. RCF	12,100 × g/14,100 × g	16,873 × g
Max. speed	13,400 rpm/14,500 rpm	14,000 rpm
Rotors available	2	1
Temperature control range		0 °C to +40 °C
Dimensions (W × D × H)	22.6 × 23.9 × 13.0 cm/22.6 × 23.9 × 12.0 cm	30 × 46 × 25 cm
Height (with open lid)	25 cm	52 cm
Weight w/o accessories	4.3 kg	22 kg
Max. power consumption	70 W/85 W	320 W
Tubes/plates capacity ^{*1}		
1.5/2.0 mL Eppendorf tubes	12	18
Eppendorf Tubes® 5.0 mL	-	-
8-place PCR strips	2	-
15 mL conical	-	-
50 mL conical	-	-
2.6 – 7 mL (13 mm)	-	-
3 – 15 mL (16 mm)	-	-
7 – 17 mL (17.5 mm)	-	-
MTP/PCR-plates	-	-
Deepwell plates (DWP)	-	-

^{*1} F = Fixed-angle rotor; S = Swing-bucket rotor

**5424 R**

The laboratory standard

24 × 1.5/2.0 mL

21,130 × *g*

15,000 rpm

4

-10 °C to +40 °C

29 × 48 × 26 cm

51 cm

21 kg

350 W

24

-

4

-

-

-

-

-

-

-

5427 R

High throughput

48 × 1.5/2.0 mL

25,000 × *g*

16,220 rpm

9

-11 °C to 40 °C

31.9 × 54.0 × 25.4 cm

30 kg

550 W

48

12

6

-

-

-

-

-

-

-

5430/5430 R

Microcentrifuge with multipurpose capabilities

48 × 1.5/2.0 mL, 6 × 50 mL, 2 × MTP

30,130 × *g*

17,500 rpm

12

-11 °C to 40 °C

33 × 42 × 25 cm/38 × 64 × 29 cm

56 cm/63 cm

29 kg/56 kg

475 W/1,050 W

48

16

8

6

6

12

12

12

2

2

DOMINIQUE DUTSCHER SAS

Centrifuge Family Overview

Eppendorf Microcentrifuges



	MiniSpin®	MiniSpin® plus	5418 R	5424 R	5427 R	5430	5430 R
Model	Entry-level		Medium capacity	The laboratory standard	High throughput	Microcentrifuge with multipurpose capabilities	
Max. Capacity	12 × 1.5/2 mL		18 × 1.5/2 mL	24 × 1.5/2 mL	48 × 1.5/2 mL	48 × 1.5/2 mL (6 × 15/50 mL, 2 × MTP)	
Max. speed (RCF)	12,100 × g	14,100 × g	16,873 × g	21,130 × g	25,000 × g	30,130 × g	
Max. rpm	13,400	14,500	14,000	15,000	16,220	17,500	
Rotor options	2		1	4	9	12	
Refrigerated	–	–	0 °C to 40 °C	–10 °C to 40 °C	–10 °C to 40 °C	–	–11 °C to 40 °C
Ordering no.	5452 000.010	5453 000.015	5401 000.010	5404 000.610	5409 000.010	5427 000.610	5428 000.610

Eppendorf Multipurpose Centrifuges



	5702	5702 R	5702 RH	5804	5804 R	5810	5810 R	5910 R	5920 R
Model	Low speed centrifuge for medium capacity needs		Active heating and cooling for high temperature accuracy	Compact high speed centrifuge for medium capacity needs. With swing-bucket rotor and deepwell plate capacity		The workhorse for medium to high-throughput laboratories: Highly versatile, high speed and high plate capacity		High throughput, high performance	Bench-top centrifuge with floor-standing capabilities
Capacity	4 × 100 mL			4 × 250 mL		4 × 750 mL		4 × 750 mL	4 × 1,000 mL
> Conical tubes	20 × 15 mL/4 × 50 mL			40 × 15 mL/16 × 50 mL		56 × 15 mL/28 × 50 mL		64 × 15 mL/ 36 × 50 mL	108 × 15 mL/ 52 × 50 mL
> Blood collection tubes	36 × 13 mm/30 × 16 mm			56 × 13 mm/52 × 16 mm		100 × 13 mm/80 × 16 mm		120 × 13 mm/ 104 × 16 mm	196 × 13 mm/ 148 × 16 mm
> Plates	–			10 × MTP/2 × DWP		16 × MTP/4 × DWP		20 × MTP/ 4 × DWP	28 × MTP/ 8 × DWP
Max. speed (RCF)	3,000 × g			20,913 × g		20,913 × g		22,132 × g	25,062 × g
Max. rpm	4,400			14,000		14,000		14,000	13,700
Rotor options	6			12		18			–
Refrigerated	–	–9 °C to 40 °C	–9 °C to 42 °C	–	–9 °C to 40 °C	–	–9 °C to 40 °C	–11 °C to 40 °C	–11 °C to 40 °C
Ordering no.	5702 000.010	5703 000.010	5704 000.010	5804 000.010	5805 000.010	5810 000.010	5811 000.015	5942 000.010	5948 000.010

Technical specifications subject to change. Errors and omissions excepted.

Your local distributor: www.eppendorf.com/contact
 Eppendorf AG · Barkhausenweg 1 · 22339 Hamburg · Germany
eppendorf@eppendorf.com · www.eppendorf.com

www.eppendorf.com/centrifugation

CryoVita® is a registered trademark of Simport Scientific, Canada. Microtainer® is a registered trademarks of Becton Dickinson and Company, USA. TÜV Nord® logo is a registered trademark of TÜV Nord AG, Germany. Reddot design award logo is a registered trademark of red dot GmbH & Co. KG, Germany. Eppendorf®, the Eppendorf Brand Design, Eppendorf Kit rotor®, OptiBowl®, Eppendorf QuickLock®, Eppendorf Safe-Lock®, Eppendorf Tubes®, Biopur®, epMotion®, epT.I.P.S.®, FastTemp pro®, MiniSpin®, MixMate® and Eppendorf ThermoMixer® are registered trademarks of Eppendorf AG. Eppendorf Quality™ is a trademark of Eppendorf AG. U.S. Design Patents are listed on www.eppendorf.com/ip. Order no.: A54X X13 020/EN4/3T/1217/NW/STEF · Carbon neutrally printed in Germany. All rights reserved, including graphics and images. Copyright © 2017 by Eppendorf AG.