

# I N S T R U C T I O N S

## Bel-Art® – SP Scienceware® Magnetic Bead Separation Racks

### WARNING!

These separation racks contain powerful Neodymium magnets. Misuse can result in injury.

- Do not use this product if you have a pacemaker or any other medical magnetized implant
- Keep away from tools and objects that could be damaged by the magnetic field
- Racks should be handled and stored away from devices subject to magnetic damage including credit cards and other products containing magnetic recording devices
- Keep away from delicate instruments, watches, electronic equipment, displays and monitors
- Magnets may attract steel or other magnetic material with high mechanical forces
- Avoid contact between two magnets
- Do not pull the magnets apart if contact has been made; twist off to prevent damage to the unit or fingers

### Cleaning and Disinfection:

Should the surfaces of your SP Scienceware acrylic products need cleaning, **DO NOT USE ANY SOLVENTS** (alcohol, acetone, etc..) or abrasives. Use of these materials **will cause cracking and/or crazing of the acrylic**. For general cleaning, wash the surfaces with a mild detergent and soft cloth, then rinse thoroughly with water.

**IMPORTANT — Do not autoclave**

Magnetic Bead Separation Rack  
for 50ml Conical Tubes  
catalog number #F19900-0000



Magnetic Bead Separation Rack  
for 5-15ml Conical Tubes  
catalog number #F19900-0001



Magnetic Bead Separation Rack  
for 1.5-2ml Centrifuge Tubes  
catalog number #F19900-0002



Magnetic Bead Separation Rack  
for 96-well Conical-Bottomed Plates  
catalog number #F19900-0003



Magnetic Bead Separation Rack  
for 96-well Flat-Bottomed Plates  
catalog number #F19900-0004



### You May Also be Interested in:

#### Bel-Art® – SP Scienceware® Micro Tube Homogenizer System

**Micro Tube Homogenizer** – Includes Cordless Motor,  
2 AA Batteries and Pestle Adapter


F65000-0000 Cordless Homogenizer Unit, –  
includes 2 AA batteries

#### Individual Pack

Catalog No.	Description	Case Qty.
F65000-0001	0.5 mL Pestle	100
F65000-0003	0.5 mL Tube	100
F65000-0005	0.5 mL Pestle and Tube	100
F65000-0002	1.5 mL Pestle	100
F65000-0004	1.5 mL Tube	100
F65000-0006	1.5 mL Pestle and Tube	100

#### Bulk Pack

Catalog No.	Description	Case Qty.
F65000-9001	0.5 mL Pestle	100
F65000-9003	0.5 mL Tube	100
F65000-9005	0.5 mL Pestle and Tube	100
F65000-9002	1.5 mL Pestle	100
F65000-9004	1.5 mL Tube	100
F65000-9006	1.5 mL Pestle and Tube	100

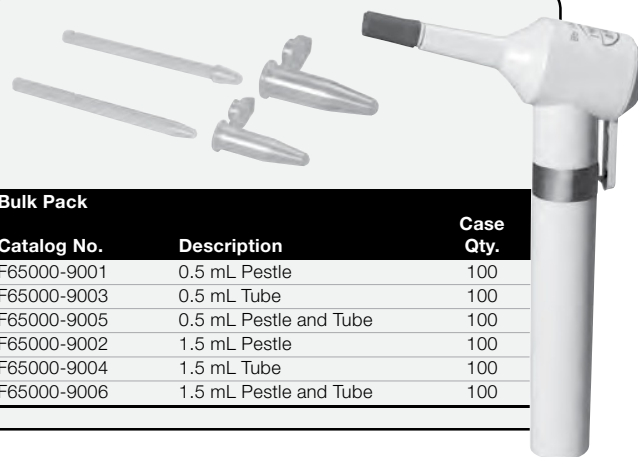

661 Route 23 South, Wayne, NJ 07470

TEL: 1-800-4BEL-ART • FAX: 973-694-7199 • www.belart.com

Bel-Art – SP Scienceware assumes no obligation or liability for any advice furnished by it, or for results obtained with respect to these products. All such advice is given and accepted at the buyer's risk.

© 2015. All Rights Reserved.

Bel-Art® – SP Scienceware® is a division of SP Industries, Inc.

Bel-Art • H-B Instrument  
**SP SCIENCEWARE™**

919900023 – 02/15

