



## **Bel-Art Products**

Wayne, New Jersey 07470

Tel: 973-694-0500 \* Fax: 973-694-7199

Internet Address: <http://www.bel-art.com>

# **Auto Secador Check List**

### **Routine Maintenance**

1. Observe if the drying module goes through its self test procedure when it is first plugged in. This should consist of 3 different stages.
2. Open door and observe if the red / green light is on inside the drying module.
3. Open door and observe if the silica gel beads are blue in color
4. Inspect door gaskets; the door gasket consists of sections of gasket material which may shift over time with the opening and closing of the door causing a gap to appear in between the sections thereby causing leaking. **Daily**
5. Door gasket may have taken a memory, this can be resolved by removing the gasket and turning over, it is important that the gasket sections meet when replacing the gasket in the door channel. **Every 3-4 Months**
6. If 1 thru 6 is okay, Suggest replacing the hygrometer.
7. If the beads do not begin to turn blue after 4 hours please call customer service for a replacement drying module or pre measured container of replacement silica gel beads. You will need the serial number in order for us to determine if you need a complete drying module or just the silica gel beads.

**PLEASE NOTE THAT THE 25% RH MENTIONED IN THE MANUAL IS BASED ON THE CABINET BEING IN AN ENVIROMENT OF 70%, ENVIRONMENTS ABOVE 70% RH MAY INCREASE THE RH LEVEL INSIDE THE CABINET. ALSO THE 25% IS BASED ON THE ENVIRONMENTS BEING APPROX. 68°F, LOWER TEMPERATURES MAY AFFECT THE ABILITY OF THE HEATING UNIT TO REGERERATE THE DESICCANT SUFFICIENTLY.**

### **Replacement Parts List**

420741001 - Shelf Vertical

420740001 - Shelf Horizontal

420701400 - Hygrometer

942070050 - Electronic Drying Module (115V & 220V)

942070053 - 115V Transformer/adapter

942070074 - 220V Transformer/adapter

420740020 – **Pre measured desiccant**

