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Counter AC-10

40345010

Counter AC-15

40345015

1 Inhalt

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40345015

11 Table of content

1.1	Important instructions for your safety	15
1.2	Unpacking the instrument.....	15
1.3	General safety warnings and instructions.....	15
2.1	Description.....	18
3	Setup of the instrument	19
3.1	Setup.....	19
3.2	Select program	19
3.3	Set countstop/stepwidth.....	20
3.4	Choose exclusive keys	20
3.5	„Count“ Mode	21
3.6	„Evaluation“ Mode	22
3.7	Resetting the device	22
3.8	Saving the counts and program parameters.....	22
3.9	PC communication.....	22
4	Cleaning and maintenance	23
5	Dismantling and disposal.....	23
5.1	Dismantling.....	23
5.2	Disposal.....	23
6	Transport and storage	23
7	Warranty.....	24
8	Technical Data	24

1 General information

The Counter is designed in accordance with Safety Class 1, and built and tested in accordance with DIN EN 60335-1; VDE 0700-1:2012-10:2012-10, safety regulations for electrical instruments. According to these regulations, the unit is designed to meet the requirements for safe and correct operations. To maintain the proper safety and operational functions of the instrument, the user should follow the instructions and safety guidelines in this manual.

1.1 Important instructions for your safety

- Every user must read and understand this manual completely before use. Failure to do so can result in serious injury or death.
- Comply with all safety and accident-prevention regulations applicable to laboratory work.
- Follow general instructions for hazard prevention and general safety instructions, e.g. wear protection clothing, eye protection and gloves.
- This operating manual is part of the product. Thus, it must always be easily accessible.
- This instruction sheet does not purport to address all of the safety problems which might result from the use of this device, chemicals, reagents, apparatus or equipment employed in any specific test or protocols. It is the responsibility of the user to consult their authorized safety advisors and establish appropriate health and safety practices and then determine the application of regulatory limitations prior to use.
- Enclose this operating manual when transferring the device to another place.
- If this manual is lost, please request another one. Please contact your dealer.

1.2 Unpacking the instrument

Unpack the outer carton containing your instrument. If there are any punctures, side walls crushed, extensive chemical stains, water marks or other physical evidence that the contents may have been damaged, notify the carrier of the potential problem and ask for specific instructions. If there is no visible damage to the carton's exterior, open it in the normal manner. When all parts are laid out, you should have:

- 1x Counter AC-10 40345010 or AC-15 40345015
- 1x external USB power adapter
- 1x instruction manual

Read the instruction manual carefully prior to operation of the instrument. Please mark points which require special attention in your field of application so they are not overlooked.

1.3 General safety warnings and instructions

Risk of explosion.

- Do not operate the device in the vicinity of highly flammable or explosive substances. The instrument is not explosion-proof.
- Do not use this device for processing any substances which could generate an explosive atmosphere.
- Do not use this device to process any explosive or highly reactive substances.
- Do not use flammable or explosive substances near the instrument.

	<p>Electric shock as a result of penetration of liquid.</p> <ul style="list-style-type: none"> ● Do not allow any liquids to penetrate the inside of the housing. ● Switch off the device and disconnect the power plug before starting cleaning or disinfection work. The On/Off Switch on the device does not disconnect the device from the power source. ● Use only closed tubes. ● Only plug the device back in if it is completely dry, both inside and outside.
---	---

Should you have any additional questions, after reading these instructions, concerning the setting up, operation or warranty terms, please contact either your distributor, or the manufacturer at the following address:

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D-97647 Sondheim/Rhön

+49 (0)9779-808-0

	<p>Electric shock due to damage to device or mains cable</p> <ul style="list-style-type: none"> ● Only connect the device to the mains supply if the device and the mains cable are undamaged. ● Only use devices that have been properly installed or repaired. ● In case of danger, disconnect the device from the mains supply by pulling the power plug from the mains socket or by using the isolating device intended for this purpose (e.g. emergency stop switch).
	<p>Lethal voltage inside the device</p> <ul style="list-style-type: none"> ● Do not open the device. ● Ensure that the housing is always closed and undamaged so that no parts inside the housing can be contacted by accident. ● The On/Off Switch on the device does not disconnect the device from the power source. Remove the plug from the AC power outlet to disconnect the instrument from the mains supply entirely. ● Do not allow any liquids to penetrate the inside of the housing. ● Repairs are only to be carried out by trained service technicians.
	<p>Damages to health due to infectious liquids and pathogenic germs</p> <ul style="list-style-type: none"> ● When handling infectious liquids and pathogenic germs, observe the national regulations, the biological security level of your laboratory, the material safety data sheets and the manufacturer's application notes.

	<ul style="list-style-type: none"> • Wear personal protective equipment. • For comprehensive regulations about handling germs or biological material of the risk group II or higher, please refer to the "Laboratory Biosafety Manual" in its respectively current valid version from the World Health Organisation.
--	--

To ensure maximum service life observe the specified ambient conditions (temperature and humidity).

Improper handling and use could lead to malfunction of the instrument or even cause hazards to the user.

	Do not use flammable or explosive substances near the instrument.
	Please comply with all safety and accident-prevention regulations applicable to laboratory work.
	<p>It is the responsibility of the user to consult and establish appropriate safety and health practices, and then determine the applicability of regulatory limitations prior to use.</p> <p>Should there be any additional questions, after reading these instructions, concerning the setup, operation or warranty, please contact either your distributor, or the manufacturer.</p>
	Use the instrument only in compliance of the intended purpose and in way that neither user nor any other persons are endangered.

2 Purpose of the Counter


Electronic count and storage instrument for statistical evaluation.

2.1 Description

The Counter AC-10 is a counting device with 10 counting keys (red), 5 program keys (green), a screen (blue), a predefined program and a freely configurable program.

Counter AC-15 is a counting device with 15 counting keys (red), 5 program keys (green), a screen (blue), 3 predefined programs and one freely configurable program.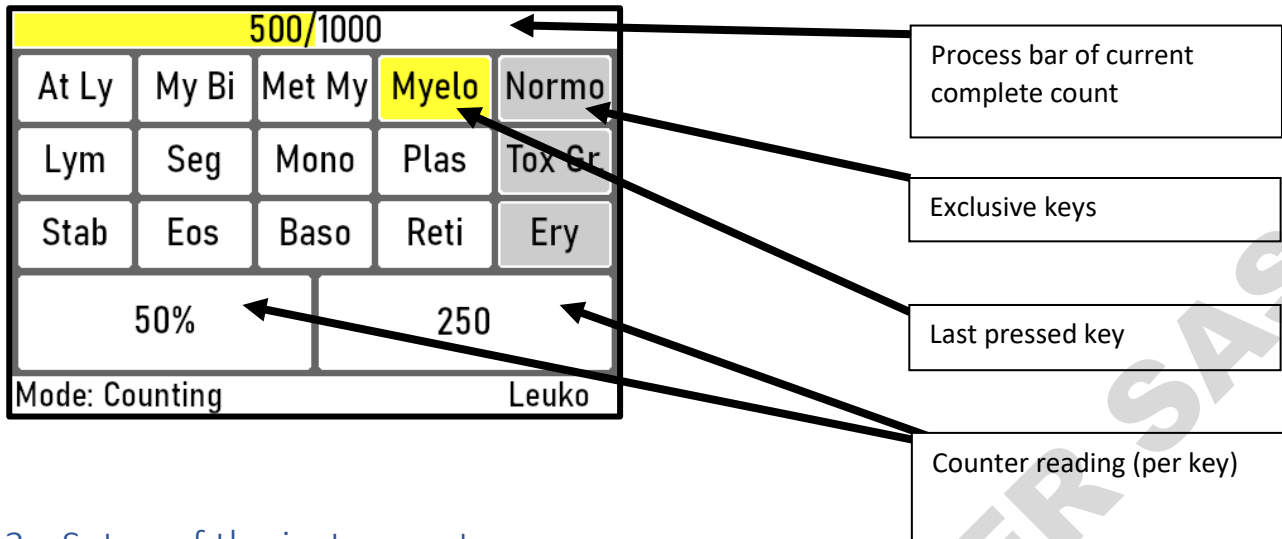

3 Setup of the instrument

After connecting to a supply voltage, the device is in standby. To start the unit, the "Standby" button must be held down for half a second. The start of the device is confirmed with a signal tone.

After the start the last selected program parameters are loaded. It is possible to skip the "Setup" mode by pressing the "Enter/Go On" key twice (three times in the "Free" program).

The device turns off via the "Standby" button.

3.1 Setup

The device is in "Setup" mode after starting. The user selects the program parameters in this mode. The parameters are set in the following steps:

1. select program
2. set end value/step size
3. select/set exclusive keys

The countstop is the value of the end counter at which the program automatically switches to the "Evaluation" mode. The end counter is incremented by the same value as the pressed key when a non-exclusive key is pressed.

In the "Reti" program, the "Div/Ery" key is incremented by the step size when pressed.

An exclusive key does not increment the end counter when pressed.

3.2 Select program


The program is selected using the "Right/+" key or "Left/-/Eval." key. The following programs can be selected.

- Leuko
 - The countstop can be set between 10 and 1000, in steps of 10 and 100.
 - The 5th column of the counting buttons is evaluated exclusively.
- Myelo
 - The countstop can be set between 10 and 1000, in steps of 10 and 100.
- Reti
 - All count buttons except the "Div/Ery" button are evaluated exclusively.

- The step size can have the following values: 1, 5, 10, 20, 30, 40, 50
- The countstop is 1000.
- Free
 - The countstop can be set between 10 and 1000, in steps of 10 and 100.
 - Exclusive keys are selected by the user.

Once the desired program is selected, it must be confirmed with the "Enter/Go On" key. The display then changes to the step size or final value setting.

Alternatively, in this phase of the "Setup" mode, the last count performed can be corrected. To do this, press any count button. If the end counter of the last created count data, is not equal to the countstop, the count will be recalled automatically so that counting can continue immediately after the unit is started.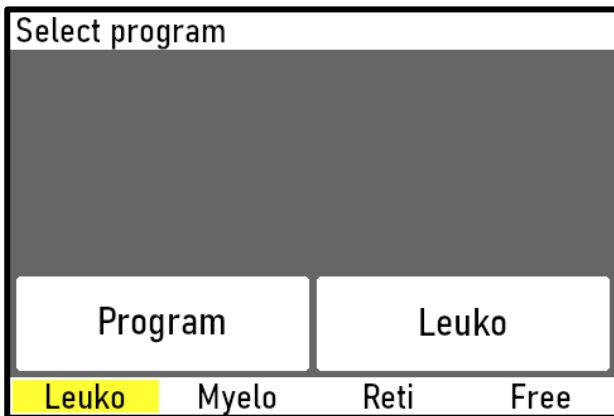


3.3 Set countstop/stepwidth

If the "Reti" program is selected, the step size is set in this step. This is done using the "Right/+" or "Left/-/Eval." keys. The step size is set after the sequence 1, 5, 10, 20, 30, 40, 50. The input of the "Left/-/Eval." key is ignored if the step size is 1. Pressing the "Right/+" key has no effect on the step size if it is 50.

Similarly, if the "Reti" program is not selected, the final value is set with the same keys. By keeping the key pressed, the final value is incremented/decremented in steps of 10. By simply pressing the keys, the final value is changed in steps of 100.

The setting of the final value or the step size is confirmed by pressing the "Enter/Go On" key.

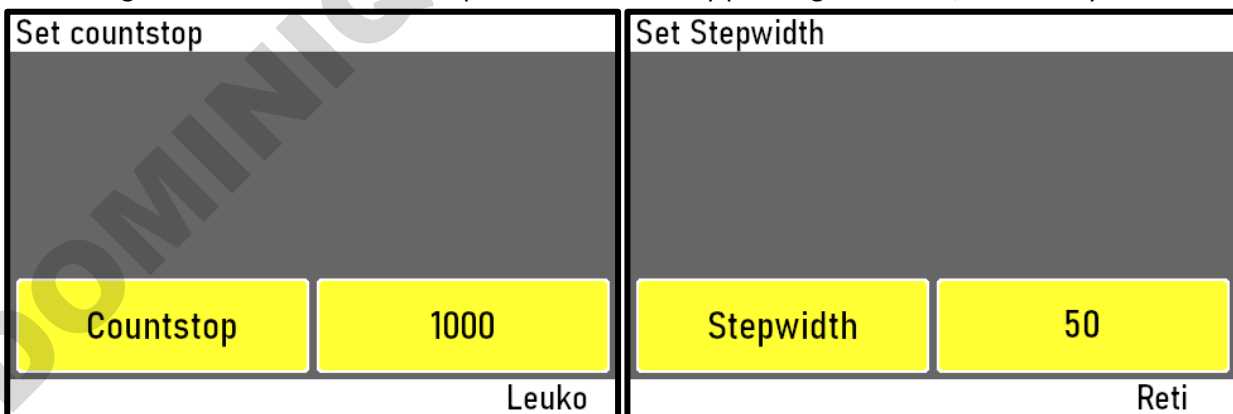3.4 Choose exclusive keys

If the "Free" program is selected, the exclusive keys are selected after setting the final value. This is done by pressing any count keys. Pressing a key that has not been exclusively evaluated marks it as an exclusive key.

If the pressed counting key has already been exclusively evaluated, it is now set as a non-exclusively evaluated key.

This setting must be confirmed using the "Enter/Go On" key. The "Count" mode then begins.

If the "Free" program is not selected, this step is skipped and the respective exclusive keys are set automatically.

Select single counts				
At Ly	My Bi	Met My	Myelo	Normo
Lym	Seg	Mono	Plas	Tox Gr.
Stab	Eos	Baso	Reti	Ery
---		---		
Free				

3.5 „Count“ Mode

After exiting the "Setup" mode, the on-screen keyboard is permanently displayed. The on-screen keyboard displays the exclusive keys and the last key pressed. Below this is the display of the current counter reading of the last key pressed. The absolute count is displayed on the right, while the count in relation to the final counter is displayed on the left. No relative count is displayed for exclusive keys.

Simply pressing the "Corr./Clear" key corrects the last count.

During the entire "Count" mode, the "Evaluation" mode can be called pressing the "Left/-/Eval." key. If the value of the final counter equals the value of the final value, the "Evaluation" mode is automatically called up.

500/1000				
At Ly	My Bi	Met My	Myelo	Normo
Lym	Seg	Mono	Plas	Tox Gr.
Stab	Eos	Baso	Reti	Ery
50%		250		
Mode: Counting		Myelo		

500/1000				
At Ly	My Bi	Met My	Myelo	Normo
Lym	Seg	Mono	Plas	Tox Gr.
Stab	Eos	Baso	Reti	Ery
50%		250		
Mode: Counting		Leuko		

500/1000				
At Ly	My Bi	Met My	Myelo	Normo
Lym	Seg	Mono	Plas	Tox Gr.
Stab	Eos	Baso	Reti	Ery
50%		250		
Mode: Counting		Free		

500/1000				
At Ly	My Bi	Met My	Myelo	Normo
Lym	Seg	Mono	Plas	Tox Gr.
Stab	Eos	Baso	Reti	Ery
---		250		
Mode: Counting		Reti		

3.6 „Evaluation“ Mode

In this mode, the counts are highlighted so that they can be easily read. The counter reading of the last count button pressed is always displayed. Pressing a count button does not cause the individual counters to increment in "Evaluation" mode.

Pressing the "Enter/Go On" key in "Evaluation" mode calls the following functions:

- The final counter is less than the final value:
Counting is continued. This switches to the "Count" mode.
- The final counter is equal to the final value:
The end value is incremented by 100 and the program switches to the "Count" mode.
- The end value is 1000:
Entering this key has no effect on AC-15. The "Evaluation" mode can only be left by holding down the Clear key.

500/1000				
At Ly	My Bi	Met My	Myelo	Normo
Lym	Seg	Mono	Plas	Tox Gr.
Stab	Eos	Baso	Reti	Ery
50%		250		
Mode: Data view			Leuko	

3.7 Resetting the device

The device is reset to the "Setup" mode by holding down the "Corr./Clear" key.

3.8 Saving the counts and program parameters

If the device is switched off correctly (holding the "Standby" button), the count data and program parameters are saved.

If an unfinished count has been saved, the unit will start in count mode immediately after switching on, using the count data of the unfinished count. Counting can continue immediately. If the stored count has been completed, it can be recalled in Level 1 by pressing any count button.

3.9 PC communication

The software is available on our website (www.assistent.eu/en) under „Downloads“

The device can transfer created counting data to the AC-15 PC Manager. For this purpose, the device must be connected to a computer using a suitable USB cable.

Reading out counting data:

- Start the software.
- Start the counting process.
- When the counting process is complete, press the "Read Out Counter Data" button in the software.
- Counter data is displayed in the software.

Saving Counter Data:

- To save the count data, press the "Save Data" button in the software.
- Fill out the form with patient data.

- Press the "Next" button to save the patient and counter data.

Load the patient information:

- Press the "Search Data" button in the software.
- Enter the data you are looking for in the search window.
- Confirm your selection.

Print the patient information:

- Click on "File" in the menu then on "Print".

Save the patient database:

- Click on "File" in the menu, then "Open Database Directory".
- Back up the database with a program of your choice.

4 Cleaning and maintenance

Surface and operating elements may be cleaned with a mild dishwashing detergent (water and a standard dishwashing detergent) and a soft, non-fuzzing moist cloth. Do not use a wet cloth. Use only a small amount of dishwashing detergent. Do not use chlorine bleach or other chlorine-based cleaning products with metallic components under any circumstances. These will damage the surface of the instrument. If you use any other cleaning method, please make sure that the intended method does not cause any damage to the instrument.

The On/Off Switch on the device does not disconnect the device from the power source. Remove the plug from the AC power outlet to disconnect the instrument from the mains supply entirely.

5 Dismantling and disposal

5.1 Dismantling

1. Switch the instrument off.
2. Disconnect the instrument from the mains.
3. Now the instrument may be removed from the working area.

5.2 Disposal

Please dispose of used instruments and defective components at your local recycling collection point. Prior to disposal, sort according to materials: metal, glass, plastic, etc. Also, be sure to dispose of the packing material in an environmental-friendly manner.

6 Transport and storage

Prior to transport:

Switch the instrument off and unplug the power supply.

Place the instrument and its parts in its original packaging or another suitable container to protect it during transport. Close the packaging with adhesive tape.

Store the instrument in a dry environment. Please observe the specified conditions of the ambient (temperature and humidity).

Do not subject the instrument to mechanical shocks or vibration during transporting it.

In case you do not use the original packaging, please mark the box with the following notes:

- Glass symbol (handle with care, fragile)
- Umbrella (keep dry)
- Content (list of content)

7 Warranty

The manufacturer agrees to correct for the original user of this product, either by repair, or at the manufacturer's discretion, by replacement, any defects in material or workmanship, which develop within 24 months after delivery of this product to the original user. In the event of replacement, the replacement unit will be warranted for the remainder of the original twenty-four (24) months period of ninety (90) days, whichever is longer.

If this product should require service, contact your distributor or manufacturer for necessary instructions.

This warranty shall not apply if the defect or malfunction was caused by accident, neglect, unreasonable use, improper service, or causes not arising out of defects in material or workmanship.

There are no warranties, expressed or implied, including, but not limited to, those of merchantability or fitness for a particular purpose, which extended beyond the description and period set forth herein.

The manufacturer's sole obligation under this warranty is limited to the repair or replacement of a defective product and the manufacturer shall not, in any event, be liable for any incidental or consequential damages of any kind resulting from use or possession of the product.

8 Technical Data

Dimensions	<ul style="list-style-type: none"> • Device: 235 x 130 x 35 mm • Packed: 300 x 230 x 90 mm
Weight	0.9 kg packed
Interfaces	USB-Mini
Network operation	USB power adapter 5VDC / 300 mA
Battery operation	External USB powerbank (not included)

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