PCRmax Alpha Cycler 1, 2 and 4

Safety Information

PCRmax rev 2.9

PCRmax

PLEASE NOTE: Before use in 120V environment

IMPORTANT NOTE:

IF USING THE AC-2 / AC-4 AT 120V, BEFORE THE UNIT IS CONNECTED TO THE MAINS AND SWITCHED ON FOR THE FIRST TIME REPLACE THE FUSES.

FOR AC4 REPLACE 8A FUSES WITH 15A ANTI-SURGE FUSES.
FOR AC2 REPLACE 5A FUSES WITH 12.5A ANTI-SURGE FUSES.

THE FUSE HOLDERS ARE LOCATED BESIDE THE POWER SWITCH OF THE UNIT.

IF THE AC-4 IS INTENDED TO BE USED IN AUSTRALIA THE UNIT MUST BE A FIXED INSTALLATION AND PERMANENTLY CONNECTED TO THE SUPPLY.

Contents

Safety Information	2
Consignes de sécurité	8
Sicherheitsinformationen	10
Informazioni sulla sicurezza	12
Información de seguridad	14
Contact Information	16
Alpha Cyler specification	17
Declaration of Conformity	19

DOMINIQUE DUTSCHER SAS

Before use

Before using the Alpha Cyclor please ensure you have read this manual carefully. If there is any doubt relating to the proper use of this equipment, please contact your local distributor or PCRmax via the website: www.PCRmax.com or email Enquiries@bibby-scientific.com.

UNPACKING

When unpacking the unit please ensure that the following have been removed from the packaging:

Alpha Cyclor with block(s).

Mains cables (UK, EU, China and US).

USB stick containing operator's manual and default protocols.

AC-4 2x 15A Fuse which must be changed in certain territories and power supplies, see below.

AC-2 2x 12.5A Fuse which must be changed in certain territories and power supplies, see below.

The user is advised to keep the original packaging in case the instrument ever needs to be returned for service or repair. PCRmax accepts no responsibility for damage incurred unless the unit is correctly packed and transported in its original packaging.

IMPORTANT NOTE:

IF USING THE AC-2 / AC-4 AT 120V, BEFORE THE UNIT IS CONNECTED TO THE MAINS AND SWITCHED ON FOR THE FIRST TIME REPLACE THE FUSES.

FOR AC4 REPLACE 8A FUSES WITH 15A ANTI-SURGE FUSES.

FOR AC2 REPLACE 5A FUSES WITH 12.5A ANTI-SURGE FUSES.

THE FUSE HOLDERS ARE LOCATED BESIDE THE POWER SWITCH OF THE UNIT.

IF THE AC-4 IS INTENDED TO BE USED IN AUSTRALIA THE UNIT MUST BE A FIXED INSTALLATION AND CONNECTED DIRECTLY TO THE MAINS.

FUSE REMOVAL AND REPLACE

The anti-surge fuses are located on the right side of the system towards the rear, beside the power switch.

- Unscrew the fuse cover, using a flat head screwdriver or similar, and remove the fuse holder.
- Pop out the two fuses that the system is shipped with and replace with AC-2 (12.5A) or AC-4 (15A).

THIS MUST BE COMPLETED BEFORE CONNECTING THE SYSTEM TO THE LIVE MAIN AND POWERING UP THE UNIT.

NOTE:

A 250mm air space should be kept free at each side of the AC-2 / AC-4 unit (left and right). If two AC-2 / AC-4 units have to be positioned next to one another, a total of 500mm should be free between both units. If the systems have to be wracked the shelves should be 600mm apart from each other vertically, this promotes the correct airflow and heat distribution from the units. As air flow with the AC-2 / AC-4 is in thought the front of the unit and out the sides, caution should be taken to not block the path of air in and out of the unit.

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Safety information

Please read all the information in this manual before using the Alpha Cyclor.

WARNING

HIGH TEMPERATURES ARE DANGEROUS: they can cause serious burns to operators and ignite combustible material. Users should be aware of the following potential hazards:

- USE CARE AND WEAR PROTECTIVE GLOVES TO PROTECT HANDS.
- DO NOT use combustible substances near hot objects.
- DO NOT operate the instrument in the vicinity of inflammable liquids or gases.
- DO NOT place any liquid directly into the instrument.
- DO NOT touch the heated lid when system is running or within 10 minutes of the system completing a run.
- DO NOT touch the block when the system is running or within 10 minutes of the system completing a run.

OPERATOR SAFETY

All operators of PCRmax equipment must have available the relevant literature needed to ensure their safety. It is important that only suitably trained personnel operate this equipment, in accordance with the instructions contained in this manual and with general safety standards and procedures. If the equipment is used in a manner not specified by Bibby Scientific Ltd. the protection provided by the equipment to the operator may be impaired.

All PCRmax instruments are designed to conform to international safety requirements and are fitted with an over-temperature cut-out. If a safety problem should be encountered, switch off the unit at the mains socket and remove the plug from the electricity supply.

INSTALLATION

The instrument should be carried using both hands. Never move or carry the instrument when in use or connected to the mains electricity supply.

The AC-2 / AC-4 have a weight of 30kgs / 47kgs respectively; it is advised that two people lift and carry the AC-2, and three people lift the AC-4.

1. All PCRmax instruments are supplied with a power cable; this may be integral or plug-in.
2.  Before connecting the instrument to the mains electricity supply, check the voltage against the rating plate (located on the back of the unit). **Please note that the unit must be earthed to ensure proper electrical safety.**
3. The units are rated to operate at: for the AC-1 100-230V, 50/60Hz and for the AC-2 / AC-4 100-230V 50/60Hz.
4. Do not allow the Alpha Cyclor to overhang the bench.
5. Ensure the bench for the AC-2 / AC-4 can withstand 30kgs / 47kg of weight.
6. Place the unit on a suitable flat and level bench or in a fume cupboard if required, ensuring that the air inlet and outlet vents on the underside and rear are free from obstruction.
7. Ensure that the feet of the instrument do not overhang the side of the bench.
8. Plug the mains cable into the socket on the back (side in AC-2 / AC-4) of the instrument.
9. Switch on the instrument using the switch located on the back (side in AC-2 / AC-4) of the unit.

REPLACEMENT CABLE

Should the mains lead need replacement, a cable of 1.5 mm² of harmonized code H05VV-F 3G connected to an IEC320 plug should be used. **IF IN DOUBT CONSULT A QUALIFIED ELECTRICIAN.**

ENVIRONMENTAL CONDITIONS

The Alpha Cyler is designed operate under the following conditions:

- Indoor use
- Ambient temperature range +5°C to +35°C
- Altitude to 2000m
- Relative humidity not exceeding 80%
- Mains supply fluctuations not exceeding 10%
- Over voltage category II IEC 60364-4-443
- Pollution degree 2

Note: The control specifications are quoted at an ambient temperature of 20°C ±2°C.

The instrument has been tested for radio frequency interference and is certified under EN61326.

WARRANTY

Bibby Scientific Ltd warrants this instrument to be free from defects in material and workmanship, when used under normal laboratory conditions for two (2) years. In the event of a justified claim, Bibby Scientific Ltd will replace any defective component or replace the unit free of charge. This warranty does not apply if damage is caused by fire, accident, misuse, neglect, incorrect adjustment or repair, damage by incorrect installation, adaption, modification, fitting of non-approved parts or repair by unauthorised personnel. Bibby Scientific Ltd liability is limited to the cost of repair or replacement of the product and excludes in particular, indirect and consequential loss, damage, costs or expenses, including but not limited to wasted time, materials and expenditure or loss of use, profit, production, revenue, expected savings or goodwill. To make a claim please contact the supplier of the instrument. This warranty is in addition to, and does not affect any statutory rights.

This manual has been prepared for the convenience of PCRmax's customers and nothing in this manual shall be taken as a warranty, condition or representation concerning the description, merchantability, fitness for purpose or otherwise of the unit or components.

Notwithstanding the description and specification(s) of the instruments contained in the operator's manual, PCRmax reserves the right to make such changes as it sees fit to the instruments or to any of the components.

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Contact Information

For technical, sales or servicing information, contact your local PCRmax distributor or PCRmax directly at:

PCRmax

Beacon Road
Stone
Staffordshire
ST15 0SA
UK

Tel: +44 (0)1785 812121

Fax: +44 (0)1785 810405

E-mail: Enquiries@bibby-scientific.com

Help: techsupport@bibby-scientific.com

Web: www.pcrmax.com

User maintenance

Before cleaning your unit, disconnect it from the power supply. The outer case of the Alpha Cycler may be cleaned with a cloth dipped in water or ethanol: hexane or 50% methanol can also be used. No part of the case or cover should be immersed in the solvents. Do not use aggressive solvents such as acetone or abrasive cleaners. The block may be wiped with water, ethanol or propan-2-ol and may be decontaminated by wiping with 2% Neutracon® or 1% bleach solution.

Before using any cleaning or decontamination method except those recommended here, the responsible person should check with PCRmax that the proposed method will not damage the equipment.

Max Strength

70% IPA

2% neutracon detergent

Alpha Cycler Specification

Temperature

Block temperature range	4°C (10°C during cycling) to 100°C
Block uniformity (at 55°C)	±0.3°C
Temperature accuracy (at 55°C)	±0.25°C
Gradient availability	Between 10°C and 100°C
Maximum gradient	29°C on a 96 well block and 19°C on a 384 well block
Minimum gradient	1°C
Column uniformity with a 10°C gradient	±0.3°C
Final store	4°C to 100°C
Temperature set point adjustment	0.1°C

Heating/Cooling Rate

Maximum heating rate	3.4°C/s
Minimum heating rate	0.1°C/s
Maximum cooling rate	1.0°C/s
Minimum cooling rate	0.1°C/s

Heated lid

Selectable heated lid temperature	35 to 115°C or off
Lid uniformity	±5°C
Pre-heat lid	Yes
Warm up time from ambient	<2 min
Over-temperature cut-out	Yes
Regulated heated lid pressure	Manually adjustable

The heated lid is only operational if the block temperature is set above 35°C. Above figures are quoted at an ambient temperature of 20°C ±2°C

Programming

Program interface	10" and 7" HD touch screen
Operation system	Android
Maximum number of programs stored	1000
Maximum number of stages per program	25
Maximum number of steps	10
Maximum number of cycles per stage	99
Programmable ramp rate	Yes, 0.1°C/s steps
Maximum hold time	4h 59m 59s
Minimum hold time	1s
Program Wizard	Yes
Active sample cooling	Yes, 4°C
Advanced reporting options	Yes
Touchdown	Yes

Incremented/decremented temperature	Yes
Incremented/decremented time	Yes
Pause facility	Yes
Program naming	Alpha numeric plus symbols
Password protection	Yes
Keyword selection	Yes
Run completion time	Yes
Auto resume on power failure	Yes, always, never or user-defined
Oligonucleotide Tm calculator	Yes, based on the Nearest-Neighbour method ¹
Software updates	Free of charge from www.PCRmax.com

Communication ports

USB port on the front of the AC-1	Connection for USB memory stick
USB port on side of the AC-2 / AC-4	Connection for USB memory stick
Type B USB port on rear of unit	For Internal use only.

Dimensions

	AC-1	AC-2	AC-4
Height (mm)	330	330	330
Width (mm)	260	535	535
Length (mm)	430	470	700
Weight (kg)	10	30	47

Power

	AC-1	AC-2	AC-4
Voltage	100-230V, 50/60Hz	100-230V, 50/60Hz	100-230V, 50/60Hz
Power	425W	900W	1,600W
Fuse ratings		220V: 8 Amp x 2 120V: 15 Amp x 2	220V: 8 Amp x 2 120V: 15 Amp x 2

¹Breslauer, K.J.; Frank, R.; Blocker, H. and Marky, L.A. (1986) Proc. Natl. Acad. Sci. USA 83, pp 3746-3750.

Declaration of Conformity

PCRmax Alpha Cycler,

Alpha Cycler 1 AC-1, Alpha Cycler 2 AC-2 and Alpha Cycler 4 AC-4

AC196 96 well
AC1384 384 well
AC296 2 x 96 well
AC2196 1 x 96 / 1 x 384 well
AC2384 2 x 384 well
AC496 4 x 96 well
AC4384 4 x 384 well
AC4296 2 x 96 / 2 x 384 well
AC4396 3 x 96 / 1 x 384 well
AC4196 1 x 96 / 3 x 384 well

These products comply with the requirements of the EU Directives listed below:

2014/30/EU **EMC Directive.**
2014/35/EU **Low Voltage Directive (LVD)**
2011/65/EC **RoHS Directive**

Compliance with the requirements of these Directives is claimed by meeting the following standards:

EN 61326-1:2013 (Electrical Equipment for Measurement, Control and Laboratory use)
EN 61010-1:2010 (3rd edition) (Safety Requirements Electrical Equipment for Measurement, Control and Laboratory use)
EN 61010-2-010:2014 (Particular Requirement for Laboratory Equipment for Heating of Materials)

The above certificates and reports are available upon request.

Signed:  (Mr S. Marriott)

Date: **20/10/2016**

Technical Director
Bibby Scientific Ltd

Bibby Scientific Ltd - Stone - Staffs - ST15 0SA - UK
Tel: +44 (0) 1785 812121 - Fax +44 (0) 1785 813748

OMINIQUE DUTSCHER SAS

Bibby Scientific Limited

Beacon Road Stone
Staffordshire ST15 0SA
United Kingdom
Tel: +44 (0)1785 812121
Fax: +44 (0)1785 813748
e-mail: enquiries@pcrmax.com
www.pcrmax.com