

HyCell™ TransFx media

Enhance transfection, increase cell viability, and optimize productivity in HEK 293 and CHO cells



HyClone®
HyCell TransFx - C™
Without L - Glutamine
Without Pluronic F - 68

0.2 um Sterile Filtered
Cat No.: SH30941.01
Lot No.: **ABC12345**
Bottle No.: 0001
Volume: 1000mL
EXP Date: MMM/YYYY
Store At: 2 to 8° C

FOR FURTHER MANUFACTURING OR RESEARCH USE. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

HyClone Laboratories, Inc. For questions or comments visit
305 West 1000 South Or call technical support at
Logan, Utah 84321 + 1435/797-8000

SH30941-01

HyClone®
HyCell TransFx - H™
Without L - Glutamine
Without Pluronic F - 68

0.2 um Sterile Filtered
Cat No.: SH30939.01
Lot No.: **ABC12345**
Bottle No.: 0001
Volume: 1000mL
EXP Date: MMM/YYYY
Store At: 2 to 8° C

FOR FURTHER MANUFACTURING OR RESEARCH USE. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.

HyClone Laboratories, Inc. For questions or comments visit
310 West 1000 South Or call technical support at
Logan, Utah 84321 + 1435/797-8000

SH30939-01

Simplified transient transfection with HyCell TransFx media

HyClone™ HyCell TransFx-H and TransFx-C media are optimized for human embryonic kidney (HEK) 293 cells and Chinese hamster ovary (CHO) cells, respectively, to reduce testing and validation times, from research to scale-up. These regulatory-friendly media were developed through the HyClone Metabolic Pathway Design process and provide an optimized, nutrient-rich formulation, minimizing the requirements for further supplementation and the need to switch medium during the process.

Enhance transfection

Versatile and efficient

HyCell TransFx-H and TransFx-C media show enhanced transient transfection efficiencies compared with evaluated competitor products.

- Work across a variety of HEK 293 and CHO cell lines
- Apply across small and large production platforms

Consistent and scalable

HyCell TransFx media have been developed to provide consistent performance and to maximize process yields.

- Increase post-transfection viability
- Support high cell densities

Optimize productivity

Improve results and lower costs

Rapidly produce recombinant proteins for structural and functional characterization or screening campaigns.

- Rapid production of recombinant proteins
- Cost savings through fewer production runs
- Meet stringent timelines

Regulatory-friendly


Traceable and consistent

In addition to component traceability and regulatory-friendly characteristics, we maintain cGMP and ISO-compliant manufacturing to provide a quality product for cell culture and bioprocessing applications.


- Animal-derived component-free (ADCF) formulation
- Hydrolysate-free



Dramatically enhanced transient transfection and production

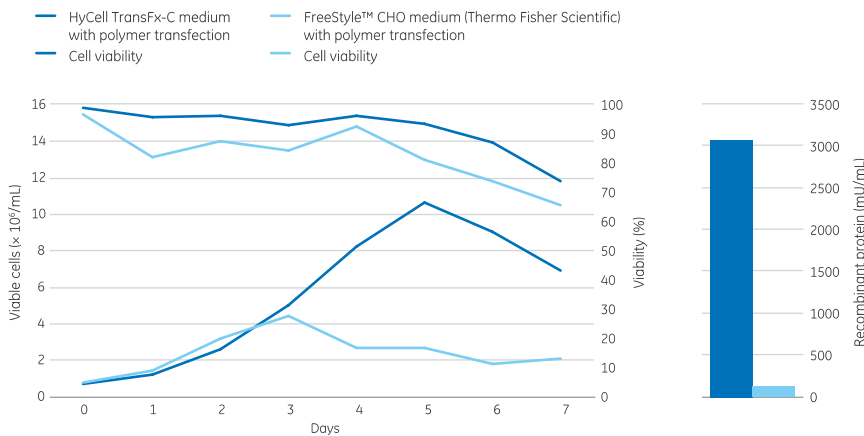

Improved growth, viability, and productivity profiles for CHO cells transiently producing recombinant protein in HyCell TransFxC medium.

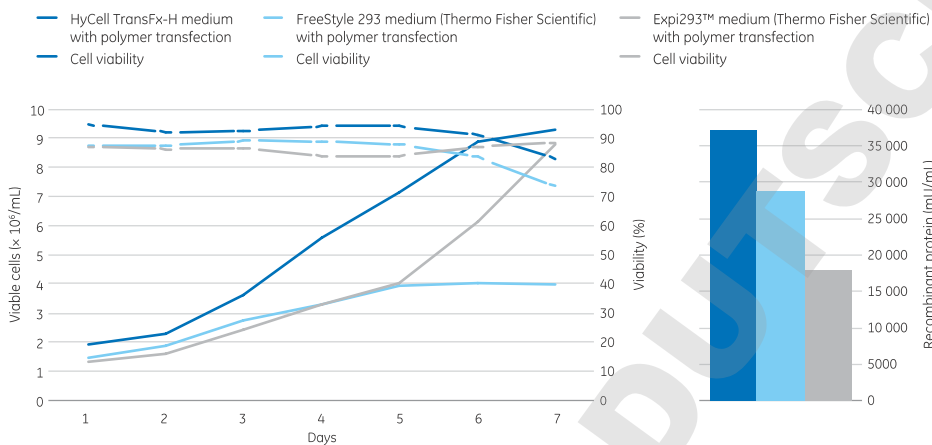

Improved growth, viability, and productivity profiles for HEK 293 cells transiently producing recombinant protein in HyCell TransFxC-H medium.

Meeting your cell viability and productivity needs

Balancing nutrient supply with metabolic waste to achieve high cell viability and productivity

HyClone Metabolic Pathway Design process is our proprietary approach to optimize and customize cell culture media. Over the last 45 years, our experts have refined the design process that analyzes the depletion of key nutrients and subsequent build-up of waste in the microenvironments of the cell culture by evaluating the distinct metabolic pathways of cells.

The Metabolic Pathway Design process enables optimization of medium nutrient consumption, while maximizing cell viability and productivity specific to your cell culture needs.

Request a FREE sample

For more information, visit www.gelifesciences.com/hyclone or contact your local sales representative today and request your FREE sample of HyCell TransFxC-H or TransFxC medium.*

*Up to 5 L medium. One request per customer please.



HyCell TransFx-C liquid medium

Without L-glutamine, without poloxamer 188

Description	Product code	
Liquid, 500mL PETE bottle	SH30941.01	○
Liquid, 1 L PETE bottle	SH30941.02	●
Liquid, 1 L BPC	SH30941.03	○
Liquid, 5 L BPC	SH30941.04	○
Liquid, 10 L BPC	SH30941.05	○
Liquid, 20 L BPC	SH30941.06	○
Liquid, 50 L BPC	SH30941.07	○
Liquid, 100 L BPC	SH30941.08	○
Liquid, 200 L BPC	SH30941.09	○
Liquid, 500 L BPC	SH30941.10	○

HyCell TransFx-H liquid medium

Without L-glutamine, without poloxamer 188

Description	Product code	
Liquid, 500 mL PETE bottle	SH30939.01	○
Liquid, 1 L PETE bottle	SH30939.02	●
Liquid, 1 L BPC	SH30939.03	○
Liquid, 5 L BPC	SH30939.04	○
Liquid, 10 L BPC	SH30939.05	○
Liquid, 20 L BPC	SH30939.06	○
Liquid, 50 L BPC	SH30939.07	○
Liquid, 100 L BPC	SH30939.08	○
Liquid, 200 L BPC	SH30939.09	○
Liquid, 500 L BPC	SH30939.10	○

HyCell TransFx-C powder medium

Without L-glutamine, without poloxamer 188

Description	Product code	
Powder, 5 L in HDPE bottle	SH30942.01	●
Powder, 10 L in HDPE bottle	SH30942.02	●
Powder, 50 L in HDPE bottle	SH30942.03	●
Powder, 100 L in HDPE bottle	SH30942.04	●
Powder, 500 L in Poly bag/pail	SH30942.05	●
Powder, 1000 L in Poly bag/pail	SH30942.06	●

HyCell TransFx-H powder medium

Without L-glutamine, without poloxamer 188

Description	Product code	
Powder, 5 L in HDPE bottle	SH30944.01	●
Powder, 10 L in HDPE bottle	SH30944.02	●
Powder, 50 L in HDPE bottle	SH30944.03	●
Powder, 100 L in HDPE bottle	SH30944.04	●
Powder, 500 L in Poly bag/pail	SH30944.05	●
Powder, 1000 L in Poly bag/pail	SH30944.06	●

● Item in stock; ○ Item is made to order. Lead times and minimum order quantities apply.

www.gelifesciences.com/hyclone

GE Healthcare Bio-Sciences AB
Björkgatan 30
751 84 Uppsala
Sweden

GE, GE monogram, HyClone, and HyCell are trademarks of General Electric Company.
Expi293 and FreeStyle are trademarks of Life Technologies Corp. All other third-party trademarks are the property of their respective owner.
© 2014–2015 General Electric Company. First published Mar. 2014
All goods and services are sold subject to the terms and conditions of sale of the company within GE Healthcare which supplies them.
A copy of these terms and conditions is available on request. Contact your local GE Healthcare representative for the most current information.
GE Healthcare UK Ltd., Amersham Place, Little Chalfont, Buckinghamshire, HP7 9NA, UK
GE Healthcare Europe GmbH, Munzinger Strasse 5, D-79111 Freiburg, Germany
GE Healthcare Bio-Sciences Corp., 100 Results Way, Marlborough, MA 01752, USA
GE Healthcare Dharmacon Inc., 2650 Crescent Dr, Lafayette, CO 80026, USA
HyClone Laboratories Inc., 925 W 1800 S, Logan, UT 84321, USA
GE Healthcare Japan Corp., Sanken Bldg., 3-25-1, Hyakunincho Shinjuku-ku, Tokyo 169-0073, Japan
For local office contact information, visit www.gelifesciences.com/contact

29105173 AC 10/2015